

Download site ALEXYS data system

Please check for updates of:

1. The ALEXYS data system
2. The USB dongle driver

On the following location: <http://www.antecleyden.com> (support section).

Never insert the dongle in the computer before installation of the driver software.

Instruction for USB dongle

The ALEXYS data system is protected by means of the supplied USB dongle. Always connect the USB dongle to a free USB port on your computer when running the ALEXYS data system. In the case no dongle is detected the ALEXYS software will start up in DEMO mode.

Under the Windows XP operating system the USB dongle is auto-detected and all necessary drivers are installed automatically. Under Windows 98 and Windows 2000 the dongle drivers should be installed manually. This should be done via the following procedure:

1. Run the program HDD32.exe
2. Restart the computer after installation
3. After restart insert the USB dongle to complete the installation.

In case you are unable to download the ALEXYS data system or dongle driver from the internet site please contact your local supplier for a request to ship the software to you on an installation CD.