


Legend LC connectors

- 250.1550 Ferrule Tefzel, 1/8", flangeless
250.1552 Nut PEEK, 1/8", flangeless, 1/4-28
- 250.1566 Ferrule CTFE, 1/8", collapsible
250.1568 Nut PEEK, 1/8", hex, 5/16-24
- 250.1558 Ferrule SS, 1/16", Valco
250.1560 Nut SS, 1/16", Valco, 10-32
- 250.1572 Fingertight fitting PEEK, hex, 10-32
- 250.1570 Fingertight fitting, PEEK, 10-32
- 250.1571 Fingertight Flex/Sencell, 10-32
- ⊗ 250.1574 Fingertight plug PEEK 1/16" 10-32
- ⊗ 250.1576 Fingertight plug PCTFE 1/4"-28

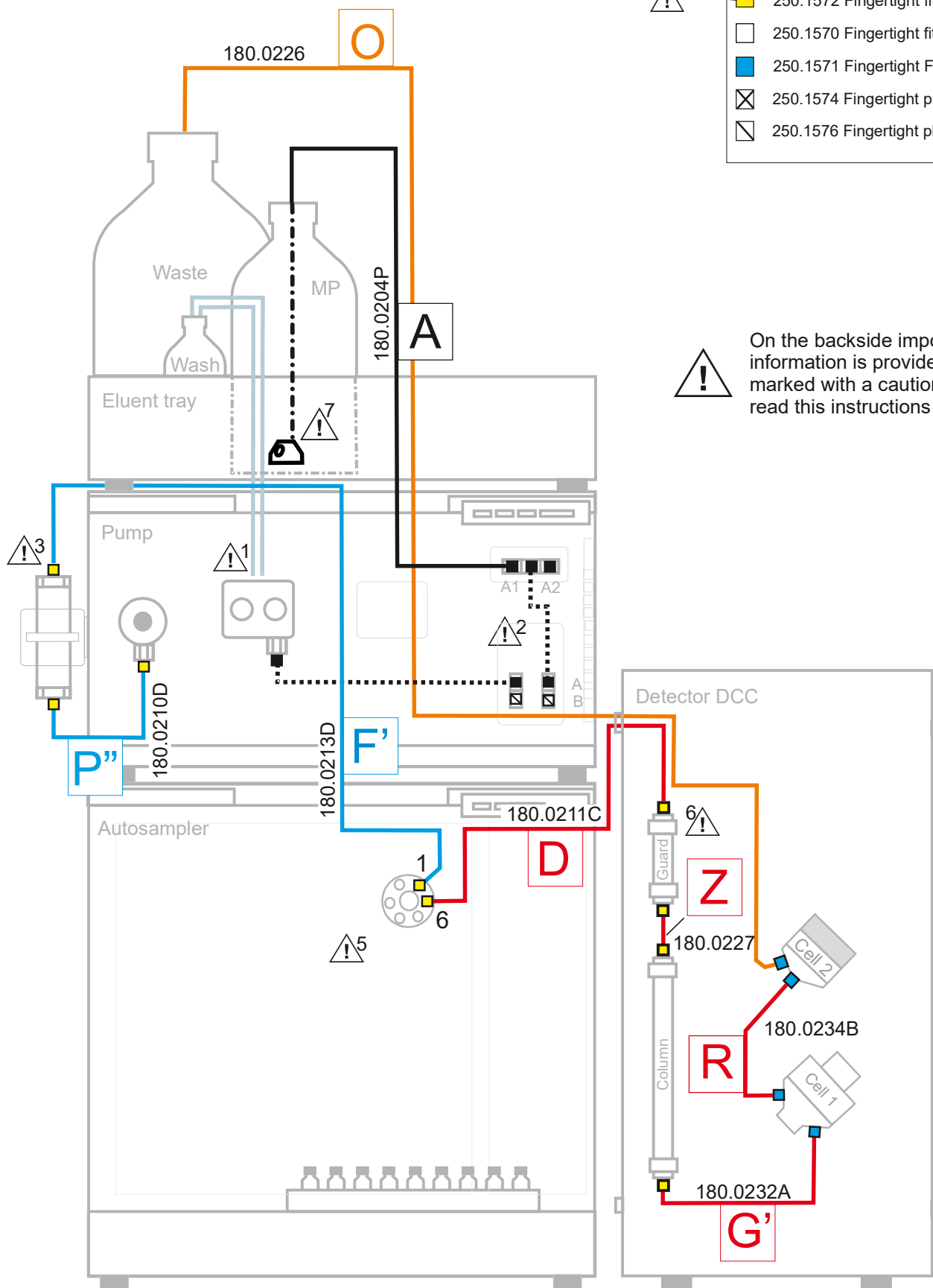

On the backside important installation information is provided for the parts marked with a caution sign. Please read this instructions before starting.





LC equipment should be used by trained laboratory personnel only. Use proper eye and skin protection when working with solvents under high pressures. Additional safety requirements or protection may be necessary depending on the chemicals used with this equipment.



IMPORTANT INSTALLATION INFORMATION

Read the LC connection kit install guide (pn 180.7001W), before installation of the LC connection kit.

1. The piston wash tubing is not part of the LC connection kit, it can be found in the P6.1L shipkit. Read instructions in manual pn 194.0010 how to install. The same applies for the AS 6.1 L waste tubing, it is part of the autosampler ship kit. Also install the supplied drainage system (funnels, hoses, nozzles) for leak management as described in the user manual of pump and autosampler.
2. In case of a P6.1L pump with integrated degasser, the tubing assembly (dotted lines) are already pre-installed. Connect the mobile phase bottle to port A1 of the solvent selection valve (SSV) using tubing assembly 180.0204 (B).
3. The pulse damper, p/n 250.AZZ00NB has to be mounted on a bracket on the side of the P6.1 L pump using the parts and instructions supplied in the shipping box of the damper.
4. All tubing connections in the high pressure part of the LC system are made with PEEK HEX head fingertight fittings (pn 250.1572). Use the tightening tool (pn 250.0094) delivered with the kit to tighten the finger tight fittings firmly. It will prevent that the tubing will slip out of the connector when the system is under pressure. Do not use a wrench, because with a wrench one might apply too much force resulting in rupture/damage of the plastic connector.
5. For optimal performance it is required to passivate all metal parts in this system using a 15% solution of HNO₃. See LC connection kit install guide for instructions.
6. In the case a guard column is used, connect the guard to the analytical column using pn 180.0227 precolumn assembly.
7. Replacement filters can be ordered as '*PEEK filter for 1/8" MP inlet, 2um*' (pn. 250.1704). When replacing, push the tubing deep into the port that is connected to the bottom inlet until it holds.

Legend LC tubing*:

 FEP 1/8" OD, 1.59 mm ID, transparent	 PEEK 1/16" OD, 0.064 mm ID, pink-striped
 Stainless Steel 1/32" OD, 0.25 mm ID, with 1/16" ends	 Silicone 3 mm OD, 1 mm ID, transparent
 PEEK 1/16" OD, 0.25 mm ID, blue-striped	 PEEK 1/16" OD, 0.50 mm ID, orange-striped
 PEEK 1/16" OD, 0.13 mm ID, red-striped	 Silicone 10 mm OD, 8 mm ID, transparent

*) Note that the color coding of the LC tubing in the drawing does not necessarily reflect the real color of the corresponding tubing.