

Clarity

Reference Guide

Clarity

ENG

Code/Rev.: M018/50A
Date: 9. 9. 2013

Phone: +420 251 013 400
Fax: +420 251 013 401
clarity@dataapex.com
www.dataapex.com

DataApex Ltd.
Petrzilkova 2583/13
158 00 Prague 5
The Czech Republic

Sections of the manual connected only to the **Clarity Full** version are marked with the  icon.

Clarity[®], DataApex[®] and [®] are trademarks of DataApex Ltd. Microsoft[®] and Windows[™] are trademarks of Microsoft Corporation.

DataApex reserves the right to make changes to manuals without prior notice. Updated manuals can be downloaded from www.dataapex.com.

Author: DM

Contents

1 Clarity	1
1.1 System	2
1.1.1 User Accounts	3
1.1.2 User Details	8
1.1.3 Configuration	9
1.1.3.1 Left section of the System Configuration dialog	10
1.1.3.2 Available Control Modules	12
1.1.3.3 A/D Converter Setup	15
1.1.3.4 Right section of System Configuration dialog	16
1.1.4 Directories	21
1.1.5 GLP Options...	23
1.1.6 Digital Outputs	25
1.1.6.1 Select Control	26
1.2 Login	27
1.3 Logout	28
1.4 Lock	29
1.5 View	30
1.5.1 Customize...	31
1.6 Help	32
1.6.1 Register	32
1.6.1.1 Register Clarity question	34
1.6.2 Check for Updates	35
1.6.3 User Code...	36
1.6.3.1 Language Selection	37
1.6.4 Send Report by E-Mail	38
1.6.5 About...	39
2 Instrument	40
2.1 Clarity Lite window specifics	41
2.2 File	42
2.2.1 Open Method	44
2.2.1.1 Open Method (Sequence, Chromatogram, Calibration)	45
2.2.2 Send Method by Email	47
2.2.3 Project	48
2.2.3.1 Error Opening Project	49
2.2.4 Archive	50
2.2.5 Restore	52
2.3 Monitor	54
2.3.1 Device Monitor	55
2.3.1.1 File	56
2.3.1.2 Control	56
2.3.1.3 View	57
2.3.1.4 Window	57
2.3.1.5 Help	57
2.4 Analysis	58

2.4.1 Single	59
2.4.1.1 Chromatogram File Name	64
2.4.1.2 Rich Text Edit	66
2.4.2 Sequence	66
2.4.3 Batch	67
2.5 Method	71
2.5.1 Event Table	73
2.5.1.1 Events used in Active Sequence	79
2.5.2 Measurement	80
2.5.3 Acquisition	82
2.5.4 Integration	84
2.5.4.1 Integration table items	85
2.5.5 Calculation	88
2.5.6 Advanced	92
2.5.7 GC (LC) Control, LC Gradient	95
2.5.8 AS	95
2.6 Setting	96
2.6.1 Export Data	96
2.6.2 Postrun...	100
2.7 View	103
2.7.1 Options...	103
2.7.1.1 General tab	104
2.7.1.2 Graph tab	107
2.7.1.3 Axes Appearance tab	110
2.7.1.4 Signals & Curves tab	111
2.7.1.5 Auxiliary Signals tab	112
2.7.1.6 Gradient tab	113
2.7.1.7 Directories tab	115
2.8 Window	116
2.9 Help	117
3 Chromatogram	118
3.1 File	122
3.1.1 Overlay Mode	126
3.1.2 Open Chromatogram	127
3.1.3 Open Chromatogram Sequentially	129
3.1.4 Open Chromatogram from Sequence	131
3.1.5 Save As...	132
3.1.6 Import Chromatogram...	133
3.1.6.1 AIA	134
3.1.6.2 Text and EZChrom ASCII	136
3.1.6.3 Import RAW...	139
3.1.7 Export Chromatogram	140
3.1.8 Sign	142
3.1.9 Verify	144
3.1.10 PostRun Setting (From Chromatogram Window)	145
3.2 Edit	148
3.2.1 Add User Column	151

3.2.2 Setup Columns...	154
3.3 Display	156
3.3.1 Preview Graph	157
3.3.2 Properties	159
3.3.2.1 Graph Properties - Graph tab	160
3.3.2.2 Graph Properties - Axes Appearance tab	161
3.3.2.3 Graph Properties - Time Axis tab	162
3.3.2.4 Graph Properties - Signal Axis tab	164
3.3.2.5 Graph Properties - Signals tab	166
3.3.2.6 Graph Properties - Auxiliary Signals tab	169
3.3.2.7 Graph Properties - Gradient tab	170
3.4 Chromatogram	171
3.4.1 Baseline	174
3.4.1.1 Lock	175
3.4.1.2 Valley	175
3.4.1.3 Together	175
3.4.1.4 Forward Horizontal	175
3.4.1.5 Backward Horizontal	175
3.4.1.6 Front Tangent	176
3.4.1.7 Tail Tangent	176
3.4.1.8 Clamp Negative	176
3.4.1.9 Cut Negative	176
3.4.1.10 Reject Negative	176
3.4.1.11 Allow Crossing	176
3.4.1.12 Spike Removal	177
3.4.2 Peak	178
3.4.2.1 Start	178
3.4.2.2 End	179
3.4.2.3 Both	179
3.4.2.4 Add Positive	179
3.4.2.5 Add Negative	180
3.4.2.6 Force Peak Name	180
3.4.2.7 Solvent Peak	181
3.4.2.8 Groups	182
3.4.3 Integration	183
3.4.3.1 Local Peak Width	184
3.4.3.2 Local Threshold	186
3.4.3.3 Integration Interval	187
3.4.3.4 Detect Negative	187
3.4.3.5 Minimal Area	188
3.4.3.6 Minimal Height	188
3.4.3.7 Minimal Half Width	188
3.4.3.8 Detector Delay	189
3.4.3.9 FFT Filter	189
3.4.3.10 Savitzky-Golay Filter	189
3.4.3.11 Moving Average Filter	189
3.4.3.12 Spike Filter	189

3.4.4 Separation	190
3.4.4.1 Valley To Valley Slope	190
3.4.4.2 Tangent Area Ratio	190
3.4.4.3 Tangent Slope Ratio	190
3.4.5 Noise & Drift	191
3.4.5.1 Noise Evaluation	192
3.4.5.2 ASTM Noise Evaluation	193
3.4.5.3 6-Sigma Noise Evaluation	193
3.4.5.4 Drift Evaluation	193
3.4.6 Overlay	194
3.4.6.1 Move	195
3.4.6.2 Scale	196
3.4.6.3 Original	196
3.4.6.4 3D View	196
3.4.6.5 Clear 3D	197
3.4.6.6 Properties... ..	197
3.4.6.7 Mathematics... (Mathematical Operations dialog)	198
3.4.7 Chromatograms	200
3.4.8 Create Label	201
3.4.9 Set Signal Names	204
3.5 Method	205
3.5.1 Calculation	207
3.5.2 Integration	212
3.5.3 Measurement	216
3.6 Results	220
3.6.1 Local Table Menu	220
3.6.2 Result Table	222
3.6.3 Summary Table	228
3.6.3.1 Summary Options	230
3.6.4 Performance Table	232
3.7 SST	235
3.8 View	236
3.9 Window	237
4 Calibration	238
4.1 Calculate By	251
4.2 File	252
4.2.1 Save As... ..	255
4.3 Edit	256
4.4 Display	257
4.4.1 Graph Properties	258
4.5 Calibration	260
4.5.1 Add All	263
4.5.2 Add Existing	264
4.5.3 Add Peak	265
4.5.3.1 Calibrate Peak	265
4.5.3.2 Recalibrate Peak	267
4.5.4 Add Group	269

4.5.4.1 Calibrate Group	269
4.5.4.2 Recalibrate Group	269
4.5.5 Options	270
4.5.5.1 Options	275
4.6 View	277
4.7 Window	278
5 Data Acquisition	279
5.1 File	281
5.2 Analysis	282
5.3 Display	285
5.3.1 Properties... ..	286
5.4 View	289
5.4.1 Axes Ranges	291
5.5 Window	293
6 Sequence	294
6.1 File	301
6.1.1 Import... ..	303
6.2 Edit	306
6.2.1 Fill Series	308
6.3 Sequence	309
6.3.1 Sequence Options	313
6.4 View	317
6.5 Window	317
7 Audit Trail	318
7.1 Station Audit Trail	319
7.1.1 File	319
7.1.2 Open Audit Trail (Append)	320
7.1.3 Export	321
7.1.4 Edit	322
7.1.5 Filter	323
7.1.6 View	323
7.1.7 Properties... ..	324
7.1.8 Window	326
7.2 Local Audit Trails	327
7.2.1 File	328
7.2.2 Edit	328
7.2.3 View	329
7.2.4 Properties... ..	330
8 Report Setup	333
8.1 List of tabs:	335
8.2 Page Setup	336
8.3 Lab. Header	338
8.4 Report Header	340
8.5 Method	342
8.6 Calibration	344
8.7 Chromatogram	346
8.8 Results	348

8.9 Sequence	349
8.10 Audit & Signatures	350
8.11 Print Preview	352
8.12 Print	354
9 Troubleshooting	355
9.1 Report error after Clarity crashed	355
9.2 How to get help	355
9 Appendices	356
9.3 List of terms	356
9.4 Shortcuts	362
10 Technical Specification	364
10.1 Declaration of Software Validation	364
10.2 Command line parameters	364
10.2.1 List of commands	365
10.2.2 Examples	386
10.2.2.1 Example 1:	386
10.2.2.2 Example 2:	386
10.2.2.3 Example 3:	386
10.2.2.4 Notes:	387
10.3 clarity.ini file	388
Index	390

To facilitate the orientation in the **Reference Guide** manual and **Clarity** chromatographystation, different fonts are used throughout the manual. Meanings of these fonts are:

Instrument (blue text) marks the name of the window, to which the text refers.

Open File (italics) describes the commands and names of fields in **Clarity**, parameters that can be entered into them or a window or dialog name (when you already are in the topic describing the window).

WORK1 (capitals) indicates the name of the file and/or directory.

ACTIVE (capital italics) marks the state of the station or its part.

The bold text is sometimes also used for important parts of the text and the name of the **Clarity** station. Moreover, there are text sections written in format other than normal text. These sections are formatted as follows:

Note: Notifies the reader of possibly interesting information.

Caution: Warns the user of possibly dangerous or very important information.

I Marks the problem statement or trouble question.

Description: Presents any closer information on the problem, describes its causes etc.

Solution: Marks the response to the question, presents a procedure how to remove it.

1 Clarity ✓ full version

The main window of the **Clarity** station contains Instrument icons as well as a menu for the configuration of the station, opening and closing Instruments etc.

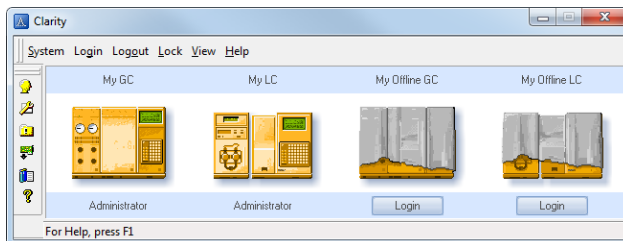


Fig 1: Main Clarity window

Clicking the Instrument icon or the *Login* button will open the **Instrument** window.

In the screenshot there are two opened Instruments (My GC and My LC), both used by the user *Administrator*. The other two instruments (My Offline GC and My Offline LC) are closed.

Above the instrument is the Name of the Instrument and below the closed instrument is the Login button and after opening it will display the Name of the currently logged user.

While trying to open **Clarity Instruments** or some other dialogs (**User Accounts**, **System Configuration** and others) or to invoke certain commands, the **Login Dialog** appears. This dialog is described in detail in the

Starting and closing Clarity

While the **Clarity** is starting or closing, the starting or closing splashscreen is visible instead of the main **Clarity** window. This might be important mainly with some of the control modules because of the quite long times when **Clarity** loads or unloads the necessary drivers.

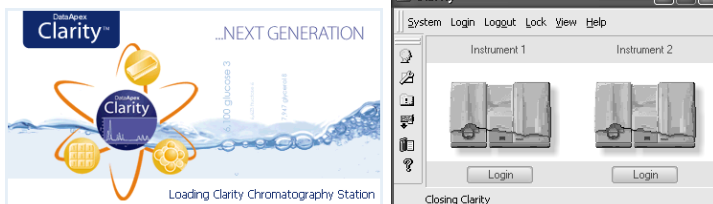


Fig 2: **Clarity** starting and closing splashscreens

Clarity Switch

It is possible to scale down the main window by double-clicking the window title bar or by the command *View - Switch*.
To restore the original size, double-click the title bar once more.

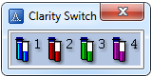


Fig 3: Main window as Clarity Switch

1.1 System

Submenu with system commands:

	User Accounts	Sets the access rights and passwords particularly for the administration of user accounts.
	User Details	Shows the details of user account settings even to the users that don't have the access to the User Accounts dialog.
	Configuration	Sets the station HW configuration.
	Directories	Sets directories for individual Instruments.
	Clarity2Go	Sets the communication with server for watching instruments in your iPhone or Android phone.
	Digital Outputs	Sets the outputs control.
	Audit Trail	Displays the records of selected station tasks and sets their logging.
	Exit	Exits the program. If the measurement is active on any Instrument, you will be notified. You will also be prompted to save changed files.

Invoking commands from the *System* menu or the *Help - User Code...* command requires all instruments to be closed. The user has to Logout all instruments. In case the *Logout All Now* button is disabled, an **Instrument** is locked. Unlock the **Instrument** manually (for more details see the chapter **Lock** on pg 29)..

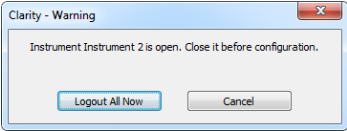


Fig 4: Logout all instruments

1.1.1 User Accounts

Administration of user accounts, access rights and passwords. Settings are saved in the CLARITY.PSW file. When the *System - User Accounts...* command or  icon is invoked, the **User Accounts** dialog will appear. Here you will be able to define all users who work with the station and set their access rights.

Fig 5: User Accounts

Note: Setting access rights is used predominantly to prevent intervention by untrained persons. This is not conceived as protection of data against theft - such protection would require setting access rights in **MS Windows**. If any file is modified outside the **Clarity** station environment, it would be identifiable due to built-in security features.

User List

Provides a complete list of previously defined users. The name of the user to whom the content of all displayed fields relates will be displayed in inverse mode. Other users are selected by using arrow keys or clicking your mouse on the desired user name.

New

Adds a new user. During this operation, the new *Username* is checked against existing usernames for possible collision.

Delete

Erases a selected user from the user list. When deleting all users the station automatically switches to “all users” (no password) mode.

Password Restrictions

Caution: Password Restrictions parameters are common to all users.

Min. Length

Refers to the minimum length of the password entered. A password shorter than the prescribed number of characters will not be accepted.

Life Time

Refers to the period of a password's validity. After its expiration the user must select a new password.

Expiration Warning

Alerts the user as to the number of days prior to a password's expiration.

Password Reuse

Relates to the period from which an already used password can be used again.

Auto Lock

If the user doesn't have any activity in Windows, all opened Instruments will be locked after specified time in minutes. All windows opened on instruments will be hidden and the user has to unlock the [Instrument](#) by clicking on its picture in the [Clarity](#) window.

User Details for:

The fields displayed in the lower part of the window always relate to the user whose username is selected from the *User List*. The selected name will be displayed in inverse mode (i.e. white on a blue background). Only a user with adequate access rights can change the settings of those access rights (selected ☒ *Open User Accounts* under *User Access Rights*).

User**User Name**

The User's name will automatically be placed in all headers of files that are created while working in the station. The username will be displayed along with the *Created by* or *Analyst* phrase when a file is selected. It is compulsory for a user to fill in this field and forbidden for two users to have the same username.

Desktop File

Refers to the name of the file (*.DSK) in which the desktop settings are saved. The file contains settings regarding size, location and visibility of the Instrument windows as well as all the amendable Instrument

parameters which are not part of system files. If this field is left empty, it will be filled in automatically according to the *User Name* field.

This setting is common to all Instruments the user has available.

Note: If you wish to share the same settings for several users, you can link them to the same desktop file.

Description

Provides a user's detailed description.

Note: *User Name* and *Description* are part of signature during electronic signing.

Other Users Can

Read & Write

Other users can view and edit your files.

Read

Other users can view your files but cannot edit or otherwise modify them.

No Access

Other users does not have any access to your files (even for reading them).

Access To

Instrument X

Defines which Instruments the user may access within a station.

User Access Rights

Open User Accounts

A user with access to user accounts can add, delete or change the access rights of other users. A selected checkbox is essentially equivalent to allocating administrator rights. It also allows the user to use the *System - GLP Options...* command to change the rules for files overwrite.

Open Configuration

Checking this field permits a specific user to change the configuration settings of the station.

Edit Method

Enables the user to create, delete or edit a method.

Note: If the method is changed by any reason (the MODIFIED inscription is shown in the [Instrument](#) window header) and the user without the *Edit Method* privilege should leave the [Instrument](#) window, all changes to the method are automatically discarded.

Edit Chromatogram

Enables the user to edit chromatograms.

Edit Calibration

Enables the user to create and edit calibrations.

Projects

Enables the user to change projects.

Import Chromatogram

Enables the user to import chromatograms.

Edit Sequence

Enables the user to create and edit sequences.

Edit Report Style

Enables the user to create and edit report styles.

Select Method

Enables the user to select a method for single analysis (Single Run).

Open Audit Trail Settings

Accesses the setting changes in station task logs.

Post Run Settings

Enables the user to change the post run settings in the [PostRun Settings](#) dialog, [Sequence](#) window and the [Batch](#) dialog.

Archive/Restore

Permits the user to archive/delete/restore files and projects in the **Clarity** software.

User Info**Password**

Notifies whether a user has already selected a password or not. Possible conditions are *Submitted* - password was already entered, *Blank* - password not entered yet.

Password Changed:

Date of last password change.

Last Login

Date of last login.

Change Password

Changes password. After the prompt a [New Password](#) dialog appears. Here you may enter a new password.

Certificate

Sets a certificate which will be available as a default option while using the electronic signature - the *Sign* function in the [Chromatogram](#) window.

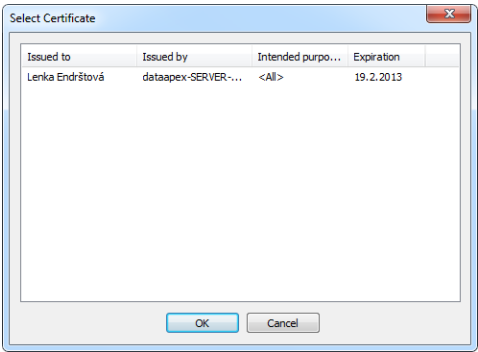


Fig 6: Select Certificate

Select Certificate

In the [Select Certificate](#) dialog selects one of the certificates installed on your computer.

Clear Certificate

Cancels an implicit certificate.

Certificate For Exporting To PDF

Invokes the [Open](#) dialog for selecting the PKCS#12 private key. Documents exported from **Clarity** are then signed automatically with this certificate.

Note: Certificates are issued by independent agencies, e.g. Thawte (www.thawte.com) or Verisign (www.verisign.com).

1.1.2 User Details

In the **User Details** dialog (accessible from **Clarity** window using the **System** menu) can users without the authorization to enter the **User Accounts** dialog view their own user account settings and enter or edit the password to their own account. Here they may also select a certificate for digital signing.

Note: If there are no users defined, the *User Details* command will not be accessible.

The description of individual fields of this dialog is identical to the corresponding fields in the **User Accounts** dialog

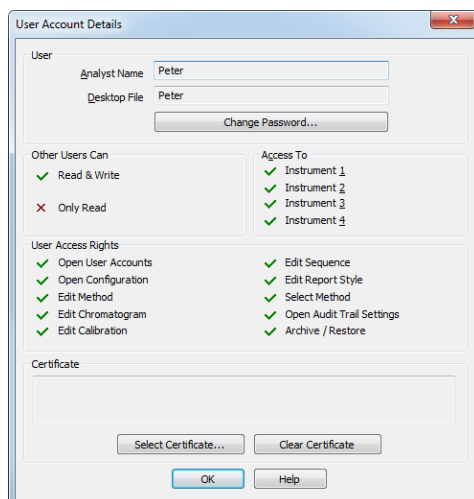


Fig 7: User Account Details

1.1.3 Configuration

Serves for the setting of the station's configuration, which is saved in CLARITY.CFG file.

After prompting the **System - Configuration...** command or pressing the  icon, the **System Configuration** dialog will be opened. Access to the **System Configuration** dialog is permitted only to users with adequate access rights settings. Access is also denied if any of the **Clarity Instruments** is not closed (you will be prompted to close opened **Instruments**).

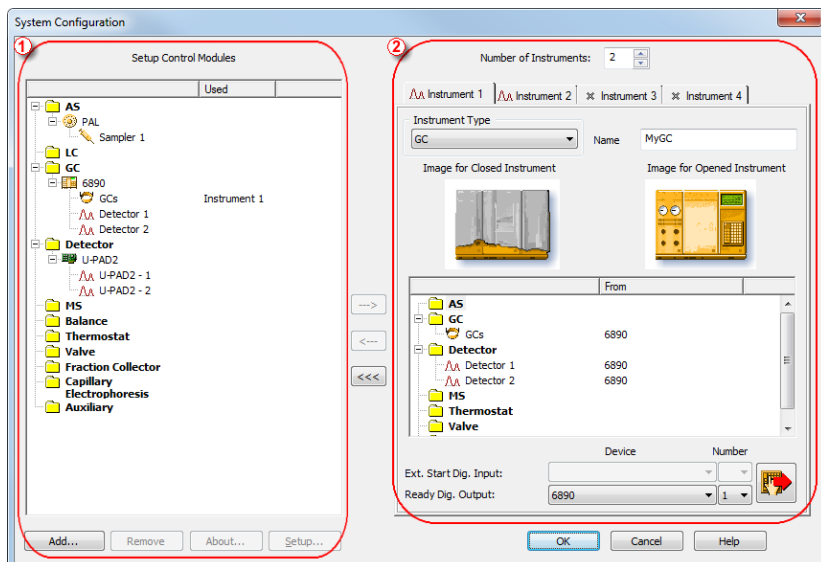


Fig 8: System Configuration

The left section of the dialog ① is used to configure equipment in the **Clarity** station, while the right section ② is where the equipment is allocated to individual Instruments.

Note: Step by step procedure of configuration is described in **Clarity Getting Started** manual. The configuration of individual Control modules can be found in corresponding **Clarity Control** manual.

1.1.3.1 Left section of the System Configuration dialog

Setup Control Modules

Lists the configured equipment (e.g. internal or external A/D converters for data acquisition, directly controlled chromatographs, autosamplers, LC pumps, etc.).

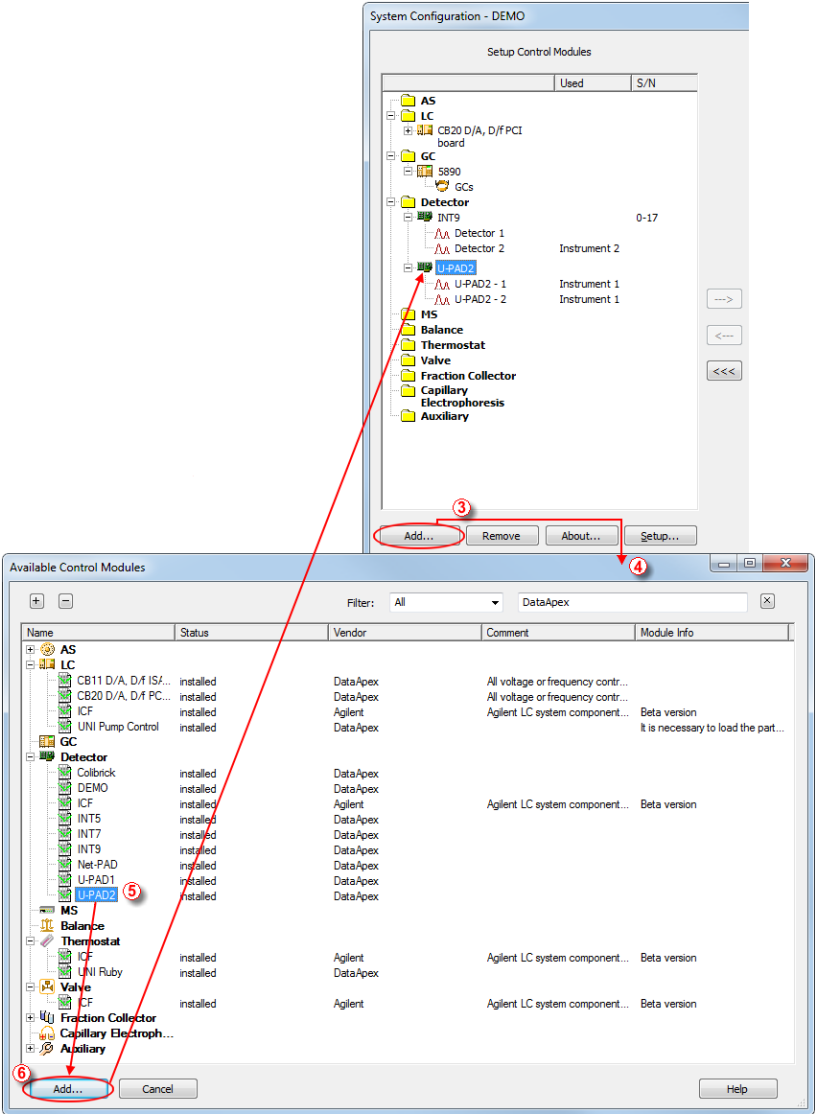


Fig 9: System Configuration - left part and Available Control Modules dialogs

Add ③

Adds and configures drivers for any available equipment. After prompting, the [Available Control Modules](#) dialog ④ will appear with a list of all available drivers for data acquisition and chromatographic equipment control (for more details see the chapter **Available Control Modules** on pg 12).

Note: If you wish to use one control module for two various Instruments, you have to add it to the **Clarity** station twice.

Select the required item (name of the driver) ⑤. By clicking the *Add* ⑥ button or double clicking on the item you can add the selected driver to the [Setup Control Modules](#) list. [X Driver Setup](#) dialog (where X stands for the device name) will appear, enabling you to fill in the exact configuration of the device (for more details see the chapter **A/D Converter Setup** on pg 15).

Note: Drivers correspond to CSW???.DLL files, which are present in the station's main station directory. The physical presence of these files in the directory is not sufficient. Files must be "registered", which is done automatically by the installation program. Files both registered and present can be checked in the station's main window using *Help - About* command. The list can be found on the [System Files](#) tab. Reinstalling the station easily resolves all file registration problems.

Remove

Removes the selected configured equipment from the **Clarity** station.

About

Displays the information regarding the selected equipment's driver.

Setup

Sets driver parameters for selected equipment in [X Driver Setup](#) dialog, where X is the equipment name, e.g. INT7, U-PAD, HP6890, etc.

Note: The [X Driver Setup](#) dialog can be also called up by double clicking on the name or icon of the device and is automatically displayed while adding the device to the **Clarity** station.

1.1.3.2 Available Control Modules

The **Available Control Modules** dialog allows to add the particular control module driver into the **System Configuration** dialog. Try to search the driver by its name using the *Filter* function or select it manually in the list of the devices section.

Note: OEM versions of various instruments can be usually set both under their original name or the OEM names.

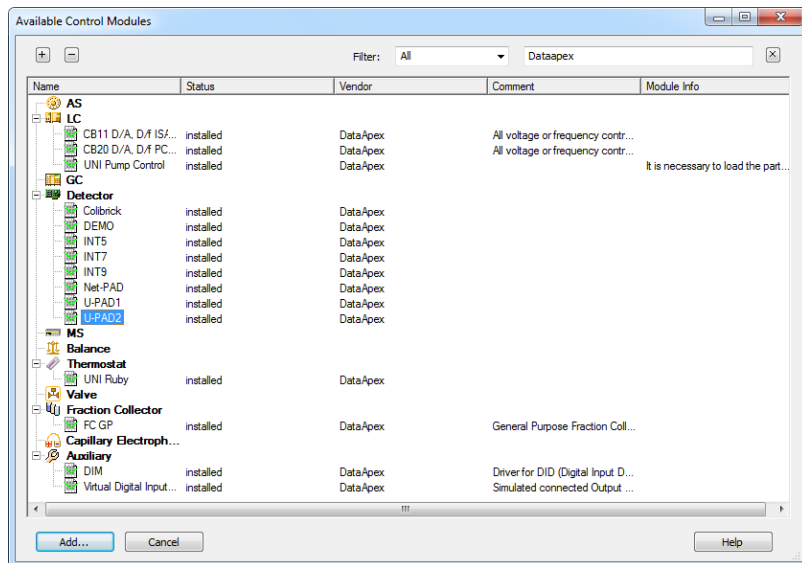


Fig 10: Available Control Modules

+ / - Buttons

Click the **+** button to unfold all sections in the **Available Control Modules** dialog or click the **-** button to fold all sections.

Filter

Serves for the simplification of the driver selection. Simply enter any text string in the second field to search all columns for this string, or restrict the search to any desired column by selecting the column name from the drop down list in the *Filter* field. The **Fig 10** on pg 12. shows all drivers containing the "DataApex" keyword filtered.

Add

Opens the setup dialog for the device selected in the list of the drivers (as the U-PAD2 is in the **Fig 10** on pg 12.). After selecting the desired driver and clicking this button the setup dialog for the driver is displayed (for more details see the manual of the appropriate control module or hardware). When correctly configured, the driver is added to the left

section of the Configuration Dialog and the [Available Control Modules](#) dialog is closed.

Cancel

Exits the dialog without any action taken.

List of the drivers:

The content of the driver's list section is dependent on the *Filter* used. All available drivers are displayed only in case no filter has been used.

Name

Contains the names of the drivers grouped in sections, sorted alphabetically.

Status

Informs the user if the control module is installed on the system or not. In the list you can see all the control modules and hardware which are controlled by **Clarity**. In typical **Clarity** Installation, not all drivers are installed on the system. By clicking on the *Add...* button an information how to install the driver will be displayed.

Vendor

Holds the information on the device vendor.

Comment

This column may hold some other type of driver description (e.g. for what sub-versions of instruments is the driver suitable and so on).

Module Info

Contains the information on the control module state (e.g. whether it is an Alpha or Beta version) and the author of the device driver. Items with the Module Info column empty are both in the Release version (had been thoroughly tested) and have been developed by the

DataApex company.

1.1.3.2.1 Control Module Is Not Installed

This dialog is displayed in case you tried to add a control module to **System Configuration** which is not installed in **Clarity** yet. You can find here an information how to install the module. Mostly it is necessary to reinstall **Clarity** and select the *Full* installation or select the specific module to install. In other cases, you must contact the control module vendor to obtain it, etc.

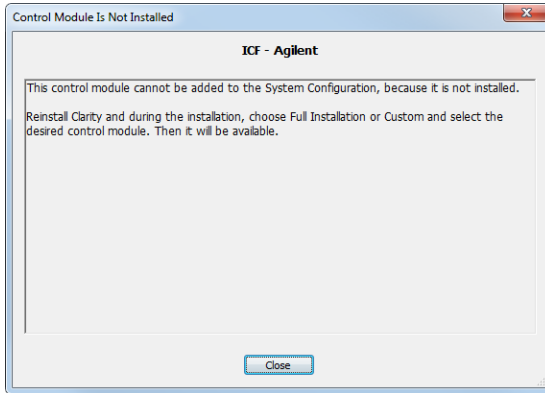


Fig 11: Control module is not installed dialog

1.1.3.3 A/D Converter Setup

Signal Names and *Units* are stored for individual signals according to the configuration of the detector. For **DataApex A/D Converters** the signal units can be configured in the **Setup** dialog accessible from **System Configuration** dialog.

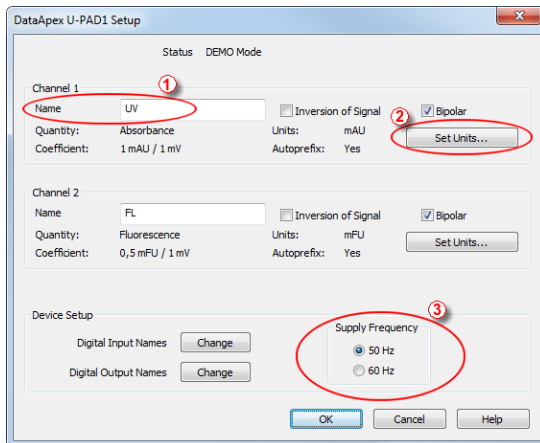


Fig 12: DataApex A/D Converter Setup dialog

For each channel provided by the A/D Converter, the name of the signal can be edited in the Name ① field of the corresponding **X Device Setup** dialog (where X is the name of the device). The Set Units... ② button can be used to change other signal parameters.

Note: In **X Device Setup** dialog of some **DataApex A/D Converter** types, desired **Supply Frequency** ③ can also be selected.

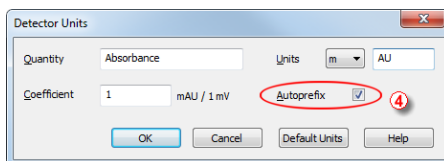


Fig 13: Detector Units

When the unit prefix in **Graph** window should be adjusted automatically, enter the units and prefix separately and check the *Autoprefix* ④ box. In tables, the units with the prefix entered will be used.

1.1.3.4 Right section of System Configuration dialog

Sets the number of displayed Instruments and allocates equipment to the relevant Instruments.

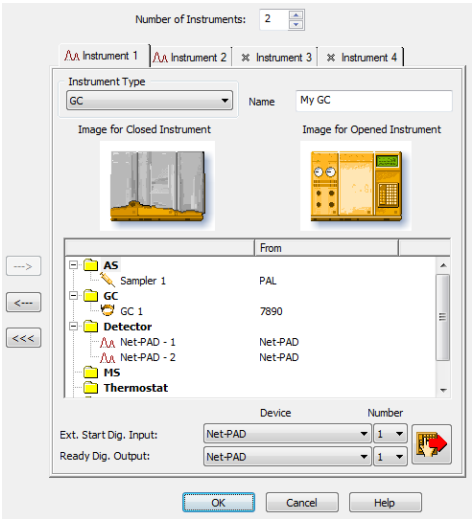


Fig 14: System Configuration - right section

Allocate the configured equipment (displayed on *Setup Control Modules* pane in the left section) to the relevant Instrument (in the right section) by “drag and drop” with the mouse or using the following buttons located in the center of the window:

Button	Description
--->	Allocates selected equipment to an active Instrument.
<---	Removes selected equipment from an active Instrument.
<<---	Removes all equipment from an active Instrument.

Number of Instruments

Provides the possibility to set the number of displayed Instruments. 1 to 4 Instruments can be displayed at a time, but you are only allowed to measure with as many Instruments as you have purchased (these are so called “on-line Instruments”). However, you can display more Instruments and use them for “off-line” preparation of sequences and evaluation of chromatograms. On-line Instruments have a red chromatogram  symbol in their header, while off-line ones have a blue  symbol. The tab(s) of any off-line Instrument(s) will have the “Offline” inscription on them. In the following image there are three Instruments set for display. The first two are on-line and the third is off-line.

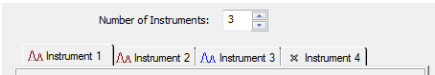


Fig 15: Tabs of individual Instruments

Tabs of Instruments 1 - 4

Tabs of individual Instruments. You can enter an Instrument name and Instrument type on each tab and connect drivers which will enable the measuring of a signal and directly control the chromatograph and other devices.

Tabs are indicated in the header with icons:

 (red chromatogram) indicates an online Instrument, which is usable for the data acquisition (as well as to evaluate chromatograms, prepare methods, etc.).

 (blue chromatogram) indicates an offline Instrument, which cannot be used to acquire data, but can be used to prepare sequences or evaluate already measured chromatograms.

 indicates an Instrument that has been switched-off. The corresponding chromatograph symbol is not displayed in the main Clarity window.

Note: The tab of a switched-off Instrument is accessible only to allow for the removal of modules previously installed on it.

Instrument Type

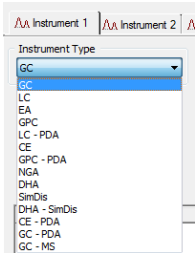


Fig 16: Selection of instrument type

Selects the type of the analysis that will be performed on this **Clarity** Instrument. In addition to standart GC and LC types, optional **Clarity Extensions** provide additional Instrument Types (EA, GPC, LC - PDA, CE and GPC - PDA).

Note: Description of these Instrument types can be found in separate **Clarity Extensions** manuals.

Depending on the selected Instrument Type, the station's main window displays an implicit symbol of a LC or GC chromatograph (if you have not set your own image) and the **List of Equipment** connected to the Instrument (see below) displays only the sections relevant to selected **Instrument Type**.

Name

Holds any arbitrary Instrument description and will be displayed in the main **Clarity** window above the picture of the chromatograph.

Instrument Images

Previews open and closed Instruments. The image is entered by clicking on it. Clicking on the symbol of a closed or opened Instrument will open the **Closed** or **Opened Instrument Image** dialog. From here you can browse for your own image using the  button.

Custom Image checkbox can be used to permit or forbid use of a user's image.

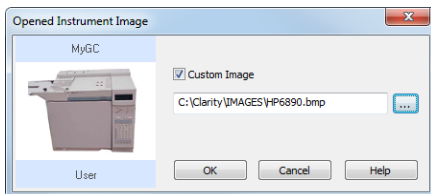


Fig 17: Opened Instrument Image dialog

If the custom image is not available, the preview will display a symbol of missing bitmap and in the main **Clarity** window the default image will be used.

List of the Equipment Connected to the Instrument

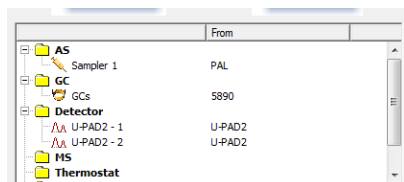


Fig 18: List of equipment connected to an Instrument

Detectors

List of data sources for the **Clarity** station.

GC/LC

Sets direct control of the used chromatograph, detector or pump.

AS

Sets direct control of the used autosampler.

Note: Other categories may appear, including Valves, Fraction Collectors, Thermostats and other miscellaneous directly controlled equipment.

Setting of digital inputs/outputs for external Start/Stop:

	Device	Number
Ext. Start Dig. Input:	Net-PAD	1
Ready Dig. Output:	Net-PAD	1

Fig 19: Digital Outputs for external Start/Stop


If you have more than one A/D converter card installed, select the appropriate card, whose digital input/output you want to use, in the *Acquisition Card* field.

User can define which digital input the selected Instrument will use to start or stop the measuring in the *Number* field of the *Ext. Start Dig. Input* line. Similarly, it can be set which digital output the selected Instrument will use to permit sampling in an active sequence in the *Number* field of the *Ready Dig. Output* line.

The option "--" (*None*) for the *Start In* and *Ready Out* fields is enabled for selected devices. It can be used in the case when the START synchronization is ensured by other means (e.g. with directly controlled autosamplers over digital line).

Instrument Method Sending

The button  invokes the [Instrument Method Sending](#) dialog.

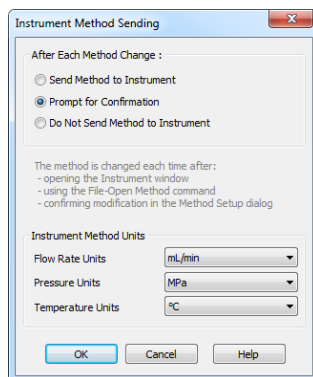

The dialog box titled "Instrument Method Sending" contains the following elements:

- After Each Method Change :**
 - ☐ Send Method to Instrument
 - ☒ Prompt for Confirmation
 - ☐ Do Not Send Method to Instrument
- The method is changed each time after :**
 - opening the Instrument window
 - using the File-Open Method command
 - confirming modification in the Method Setup dialog
- Instrument Method Units**
 - Flow Rate Units: mL/min
 - Pressure Units: MPa
 - Temperature Units: °C
- Buttons: OK, Cancel, Help

Fig 20: Instrument Method Sending

This dialog sets whether the LC/GC/AS method has to be sent to the instrument (device) at each change of the template method and allows the user to select the units for *LC Flow Rate*, *LC Pressure* and *Temperature Units*.

After Each Method Change

The method change occurs when:

- Opening the [Instrument](#) window (which automatically loads template method).

- Loading new template method to the **Instrument** window using the *File - Open* command.
- Changing the currently opened template method (i.e. after editing the **Method Setup** dialog).

There are three possible regimes available:

Send Method to Instrument - The LC/GC/AS method will be sent to the instrument (device) automatically without prompt.

Prompt for Confirmation - At each method change, the **Clarity** station will display the **Send Method Confirmation** dialog in which the sending of the method may be confirmed or suppressed.

Do Not Send Method to Instrument - **Clarity** will not send LC/GC/AS method to instrument (device) on every change.

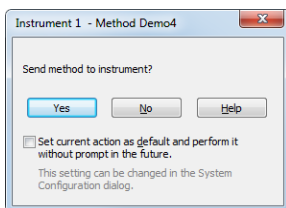


Fig 21: Send Method Confirmation

This setting is individual for each **Clarity** Instrument, but common to all instruments (GC/LC/AS devices) attached to the same **Clarity Instrument**. Regardless of the setting, the LC/GC/AS method will be sent:

- From the running **Active Sequence** according to the **Sequence Table**.
- Manually using the **Send Method** button from the **Single Analysis** dialog.

Instrument Method Units

These settings affect the Gradient Table, tabs of auxiliary pumps, Device Monitor, monitors of auxiliary pumps and tabs of other controlled devices in **Method Setup** and **Chromatogram** windows.

1.1.4 Directories

Sets the working directory ① that contains PROJECTS subdirectory ② with the list of available projects ③. Each project specifies by its name the subdirectory for saving data ④. A working directory is the directory where existing projects are sought and new projects established; it can be set independently for each Instrument. The *Directories...* command or the  icon opens **Instrument Directories for Projects** dialog containing four **Instrument 1 to Instrument 4** fields.

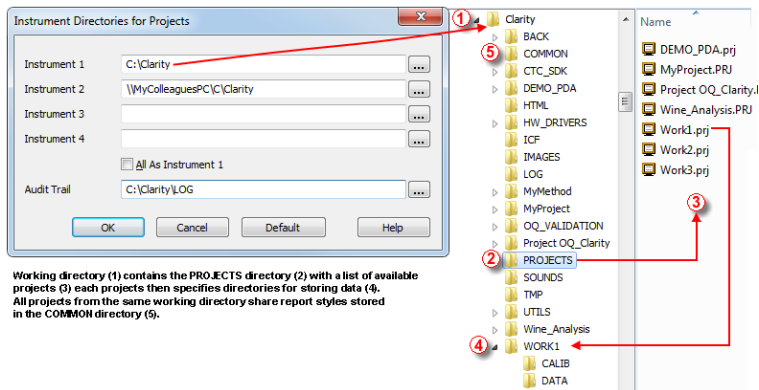


Fig 22: Instrument Directories

Note: In the above image the **Instrument 2** shares projects with another **Clarity** station that has been installed on a networked computer named **MyColleaguesPC**.

Each field specifies the working directory for the corresponding Instrument. If a field is left blank, the working directory will be set to the directory from which the CLARITY.EXE executable file was run (C:\CLARITY by default). This directory may be set any time by the *Default* button.

Use the  button to find the required directory. The *All As Instrument 1* checkbox sets a single directory for all Instruments.

Note: When defining a new directory, user is asked whether he wants to create the directory (nothing with the same name exists) or whether he wants to create the COMMON sub-directory too and fill it with the default files.

The *Audit Trail* line specifies the directory to which the files with all logged activities of the station are stored. If the cell is left blank, the log files will be stored by default in the LOG subdirectory of the main station directory. Pre-set directories are stored in the CLARITY.CFG configuration file, thus being common for all users, even in the protected mode.

Opening of a project on [Instrument](#) is described in the chapter "**Project**" on pg **48**.

In case there is an error during opening a project, [Error Opening Project](#) dialog is displayed.

Caution: If you reinstall or update the **Clarity** station, you should consider that the installer updates files in the COMMON directory located in the **Clarity** installation directory only. It is necessary to update the content of COMMON directories in your custom project directories manually.

1.1.5 GLP Options...

This dialog is used to set the global regulated-environment connected options. After prompting the *System - GLP Options...* command, the user is asked to enter the password for the **Clarity** station. Only users allowed to modify the User Accounts (for more details see the chapter **User Accounts** on pg 3). are allowed to enter this dialog. All changes performed in this dialog are logged into the station audit trail.

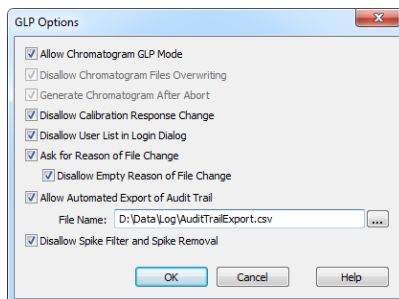


Fig 23: GLP Options dialog

Allow Chromatogram GLP Mode

This checkbox ensures that settings preventing the possibility of losing the data from chromatograms are correctly set. If checked, both the *Disallow Chromatogram Files Overwriting* and *Generate Chromatogram After Abort* checkboxes will be checked.

Disallow Chromatogram Files Overwriting

This checkbox cancels the possibility to overwrite the chromatogram if a file with the same name already exists in the selected directory. The behavior differs according to the action that should cause the chromatogram to be overwritten - when it is a user action (using the *Save As* command for example), the user will be prompted to change the file name to any still unused file name. If the overwriting should be caused by the **Clarity** data acquisition (e.g. creating new chromatograms caused by **Single Run** or **Sequence** measurement), a unique 4- to 6- digit number is added at the end of the original filename. In the improbable case that all of these combinations are already taken, the **Clarity** will ask the user to change the file name personally so it is unique.

Generate Chromatogram After Abort

This feature will cause the generation of the chromatogram in case the analysis is aborted by any reason.

Disallow Calibration Response Change

When checked, the manual change of the response in the [Calibration](#) window is disabled.

Note: The manual change, if allowed, is still marked in the [Calibration Audit Trail](#), but it breaks the link between the calibration standard and the calibration itself and is not recognizable in the chromatogram linked to the calibration file.

Disallow User List in Login Dialog

This checkbox disables the display of all available **User Names** in the [Login](#) dialog. The user is then required to enter two unique identification components to successfully log in.

Ask for Reason of File Change

The *Ask for Reason of File Change* checkbox sets the possibility to fill in the reason for a change. When saving a chromatogram, method, sequence, calibration or GPC calibration file, the [Reason for Chromatogram \(Method, Sequence, ...\) Change](#) dialog will appear. The user can (but is not obliged to) fill in the reason for the change.

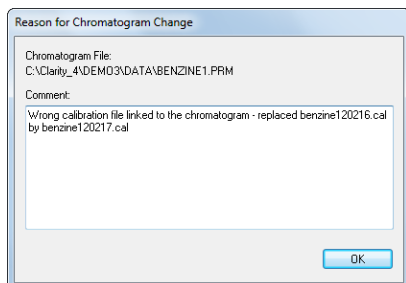


Fig 24: Reason for Chromatogram Change

Note: The reasons of changes are displayed in the **Audit Trails**, separated by a dash from the „file has been saved“ event.

Disallow Empty Reason of File Change

By checking this checkbox the user becomes obliged to fill in the reason of the change. No empty reason field nor reasons solely composed from spaces are now allowed.

Allow Automated Export of Audit Trail

By checking this checkbox the station audit trail file will be automatically exported to the file and location specified in the *File Name* field. To set the appropriate location, use the  button to browse to the desired directory. Changes to this export file will be made continuously and the exported file will be marked as read-only for as long as the **Clarity** is opened.

Disallow Spike Filter and Spike Removal

By checking this checkbox, the *Spike Filter* and *Spike Removal* functions from the [Chromatogram](#) window will be disabled. These are potentially dangerous operations to work with in respect to the GLP practice. The operations are only available in **Clarity** version 4.0 or higher.

1.1.6 Digital Outputs

This dialog, accessed through the *System - Digital Outputs...* command or by  icon, is used to control the digital outputs of the installed A/D boards. These TTL outputs can control, via the **Event Table** (for more details see the chapter **Event Table** on pg 73)., the external measuring devices during the analysis. The command enables one to set up the initial and current state of these outputs by means of the displayed **Digital Output Control** dialog. Any changes performed in the dialog will be saved at the moment of the dialog closing.

Note: If more than one board has been installed, the **Select Control** dialog containing a list of all boards will open first. If a converter has not been installed or is not functional, the *Digital Outputs...* command will not be active (dimmed).

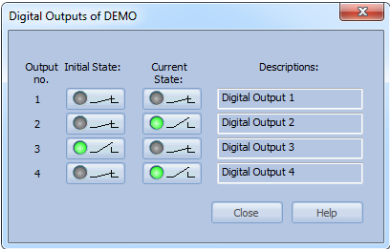


Fig 25: Digital Output Control

Output no.

Lists the numbers of individual outputs.

Initial State

Selects the **Initial** status of the digital output when the station starts. The status is indicated in two halves: first shows the digital output (*Output*) and second its usability as a relay (*Relay*). Where no relay contact can be used, only the state of the output is shown.

Current State

Displays and enables modification of the current status of outputs. The status is indicated in two halves: first shows the digital output (*Output*) and second its usability as a relay (*Relay*). Where no relay contact can be used, only the state of the output is shown.

Change in these columns takes effect immediately. To change the status click the corresponding field with the inscription *HIGH* or *LOW* indicating the output value. The output status is also indicated by color, where gray corresponds to logical zero (*LOW*) and green to logical one (*HIGH*).

Descriptions

Outline description indicating the meaning of individual outputs. The names of individual digital outputs can be edited in the **System**

Configuration dialog for the actual device providing the outputs.

Note: It is not possible to access the **Digital Outputs** dialog of the device that is configured solely on the Instrument you do not have the access to. The access to particular Instruments for different users may be set in the **User Accounts** dialog.

1.1.6.1 Select Control

If more than one board is installed the **Select Control** dialog containing a list of all boards opens first.

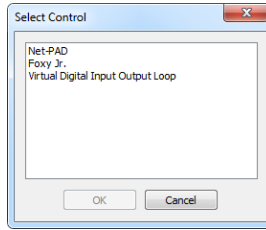


Fig 26: Select Control

Note: If a converter has not been installed or is not functional, the **Digital Outputs** command will not be active (dimmed).

1.2 Login

Submenu for opening an Instrument, which contains a list of all configured Instruments. Opened Instruments are checked.

The **Instrument** window can also be opened by clicking directly on the Instrument's picture in the main **Clarity** window or using a **Ctrl** + **1** shortcut (2, 3, 4 for further Instruments).

If you have an user account, the **Login Dialog** will be displayed after the prompt. Here you select your username from the drop-down list. Choose the user name and fill in the password to log to the **Clarity Instrument**. For heightened security, asterisks ("*") are displayed on your screen when entering a password instead of characters.

Note: After the password has been entered erroneously three times, it is necessary to restart the **Clarity** station.

If the station is in "not restricted" (no password) mode, and therefore no user account exists, the dialog for entering your name will be displayed. The entered name will be saved in all files which you create during your session. You can quickly change the current project or create a new project in *Select Project* combo box. Project offered by default is the last one used by selected user (as saved in user's desktop file) or the last one used in default desktop file (where no user's are created). In case Clarity's command line parameters are used, it is possible to choose project offered through these parameters, either by directly selecting appropriate project or by selecting user with particular project in his desktop file.

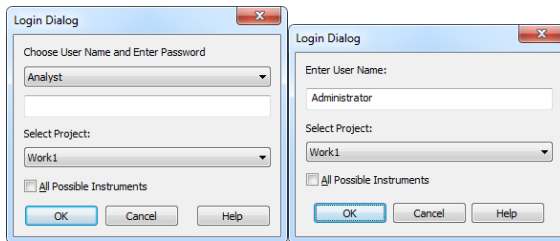


Fig 27: Login dialog and Username dialog

All Possible Instruments

If you have checked *All Possible Instruments* box, then all available Instruments that you have access to will be opened.

Note: A modification of the **Login Dialog** also appears when configuring, setting user accounts or creating electronic signatures.

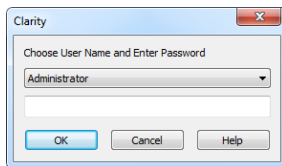


Fig 28: Login dialog when opening System Configuration or User Accounts

Initial password

After receiving your username from the administrator and logging in to the station for the first time, you will receive a verification notice and be prompted to enter your password.

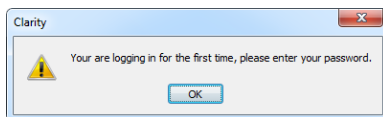


Fig 29: Initial password warning

To prevent keying mistakes during the initial login process, you will be prompted to enter your password two times.

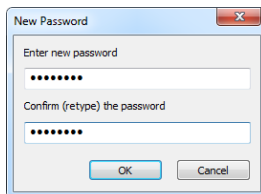


Fig 30: New Password dialog

The administrator can set and control the conditions of a password, allowing for a minimum number of characters and/or the time limit of a password's validity. After a password expires, you will be prompted to enter a new password.

1.3 Logout

This submenu contains a list of all configured **Clarity** Instruments, which can be closed from here. The names of Instruments that are already closed will be dimmed. The user will be notified of an analysis or sequence in process as well as of all modified files that have not yet been saved. Modifications not yet saved will always be marked after the filename by the *MODIFIED* inscription.

Logout All

Closes all opened Instruments at once.

1.4 Lock

This is a submenu used for locking/unlocking open Instruments in the protected mode. This enables the user to guard against an Instrument's unauthorized intervention during his/her absence. Locked Instruments are denoted with a check mark next to its name in the displayed list and by the "lock" symbol before the user's name in the main **Clarity** window. Locking an Instrument affects neither the proceeding data acquisition nor a running sequence. An Instrument can be unlocked by entering the same password used to login to the Instrument in the first place.

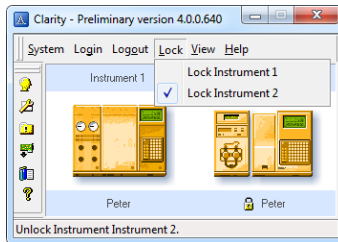


Fig 31: Opened "MyLC" and locked "My GC" Instruments

Note: If the Instrument has been locked by a user, it can be also unlocked by the **Clarity Administrator** (for more details see the chapter **User Accounts** on pg 3). The record of this action is saved into the audit trail and the person that unlocked the Instrument is now logged on it.

1.5 View

This is a submenu for setting up the appearance of the main **Clarity** window. Similar ones can be found in all major **Clarity** windows.

Note: If the main **Clarity** window covers any of the dialogs, the workstation will look as if it has not reacted to any of your commands. In this case it is important to shift the main window aside and react to the covered dialog report.

Switch Scales down the main **Clarity** window. The same effect is achieved by double-clicking on the title bar. To restore the original size, double-click on the title bar once more.

Note: It is advisable to combine this command with the *Always on Top* command.

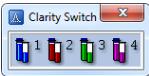


Fig 32: Main window as Clarity Switch

Hide All Minimizes (hides) all open windows.

Show All Displays all open and minimized (hidden) windows.

Toolbar Hides/shows all toolbars with the command icons set by *Customize* command. By using the mouse, it is possible to move or hide toolbars and move the menu bar on the screen.

Note: In case you accidentally move the menu bar out of the screen, you can return it back by using the shortcut **Alt + V**.

Customize... User set toolbars.

Reset All Sets all toolbars and menus back to their initial state - that which the **Clarity** workstation was in when you installed it.

1.5.1 Customize...

Customizes the toolbars.
The **Toolbars** tab lists all created toolbars. To display/hide a toolbar check/uncheck the box next to its name. The *New* command will create a new, empty toolbar; the *Reset* command restores the original setting.

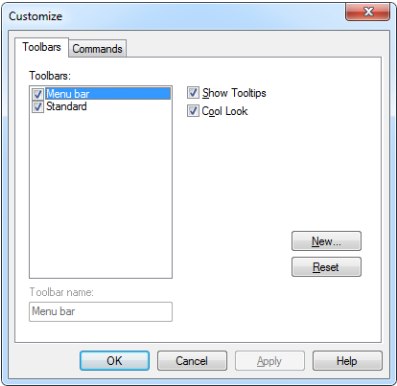


Fig 33: Customize - Toolbars

The **Commands** dialog enables the user to arrange, at will, available icons of individual toolbars. Available icons will be arranged in categories that roughly correspond to menu items. Add/remove an icon to/from the toolbar by dragging it with the mouse.

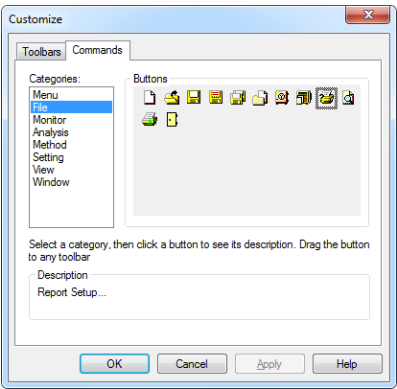


Fig 34: Customize - Commands

Note: Toolbars are set separately for each window and for each user (except the main **Clarity** window). The set of available icons differs from window to window.

1.6 Help

This is an entry point to a submenu of commands that displays fundamental information about the **Clarity** chromatographic station. An analogous menu is available in all basic station windows.


Help	Displays the introductory help item concerning the main Clarity window.
Register...	Allows to register your workstation online.
Check for Updates...	Automatically determines the presence of new versions of Clarity from the Internet.
Languages	Allows you to select the language. You will have to restart Clarity for the change to have an effect.
User Code...	Provides the possibility to enter a user code.
Trial Prolongation...	It opens a dialog that allows you to extend the trial period.
Send Report by Email	Sends information regarding the state of Clarity to the DataApex user help.
Clarity On-Line	Invoking this command opens your default web browser and connects to the DataApex website. Your computer must be connected to the Internet to be able to do so.
About...	Displays information about Clarity .

1.6.1 Register

Online registration of your **Clarity** station. Registration is voluntary and it automatically allows you to download updates from the **DataApex** website directly from the **Clarity** station environment using the *Check for Updates* function.

The selected username and password is also used to gain access to our protected website sections at www.dataapex.com, which are reserved for registered users only. If you have already registered on our website, use this same username and password in the **Clarity Registration** dialog.

In case your computer is not connected to Internet, after pressing *Register...* menu item an off-line registration form will be opened. You have to fill it, print it and send it to us by mail or fax.

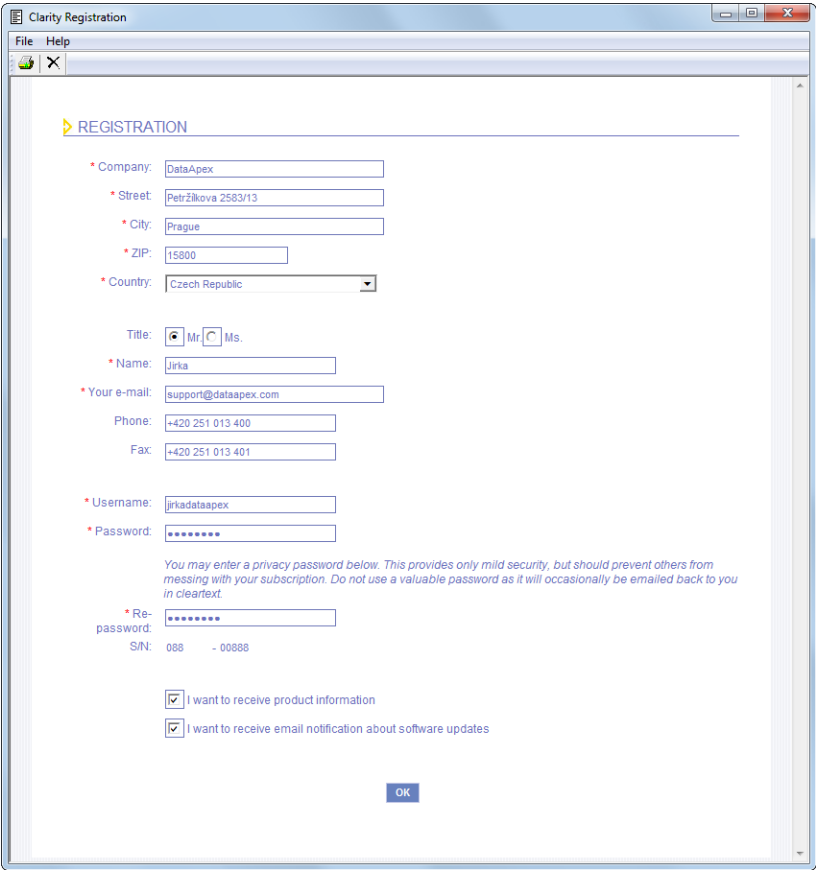

The image shows a 'Clarity Registration' dialog box with a standard Windows-style title bar and menu bar (File, Help). The main area is titled 'REGISTRATION' and contains a form with the following fields and options:

- * Company: DataApex
- * Street: Petržilkova 2583/13
- * City: Prague
- * ZIP: 15800
- * Country: Czech Republic (dropdown menu)
- Title: ☒ Mr. ☐ Ms.
- * Name: Jirka
- * Your e-mail: support@dataapex.com
- Phone: +420 251 013 400
- Fax: +420 251 013 401
- * Username: jirkadataapex
- * Password: (masked with dots)
- You may enter a privacy password below. This provides only mild security, but should prevent others from messing with your subscription. Do not use a valuable password as it will occasionally be emailed back to you in cleartext.*
- * Re-password: (masked with dots)
- SIN: 088 - 00888
- ☒ I want to receive product information
- ☒ I want to receive email notification about software updates
- OK button

Fig 35: Clarity Registration dialog

OK

Sends the registration data to **DataApex** through the Internet. Your computer must be connected to the Internet in order to register successfully.

Note: The next screen allows you to check and correct all the information before pressing the *Send* button.

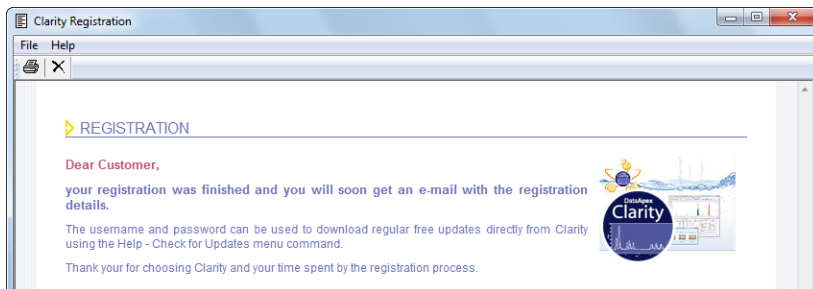


Fig 36: Successful registration

Print

Prints the registration data. Printed registration can be sent by mail or fax.

1.6.1.1 Register Clarity question

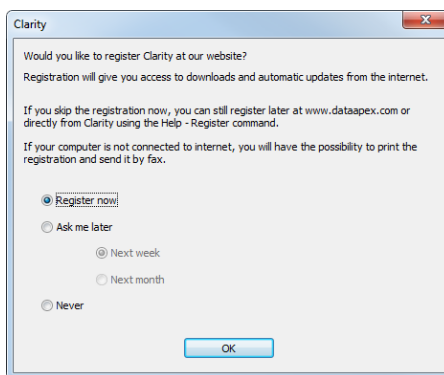


Fig 37: Clarity registration message

This dialog allows you to register your **Clarity** station online at the **DataApex**. Registration is voluntary and it automatically allows you to download updates from the **DataApex** website directly from the **Clarity** station environment using the *Check for Updates* function. Select the *Register now* and click the **OK** button to open the **Clarity Registration** dialog (and proceed with the registration) or skip the registration for a time set by the *Ask me later* option. You will be notified of the registration once the set time has passed. You can select the *Never* option and click the **OK** button to skip the registration whatsoever. The registration may be performed anytime by invoking the *Help - Register* command from the main **Clarity** window. If you have already registered on our website, use this same username and password in the **Clarity Registration** dialog.

1.6.2 Check for Updates

Automatic survey of the newest available updates from the Internet. If your computer is connected to the Internet, the **Clarity** workstation connects to an updating server. The **Check for Updates** dialog opens and displays the phrase “*Connecting to server*”. When a successful connection is made, the program for updating compares the number of your version to the last available version on the server. If you have the newest available version, the message “*No updates available at this time*” will be displayed. If you get the *Failed to connect to server* message, please check your internet connection, for example by opening any internet page in your browser.

Note: Automatic update feature can be used only if you have registered your **Clarity** workstation through the *Register* command.

When the update program determines that your version is older than that available on the server, it will offer you an automatic download of newer version from the Internet.

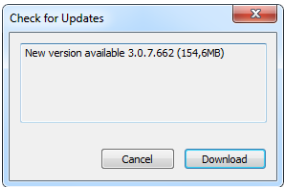


Fig 38: Check for Updates dialog

Confirm the update with the *Download* button.

Note: If the user who performs the download does not posses appropriate rights to install applications in Windows system, a warning message will be displayed. It will offer the possibility to download the update and save it on the local hard disk.

Afterwards, the **AutoUpdate** dialog will appear. Here you can see the rate and progress of downloading of the update file from the Internet.

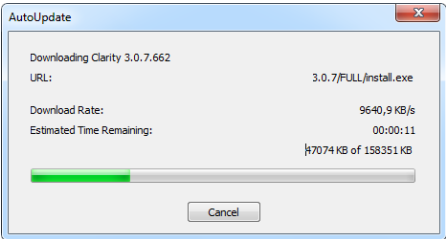


Fig 39: AutoUpdate dialog

While downloading the update from the Internet, you will not be able to use the **Clarity** workstation for measuring and processing of analyses. After completion of the download from the updating program, the **Clarity** workstation automatically switches off and runs an Installation program. This will take you step-by-step through the update of the workstation.

1.6.3 User Code...

Enters, displays and edits the **User Code**. The user code is the 16-character encoded form of the Serial Number of your **Clarity** station.

Note: The **Serial Number**(S/N) is a unique number with the basic structure AAA-BBBBB+CCC/DD, which must correspond to the number in the protective hardware key. It can be found on the installation DVD-ROM or in the [About](#) dialog in **Clarity**.

The **User Code** is saved in the CLARITY.SNO text file in the root folder of the **Clarity** workstation (implicitly C:\CLARITY).

The **User Code** also defines the number of **Clarity Instruments** available for measurement and which **Control Modules** you are allowed to use.

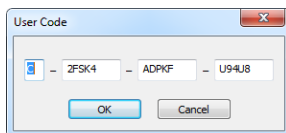


Fig 40: User Code dialog

Trial Mode

If you do not enter the **User Code** (accomplished by hitting *Skip* in the [Registration](#) dialog during the installation), **Clarity** will be able to run 30 days or 100 starts (whatever comes first) in **Trial Mode**. **Trial Mode** is indicated in the header of the main [Clarity](#) window.

1.6.3.1 Language Selection

This dialog is displayed in case the **Clarity** is set to a language which is not available for your licence (User Code). Please, contact **DataApex** to consult the languages availability.

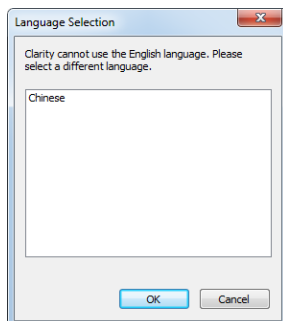


Fig 41: Language Selection dialog

After you select one of the listed languages, restart of the **Clarity** Chromatography Station is performed to change the environment into selected language.

1.6.4 Send Report by E-Mail

Sends the information on the state of the **Clarity** workstation to the user help desk of **DataApex**. Use this command in the case that you have problems with the functionality of the **Clarity** workstation. The command creates a new e-mail message addressed to our technical support.

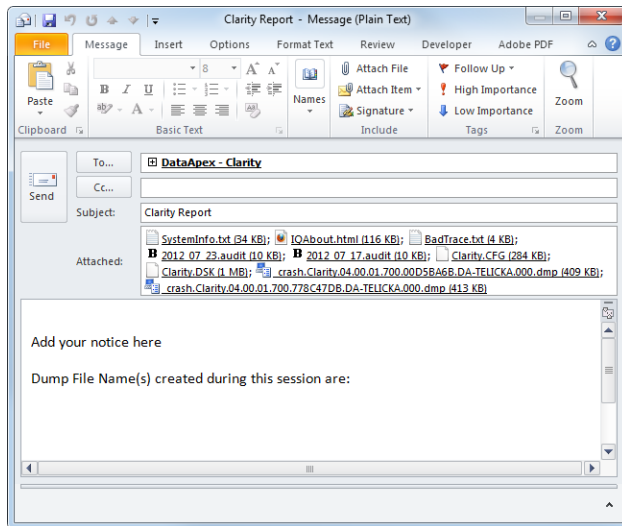


Fig 42: E-mail message with information about Clarity

In the appendices are the SYSTEMINFO.TXT, BADTRACE.TXT files, current desktop, current configuration and two last files of the station audit trail. All these files contain information describing the state of the **Clarity** workstation at the time when the problem(s) occurred.

Note: The SYSTEMINFO.TXT and BADTRACE.TXT files can be found in the root folder of the workstation (implicitly C:\CLARITY). The content of the SYSTEMINFO.TXT file is also displayed in the [About - System](#) dialog. You may read through the files of the log in the [Audit Trail](#) window by using the [Open](#) command (for more details see also chapter [Audit Trail](#) on pg 318).

Should the work with the **Clarity** workstation have been terminated improperly, an offer to send an e-mail with the error description will automatically be displayed on the next start up. The data will be send under no circumstances until you approve it.

Note: You can deactivate the automatic offer to send the error description by checking the *Send Reports About Unsuccessfully Finished Sessions* checkbox on the [General](#) tab of the [User Options](#) dialog described on pg 104.

1.6.5 About...

Displays the version number and date of the release, serial number and other system information.

Clarity tab

Displays the name of the release followed by its topical number in the format W.X.Y.Z. The compilation date will be prefaced by *Build*. The *S/N* abbreviation prefaced the serial number of your station in the format XXX-YYYYY+ZZZ/NN.

What's New tab

The tab shows the contents of the WHATSNEW.TXT text file from the main directory of the **Clarity** station. The file contains a chronological list of all changes made in each new version of the station.

Readme tab

Displays a brief description of the station.

System Files tab

The tab shows the contents of the IQ.HTML text file from the main directory of the **Clarity** station. This file presents fundamental information about both the station and the computer and is updated each time the station is run. If you encounter any problems, send the file to your vendor or directly to **DataApex**. This tab also displays the number of Instruments allowed and lists available control modules.

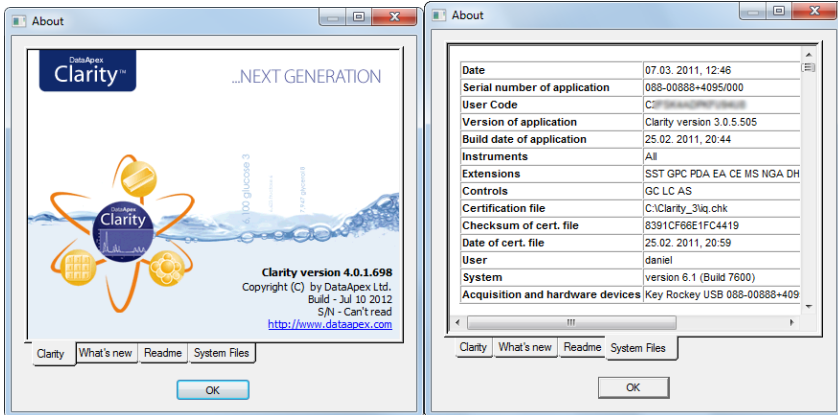


Fig 43: About - Clarity & System Files

2 Instrument

Caution: In **Clarity Lite**, this window is also called the (Main) **Clarity Lite** window and supplies some of the properties of **Clarity** window. Any differences between **Clarity Instrument** window and **Clarity Lite** window will be discussed later, most notably on pgs. 41., 42. and 117.

This chapter describes the commands of the **Instrument** window - the fundamental window for chromatogram measurement and evaluation. The station allows for up to four **Instrument** windows to be displayed at one time. The number of displayed Instruments may exceed the number of Instruments actually purchased, but the excess Instruments may not be used for data acquisition.

Note: Users commonly utilize these so called **Offline Instruments** for method development or evaluation of already measured analyses.

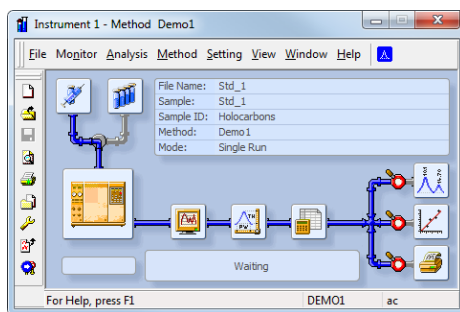


Fig 44: Instrument

The **Instrument** window displays a logical sequence of operations and commands necessary for the comprehensive processing of an analysis with icons. This **Scheme of Data Processing** starts from a sample injection taken and finishes with a printout of the results. Each Instrument displays the scheme in a dedicated color. Each icon corresponds to one menu command. In addition to the menu and the scheme of data processing the window contains the **Information Table** providing information about the measured chromatogram and **Status Line** providing information about elapsed time and current status of the measurement.

Note: If the instrument is in the **Control Time**, the time is displayed as a negative number, counting up to zero where the **Control Time** will end and a new analysis (if set) will begin.

The lower end of the **Instrument** window then holds the information on how to invoke the context help and the information on the project opened and the name of the **User** who is logged to the **Instrument**.

To open or display an Instrument, invoke the command from the *Login* submenu or click on its icon in the main **Clarity** window. The Instrument name is taken from the configuration, where you can set it (for more details see the chapter **Right section of System Configuration dialog** on pg 17).

Scheme of Data Processing

The icons from the scheme serve for easy access to all dialogs necessary for the management of an analysis.

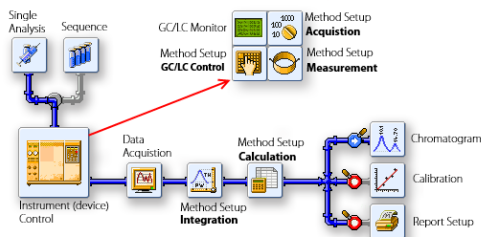


Fig 45: Scheme of Data Processing

When the sequence is running, the **PostRun** checkboxes (, ) will be locked to the state corresponding to the setting on the currently running row (for more details see the chapter **Postrun...** on pg 100).

2.1 Clarity Lite window specifics

In the **Clarity Lite** version, after starting the software, the dialog for entering your name will be displayed. The entered name will be saved in all files which you create during your session.

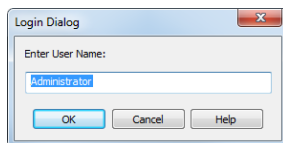


Fig 46: Username dialog

Then the **Clarity Lite** window (in all important respects same as **Instrument** window) will open.

2.2 File

An entry point to a submenu of commands for working with system files, setting up the printer and exiting an Instrument.

	New Method	Creates a new method. When invoked a Save As dialog is invoked and the user has to enter the file name of the new method. User will be notified if the previously opened method was modified and not yet saved. The modification of the method is indicated in the title of the Instrument window next to the filename by the "(MODIFIED)" inscription. No file may be saved under the name NONAME.
	Open Method	Selects and open an existing method.
	Save Method	Saves the method. If a new file is to be saved, this command has no effect. In this case use <i>Save Method As</i> command instead. The default name NONAME is reserved for new, as yet not saved files.
	Save Method As	Saves a method under a selected name and/or to another directory. The command opens the Save As dialog (similar to the Open Method dialog). Here a new name can be entered or a new directory selected. If a file of a particular name already exists and has been read in some other Instrument or window, the name cannot be used. Moreover, in the protected mode (full version  only) a user may not overwrite a file he/she is not entitled to modify (see the <i>User Account</i> command).
	Send Method by Email	Sends the displayed method file as an email attachment.
	Project...	Selects project directory.
	Archive... 	Backs up files and projects of Clarity station.
	Restore... 	Restores files and projects of Clarity station.
	Report Setup...	Selects a report style. This command opens the Report Setup dialog (for more details see also chapter Report Setup on pg 333).

 Print Preview	Displays a preview of the printout prior to actual printing. This command opens the Print Preview dialog (for more details see also chapter Print Preview on pg 352).
Print to PDF	Prints report to PDF file.
Send Printed PDF by Email	Opens new email message in default email client, prints to PDF and adds created PDF file as an email attachment.
 Print...	Selects a printer and start printing. Invoking this command opens the Print dialog (for more details see also chapter Print on pg 354).
-----	Lists recently opened files so that you can easily access method files you have been currently working with. The number of files in the list is determined by the <i>Recent Files</i> parameter in the User Options - General dialog available from the Instrument window through the <i>View - Options</i> command.
Exit	Terminates work with an Instrument and closes all its windows. A warning is issued if an analysis is in progress. The user is also notified of modified files that are unsaved. In the full version , an Instrument can also be closed by using the <i>Logout</i> command from the main Clarity window or gained by right-clicking on its icon.

2.2.1 Open Method

Selects and opens a method. When invoked the opened [Open Method](#) dialog enables the user to select the desired method. The method will then be loaded and its name displayed in the **Information Table** and in the [Instrument](#) window title bar. The user is notified if a previous modified method has yet to be saved. Unsaved modifications are indicated by the inscription MODIFIED next to the filename. To select another method click the  icon.

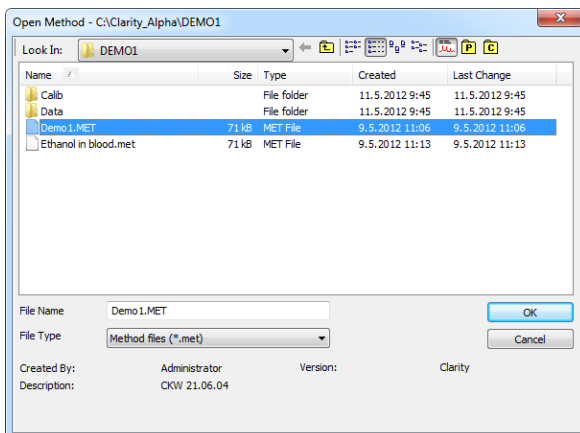


Fig 47: Open Method

Note: When the mouse cursor is placed over the title of the template method in the **Information Table**, the yellow box (tooltip) will open, displaying information about whether a measurement is currently taking place according to this method.

The template method name is displayed both in the **Information Table** and in the window title bar. The station enables measurement according to template method while simultaneously editing another template method (current instrument method). The information table shows the name of the template method used in the proceeding analysis, while the [Instrument](#) window title bar indicates the method currently being displayed and, accordingly, editable.

From this description it follows that until a measurement begins, the current Instrument method and template method will be the same.

Note: To display the method used in the measurement in progress, invoke the *Method* command from the [Single Analysis](#) dialog or the *Sequence - Edit Current Method* command from the [Sequence](#) window.

Note: If a measurement is not taking place, you may open the method by clicking on its name in the **Information Table**.

2.2.1.1 Open Method (Sequence, Chromatogram, Calibration)

This is a universal dialog for selecting files. As it differs somewhat in appearance from the standard window for file selection of the **Windows** operating system, it is described in more detail below:

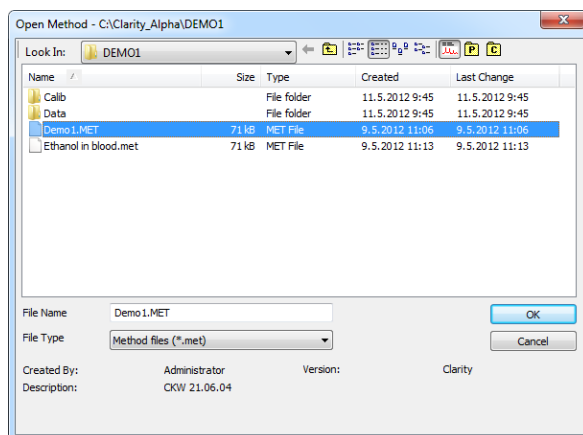


Fig 48: Open Method

List of files

Displays all files of the selected *File Type* from the previously opened directory.

The icon displays files from the current project. The icon displays files from the COMMON directory.

The icon returns you to the previous directory.

The icon transfers you up to the parent directory.

With the icon found in the *Look In* combo box you may unroll the tree structure of the disks of your computer and quickly select any folder in that structure.

Note: When selecting a chromatogram the (yellow) and (red color) icons enable one to switch to the data or calibration subdirectory of the current project.

Files are sorted and displayed in conformity to their previous setting. The layout can be changed by using the button to access the *Detailed View* menu. In the *Detailed View* layout you can change the sorting order by clicking on the heading of the desired column. A second click reverses the order. The type and method of file sorting is indicated by a small arrow next to the column name.

The mode in which files are displayed and the ordering method are automatically saved in the current desktop file (*.DSK).

Note: In the non-protected mode all settings are the same for the entire station. However, in the protected mode each user may have a dedicated *.DSK file.

File Name

This field contains the name of the selected file. The name can be entered directly or selected from the list of filenames. When the name is being entered directly, only files whose names begin with the characters already entered will be displayed. The files and directories that match the so-far entered letters in the given actual directory will be shown in the listbox on the *File Name* row.

Note: If you search for files containing only some characters, enter an asterisk (*) - which stands for any number of characters) or several question marks (?) - substitutes for a single character) before those characters.

Use the  button to display/hide the following **detailed information** about the selected file:

Created by

Indicates the name of the user who created the file. Each file contains the name of the user who created it by use of either the *New* or *Save As* commands.

Description

A supplementary file description that enables to fill in or update the information in the corresponding editing window of the system file (this field serves for information only).

Version

Displays the version of the station in the format from which the file was last saved.

2.2.2 Send Method by Email

Sends the displayed method file as an email attachment. The command will create a new email message in your default e-mail client.

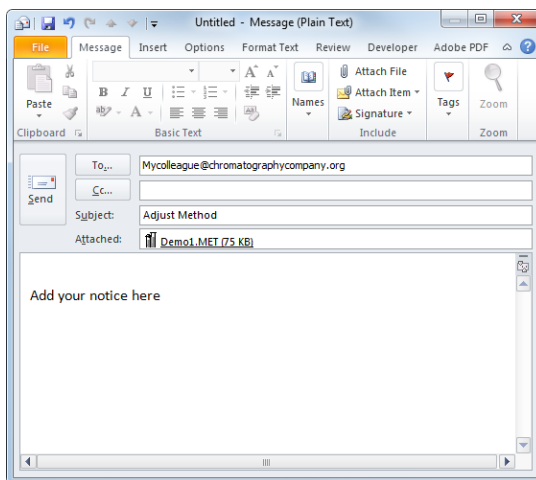


Fig 49: Send method by email

Caution: The proper function of this command is conditional upon having a working email account.

The command is used in all major **Clarity** windows and serves for sending the respective files handled in these windows (e.g. *Send Calibration by Email*, *Send Chromatogram by Email*, etc.)

2.2.3 Project

Use the *Project...* command (icon ) when working with project directories.

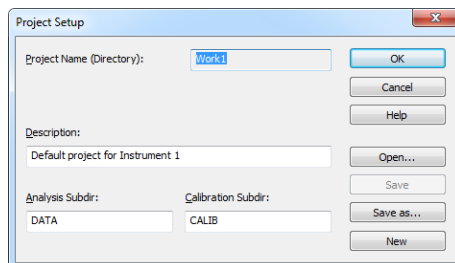


Fig 50: Project Setup

Invoking this command will open the **Project Setup** dialog. Here the setup of the current project can be updated or another project opened or created.

Project Name (Directory)

Contains both the project name and the project directory. Only methods and sequences will be stored in the project directory (chromatograms and calibration files are stored in the corresponding subdirectories).

Description

Description of the project (this field serves for information only).

Analysis Subdir

Subdirectory for storing data (chromatograms).

Calibration Subdir

Stores calibration standards and calibration files.

Note: More information on the project directory structure can be found in the chapter "**Directories**" on pg 21.

Open...

Selects a new project from a list of existing ones. In the full version , the same project can be opened in two different Instruments. This configuration can be useful, for example, when preparing a new sequence during measurement.

Caution: Although the same **project** can be opened in two **Clarity** Instruments within one station, the same **file** cannot. However, with multiple **Clarity** (Lite) stations running in a PC network, this limitation does not apply and the same file can be displayed on any number of stations.

Note: The list of existing projects corresponds to the list of *.PRJ files in the PROJECTS directory.

Save

Saves the displayed project parameters.

Save as

Copies the current project and saves it under another name. The project description and the name of data subdirectories are copied automatically. Copies of all files are created only after confirmation.

New

Creates a new project. Invoking the command displays a list of existing projects where the name of the new project can be entered. Two data subdirectories are offered with the default names DATA and CALIB.

2.2.3.1 Error Opening Project

This error dialog is displayed if there is an error during opening a project, it also allows you to decide, what will **Clarity** do in this type of the situation.

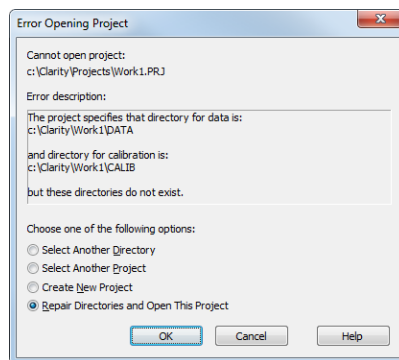


Fig 51: Error Opening Project

Select one of the option below and click the **OK** to perform the action and continue.

Select Another Directory

Set this option in case the directory with your projects has moved.

Select Another Project

Allows to select another project within the projects directory.

Create New Project

Allows to create new project in existing the projects directory.

Repair Directories and Open This Project

Allows to create required subdirectories in the project's directory.

2.2.4 Archive ✓ full version

Archives or erases selected files or entire projects in the **Clarity** station. The command opens the **Backup - Create Archive** dialog. Files from the current project can be backed up to any directory on any disk either as simple copies of individual files or as a single compressed *.DGZ file.

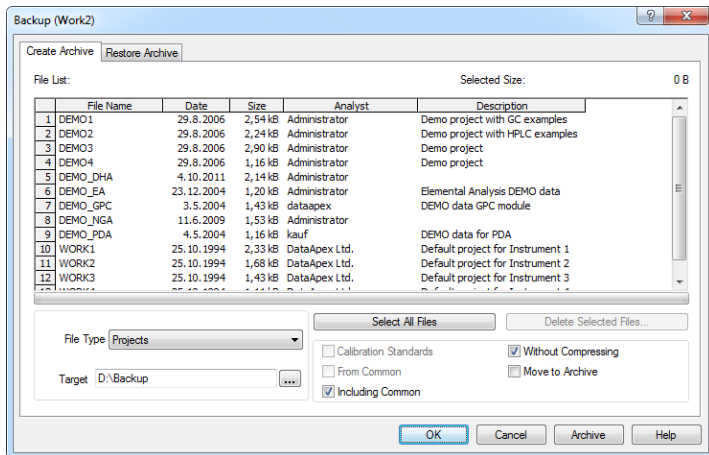


Fig 52: Backup - Create Archive

File list

Lists all selected files from the current project. Click on files to select them, click again to deselect the file. Any number of the displayed files can thus be selected. In addition to the filenames, the list contains the date when it was last saved, file size, and the name of the user who created the file. *Sample ID* and *Sample* Columns will be shown for chromatograms. *Description* will be displayed for the other files and projects.

Selected Size

Indicates the size of the selected files (in kB).

File Type

Selects the type of file: *Chromatograms*, *Sequence files*, *Method Templates*, *Calibration Files*, *Report Styles*, *SST Method files*, *GPC Calibration files*, *PDA Library files*, *Projects* (meaning all files of a given project), *Audit Trail Files*, *PetroCal Files* and *NGA Method Files*.

Note: Archiving the whole project will also archive all files from the common directory (C:\COMMON by default) that contains for example report styles.

Select All Files

Rapidly selects all files in a list.

Delete Selected Files

Deletes selected files. Once deleted, files cannot be restored from within the **Clarity** station!

Calibration Standards

Determines whether chromatograms from the DATA or from the CALIB directory will be listed.

Without compressing

If checked, the selected files will be copied or moved, otherwise a single compressed *.DGZ file will be created.

Note: Previously existing files of identical names will be rewritten and not extended!

From Common

Displays system files only from the common directory COMMON. This checkbox is active only for system files.

Move to Archive

After a successful backup, the files will be deleted from the current directory.

Including Common

While this checkbox is checked, the COMMON subdirectory will be also archived with other files of the archived project. This option can only be checked while archiving whole projects.

Target

Defines the target directory for copying the files.

OK

Saves any modifications made in the archive settings and closes the [Backup](#) dialog.

Cancel

Closes the [Backup](#) dialog without saving any of the modifications made in the archive settings.

Archive

The proper command for backing up (copying or moving) selected files. A warning is given if files with identical names exist in the target directory.

2.2.5 Restore ✓ full version

Restores or deletes files and projects in the **Clarity** station. This command invokes the **Backup - Restore Archive** dialog. The *Restore* command restores files from any directory to the current project.

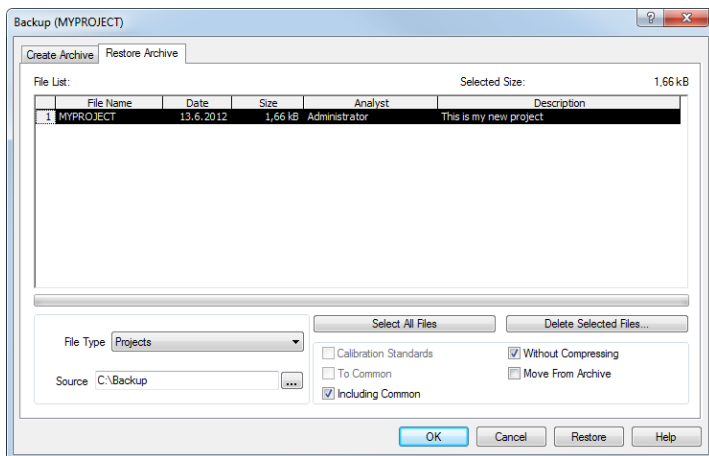


Fig 53: Backup - Restore Archive

File list

All files of the selected type from the directory selected by *Source* command. Files are selected by clicking once and deselected by clicking a second time allowing for any number of displayed files to be selected this way. The list that is displayed contains file names, size, the date on which they were last saved, and the name of the user who created the file. The *Sample ID* and *Sample* columns will be shown for chromatograms. The *Description* column will be displayed for other files and projects.

Selected Size

Indicates the size of the selected files (in kB).

File Type

Selects the type of file: *Chromatograms*, *Sequence files*, *Method Templates*, *Calibration Files*, *Report Styles*, *SST Method files*, *GPC Calibration files*, *PDA Library files*, *Projects* (meaning all files of a given project) and *Audit Trail Files*.

Note: Restoring the whole project will also restore all files to the common directory (C:\COMMON by default) that contains for example report styles.

Select All Files

Rapidly selects all files in a list.

Delete Selected Files

Deletes selected files. Once deleted, files cannot be restored from within the **Clarity** station!

Calibration Standards

Determines whether the chromatograms will be restored to the DATA or CALIB directory.

Without compressing

If checked, the selected files will be simply copied or moved, otherwise the content of a single compressed *.DGZ file will be displayed.

To Common

Restores system files to the COMMON directory. This checkbox is active only for system files.

Move from Archive

Files successfully restored will be deleted (the *Restore* command functions similarly to the *Move* function).

Including Common

While this checkbox is checked, the project will be restored including the COMMON subdirectory (if it is included in it). The option is only available for restoring whole projects. Note that files restored from the project will try to overwrite the actual content of the COMMON subdirectory.

Source

Identifies the source directory (of the *.DGZ file) where selected files will be sought.

OK

Saves any modifications made in the settings of the **Backup - Restore** dialog and closes the **Backup** dialog.

Cancel

Closes the **Backup** dialog without saving any of the modifications made in the settings of the **Backup - Restore** dialog.

Restore

The proper command for restoring (copying or moving) the selected files. A warning dialog is shown if files with identical names exist in the target directory.

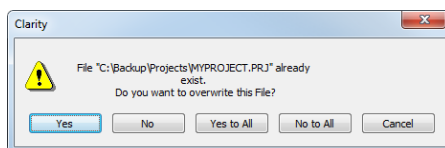


Fig 54: Warning message when the filename already exists

2.3 Monitor

An entry point to a submenu of commands for displaying the signal from connected detectors or monitoring the status of a directly controlled chromatograph.



Data Acquisition

Displays the signal of the connected detectors. Displays the signals of connected detectors. After selection, the [Data Acquisition](#) dialog described on pg 279. appears. You may also select this dialog by using the  icon located in the [Instrument](#) window.



Device Monitor

Monitors chromatograph status.

2.3.1 Device Monitor

Displays the current status of a directly controlled chromatograph and other controlled modules. Invoking this command opens the [Device Monitor](#) window, which accommodates monitors from multiple devices.

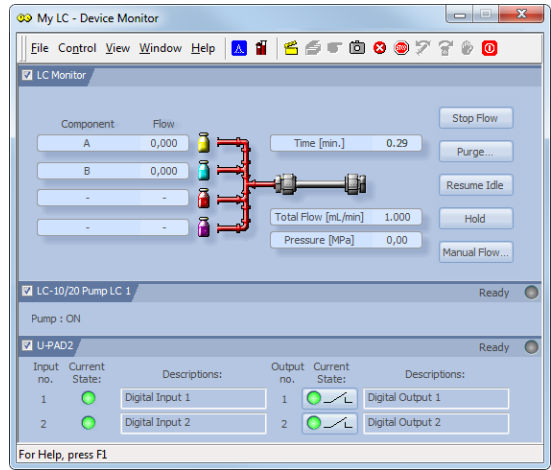


Fig 55: Device Monitor

Menu

File	Submenu for controlling the Device Monitor window.
Control	Submenu for controlling the device.
View	Submenu of commands for setting up the appearance of the Device Monitor window.
Window	Submenu containing commands governing the appearance of base windows.
Help	Submenu of commands that displays help.

Toolbar

The toolbar includes buttons that can be used for controlling the analysis process (*Run Single*, *Run Sequence*, *Stop*, *Abort*, *Shutdown*, etc.) (for the description of these functions see the chapter "**Single**" on pg 59. and the chapter "**Sequence**" on pg 294.). Additionally the (*Stop Flow*, *Purge...*, *Resume Idle*, *Hold* and *Manual Flow...*) functional buttons will be available in selected monitors after the appropriate control module has been installed and configured in the [System Configuration](#) window. A

detailed description of the particular sections of the [Device Monitor](#) window can be found in corresponding manuals for Control modules.

Individual panes - monitors

Context menu in [Device Monitor](#) window (right mouseclick) enables to change the order of individual device monitor panes.

Individual monitor panes can be collapsed/expanded using the checkbox in the header of the pane (left to the Device Name). In the collapsed form only the state is displayed.

Note: Availability of this command depends on the capabilities of pumps used in the gradient. The command is only available during the analysis run.

Note: Availability of this command depends on the capabilities of pumps used in the gradient. The command is only available during the analysis run.

2.3.1.1 File

Close	Closes the Device Monitor window.
--------------	---

2.3.1.2 Control

Shut Down	Function is individual for each device. Detailed description is in a separate manuals for Control modules.
Stop Flow	Function is individual for each device. Detailed description is in a separate manuals for Control modules.
Purge...	Command for flushing/purging the pump outside a run.
Resume Idle	Resumes back to <i>IDLE</i> state after flushing/purging the pump using the <i>Purge</i> command.
Hold	Pauses the execution of the Gradient Table at the actual flow and mobile phase composition. The same flow and composition will be maintained until the <i>Resume</i> button is pressed.
Resume	Resumes the execution of the Gradient Table after the <i>Hold</i> button has been pressed.
Manual Flow...	Opens the LC Control Manual Flow dialog allowing to set custom flow and mobile phase composition, disregarding the Gradient Table set in the method.

2.3.1.3 View

Fit to Content	Resizes the window according to its content.
Toolbar	Displays/hides all toolbars. The command is analogous to that shown in the Clarity window (for more details see also chapter View on pg 30).
Customize...	User customized toolbars. The command is analogous to that shown in the Clarity window (for more details see also chapter Customize... on pg 31).
Component Flow in %	Display the component Flow in % of the <i>Total flow</i> instead of the volume units.
Reset All	Returns all menus and toolbars in this window into their initial state.

2.3.1.4 Window

Main	Displays the main Clarity window (for more details see also chapter Clarity on pg 1).
Instrument	Displays the Instrument window (for more details see also chapter Instrument on pg 40).

2.3.1.5 Help

Help	Displays the help item concerning the Device Monitor window.
-------------	--

2.4 Analysis

This is a submenu containing commands for measurement control:

	Single	Prepares and measures a single analysis.
	Sequence	Prepares and measures a series of analyses.
	Batch ✓ Full version	Processes previously performed analyses.

2.4.1 Single

Prepares and runs a single analysis. Invoking this command opens the [Single Analysis](#) dialog, which can also be opened using the  icon in the [Instrument](#) window. The dialog provides commands for controlling a measurement and its fundamental description. The name of the current method is displayed in the title bar.

Note: If the [Single Analysis](#) dialog contains other fields than described here, then you are probably using one of the **Extensions** (such as **GPC** or **EA**), refer to the corresponding manual.

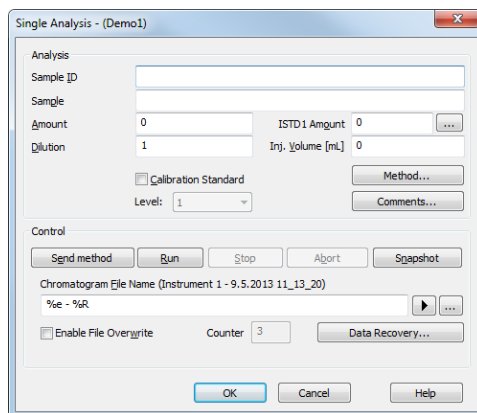


Fig 56: Single Analysis

Analysis

Comprises the description and parameters of an analysis. All entered values will be transferred to the chromatogram header after completion of the analysis.

Sample ID

Sample identification field used for information purposes only. Maximum number of characters in this field is limited to 64. Variables that can be used are described in the *Chromatogram File Name* field (for more details see the chapter **Chromatogram File Name** on pg 64).

Sample

Sample description used for information purposes only. Maximum number of characters in this field is limited to 64. Variables that can be used are described in the *Chromatogram File Name* field (for more details see the chapter **Chromatogram File Name** on pg 64).

Amount

Sample amount. The value is used in calibration calculations of individual components percentage in the *ESTD* and *ISTD* methods. In *ISTD* enter the

sample amount without the addition. Enter values using the same units that were used in the calibration file (*Units - Compound* parameter in [Calibration Options](#) window). The default value is 0.

ISTD Amount

An added amount of the internal standards for the ISTD method - the amount of the ISTD1 can be directly edited in the [Single Analysis](#) dialog, the amounts for ISTD2 - ISTD5 can be set using the [Edit ISTD Amounts](#) dialog accessible through the  button next to it. Always state this parameter in the same units that are used in the calibration file (*Units - Compound* parameter in [Calibration Options](#) window). The default value is 0.

Dilution

A sample dilution factor. This value is used to multiply each calibrated amount in the *Amount* column of the *Result Table*. The default value is 1.

Inj. Volume

An injected volume of the sample. The responses of all compounds will be corrected by the ratio of this value and the *Default Injected Volume* parameter from the calibration file in recalibration or calibration calculations. The default value is 0.

Calibration Standard

Marks the sample being analyzed as a calibration standard. The resulting chromatograms will be stored in the calibration subdirectory. This option cannot be changed during measurement.

Level

Sets the calibration level for calibration or recalibration. This field is active only when *Calibration Standard* checkbox is selected.

Note: If you select "None" then resulting chromatogram will not be used for calibration (it will only be copied to calibration subdirectory).

Method...

Used to display and edit the current method. Unless an analysis is in progress the method set in the [Instrument](#) window will be displayed. During analysis another method may be read into the [Instrument](#) window. Parameters of the current analysis can only be changed here.

Note: If you edit the current method or open a new one on the [Instrument](#) while the **Single Run** analysis is running, you will be prompted to send the newly opened method to the controlled devices after the run is complete.

Comments...

Opens the [Rich Text Edit](#) dialog, which serves for adding some information about the chromatogram created by current method.

Control

Section with commands and parameters for controlling an analysis.

Send Method

Sends actual method parameters to controlled modules, regardless on the setting of the LC/GC/AS method.

Note: The *Send Method* button is visible only when the instrument is configured with a control module.

Run

Runs an analysis. Analysis in progress is indicated by the inscription **RUNNING** in the **Status Line** below the  icon. The parameters of the current measurement method in the [Method Setup - Acquisition](#) dialog cannot be changed while an analysis is in progress.

Invoking the command during analysis is possible only when *External Start* setting in [Method Setup - Measurement](#) dialog is set to *Start-Restart*. Invoking the command terminates running analysis and at the same time starts another one - this operation procedure is often called continuous measurement.

If you enter the *Chromatogram File Name* including a subfolder, for example *%B_%Y%e - %R%e - %R*, then the [Clarity - Directories](#) dialog will be displayed.

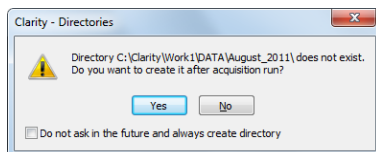


Fig 57: Clarity - Directories dialog when the subdirectory doesn't exist

Click **Yes** if you want to allow Clarity to create the subfolder after the acquisition is finished. If you press **No**, then you must delete the directory and the backslash from the *Chromatogram File Name*. In the dialog you can see that it is also possible to use the parameters in the subfolder name as in the chromatogram name.

Note: Invoking the command during analysis is possible only when there are only analog detectors (INT7, Net-PAD,...) configured on Instrument. If there is one or more digital detectors on Instrument, invoking the command during analysis is not possible.

Stop

Terminates the running analysis and processes the results - a chromatogram is created with the name entered in the *Chromatogram File Name* field. The method used in its processing is the measurement method (its name is found in the **Information Table**). Events following the completion of an analysis are specified by items from the [PostRun Setting](#)

dialog (some of them correspond to the  and  icons in the [Instrument](#) window). The Instrument returns to the *STOP* or *WAITING* state. While the program controlling the chromatograph is running, *CONTROL* is displayed.

Abort

Aborts an analysis without storing or processing the data. The instrument will return to either the *STOP* or *WAITING* state, the chromatograph controlling program, if any, is aborted.

Snapshot

Processes the existing part of the analysis without its termination. A chromatogram is created using the name specified in *Chromatogram File Name* field. The processing mode is the same as with the *Stop* command, the Instrument, however, remains *RUNNING*. If you are about to replace an already existing file, you will receive a warning, but only upon the first use.

Chromatogram File Name

Sample name for created chromatograms. The name can compose of fixed expressions, as well as variable parameters (e.g. adding time, date and/or sample ID and many other parameters). Fixed part may be combined with the variable part in a single chromatogram name.

After clicking the  button a list of available parameters and a brief description of their functions is displayed. You may stepwise compose the title of the file by selecting the required variables (for more details see the chapter **Chromatogram File Name** on pg 64).

It is possible to include also the name of a subfolder in the name to store chromatograms into subfolders of the [Data Subdirectory](#). E.g. the *MySubProjectMyFileName* in the DEMO1 project will result in the MYFILENAME.PRM stored in the C:\CLARITYDEMO1\DATA\MYSUBPROJECT folder. If the subfolder doesn't exist, then it will be created when the chromatogram will be stored.

Note: *Chromatogram File Name* field content can be changed when the analysis is running and will influence the name of its resulting chromatogram, when the change is accepted by pressing the OK button afterwards.



Selects a chromatogram location that is different from the current project. After clicking on this button the [Save As](#) dialog will appear. Here you may select a folder for saving the chromatogram.

Note: It is not particularly suitable to save chromatograms outside of their project folder.

This button is most often used for determining the titles of previously existing files in the folder and helps to avoid the use of identical titles. It is also possible to create a new file on the basis of an already existing title.

Enable File Overwrite

Enables the overwrite of existing chromatograms without giving notice. It is necessary to close the [Single Analysis](#) dialog and open it again for this setting to take effect.

In case the file overwriting is not allowed and the chromatogram with the same name already exists, a chromatogram with the name 00000000.PRM is created instead. The number for such chromatograms is then incremented in case the necessary file name is already taken.

Counter

Shows the current value of the sample counter. If automatic numbering is used for naming chromatograms in the %n format, the current value may be changed here.

Data Recovery...

Rescues data from an interrupted analysis. If an analysis has been interrupted for any reason (power failure, computer breakdown, invocation of the *Abort* command, etc.) a portion of the measured chromatogram may still be rescued.

Invoking the command opens the [Rescue Data](#) dialog where you select from the *LAST.RAW and *RUN.RAW files. If no data can be rescued, the items will be dimmed.

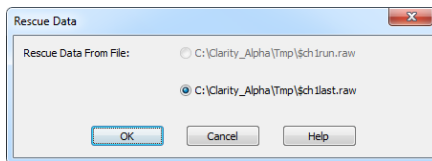


Fig 58: Rescue Data

The *LAST.RAW and *RUN.RAW files are stored in the TMP subdirectory of the main station directory (C:\CLARITY\TMP by default). The *RUN.RAW file will be created only when **Clarity** station has crashed while measuring data (e.g. after a power failure). Otherwise the last measured chromatogram will always be saved to the *LAST.RAW file. Pressing the [OK](#) button invokes the [Save As](#) dialog. During an analysis data is saved to the hard disk every minute.

2.4.1.1 Chromatogram File Name

This field contains a template for creating chromatogram names. Up to 254 characters may be entered including spaces and international characters, the following characters are, however, forbidden: , , , , , , , , .

The automatic filename composition option is a powerful tool. The station uses a set of variables preceded by the "%" character as described below. After an analysis is terminated each variable is replaced by its current value. Variables may be combined at will. The current value is displayed in parentheses above the editing field and in the **Information table** of the [Instrument](#) window.

Tab 1: List of common file name variables

Variable	Meaning	Example of use
%n	Sample serial number. The value may be modified in the <i>Counter</i> field. The value may be set automatically to 0 or 1 each time the Instrument is opened by the <i>Reset when Open Instrument</i> checkbox from the <i>Counter</i> section of the User Options - General dialog. For more details see the chapter General tab on pg 104.	
%xn	Selects fixed number of places (e.g. for x=3:).	%3n = 001-999
%n(x)	Selects step, (e.g. for x=5:)	%n(5)=1, 6, 11
%P	Name of the project	Pollutants
%J	Name of the method (without .MET suffix)	DDTInWater
%s	Name of the sequence (without .SEQ suffix)	BalticSea
%v	Vial number (Only for the sequence table).	
%i	The number of the injection taken from a given vial (Only for the sequence table).	
%L, %l	The Sequence Line Number is a number of the row in the Sequence Table (Only for the sequence table).	
%c	Number of an Instrument.	
%e	Name of an Instrument.	
%g	Analyst name/ User name.	
%Q	Sample (contents of the Sample field).	
%q	Sample ID (contents of the Sample ID field).	
%f	Vial barcode.	BarCode-00001
%%	The sign for percentage.	%
%T	Time of the analysis start (in the format hh-mm).	%T=13-41
%D	Date (in the format dd_mm_yyyy).	%D=24_02_2009

Advanced date and time formatting

Submenu with advanced date and time formatting functions.

Tab 2: List of advanced file name variables

Variable	Meaning	Example of use
%H	Hours (24-hour format).	00 - 24
%h	Hours in the (12-hour format).	00 - 12
%p	Designation a.m./p.m.	AM/PM
%M	Minutes.	0 - 59
%S	Seconds.	0 - 59
%a	Day of the week (abbreviated).	Mo, Thu, ...
%A	Day of the week.	Monday, ...
%b	Month (abbreviated).	Jan, Feb, ...
%B	Month name.	January, ...
%R, %r	Date and time of the analysis start as set in the current Windows format.	
%o	Day of month.	01 - 31
%j	Day in calendar year.	1 - 366
%m	Month.	01 - 12
%U	Week of calendar year beginning on Sunday.	00 - 53
%u	Day of week beginning on Sunday.	1 - 7
%W	Week of calendar year beginning on Monday.	00 - 53
%w	Day of week beginning on Monday.	1 - 7
%x	Date as set in the current Windows format.	
%X	Time as set in the current Windows format.	
%y	Last two digits of calendar year.	00 - 99
%Y	Full calendar year.	2001
%Z	Name of the time zone.	

The above variables may also be used in *Sample* and *Sample ID* fields both for individual injections and in a sequence table. In order to use the variables in *Sample* and *Sample ID* fields you must also use it in the filename.

Note: Some parameters (e.g. %p) may depend on the settings in the **Windows** system and may thus be shown in different format than in the table above.

2.4.1.2 Rich Text Edit

The **Rich Text Edit** dialog allows the user to enter any description of the method. The description can be formatted using the classical functions of text processors (Bold, Italics, Underlined text, text alignment, text color, text font, text size, bulleted text). This dialog is accessible (and the method description can be edited) prior to the analysis from the **Single Analysis** dialog using the *Comments...* button or after the analysis has ended from the **Chromatogram** window using the *Edit...* button in the **Chromatogram** section of the **Measurement Conditions - Instrument** tab.

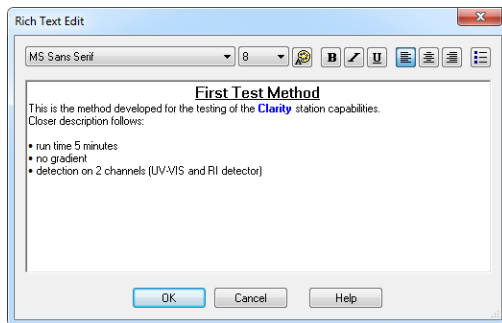


Fig 59: Rich Text Edit

2.4.2 Sequence

Command for the preparation and measurement of a series of analyses using sequence file. The **Sequence** window is displayed at the prompt. The major component is a **Sequence Table**, where each line describes the method of processing for one or more samples. Detailed information on **Sequence** window can be found in the chapter "**Sequence**" on pg 294.

2.4.3 Batch Full version

Batch processes single chromatograms or automatically and simultaneously executes selected actions with numerous chromatograms. When invoked, the **Batch** dialog will be displayed.

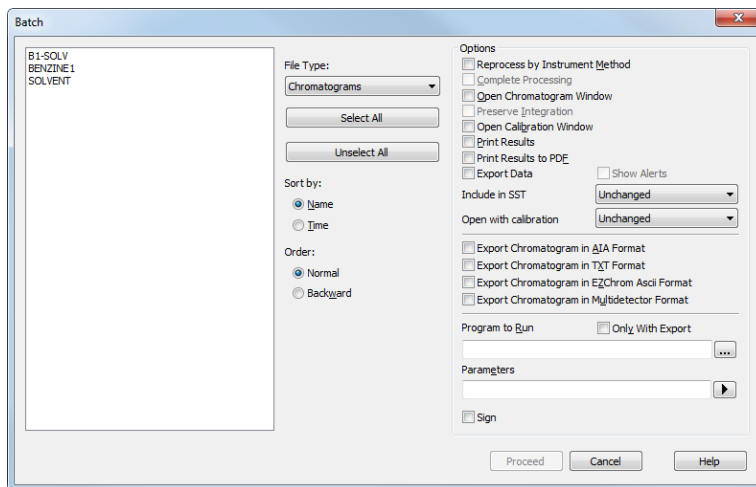


Fig 60: Batch

List of files

Lists all chromatograms or sequences from the current project according to the settings in the *File Type* field. Left click the mouse while holding the **Ctrl** key to select and deselect multiple files. This will enable you to progressively mark any arbitrary number of displayed files.

File Type

Selects the type of files to be batch processed.

Chromatograms - The list will display chromatograms from the data subdirectory.

Calibration Standards - The list will display chromatograms from the calibration subdirectory (i.e. calibration standards).

Sequence Files - Sequences from the current project will be displayed. All chromatograms which were measured in the last run of the marked sequences will be processed.

Note: Each sequence file remembers chromatograms, which were measured last and, therefore, it is not necessary to update the *File Name* column. Chromatograms not found are ignored.

Caution: When a sequence is run repeatedly, the list of measured files in that sequence will be reset and the original files cannot be processed using the *Batch* command.

Select All/ Unselect All

Quickly selects or deselects all files in the list.

Sort by Name/Time

Sorts files by name or by the time upon which they were last saved.

Sort - Normal/Backward

Sorts files normally or in reversed order. As files are processed in the order in which they appear in the list, the order of processing can be changed this way.

Options section**Reprocess by Instrument Method**

Reintegrates all selected chromatograms according to the current method.

Note: In this case, chromatograms that have been processed according to the sequence file, will merely be reintegrated. For complete processing, including updating according to the current sequence table contents (header update, recalibration), select *Complete Processing* parameter.

Complete Processing

This command is active only for processing chromatograms according to the sequence file (the *File Type* field must be set to *Sequence Files*). In this case, all chromatograms will be processed according to the current sequence table contents and methods used, including the eventual recalibration. This means that you can batch change headers of already measured chromatograms and reintegrate them, if you modify the integration parameters in used methods, etc.

Note: All other checkboxes except the *Sign* and *Preserve Integration* will be inactivated (dimmed).

Caution: Only the last set of measured chromatograms will be processed (if you used the same sequence more than once).

Open Chromatogram Window

Each chromatogram will be loaded into the [Chromatogram](#) window. Opening the chromatogram in [Chromatogram](#) window is necessary in case you want to include the chromatogram graph in the printed [Report](#).

Preserve Integration

Chromatograms will be proceeded without reintegration, original integration is preserved including hand-made modifications.

Open Calibration Window

Each chromatogram will be loaded into the [Calibration](#) window.

Print Results

Prints a report according to the report setup which is adjustable in the [Instrument](#) window using the *File - Report Setup* command or the  icon. The default print style is *Instrument*.

Print Results to PDF

Prints a report to the *.PDF file according to the report setup.

Export Data

Each chromatogram will be exported according to the settings in the [Export Data](#) dialog (for more details see also chapter **Export Data** on pg 96).

Show Alerts

If you do not wish to show system alerts, do not select this box. In this case system alerts are not displayed, but only saved in the log.

Include in SST

Selects whether the batch-processed chromatograms will be included in SST calculations. Available options are *Exclude*, *Include* and *Unchanged*.

Open with calibration

Selects the type of calibration connected to the chromatogram after the reprocess. Available options are *Linked*, *Stored* or *Unchanged*.

Export Chromatogram in AIA Format

Exports a chromatogram in AIA format. Corresponds to the *File - Export Chromatogram* command in the [Chromatogram](#) window (for more details see also chapter **Export Chromatogram** on pg 140).

The CDF format file will have the same name as the chromatogram and will be saved in the same (data or calibration) subdirectory of the current project as the exported chromatogram.

Export Chromatogram in TXT Format

Exports a chromatogram in TXT format. Corresponds to the *File - Export Chromatogram* command in the [Chromatogram](#) window (for more details see also chapter **Export Chromatogram** on pg 140).

The TXT format file will have the same name as the chromatogram and will be saved in the same (data or calibration) subdirectory of the current project as the exported chromatogram.

Export Chromatogram in EZChrom ASCII Format

Exports a chromatogram in ASC text format, as it is used by the EZ-Chrom program. Corresponds to the *File - Export Chromatogram* command in the [Chromatogram](#) window (for more details see also chapter **Export Chromatogram** on pg 140).

A file with the ASC suffix will have the same name as the chromatogram and will be saved in the same (data or calibration) subdirectory of the current project as the exported chromatogram.

Export Chromatogram in Multidetector Format

Exports a chromatogram in CHR text format so that all signals of multidetector chromatogram are exported to a single file (for more details see also chapter **Export Chromatogram** on pg 140).

A file with the CHR suffix will have the same name as the chromatogram and will be saved in the same (data or calibration) subdirectory of the current project as the exported chromatogram.

Program to Run

Names the program which is to be run after each chromatogram is processed (e.g. EXCEL.EXE).

Only with Export

When invoked, the program can only be run if the *Automatic Export* has also been selected.



Searches for the program to be run. Standard **MS Windows** window will be displayed for program search.

Parameters

Sets the parameters of an initiated program; these parameters are generally entered in the command line after the program name (e.g. name of loaded file, etc.).

The following variables can be used in the *Parameters* field:

- %f will be replaced by the name of the chromatogram after export.
- %e will be replaced by the name of the file after export. This does not apply when exporting multidetector chromatograms in the .TXT or .AIA formats.

Sign

Batch signs selected chromatograms. The [Sign](#) dialog will be displayed after running a batch, and you will be asked to enter a username, password and reason for signing.

Proceed

Runs batch processing.

2.5 Method

Displays and updates the measuring method through a submenu of commands. Every added Control module will have its own submenu command here (e.g *Thermostat Control*, *Valve Control*, *Fraction Collector Control*, ...). The method is divided into the following parts, each on a separate tab:

Tab 3: List of common tabs in the Method Setup dialog

Tabs	Contents of the tab
Event Table	Enables to trigger actions during analysis.
Measurement	Parameters that specify the measurement conditions.
Acquisition	Parameters that decide the method used to measure the detector signal.
Integration	Parameters that define the method of integration and baseline modification.
Calculation	Parameters that define the method of results calculation.
Advanced ✓ Full version	Contains <i>Subtraction Chromatogram</i> function and <i>Column Calculation</i> parameters.
GC (LC) Control ✓ Full version	Parameters that define direct control of the chromatograph.
AS Control ✓ Full version	Parameters that govern the method of an injection when using direct autosampler control.

Aside of the particular tabs the [Method Setup](#) dialog also contains a series of buttons. These buttons are present on every tab of the [Method Setup](#) dialog, in truth the tabs are only laid into the main frame which holds the buttons.

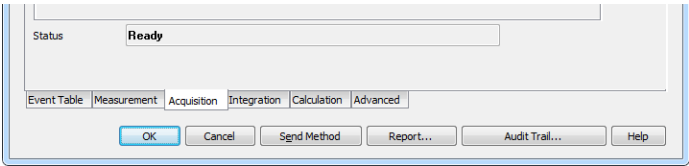


Fig 61: Method Setup - buttons

These buttons, when invoked, perform the following actions:

OK

Applies the changes performed in the method setup dialog to the opened version of the method and, if possible (acquisition is not running), sends or asks to send the method to all connected control modules/hardware (according to the settings in the [Instrument Method Sending](#) dialog). If the

acquisition is running, only the change in the *Autostop* and *Run Time* parameters from the [Method Setup - Measurement](#) tab are sent to the hardware.

Cancel

Discards all changes performed in the [Method Setup](#) dialog and leaves the dialog.

Send Method

When invoked, sends the method to all connected control modules/hardware. This button is not available during the data acquisition (when the hardware is working).

Report

Opens the [Report Setup](#) dialog on the [Method](#) tab. When there are any changes performed in the [Method Setup](#) dialog prior to using the *Report* button, the user is asked to either apply the changes or not. The report will then be made according to changed or unchanged method.

Audit Trail

Opens the [Method Audit Trail](#) window, a window holding an audit trail of the current method. When there are any changes performed in the [Method Setup](#) dialog prior to using the *Audit Trail* button, the user is asked to either apply the changes or not.

Note: The [Method Audit Trail](#) is, unlike other local audit trail windows, modal. That means that it is necessary to close it in order to be able to work with any other **Clarity** windows and dialogs.

Help

This button calls the online context help for the tab being currently opened.

2.5.1 Event Table

The *Event Table* enables to trigger action with specified Output based on events from selected Input. [Event Table](#) tab is common to all devices on the same **Clarity** Instrument.

Inputs:

- Acquisition and Idle events - Time, signal level, digital input state
- Events from Control module - Auxiliary signal level
- System events - Begin and end of sequence, acquisition etc.

Outputs:

- Digital Output control - **DataApex** A/D Converters , selected Control modules (for more details see the chapter **Events used in Active Sequence** on pg 79).
- Acquisition and Sequence control commands (*Stop, Abort, Skip Vial*, etc.).
- Run program - Launching an external program, with/without start parameters. Basic set of external tools is supplied with **Clarity**.
- None - Nothing will happen, suitable for development and testing of the method.

Each row in the table represents an independent event. When several events controlling the same output collide, the subsequent event masks the preceding one. When two events occur simultaneously, the event listed as last in the **Event Table** prevails.

Some events occurs in *RUNNING* or *IDLE* state only, these are coined with **Run** or **Idle** words in the event name.

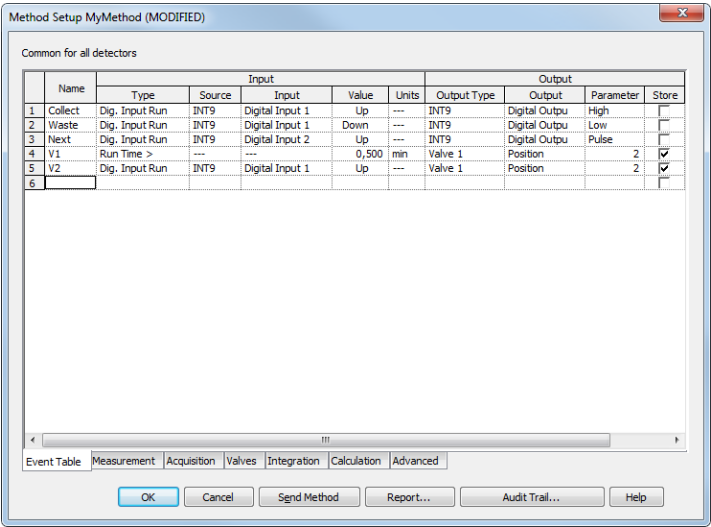


Fig 62: Method Setup - Event table

Individual lines or the entire table of events may be erased by right clicking the mouse on the available context menu above the table and using *Delete Row* or *Delete All Rows* commands. When a line becomes invalid, it will be highlighted with red background. Such row may be edited and must be repaired prior to saving the method. It is possible to select and copy the whole rows, paste them in the table and modify them. This is useful for creating sets of similar events.

Name

Arbitrary name to describe the event.

Input

The set of columns describing the Input that will trigger the action specified in *Output* columns.

Type

Sets the type of the input. Available options are shown in the **Tab 4** on pg **75**..

Tab 4: List of events available in the Event Table

Name of the event	Condition on when the event is triggered
<i>Run Time ></i>	Duration of acquisition (Instrument in <i>RUNNING</i> state) exceeds adjusted value (reset by every start of acquisition).
<i>Input Run ></i>	Measured quantity (e.g. input voltage or AU) exceeds adjusted value (works in <i>RUNNING</i> state only).
<i>Input Run <</i>	Measured quantity (e.g. input voltage or AU) drop below adjusted value (works in <i>RUNNING</i> state only).
<i>InpRel Run ></i>	Measured quantity (e.g. input voltage or AU) exceeds adjusted value+offset (works in <i>RUNNING</i> state only, offset is locked in the moment of acquisition start).
<i>InpRel Run <</i>	Measured quantity (e.g. input voltage or AU) drop below adjusted value+offset (works in <i>RUNNING</i> state only, offset is locked in the moment of acquisition start).
<i>Time Idle ></i>	Duration of idle time (Instrument in <i>WAITING</i> state) exceeds adjusted value (reset by every sending of method and by the end of the acquisition).
<i>Input Idle ></i>	Measured quantity (e.g. input voltage or AU) exceeds adjusted value (works except <i>RUNNING</i> state only).
<i>Input Idle <</i>	Measured quantity (e.g. input voltage or AU) drop below adjusted value (works except <i>RUNNING</i> state only).
<i>InpRel Idle ></i>	Measured quantity (e.g. input voltage or AU) exceeds adjusted value+offset (works except <i>RUNNING</i> state only, offset is locked in the moment of method sending).
<i>InpRel Idle <</i>	Measured quantity (e.g. input voltage or AU) drop below adjusted value+offset (works except <i>RUNNING</i> state only, offset is locked in the moment of method sending).
<i>Dig. Input Run</i>	Digital input changes its state (works in <i>RUNNING</i> state only).
<i>Dig. Input Idle</i>	Digital input changes its state (works except <i>RUNNING</i> state only).
<i>Aux Run></i>	Auxiliary input level exceeds adjusted value (works in <i>RUNNING</i> state only)
<i>Aux Run<</i>	Auxiliary input level drop below adjusted value (works in <i>RUNNING</i> state only)
<i>Aux Idle></i>	Auxiliary input level exceeds adjusted value (works except <i>RUNNING</i> state only)
<i>Aux Idle<</i>	Auxiliary input level drop below adjusted value (works except <i>RUNNING</i> state only)
<i>Ready</i>	All connected hardware becomes ready.
<i>Run Begin</i>	Run has begun, i.e. start of acquisition, start of the sequence row (before injection and after injection), but not the restart.

Name of the event	Condition on when the event is triggered
<i>Injection</i>	Start of the sequence row, Instrument is waiting for carrying out the injection.
<i>Bypass Injection</i>	Start of the bypassed sequence row (row with <i>Std = Bypass</i> value set).
<i>Acq Begin</i>	Acquisition has begun, i.e. start of acquisition, start of the sequence row (after injection or bypass), but not the restart.
<i>Restart Acq</i>	Restart of acquisition (by wire or manually by button).
<i>Acq End</i>	Correct completion of acquisition (<i>Autostop</i> or <i>Stop</i> button).
<i>Run End</i>	Completion of acquisition (<i>Autostop</i> , <i>Stop</i> button, <i>Abort</i> button).
<i>Control Elapsed</i>	Completion of acquisition and all programs on controlled devices (e.g. gradient program on pump or temperature program on column oven).
<i>HW Error</i>	Hardware error occurred (e.g. error on communication, device firmware error or physical failure).
<i>User Abort</i>	Abortion of acquisition by <i>Abort</i> button.
<i>User Shut Down</i>	System was shut down by user (using button on Device Monitor , command is available with controlled device connected only).
<i>Close Instrument</i>	Instrument was closed.
<i>SST OK</i>	On the end of acquisition, when SST resulted with OK.
<i>SST Unknown</i>	On the end of acquisition, when SST resulted with Unknown.
<i>SST Failure</i>	On the end of acquisition, when SST resulted with Failed.

Notes: *WAITING FOR INJECTION* state is considered as *IDLE* from event table's point of view, so for correct launching of e.g. *Stop Seq* command on any circumstances is necessary to create two events with the same command, one for *RUN* and one for *IDLE* state.

Notes: SST events occurs only when acquisition is finished.

Notes: User Shut Down event is available with controlled device connected only and after removing corresponding device from configuration particular event table row must be removed too.

Notes: In the event table, controlled devices are listed under names entered in their configuration dialogs, after change of name of the device, particular event table row must be corrected or removed.

Source

Sets the device which should be used as Input (e.g. detector, A/D converter, ...).

Input

Applies only for Digital Inputs, where it sets the input number.

Value

Indicates the value that is to be compared during the acquisition with current values (time, voltage on detector, digital input state etc.).

Note: When using Events evaluating Input values (Input Run, InpRel Run, Input Idle, InpRel Idle), please note that event will be triggered only when the change of input signal will be greater than EventThreshold value set in [Clarity.ini](#) file.

Units

Units of the *Value* (if applicable). For detectors, these are the units set in the [Device Setup](#) dialog of the [System Configuration](#) dialog. This field is not editable.

Output

The set of columns describing **Output** triggered by the action specified in *Input* field.

Output Type

Selects from the available output options:

- A/D converter (INT7, INT9, U-PAD, Net-PAD, etc.) - digital outputs
- Controlled device (Agilent 6890, etc.) - digital outputs
- Run Program - runs external program, command line parameters can be set.
- None - Nothing will happen, suitable for development and testing of the method. (e.g. when it is needed to log event in the chromatogram)

Note: There are two auxiliary software utilities in the UTILS subdirectory. SENDCOM.EXE for sending simple commands over serial line and BEEP.EXE for acoustic signaling, check the text files in the UTILS subdirectory for detailed description.

- Command - commands for internal controlling of the software run (for more details see the chapter **Sequence** on pg 310).

Output

Meaning of this field varies according to the *Output Type* used:

- A/D converter - sets the number of the controlling output to be set to the value indicated in the *Parameter* field.
- Controlled device - sets the name/number of the controlling output to be set to the value indicated in the *Parameter* field.
- Run Program - sets the filename of the software to be run (including the full path).
- Command - sets the command for internal controlling of the software run:

Tab 5: List of possible outputs for the Command output type

Name of the command	Meaning of the command
<i>Start Seq</i>	Start of the Sequence . Sequence file must be properly filled and saved in order to execute this command.
<i>Start Acq</i>	Start of the acquisition. When launched in <i>RUNNING</i> state, restarts the acquisition only with <i>Start-Restart</i> option checked (for more details see the chapter Measurement on pg 80). .
<i>Stop Acq</i>	Stop of the acquisition. In sequence run, current acquisition is halted, the Sequence continues on the next row.
<i>Stop Seq</i>	Stop of the Sequence . Current acquisition will be finished, then the sequence will stop.
<i>Repeat Injection</i>	Acquisition of current row in Sequence is halted immediately, the same sample is injected and analysed again.
<i>Skip Vial</i>	Acquisition of current row in Sequence is halted immediately, the Sequence continues on the next row.
<i>Abort</i>	Aborting of Single Run or the Sequence (chromatogram is always saved, regardless of setting in GLP options).
<i>Shutdown</i>	Shutting down all the devices connected (command is available with controlled device connected only).

Note: For more details see the chapter **Sequence** on pg 294.

Parameter

Its meaning varies according to the **Output Type** used:

- A/D Converter: The level to be set on the corresponding output specified in the *Output* field:

High - High level (output voltage 3 to 5 V) - corresponds to an opened contact.

Low - Low level (output voltage 0 to 1 V) - corresponds to a closed contact.

Pulse - Inverts the current level of the output for a given time and then returns back. Implicit length of the pulse is 1 sec.

Note: The duration of the pulse can be adjusted in the CLARITY.INI file (for more details see the chapter **clarity.ini file** on pg 388).

- Run Program: Adds command line parameters to be applied to the program specified in the *Output* field.

Store

Sets whether the performed event should be stored within the chromatogram file. Stored events can be displayed in the [Chromatogram](#) window.

2.5.1.1 Events used in Active Sequence

The control outputs are also used in *Active Sequence* to enable the autosampler to inject the next sample. In full version , these inputs are assigned in the [System Configuration](#) dialog.

Caution: Do not use these inputs and outputs in the **Event Table**.

Default assignment of the *Ext. Start Dig. Input* and *Ready Dig. Output* functions can be found in the manual for the corresponding hardware.

2.5.2 Measurement

The **Method Setup - Measurement** tab describes in detail the measuring conditions and determines the duration of an analysis.

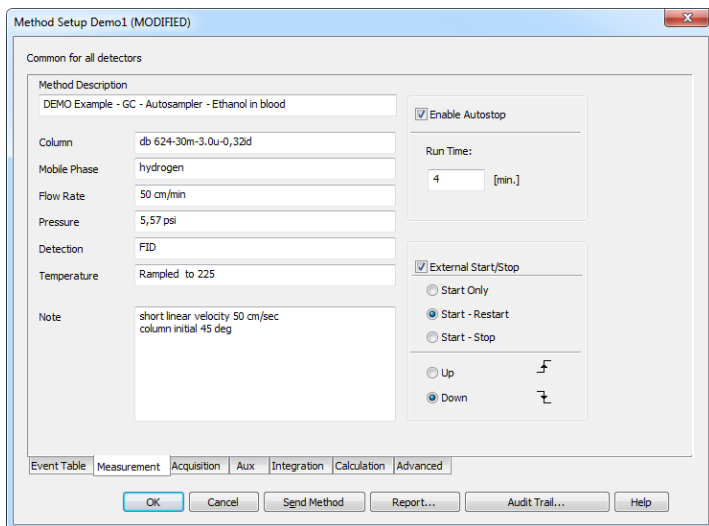


Fig 63: Method Setup - Measurement

The **Method Setup - Measurement** dialog will open with the method name displayed in the title bar when this command,  (GC) or  (LC) is invoked. While performing a multi-detector measurement, this tab is common for all signals.

Method Description

Contains the description of the method. When opening the method, the content of this field is displayed in the *Description* field of the **Open Method** dialog after the method is selected.

Column

Describes (informational purposes only) the column being used.

Mobile Phase

Describes (informational purposes only) the mobile phase used.

Flow Rate

Indicates (informational purposes only) the mobile phase flowrate.

Pressure

Indicates (informational purposes only) the mobile phase pressure.

Detection

Contains (informational purposes only) the detector characteristics.

Temperature

Describes (informational purposes only) temperature parameters.

Enable Autostop

Sets automatic termination of an analysis after the time set in the *Run Time* field.

Run Time

Specifies the duration of analysis. After the appointed run time has been reached, the analysis will automatically be terminated and processed. The possible time interval is between 0.2 - 9999 minutes. *Enable Autostop* option must be checked for this field to be editable and functional.

Note: The *Run Time* field must be set for all methods used in an active sequence since the time determines the cycle duration.

Note

An optional note (for informational purposes only).

External Start/Stop

Enables external start synchronization using an external signal from Digital Input on the A/D converter. Three types of the *External Start/Stop* operation are possible:

- *Start Only* - The external signal only starts analysis. Any external signals coming during the analysis will be ignored.
- *Start - Restart* - The external signal terminates the analysis in progress and at the same time starts a new one. The first signal starts the analysis, the second signal terminates it and at the same time starts a new one. This is known as “the continuous measurement”.

Note: The *Start/Restart* operation mode can not be performed for majority of directly-controlled instruments.

- *Start - Stop* - The external signal starts and stops an analysis. The first signal starts the analysis, the second signal stops it, third signals starts a new analysis etc.

The externally synchronized instrument may react on contact closure or contact opening. **Clarity** allows both options.

Up

The station reacts to a voltage change in the relevant controlling input from low (< 0.7 V) to high (> 3 V), or in other words, to an opened contact.

Down

The station reacts to a voltage change in the relevant controlling input from high (> 3 V) to low (< 0.7 V), or in other words, to a closed contact.

2.5.3 Acquisition

Sets the measuring conditions. When the menu command or  icon is invoked, the **Method Setup - Acquisition** dialog will be opened. The type of items displayed depends, in part, on the type of A/D converter that has been installed (or the type of chromatograph used while using direct digital acquisition). Be aware that parameters cannot be changed during analysis! The type of the A/D converter along with the name of the detector can be seen in the upper part of the tab.

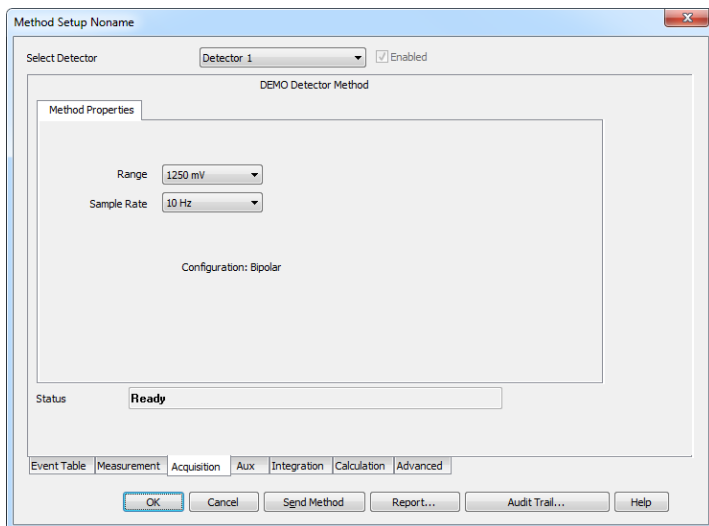


Fig 64: Method Setup - Acquisition

Note: At the beginning of the acquisition **Clarity** performs calibration of the A/D converter. Depending on the type of the converter (INT9, U-PAD2, Colibrick, etc.) and *Sample Rate* the first 20 - 200 ms of the chromatogram will not contain valid data.

Select Detector

Selects between the individual detectors in case several detectors are configured on the same Instrument (multi-detector configuration).

Caution: When using a multi-detector configuration, it is necessary to set the parameters individually for each detector.

Enabled

Enables/disables signal selected in the *Select Detector* field.

Range

Selects the voltage input range (in mV) and defines the maximum output voltage of the detector that the A/D converter can process.

Sample Rate

Specifies the sample rate (ranging from 6,3 to 480 samples per second, depending on the converter used) of a chromatogram. The selected sample rate will affect the ability to measure very narrow peaks. Sufficient number of samples (about 20) is a prerequisite for successful peak detection. This parameter decides on the size of the resulting chromatogram and the speed of its processing.

Note: Available sampling speeds may change with the acquisition equipment used (INT7, INT9, U-PAD, U-PAD2, Colibrick, Net-PAD, digital collection directly from the chromatograph, etc.).

2.5.4 Integration

Sets the integration parameters. When invoked by the menu command or by the  icon, the **Method Setup - Integration** dialog will be opened.

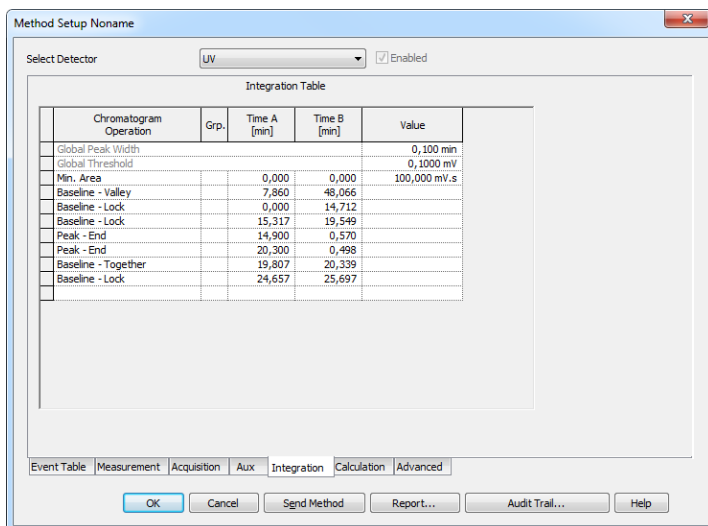


Fig 65: Method Setup - Integration

Caution: When using a multi-detector configuration, each detector will have its own integration table.

Select Detector

Selects between integration tables for the individual detectors, when using a multi-detector configuration.

Integration Table

The integration table defines the sequence of baseline operations to be conducted with a chromatogram that will be measured according to the displayed method.

Chromatogram Operation

Selects the operations to be performed on the chromatogram. Clicking the right border of the field displays a list of all possible operations. (This is identical to the list of commands in the *Chromatogram - Baseline*, *Peak*, *Integration*, *Separation* and other sub-menus of the *Chromatogram* window).

Grp.

Group identifier for the *Add Group* and *Delete Group* commands.

Time A

a) Indicates the starting time of the operation (for interval commands).

- b) Indicates the peak retention time (for commands that modify the peak start, peak end, etc.).

Time B

- a) Indicates the final time of the operation (for interval commands).
b) Indicates the position of a new start, end, etc. relative to the peak apex (for commands that modify the start, end, etc.).

Note: Zero entered in the *Time B* will set the end of the interval to the end of the chromatogram.

Note: Indicating the relative position (i.e. time) is more suitable than indicating a mere absolute value, since in using the **Integration Table** the peak with the most similar retention time is then found and the peak start/end is shifted to the point calculated from the relative shift on the actual retention time. This substantially diminishes the differences in actual retention time of the same peak in different chromatograms.

Value

Value (if relevant for the operation) with appropriate units.

2.5.4.1 Integration table items

The first two lines contain a default setting of two basic integration parameters - *Global Peak Width* and *Global Threshold*. This setting applies everywhere, where no values for *Local Peak Width* or *Local Threshold* have been entered.

Global Peak Width / Local Peak Width [min]

This parameter defines the smallest peak width, which is to be detected. The value range is between 0,001 and 10 minutes.

The value is not critical, so even narrower peaks than the entered value will be detected, but if set incorrectly, the beginning and end of a peak may be inaccurately determined and the baseline could be incorrectly inset.

The *Global Peak Width* parameter is always displayed in the first row of the integration table and is applied to the whole chromatogram.

It is possible to set a different peak width for individual sections by adding the *Local Peak Width* parameters into further rows of the integration table which then override the *Global Peak Width* settings in the given area of chromatogram.

Global Threshold / Local Threshold [mV]

Determines the sensitivity the integration algorithm has to noise. The height of all peaks (measured from the baseline to the top) that should be detected has to be at least double this parameter. The value range must be 1 μ V - 10 V.

The *Global Threshold* parameter is always displayed in the second row of the integration table and is applied to the whole chromatogram.

It is possible to set a different noise threshold for individual sections by adding the *Local Threshold* parameters into further rows of the integration table which then override the *Global Threshold* parameter in the given area of the chromatogram.

Integration Interval

Limits the time period over which the chromatogram will be integrated. This means that the chromatogram profile will be ignored beyond the set time interval. This is the main difference compared to *Lock* command, which cancels already integrated peaks only in defined sections and does not reintegrate the whole chromatogram. If the parameter is not defined, integration is done on the whole chromatogram.

Rejection section

Consists of several integration parameters which specify the conditions for including/excluding peaks in/from the integration. The peak parameters must be larger than the specified values, otherwise the peak will not be displayed.

Min. Area

This parameter sets the minimum peak area in mV.s. Peaks whose area is smaller than the specified value are not displayed.

Min. Height

This parameter gives the minimum peak height in mV. Peaks whose height is smaller than the specified value are not displayed.

Half Width [min]

This parameter gives the minimum peak half-width in minutes. Peaks whose *W05* is smaller or equal to the specified value are not integrated.

Separation section

Several parameters defining the integration of non-separated peaks.

ValleyToValley Slope

Specifies the maximum slope of the baseline. If the baseline slope is smaller than the specified value, the separating perpendicular line is canceled and the baseline passes through the valley.

The default value of 0 prevents the baseline from passing through the valley.

Tangent Area Ratio

Constitutes the first condition imposed on tangential separation. The condition is satisfied if the ratio of areas of the main and the separated peak exceeds the specified value.

The default value of 0 prevents tangential separation. If there is no *Tangent Slope Ratio* item defined in the **Integration table**, its value is considered to be 1.5.

Tangent Slope Ratio

This is the second condition imposed on tangential separation. The condition is satisfied if the ratio of slopes of the second and the first peak exceeds the specified value. The slope ratio is defined as the tangent to the first peak at its end point divided into the slope of the line connecting the beginning and end of the second peak.

The default value of 0 prevents tangential separation. If there is no *Tangent Area Ratio* item defined in the **Integration table**, its value is considered to be 0.1.

The evaluation process

The evaluation process does not proceed exactly in the same order as the parameters in the integration table are shown. The process is:

- Primarily initiated with the *Global (Local) Peak Width*, *Global (Local) Threshold* and *Integration Interval* parameters.
- Secondly the *ValleyToValley Slope* parameter will be applied. After all groups of non-resolved peaks have been tested by this parameter the remaining two parameters *Tangent Area Ratio* and *Tangent Slope Ratio* are applied to all remaining non-resolved groups. Tangential separation is used whenever the relevant criteria are met.
- Only then is the **Integration Table** used, followed by the *Rejections* parameters.

It thus follows that the separation parameters are not applied (e.g. to manually added peaks) and can be suppressed by the *Together* and *Valley* commands included in the **Integration table**.

2.5.5 Calculation

Sets the method of calculation display. When invoked by the menu command or via the  icon, the **Method Setup - Calculation** dialog will open. When utilizing a multi-detector chromatogram, the tab will be common to all signals.

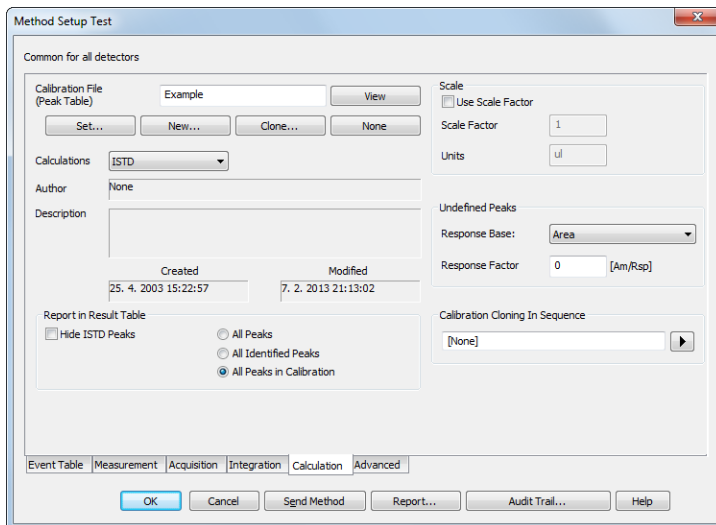


Fig 66: Method Setup - Calculation

Calibration File (Peak Table)

Name of the calibration file. Any chromatogram measured according to this method will be calibrated with it.

View

Displays calibration file in the **Calibration** window.

Note: The **View** command is contained in what is known as the modal window and, therefore, the **Calibration** window is displayed only under the **Method Setup** dialog. For a full display, you must close the **Method Setup** dialog using **OK** or **Cancel** button.

Set...

Selects a calibration file.

New...

Creates a new, empty calibration. By using this button the **Save As** dialog opens. It is possible to select the name and location of the calibration file there.

Clone...

Creates a copy of a current calibration file. By using the [Save As](#) dialog you can select the name and location of the calibration file. The copy will then be linked to the method.

Note: If the new calibration should not contain responses from the original file, they can be deleted by using the *Clear Responses* checkbox in the [Save As](#) dialog.

None

Closes a calibration file. The calculation type is changed to uncalibrated - *Uncal*.

Calculations

Selects a calculation type:

Uncal

Uncalibrated calculation not using a calibration file. The integration result is a percentage share of individual compounds in the chromatogram.

ESTD

External standard method. The result is a percentage and absolute share of individual compounds and their names.

ISTD

Internal standard. The result is a percentage and absolute share of individual compounds and their names.

NORM

This calculation type performs the calculations identically to the **ESTD** calculations according to the actual calibrations settings. Additionally it checks the condition, that all identified peaks are calibrated and all peaks in calibration were identified in the chromatogram. No results and warning message in *Result Table* header are given in opposite case. It means that *Peak Type* would be *ERROR* and *Amount%* zero.

STDADD

Standard Addition. The result is a percentage and absolute share of individual compounds and their names.

Caution: Each new unknown sample will need a new calibration file - use of the *Calibration Cloning In Sequence* function is highly recommended when the *STDADD* calculation is performed.

Integration

Selects the integration algorithm to be applied to new chromatograms measured according to this method or reprocessed according to this method. Some functions of the integration algorithm may be only present in particular versions of integration algorithm.

Author

Name of the user who created the calibration.

Description

Informative description of calibration.

Created

Date of calibration creation in the respective directory.

Caution: When copying or moving an arbitrary file, the **Created** field contains the date of this copy or move.

Modified

Date of last modification.

Report in Result Table

Determines which peaks will be displayed (and their values summed in the Total line) in the result table when calibrated calculations are in use.

Note: When the *Calculation* item is set to *Uncal*, this section will be inactive (dimmed).

Hide ISTD Peaks

The peaks identified as an internal standards will not be shown in the results table.

All Peaks

All chromatogram peaks will be displayed.

All Identified Peaks

Only peaks identified in the calibration file will be displayed.

All Peaks in Calibration

All peaks from the calibration file will be displayed regardless of how many peaks the chromatogram contains and how many were identified. This is known as the fixed table format as is made by the number of peaks in a calibrated file.

Scale

Auxiliary calculation parameters. To utilize these, do not forget to select the *Use Scale Factor* checkbox.

Note: When the *Calculation* item is set to *Uncal*, this section will be inactive (dimmed).

Use Scale Factor

Permits the use of *Scale Factor Units* and *Units After Scaling* parameters.

Scale Factor

All values in the *Amount [x]* column in the result table will be multiplied by this value.

Units After Scaling

Units for scale application. In the result table the original units adopted from the calibration file are replaced by units defined here.

Unidentified Peaks

Handles the behavior of unidentified peaks present in the chromatogram.

Note: When the *Calculation* item is set to *Uncal*, this section will be inactive (dimmed).

Uncal. Base

Selects the calculation base (*Area/Height*) for unidentified substances.

Uncal. Response

Coefficient for calculating the quantity of unidentified substances. The quantity is determined as the area (height - depending on the *Uncal. Base* item) of a peak multiplied by this coefficient. This is basically an artificial single-point calibration.

Calibration Cloning In Sequence

Sets the name for cloned calibration files used in the Standard Addition calibration method and during the Calibration Bracketing. The field enables the use of any text, however it is recommended to use variables (added using the  button) appropriate to the calibration method. For more details on the variable types and functions, see the chapter "**Chromatogram File Name**" on pg 64.

Note: It is advantageous to use the *Sample ID* (%q) as a part of the variable for the Standard Addition measurements, where all lines belonging to one sample (*Unknown*, possible *Blank*, all *Standards* on other levels of the same sample) share the same *Sample ID* in the [Sequence](#) window. For the Calibration Bracketing, the use of *Date* and *Time* (in any format) as a part of the *Calibration Cloning In Sequence* field may be advantageous. For Safe Calibration Usage, the use of *Date* (%D) variable as a part of the cloned calibration name is advantageous.

2.5.6 Advanced ✓ full version

Sets other parameters of the method, most notably the *Subtraction* and *Auxiliary Signals*. When invoked by the menu command, the **Method Setup - Advanced** dialog will open. The tab is common to all signals while performing multidetector measurement.

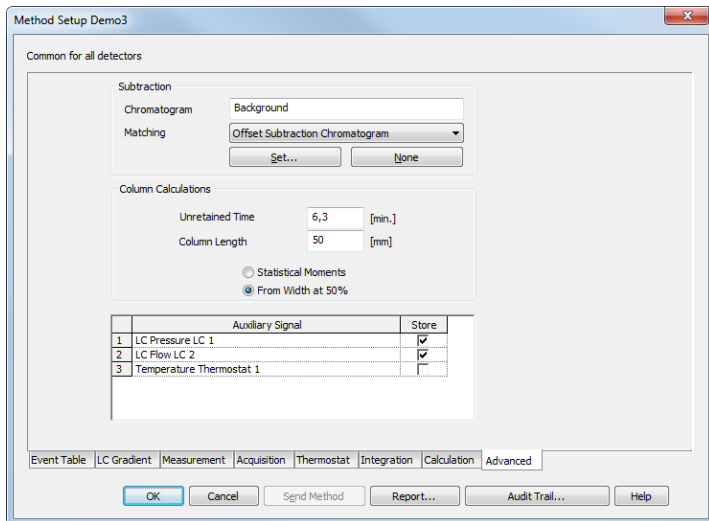


Fig 67: Method Setup - Advanced

Subtraction

Sets an automatic subtraction of a previously measured chromatogram from the currently measured chromatogram. The subtraction will be indicated in the printout with the “*Mathematically reprocessed*” inscription. Use of automatic subtraction is also shown in the Chromatogram Audit Trail.

The subtracted chromatogram is saved into the measured chromatogram, it is not simply linked to it. The changes occurring with the subtraction chromatogram thus do not influence the measured chromatogram from which the background is subtracted. The subtraction, however, does not hold its history - only last subtracted chromatogram's data is saved when it is changed, all older subtraction chromatograms are discarded.

Note: The subtracted chromatogram may be displayed during a measurement in the **Data Acquisition** window. It will appear in the background as a grey chromatogram.

Chromatogram

A dialog for specifying the subtracted chromatogram.

Matching

Sets the method of the background chromatogram subtraction.

No change - The subtraction chromatogram will be subtracted from the measured chromatogram without any additional amendments.

Offset Subtraction Chromatogram - The **Clarity** workstation attempts to “offset” the measured chromatogram with the subtraction chromatogram by shifting it on the axis X (time axis) before the actual subtracting.

Scale Subtraction Chromatogram - The **Clarity** workstation attempts to “offset” the measured chromatogram before subtracting by scaling the subtraction chromatogram on the time axis.

Note: Offsetting of chromatograms is designed according to the largest peaks in the chromatogram (i.e. either the highest peak or the peak with the greatest area).

Set...

Opens the [Open Chromatogram](#) dialog for selecting the file to be subtracted.

None

Cancels subtraction of the chromatogram.

Note: Should you wish to display e.g. the solvent only on the background of the [Data Acquisition](#) window without automatic subtraction, use the *File - Set Background Chromatogram...* command in the [Data Acquisition](#) window.

Column Calculations

Parameters affecting calculations on the [Performance](#) tab in the [Chromatogram](#) window.

Unretained Time

Determines the unretained peak time of used column in minutes. This parameter is used to calculate the capacity ratio.

Column Length

Determines the column length in millimeters. This parameter is used to calculate relative column efficiency.

Statistical Moments

Calculation runs based on statistical moments.

From Width at 50%

Calculation runs based on peak width at half the height and corresponding half-lengths.

Auxiliary Signals

This section of the [Method Setup - Advanced](#) tab is used for setting the auxiliary signals storing and possible display.

Note: The presence of auxiliary signals is dependent on the directly controlled device. Some of these have auxiliary signals, some do not.

Auxiliary signal

List of available auxiliary signals.

Store

Check the checkbox to save the particular auxiliary signal to the measured chromatogram.

Note: To display the auxiliary signals in the graph it is necessary to set their display in the [Graph Properties](#) dialog of the particular graph. It is, however, necessary to have the particular signal stored in order to be able to display it.

2.5.7 GC (LC) Control, LC Gradient

Direct instrument control pages. When prompted, the **GC** (**LC**, **LC Gradient** or other) tab of the **Method Setup** dialog will be displayed. The command is functional only if the respective control module is installed and some type of such directly controlled instrument is configured on the **Clarity** Instrument. Detailed information is included in corresponding **Clarity** Controls manual.

2.5.8 AS

Selection of autosampler sampling program. When prompted, the **AS** tab of the **Method Setup** dialog will be displayed. The command will only be functional if the respective control module has been installed and a directly controlled autosampler is configured on the given Instrument. Detailed information on **AS** tab can be found in the manual of respective control module.

2.6 Setting

Submenu with commands for setting export and post-processing of chromatograms.

Export Data

Opens the [Export Data](#) dialog which sets the method of automatic results export.



Postrun

Opens the [PostRun Setting](#) dialog which selects post-processing of a measured analyses.

2.6.1 Export Data

Sets the automatic results and chromatogram export parameters. When prompted, the [Export Data](#) dialog will be displayed. The same dialog can be called up using the *File - Export - Export Data* command from the [Chromatogram](#) window, where the command is intended for manual export of the active chromatogram and its results. Remember that this is the same window, so changes made in one place also apply to the other window. Selected data can be exported to a file or clipboard. Automatic export can be activated in the [PostRun Setting](#) window.

Note: User columns may only be exported from the **Result Table**.

Using the **Ctrl + E** shortcut it is possible to export data directly from the [Chromatogram](#) window without opening the [Export Data](#) dialog.

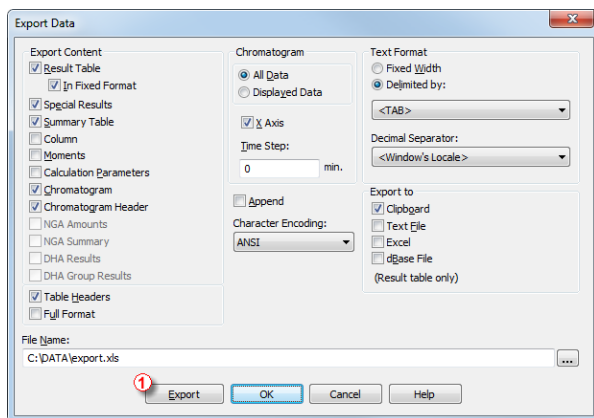


Fig 68: Export Data

Note: Besides exporting using *Export Data* command, a chromatogram can also be exported by *Export Chromatogram in AIA (TXT, EZChrom) Format* command (for more details see the chapter **Export Chromatogram** on pg 140).

Export Content

Determines the type of exported data.

Result Table

Exports integration results table.

Caution: Hidden columns will not be exported.

In Fixed Format

When checked the data will be exported always in the same format regardless of the setting on the screen. This means it will not be WYSIWYG format.

It is useful to check this option when the exported data should be automatically processed by other program. However, user columns will not be exported while the *In Fixed Format* checkbox is checked.

Special Results

Exports any data given by control modules (e.g. fraction table in case of fraction collector, ...).

Summary Table

Exports summary table.

Column

Exports *Column Performance* table counted from the half height. This table is exported in fixed format.

Moments

Exports *Column Performance* table counted from statistical moments. This table is exported in fixed format.

Calculation Parameters

Exports calculation parameters (parameters from the [Method Setup - Calculation](#) tab).

Chromatogram

Exports actual chromatogram according to specifications in the *Chromatogram* section.

Chromatogram Header

Exports chromatogram header.

NGA Result Table

Exports the Result Table from the *NGA Results* tab in the **NGA** Extension. Available only if the *Instrument Type* is NGA.

NGA Summary Table

Exports the Summary Table from the *NGA Summary* tab in the **NGA** Extension. Available only if the *Instrument Type* is NGA.

Table Headers

Exports all table headers with exported tables. (In export to Excel, Summary Table always contains table header.)

Full Format

Exports full format of all tables. Each line is extended by the filename, date and time of creation.

Chromatogram

Section determining the format of exported chromatogram. It will be active only when the *Chromatogram* checkbox in the Export Content group is checked.

All Data

Exports whole chromatogram regardless of selection.

Displayed Data

Only the displayed time period of chromatogram will be exported.

X Axis

Determines, whether or not will the data contain time in the first column.

Note: The time axis is exported with 5 decimal places precision (other data with 4 decimal places).

Time Step

Time step in minutes. Determines the length of time period from which will each exported value be calculated (using arithmetic mean). The lowest purposeful time step is a thirtieth of the *Peak Width* parameter for a specific chromatogram. If a smaller or zero value is entered, output data is exported at this, the lowest step. If you enter a time step larger than the chromatogram length, you will get a single value corresponding to the mean voltage value of the whole chromatogram.

Append

When the *File Name* field in [Export Data](#) dialog is set, then *Append* checkbox will:

- append data to existing file or
- create new file if it did not exist.

Note: If the *File Name* field is empty, it will always create separate file for each chromatogram.

Caution: Excel files cannot be appended, a new file is always created instead.

Character Encoding

Selects the encoding type of the chromatogram while exporting. *ANSI* or *Unicode (UTF-8)* encoding format is available.

Text Format

Output format selection:

Fixed Width

Items will be separated by an adequate number of spaces to maintain alignment (this occurs only for disproportional letter types).

Delimited By:

Items will be separated by: *Comma*, *Semicolon* or *Tabulator*.

Decimal Separator:

Chose decimal separator: *Window's Locale*, *Dot* or *Comma*.

Automatic Export To**Clipboard**

Automatically exports to clipboard.

Text File

Automatically exports to text file.

Excel

Automatically exports to Excel file.

Note: Exporting to Excel is not supported in any **Starter** edition of MS Office.

dBase File

Automatically exports to database file of dBase format. Only the results table is exported in database format.

If you select export only to the *dBase File*, the entire window will be grayed and *Results* and *Full Format* checkboxes will be checked.

File Name

Filename for export to a file. If this field is not filled, a file with the same name as the exported chromatogram will be created with the TXT suffix, and located in the same directory.



Searches for a file or directory that is to be entered to the *File Name* field.

Export

Performs the export to destinations selected in *Automatic Export To* area.

Note: The *Export* button  is only available if you open this dialog from the [Chromatogram](#) window.

OK

Saves the actual settings.

Cancel

Leaves the dialog without saving any changes.

2.6.2 Postrun...

Post-processing of a chromatogram. Selected actions are always executed following the completion of each **Single Analysis** measurement. When prompted, the **Postrun Setting** dialog will be displayed.

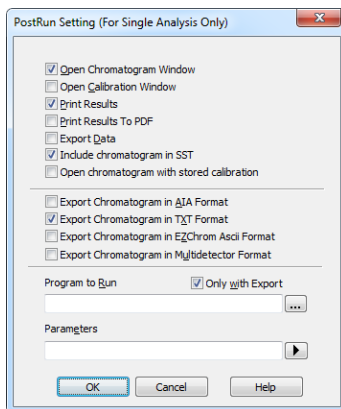


Fig 69: Postrun Setting

Note: Sequence measurement uses its own setting from the **Sequence Table**.

Open Chromatogram Window

The created chromatogram will be automatically loaded into the

Chromatogram window. A selected checkbox corresponds to the  symbol found next to the  icon in the **Instrument** window.

The minimized **Chromatogram** window is not shown up front at chromatogram upload when the *Open Chromatogram* command in the **Postrun Setting** dialog has been activated.

Open Calibration Window

The created chromatogram will be automatically loaded into the

Chromatogram window. A selected checkbox corresponds to the  symbol found next to the  icon in the **Instrument** window.

Print Results

A report will be printed according to the print style set in the **Instrument** window by using the *File - Report Setup* command. The default style is INSTRUMENT. A selected checkbox corresponds to the  symbol next to the  icon in the **Instrument** window.

Print Results To PDF

A report will be printed to *.PDF file according to the print style set in the **Instrument** window by using the *File - Report Setup* command. The

default style is INSTRUMENT.

Export Data

The created chromatogram will be exported according to settings made in the [Export Data](#) dialog (for more details see also chapter **Export Data** on pg 96).

Include Chromatogram in SST

Sets whether the created chromatogram will be included in the SST calculations or not.

Open Chromatogram with Stored Calibration

Saves and opens the chromatogram with the stored version of the calibration file, as opposed to the linked calibration when not checked.

Export Chromatogram in AIA Format

Exports a chromatogram in AIA format. This corresponds to the *File - Export Chromatogram* command in the [Chromatogram](#) window. The exported chromatogram will be saved under the same name as the source chromatogram (with the *.CDF suffix) and located in the same directory (for more details see also chapter **Export Chromatogram** on pg 140).

Export Chromatogram in TXT Format

Exports a chromatogram in the TXT format. Corresponds to the *File - Export Chromatogram* command in the [Chromatogram](#) window (for more details see also chapter **Export Chromatogram** on pg 140).

The exported chromatogram will be saved under the same name as the source chromatogram (with the *.TXT suffix) and located in the same directory.

Export Chromatogram in EZChrom ASCII Format

Exports a chromatogram in the ASC text format, which is used by the EZ CHROM program. This corresponds to *File - Export Chromatogram* command in the [Chromatogram](#) window (for more details see also chapter **Export Chromatogram** on pg 140).

The exported chromatogram will be saved under the same name as the source chromatogram (with the *.ASC suffix) and located in the same directory.

Export Chromatogram in Multidetector Format

Exports a chromatogram in the TXT format. Multiple signals will be stored in separate columns of the resulting file. Corresponds to the *File - Export Chromatogram* command in the [Chromatogram](#) window (for more details see also chapter **Export Chromatogram** on pg 140).

The exported chromatogram will be saved under the same name as the source chromatogram (with the *.TXT suffix) and located in the same directory.

Program to Run

Sets the program which will be run after the completion of the measurement.



Searches for the program which should be entered in the *Program to Run* field.

Only with Export

This item has two functions:

- The program specified in the *Program to Run* field will only be run if the *Automatic Export* checkbox has been selected.
- While manually exporting from the [Chromatogram](#) window using **Ctrl + E** shortcut this item enables the simultaneous running of the program specified in the *Program to Run* field.

Parameters

Parameters of the run program, which are normally entered in the command line behind the program name (e.g. name of loaded file, macro, etc.).

The following variables can be used in the *Parameters* field:

- Variable %f will be substituted by the name of the chromatogram at the moment of processing.
- Variable %e will be substituted by the name of the file to be exported at the moment of processing. (This does not apply when exporting multidetector chromatograms in TXT or AIA formats.)

Click the  button next to the *Parameters* field to enter the variables automatically.

For facilitating of exported chromatogram files copying you can use COPYFILES.EXE utility, located in UTILS subfolder. It allows you to copy all the exports from one chromatogram (.TXT, .PDF) at once. An utility is thoroughly described in COPYFILES.TXT file in the same folder.

2.7 View

This is an entry point to a submenu of commands for setting up the appearance of the [Instrument](#) window.

Toolbar	Displays/hides all toolbars. The command is analogous to that shown in the main Clarity window (for more details see also chapter View on pg 30).
Customize...	User customized toolbars. The command is analogous to that shown in the main Clarity window (for more details see also chapter View on pg 30).
Reset All	Returns all menus and toolbars in the Instrument window into their initial state.
Options...	Sets up Instrument properties.

2.7.1 Options...

Sets [Instrument](#) window properties. When prompted, the [User Options](#) dialog will be displayed.

Settings are saved in the appearance settings file (*.DSK). This means that when "all user" mode is being used (no password entered) these settings are common to all users. These settings are common to all [Instrument](#) windows in the **Clarity** station.

 If the station is used in "login" mode (i.e. all users have a password), settings are derived from the settings file of the respective user.

Note: Location of the *.DSK file can be set in the [User Accounts](#) dialog in the *Desktop File* field (for more details see also chapter **User Accounts** on pg 3).

Currently logged user and his desktop file name is shown in the title bar of the dialog. In the **Fig 70** on pg **103** is example of "all user" mode configuration, when *Administrator* is the default user and *CLARITY* is the default desktop file name.

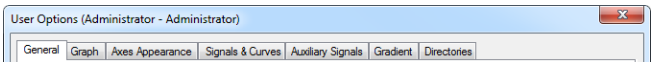


Fig 70: Title bar of the User Options dialog

2.7.1.1 General tab

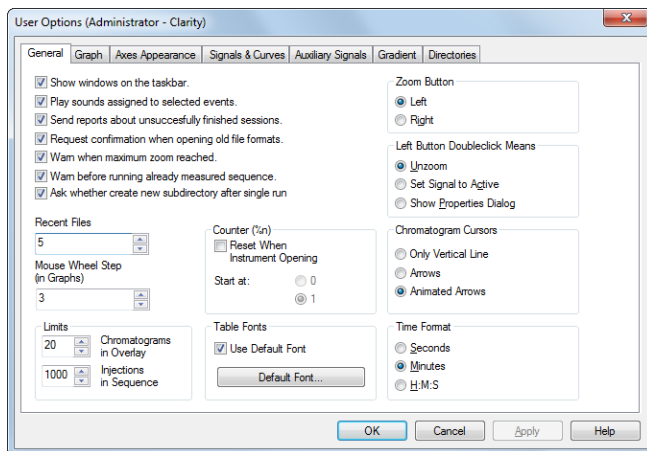


Fig 71: User Options - General

Show windows on the taskbar

All open windows can be displayed as icons on the taskbar or in the reverse case, it is possible to choose to have only the opened **Instrument** and the station's main **Clarity** window displayed on the taskbar.

Play sounds assigned to selected events

Permits sounds to be played with selected actions: *Start, Stop, Abort, Snapshot, Open Chromatogram, Close Chromatogram*. Other sounds can be selected on the **Sounds** tab of the **Control Panels** Windows dialog in the *DataApex - Clarity* section.

Send reports about unsuccessfully finished sessions

Permits the automatic delivery of a report that describes the problem to **DataApex** tech support. When this checkbox has been selected, a message box will automatically appear on the very next start-up after an error has occurred causing an improper shutdown of the system. This message box will enquire whether or not you would like to e-mail error information directly to the **DataApex** tech support.

Request confirmation when opening old file formats

When this checkbox is selected, the warning appears while opening any file saved in a station's previous version's format informing the user of the existence of this older format. You will then be advised how to resave the file, either in the original format or in the new version's format. If this option is unselected, the station will automatically convert all open files to the new version's format.

Warn when maximum zoom reached

Activates/deactivates the warning message display when attempting to exceed the maximum zoom of the graph.

Warn before running already measured sequence

Activates/deactivates the warning message display when attempting to run already measured sequence.

Note: Only the data from the last measurement of a given sequence may be reprocessed by the *Batch* command with *Sequence Files* parameter.

Ask whether create new subdirectory after single run

Enables/disables the confirmation message for creation of subdirectories after the single run measurement, if the subdirectory defined in the file name is not present. Similar feature is present in the Sequence measurements too, but the creation of the subdirectories is always silent.

Recent Files

Sets the number of recently opened files that will be displayed in the *File* menu. This setting is common to the [Instrument](#), [Sequence](#), [Chromatogram](#) and [Calibration](#) windows.

Mouse Wheel Step (in Graphs)

Sets the mouse wheel sensitivity on zooming the graphs.

Limits

Sets the general limits in **Clarity**.

"Chromatograms in Overlay"

Limits the number of chromatograms that can be simultaneously displayed in the [Chromatogram](#) window. The chromatogram displayed longest will be closed when the newer chromatogram is opened. 10,000 is the maximum number, the preset default is 20.

Injections in Sequence

Limits the maximum number of analyses measured from a single sequence. When the number of measurements reaches the set limit, the check of the sequence prevents from running it. This limit is set to prevent the lowering of the **Clarity** performance - sequence check may be unnecessarily long when there are several thousands of injections there.

Whole chromatograms are always counted regardless of the number of signals in the multi-detector chromatogram(s).

Counter (%n)

Allows for automatic reset of the %n counter used in the [Single Analysis](#) dialog. After checking, the counter will reset to 0 or 1 (according to further setting) at each [Instrument](#) window start.

The setting is common for one user, but the counter of each Instrument increases its value independently.

Note: Similar setting for sequential measurement can be found in the [Sequence Options](#) dialog (for more details see also chapter **Sequence Options** on pg 313).

Table Fonts

Selects default fonts for all **Clarity** station tables.

Use Default Font

When selected, the station will use the font set by the *Default Font...* button in all tables. If the box is checked and filled with gray color (instead of black), it means that at least one column in one table contains a different font.

Default Font...

After clicking this button the [Font](#) dialog will appear where you may choose the font, style, color and size of the table text font.

Note: If you wish to use a different font in any table, right click the mouse on the table and select the *Setup Columns...* command. This will open the [Setup Columns](#) dialog where it is necessary to deselect the *Use Default Font* checkbox in favor of desired font for specific column by using *Font* button.

Zoom Button

Select whether the selections from a graph should be made using the right mouse button (as in station **CSW17**) or the left mouse button (which is the new standard). If you use the left mouse button, the right button will be used to open a local menu.

Left Button Doubleclick Means

Defines what will occur when doubleclicking the left button in graph:

Unzoom - The graph returns to the appearance before zooming.

Set Signal to Active - If more chromatograms than one are displayed, doubleclicking over a specific chromatogram will set it as active.

Show Properties Dialog - Double-clicking on the graph opens the [Graph Properties](#) dialog for setting graph and chromatogram properties.

Chromatogram Cursors

Sets the cursor style for chromatogram modification. Cursor is shown as:

<i>Only Vertical Line</i>	A vertical line throughout the whole graph.
<i>Arrows</i>	Vertical line with two arrows at the point where this line intersects the chromatogram curve.
<i>Animated Arrows</i>	Animated arrows that function as those listed above.

Time Format

Determines the format of the graph's time axis in the [Data Acquisition](#), [Chromatogram](#) and [Calibration](#) windows.

Note: Times in all tables are always displayed in minutes!

2.7.1.2 Graph tab

Sets the appearance of the default graph. This selection applies to all graphs when the local setting in the [Graph Properties - Graph](#) dialog has the *Use User Options* checkbox selected.

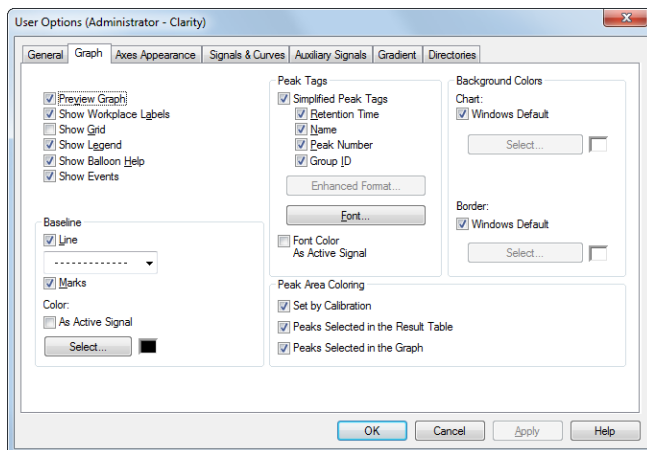


Fig 72: User Options - Graph

Preview Graph

Shows/hides a preview graph.

Show Workplace Labels

Shows/hides all labels and lines created using the *Create Label* command.

Show Grid

Shows/hides a grid for simplified graph reading.

Show Legend

Shows/hides a legend that displays chromatogram names and their signals.

Show Balloon Help

Displays balloon help next to the cursor lines when user performs interactive modifications of the chromatogram. This feature can also be switched on/off using the  icon on the toolbar in the [Chromatogram](#) window.

Show Events

Permits the display of events from the [Event Table](#) tab of the [Method Setup](#) dialog.

Baseline

Line - Shows/hides the baseline.

Marks - Shows/hides auxiliary marks of peak starts and ends.

Color

As Active Signal - Determines whether baseline color will be the same as the chromatogram color. If not, it will be the color set by the *Select* button.

Select - Button for color selection.

Peak Tags

Allows to shows, hide peak tags or set enhanced format of tags. The format of fields and values is defined by the format set for the **Result Table**.

Simplified Peaks Tags

Enables/disables checkboxes concerning simplified peaks tags. Simplified tags follows the changes of retention time caused by *Scale* and *Offset* parameters of [Time Axis](#) settings.

Retention Time - Retention time.

Name - Peak names made according to calibration file.

Peak Number - Peak serial number.

Group ID - Name of group to which a peak pertains.

Enhanced Format...

Displays the [Peak Tags Format](#) dialog which allows to set custom format of peak tags. Available only if the *Simplified Peaks Tag* checkbox is unchecked. Enhanced peak tags are derived from the result table, thus changes on [Time Axis](#) does not influence them.

Font...

Sets font for peak tags. When prompted, the standard [Font](#) dialog will be displayed.

Font Color As Active Signal

Determines whether the peak tag color will be the same as the chromatogram color. This overrides the color set by the *Font...* button.

Peak Area Coloring

Set by Calibration - The area of peaks will be colored based on the color specified in **Calibration Summary Table** (*Peak Color* column).

Peaks Selected in the Result Table - Peaks selected in the *Result Table* will be highlighted in the chromatogram.

Peaks Selected in the Graph - Peaks selected in the chromatogram graph will be highlighted in the [Result Table](#).

Background Colors

Selects the background color - *Chart* section - and border color - *Border* section - of a graph.

Windows Default - Color will automatically be adopted from the **MS Windows** appearance scheme. If deselected, the color will be set according to the *Select* button.

Select - Color selection button.

2.7.1.2.1 Peak Tags Format

Peak Tags Format dialog allows custom setting of the peak tag. The values are formatted according to the **Result Table**.

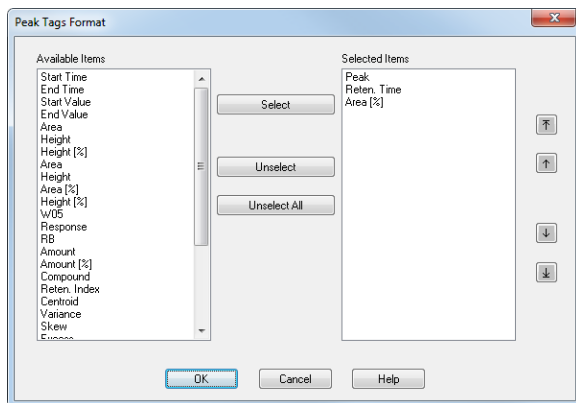


Fig 73: Peak tags format dialog

Available Items

Allows to select an item from the list.

Select

Moves the selected item from the *Available Items* list to the *Selected Items* list.

Unselect

Moves the selected item from the *Selected Items* list to the *Available Items* list.

Unselect All

Moves all items from the *Selected Items* list to the *Available Items* list.

Selected Items

List the items those will be displayed in the peak tag. To change their order, use the arrows on the right side of the dialog.

2.7.1.3 Axes Appearance tab

Axes appearance in the graph.

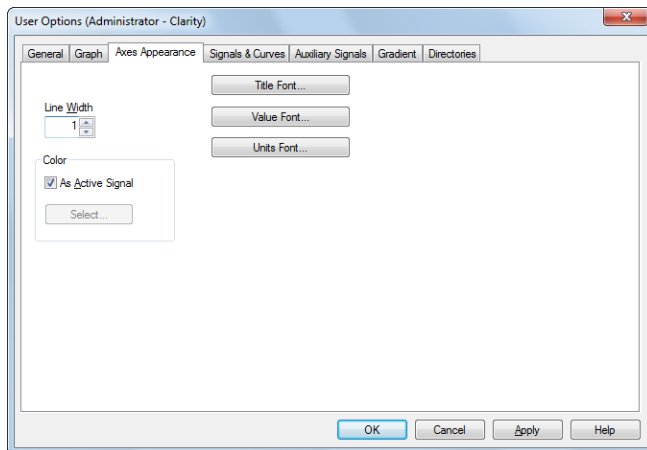


Fig 74: User Options - Axes Appearance

Line Width - Axis width in pixels.

Color - Axes color selection.

As Active Signal - Color will be the same as the color of the active signal.

Select... - Color selection button.

Title Font... - Selects font for axes graph titles.

Value Font... - Selects font for axes numerical values.

Units Font... - Selects font for axes units.

2.7.1.4 Signals & Curves tab

Sets the default curve properties.

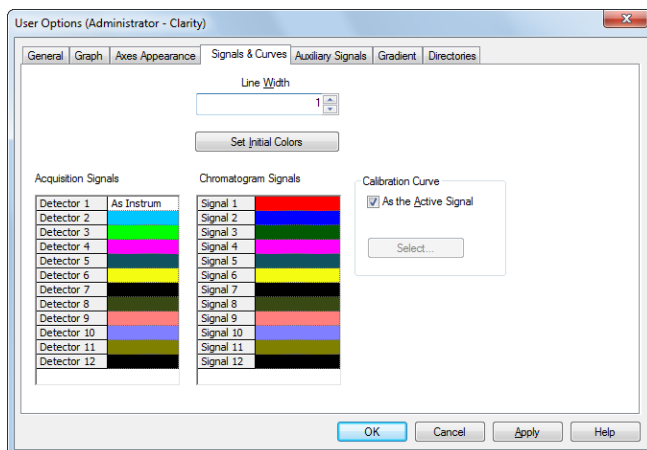


Fig 75: User Options - Signals & Curves

Line Width

Selects the default curve width.

Set Initial Colors

Resets to initial colors (the default setting returns colors to the state they were in after the initial installation of the **Clarity** station).

Acquisition Signals and Chromatogram Signals

Actual settings of colors for particular signals can be seen in these sections. Clicking on the appropriate button sets the color from a pre-set palette or will allow the user to create a newly defined color.

Calibration Curve

Section for changing the calibration curve color.

As the Active Signal

Sets the color of the calibration curve to the color of the active signal.

Select...

Button for the color selection.

2.7.1.5 Auxiliary Signals tab

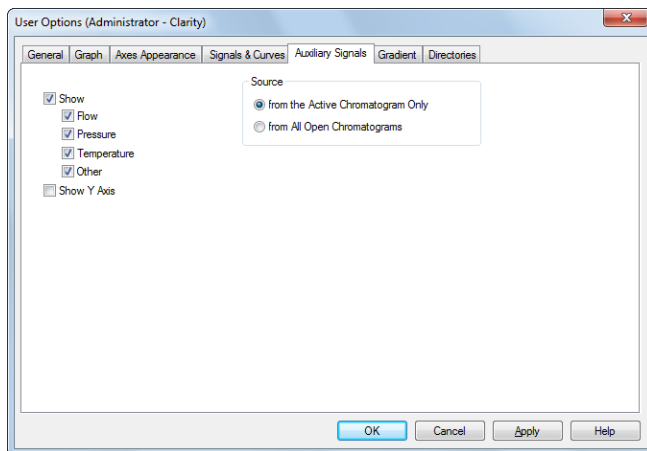


Fig 76: User Options - Auxiliary Signals

This tab determines which auxiliary signals will be displayed in the graph. Each auxiliary signal is included in one of following categories: *Flow*, *Pressure*, *Temperature*, *Other*. The display of all signals in each category is allowed together.

Show

If this field is left unchecked, no auxiliary signal will be displayed.

Note: To display any auxiliary signals it is also necessary to set the storing of the particular signal in the [Method Setup - Advanced](#) dialog (for more details see also chapter **Advanced** on pg 92).

Show Y Axis

Determines whether the secondary signal axis (displayed on the right of graphs, context to the auxiliary signal selected at the moment) will be shown.

Source

This section governs which auxiliary signals will be shown.

from the Active Chromatogram Only

Selected types of auxiliary signals will be displayed only in the active chromatogram.

from All Open Chromatograms

Selected types of auxiliary signals will be displayed in all opened chromatograms.

2.7.1.6 Gradient tab Full version

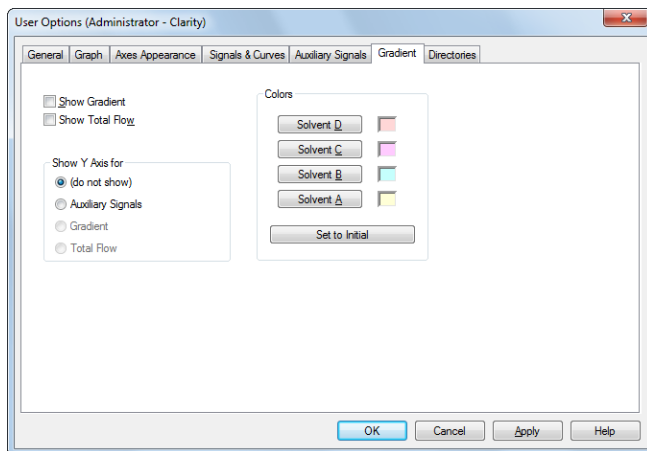


Fig 77: User Options - Gradient

Selects the gradient graph display (in the [Chromatogram](#) and [Data Acquisition](#) windows) when utilizing direct control of HPLC pumps.

Show Gradient

Switches the display of the gradient on/off and activates the possibility to select the gradient as the right Y axis.

Show Total Flow

Switches the display of the total flow on/off and activates the possibility to select the total flow as Y axis.

Show Total Flow

Switches the display of the total flow on/off and activates the possibility to select the total flow as Y axis.

Show Y Axis for

Determines the appearance of the right (second) Y axis. Available options are:

(do not show) - Not visible.

Auxiliary Signals - Shows the axis for auxiliary signals. To actually display the axis, it is necessary to select the particular auxiliary signal in the graph. In case that several auxiliary signals are in the same chromatogram, the axis of auxiliary signal that was active last will be displayed.

Gradient - The percentage share of individual solvents.

Total Flow - Flow rate in units set in the [Instrument Method Sending](#) dialog accessible from the [System Configuration](#) dialog.

Colors

Selects colors for the display of individual gradient components.

Set to Initial

Returns to default colors.

Solvent D-A

Selects the colors of each of four possible gradient components. Buttons are arranged in the same order as in the graph (i.e. from the bottom).

2.7.1.7 Directories tab

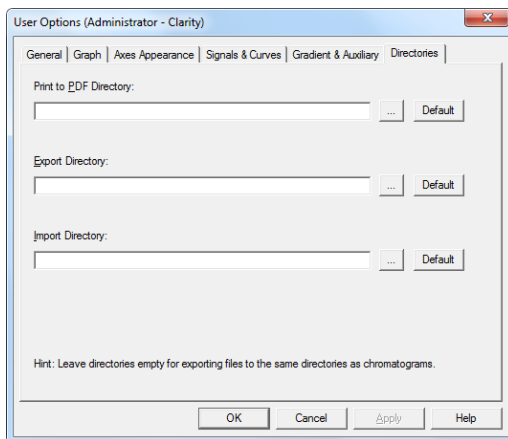


Fig 78: User Options - Directories

Sets the directories that will be used for the automatically created *.PDF files (Print to PDF Directory) and exported files (Export directory) and where **Clarity** will look for files to import (Import directory). If the fields are left empty, **Clarity** will use the DATA subdirectory of the current project as default directories for the respective actions. The *Export Directory* can be overridden when the *File Name* field in the [Export Data](#) dialog is specified.

2.8 Window

The commands for working with windows of given Instrument.

Hide All	Hides all open windows of the Instrument.
<i>Note:</i> When the <i>Show windows on the taskbar</i> checkbox in the User Options - General (described in the chapter "General tab" on pg 104.) is left unchecked, the windows will be hidden.	
Show All	Displays all hidden windows.
Main ✓ Full version	Displays the main Clarity window described on pg 1.
Chromatogram	Displays the Chromatogram window described on pg 118.
Calibration	Displays the Calibration window described on pg 238.
Single Analysis	Displays the Single Analysis dialog described on pg 59.
Sequence	Displays the Sequence window described on pg 294.
Station Audit Trail ✓ Full version	Displays the Audit Trail window described on pg 319.
Data Acquisition	Displays the Data Acquisition window for online monitoring of a signal from detector(s) described on pg 279.

2.9 Help

In the **Clarity** version, this submenu substitutes the missing *Help* submenu from the main [Clarity](#) window of the full version. All the submenus are the same as described in the chapter "**Help** " on pg **32**. In the full **Clarity** version, only the *Help* command is accessible from the [InstrumentHelp](#) submenu.

3 Chromatogram

This chapter is dedicated to describing the **Chromatogram** window, a window which is designed for displaying, adjusting and evaluating chromatograms. The window can be displayed using the *Window -*

Chromatogram command or the  icon from the **Instrument**, **Calibration** or **Sequence** window. The **Chromatogram** window may also be opened automatically when a new chromatogram is created, if the symbol next to the  icon in the **Instrument** window is set to  (in case of the single analysis) or  (sequence measurement).


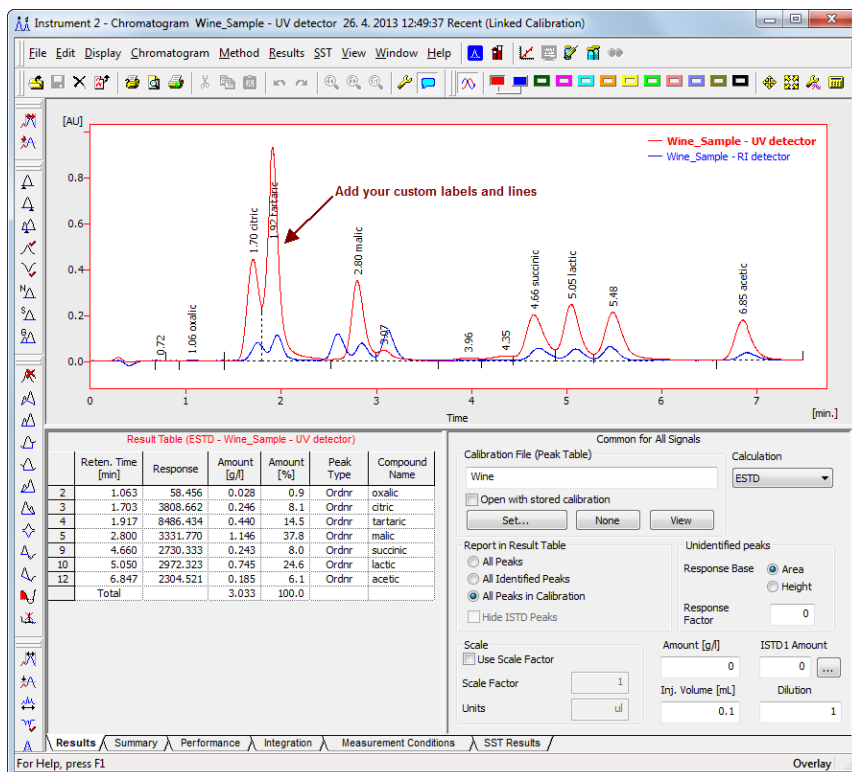


Fig 79: Chromatogram

Window header

Displays the name of the active chromatogram. The name of the active signal is displayed next to the name of the active chromatogram when utilizing a multi-detector chromatogram. When utilizing the display from any previous state of the chromatogram (using the *Method* field from the

Open Chromatogram dialog), both the date of this version and a flag reporting whether the used calibration is linked or saved in the chromatogram are also displayed next to the name of the chromatogram. The window is divided into an upper panel with a graph and two lower panels with tabs of the results and the method of the active chromatogram. The border between the panels may be freely shifted using the splitter. Double-clicking around the border (the cursor changes to ) enlarges the panel across the whole window, double-clicking again returns the panel to its original state.

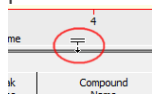

Orientation graph

An auxiliary graph where the chromatogram is always displayed in its original size. The graph is designated as an illustration of the cut-out performed in the main graph. Cut-outs may also be performed here, which is particularly suitable for shifting and zooming. The orientation graph is hidden by default, but can be displayed or hidden again by using the *Display - Preview Graph* command.

Note: After reducing the area of the graph below a certain level (e.g. by shifting the border "splitter" upwards) or reducing the overall size of the window, the small overview window (if opened) will hide and to be displayed again after the enlarging the graph, the command *Display - Preview Graph* have to be invoked once more.

Main graph

Displays the chromatogram including the baseline, thus enabling interactive adjustment of the baseline and many other functions. Method of graph display is set by *Display - Properties* command.

It is possible to select cut from chromatogram, simply by dragging the cursor over desired area with left mouse button pressed. The cut can moved along the whole chromatogram, by dragging with mouse (with Ctrl + left mouse button pressed), or by using keyboard arrows. See also the chapter "**Display**" on pg 156.

Orientation in peaks and their corresponding rows in [Result Table](#) is accomplished by filling the peak with a color after its selecting by clicking the cursor into its area. Corresponding row in [Result Table](#) is also highlighted. Selection is abandoned by clicking the cursor out of the peak area. It is also possible to select the row in [Result Table](#) by clicking into its first cell on the left (the one with number of peak displayed), corresponding peak in graph will be filled in. Multiple peaks can be selected at a time by clicking into their areas while holding **Ctrl** key or by clicking into the first and last one while holding **Shift** key. By default the same color is used for all the peaks (the same as the color of signal curve itself), custom colors can be set for every compound in Calibration (*Peak*

Color column). This feature can be further set in the [Graph Properties - Graph tab](#).

Current co-ordinates of cursor

The current position of the cursor is displayed in the right part window in the space between the orientation and main graphs or above the main graph, if the orientation graph is not displayed.

Graph legend

A legend describing the individual chromatograms (signals) is displayed in the upper right-hand corner of the graph. If you have more chromatograms opened than can fit on the screen, **▼▼** signs for scrolling the text will appear above and/or below the legend. The currently displayed chromatogram (signal) is highlighted in bold letters.

Double click the Chromatogram/Signal description in the legend to make it active.

Signals (or chromatograms in **Overlay** mode) can be hidden from view by **Ctrl** + left mouse click on the relevant colored  icon in the **Overlay** toolbar or **Ctrl** + double-click on its description in the legend.

Hidden Signal will be indicated in the **Overlay** toolbar by a cross .

Note: The legend is displayed or hidden by (un)checking the *Show Legend* checkbox in the [Graph Properties - Graph](#) dialog.

Balloon Help

When user performs interactive modifications of the chromatogram (adding/removing peaks, shifting peak start, etc.), **Clarity** displays balloon help tips next to the cursor lines.

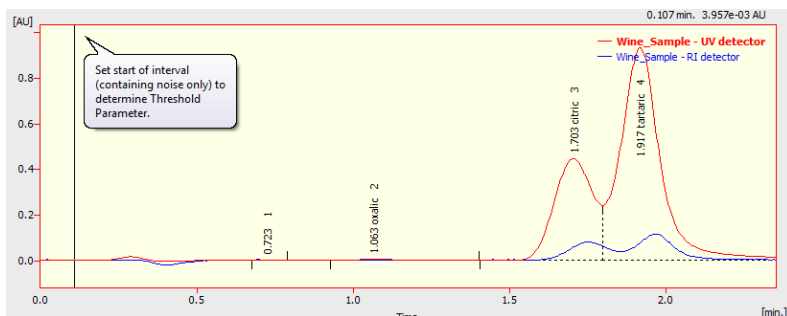


Fig 80: Balloon Help

This feature can be switched On/Off using the  icon on the **Standard** toolbar in the [Chromatogram](#) window.

Results panel

Displays a number of tabs in the lower part of the window, which contain result tables and other parameters of the chromatogram. They also enable the user to make adjustments to evaluations or display a description of the chromatogram.

Some tabs may not be visible following the station start-up, in which case the settings can be checked using the *View - Show Tables* command from the [Chromatogram](#) window menu.

Results

Contains result tables and parameters from the [Calculation](#) section of the method from which the last measurement was taken (for more details see the chapter **Result Table** on pg 222).

Summary

Contains a summary table of results for displaying the results of a number of chromatograms (signals) all at the same time (for more details see the chapter **Summary Table** on pg 228).

Performance ✓ full version

Contains a table of chromatogram parameters that define the quality of the system as well as relevant parameters from the [Calculation](#) section of the method according to which measurement was taken (for more details see the chapter **Performance Table** on pg 232).

Integration

Contains an **Integration Table** from the [Integration](#) section of the method according to which measurement was taken (for more details see the chapter **Integration** on pg 212).

Measurement Conditions

This tab has sub-tabs containing parameters from the sections [Measurement](#), [Event Table](#) and possibly also [GC,LC](#), [LC Gradient](#) or [AS](#) sections, according to which measurement was taken (for more details see the chapter **Measurement** on pg 216).

SST Result ✓ full version

Stands for “**System Suitability Test**” and is used for monitoring the quality of a measurement. This tab will be displayed only when a licence for the **SST** module has been purchased. When selected, the [SST Result](#) tab will be displayed. A detailed description of this item is included in a separate manual.

3.1 File

This is a sub-menu with commands for working with chromatograms and other specific system commands.

	Overlay Mode	Displays a number of chromatograms simultaneously.
	Open Chromatogram...	Selects and displays chromatogram (Ctrl + O).
	Open Chromatogram Sequentially	Opens and browses chromatograms from the last active folder. In <i>Overlay</i> mode browses through currently opened chromatograms.
	Open Chromatogram From Sequence	Displays all chromatograms measured according to the selected sequence.
	Close	Closes the active chromatogram (Ctrl + W). If the chromatogram has been changed and not yet saved, you will be notified accordingly. Modifications not yet saved are always marked by the inscription MODIFIED after the filename.
	Close All	Used to close all displayed chromatograms (Ctrl + Shift + W). If any chromatogram has been changed and not yet saved, you will be notified accordingly. Modifications in not saved chromatograms are always marked by the inscription MODIFIED after the filename.
	Save	The command saves the active chromatogram (Ctrl + S). When a mathematical curve created in the <i>OVERLAY</i> mode by the <i>Chromatogram - Overlay - Mathematics</i> command is to be saved, a chromatogram will be created and saved according to the current method shown in the Instrument window.
	Save As...	Saves active chromatogram under selected name (Ctrl + Shift + S).

Send Chromatogram by Email	Sends the displayed active chromatogram as an email attachment. In the Overlay mode all displayed chromatograms will be added to the email. The command is analogous to the <i>Send Method by Email</i> command described in the chapter "Send Method by Email" on pg 47.
Import Chromatogram...	Imports chromatograms in the text, AIA or EZChrom ASCII format.
Export	Initial menu for other Export commands:
Export Data...	Exports results and data of the active chromatogram. After selection, the Export Data dialog described on pg 96. will be displayed. By using the Ctrl + E shortcut it is possible to export the data without opening the Export Data dialog.
Caution: User columns are exported only from the Result Table, when the <i>In Fixed Format</i> checkbox is not checked.	
Export Chromatogram...	Exports chromatographic curve into AIA, text or EZChrom ASCII format.
Export Summary Table...	Exports <i>Summary Table</i> of results into a text file. After selecting the command the Export Summary Table To dialog will be displayed for selecting the name of the file and directory. Individual items of the summary table are divided in the file by a tabulator.
Caution: Chromatograms containing unsaved modifications will be saved before exporting. Export is logged in the Audit Trail.	
Caution: The <i>Summary Table</i> is exported in "as on screen" format.	
Export as Picture to Clipboard	Copies currently displayed chromatogram into clipboard as a vector picture.

Export as Picture to File...	Exports chromatogram as a picture in vector format. All chromatograms from the main graph including all labels are always exported in the format EMF format (Enhanced Metafile). The exported graph is saved into the file. After selecting the command, the Save As dialog will be displayed for selecting the name of the file and directory.
Sign... 	Electronically signs the active chromatogram.
Verify... 	Displays and verifies electronic signature.
 Report Setup...	Selects print configuration (Ctrl + Alt + P). After selection, the Report Setup dialog described on pg 333 . will be displayed.
 Print Preview	Previews selected print configuration before printing (Ctrl + Shift + P). After selection, the Print Preview dialog described on pg 352 . will be displayed.
 Print to PDF	Prints report to PDF file.
 Send Printed PDF by Email	Opens new email message in default email client, prints report to PDF and adds created PDF file as an email attachment.
 Print...	Command for selecting printer and actual printing (Ctrl + P). After selection, the Print dialog described on pg 354 . will be displayed.
 Perform Postrun...	Command for performing the postrun operations from the Chromatogram window. After selection, the PostRun Setting (From Chromatogram Window) dialog described on pg 145 . will be displayed.
-----	In this place the <i>File</i> menu contains a list of the last open chromatograms. This list serves for quick access to the currently used chromatograms.
Note:	The number of displayed files is set using the <i>Recent Files</i> field in the User Options - General dialog, accessible from the Instrument window using the <i>View - Options</i> command.

Exit

Closes the [Chromatogram](#) window. If any chromatogram has been changed and not yet saved, you will be notified accordingly. Modifications not yet saved are always marked by the inscription MODIFIED after the filename.

3.1.1 Overlay Mode

The user can work with multiple chromatograms in the *OVERLAY* mode. The activation of this mode is indicated by a checkmark next to the *Overlay Mode* command, an activated  icon in the *Overlay* toolbar and the display of the *OVERLAY* inscription in the status bar of the *Chromatogram* window (located in the lower right-hand corner of the window). In the *OVERLAY* mode it is possible to display a larger number of chromatograms at the same time by simultaneously or consecutively using the *Open* command or the  icon.

When a fixed range of display has not been set, the main graph will re-plot to display all chromatograms in their entirety even with the selection of variably sized chromatograms. Only the active chromatogram will display the descriptions of peaks and indicate the course of the baseline. The active chromatogram (signal) will have the relevant colored  icon depressed in the *Overlay* toolbar. Its name will be stated in the window header and a bullet will be displayed next to its name in the list of open chromatograms (signals) in the *Chromatogram* menu.

Note: Be careful not to confuse this list of currently open chromatograms with the list of last open chromatograms in the *File* menu.

The active chromatogram is also stated in bold letters in the legend on the right-hand side of the main graph. If no fixed color of the graph axes is selected, these axes always have the color of the active chromatogram.

Another active chromatogram is selected by:

- Clicking on the icon  with the color of the required chromatogram.
- Clicking on its name in the list of open chromatograms (signals) in the *Chromatogram* menu.
- Double-clicking on its name in the graph legend.

When using a multi-detector chromatogram, the icons of signals belonging to one chromatogram will be connected by the  link. After canceling the *OVERLAY* mode, only the active chromatogram will remain displayed.

3.1.2 Open Chromatogram

Selects and displays a chromatogram (**Ctrl** + **O**). After selection, the **Open Chromatogram** dialog (similar to the **Open Method** dialog for selection of method) will be displayed. The action of opening of the chromatogram is described in the chapter "**Open Method (Sequence, Chromatogram, Calibration)**" on pg 45.

Only the fields not described in the chapter Open Method are listed here.

Signals

When a multi-detector chromatogram is being used, this command enables a display selection of certain signals only.

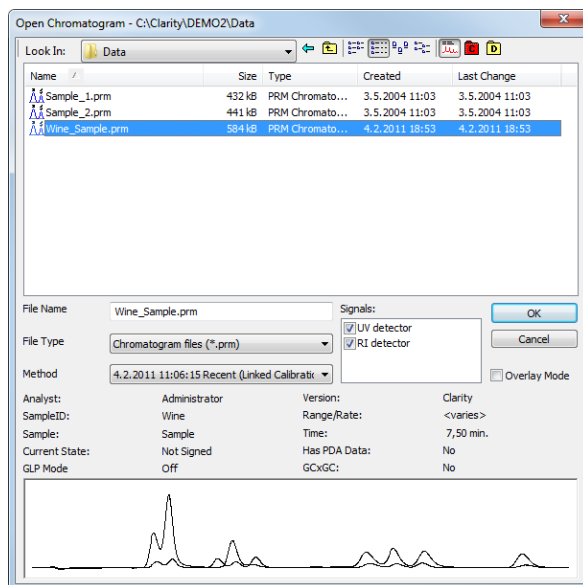


Fig 81: Open Chromatogram

Method

Enables the option of opening a previous version of a chromatogram. The chromatogram file maintains a history of adjustments as a list of methods. Current version of the method is saved, along with a copy of the currently attached calibration, into this list whenever the chromatogram is saved.

Note: In addition to manually saving a chromatogram using the Save and Save As commands, the station will automatically save each version when the chromatogram is printed, signed electronically or exported.

The calibrations stored in the chromatogram are stored without their own history and without calibration log. To conform with GLP it is

recommended to preserve the linked calibrations containing complete history of the calibration.

Detail Info

The *Detail Info* pane switched on/off using the  icon contains the detailed description of the chromatogram (information from the Chromatogram header, information about used Signature, presence of PDA data, software used for acquisition) and a preview of the curves.

Method (stored x linked)

By design, the chromatogram will be opened according to the method and calibration reference in which it was last saved. The first row of the *Method* field always ends with the *Recent (Linked Calibration)* inscription.

To display the chromatogram along with the last saved method including the saved calibration, select the second row of the **Method** field ending with the *Recent (Stored Calibration)* inscription. To display another version, select the row with the relevant date in the *Method* field.

When **Clarity** cannot find the linked calibration while opening a chromatogram, it will invoke a message offering to use calibration stored in the chromatogram.

Overlay Mode

This option is preset according to *OVERLAY MODE* option set by the *File - Overlay Mode* command in the [Chromatogram](#) window. If checked, more than one chromatogram may be selected simultaneously using the **Ctrl** or **Shift** shortcut.

Note: Continuous selection can be performed by clicking on the first file, pressing and holding down the **Shift** key, and then clicking on the last file in the row. All chromatograms will then be highlighted. Click on the **OK** button or hit the **Enter** key to then open these files.

Note: Non-continuous selection can be performed by using the **Ctrl** key. Hold this key down and click on the name of each desired chromatogram to highlight it. Click on the **OK** button or hit the **Enter** key to then open these files.

When attempting to open a chromatogram in an *OVERLAY MODE* which has already been opened and contains unsaved changes, a message will appear asking whether you wish to return to the original version of the chromatogram and thus discard all the unsaved changes.

Likewise, if the *OVERLAY MODE* is not switched on yet and there is an open chromatogram containing changes, a message will appear asking whether you wish to return to the original version of the chromatogram and thus discard all the unsaved changes.

Overlay Mode is not available when opening a chromatogram of the standard in the [Calibration](#) window.

3.1.3 Open Chromatogram Sequentially

Opens chromatograms from the last folder opened using *Open Chromatogram* command. In *Overlay* mode serves for browsing through currently opened chromatograms.

Caution: The **Sequential Opening** will be active only if the *Overlay* mode is **OFF**. With *Overlay* mode **ON** you can browse through previously opened chromatograms only.

◀ First	Opens the first chromatogram from current folder or currently overlaid files(Ctrl + F6)
◀ Previous	Opens the previous chromatogram from current folder or currently overlaid files(F6)
▶ Next	Opens the next chromatogram from current folder or currently overlaid files(F8)
▶ Last	Opens the last chromatogram from current folder or currently overlaid files(Ctrl + F8)

Filters and **sorting** applied in the *Open Chromatogram* dialog can limit the list of chromatograms for the **Sequential Opening** that will be performed later using the icons in the *Chromatogram* window.

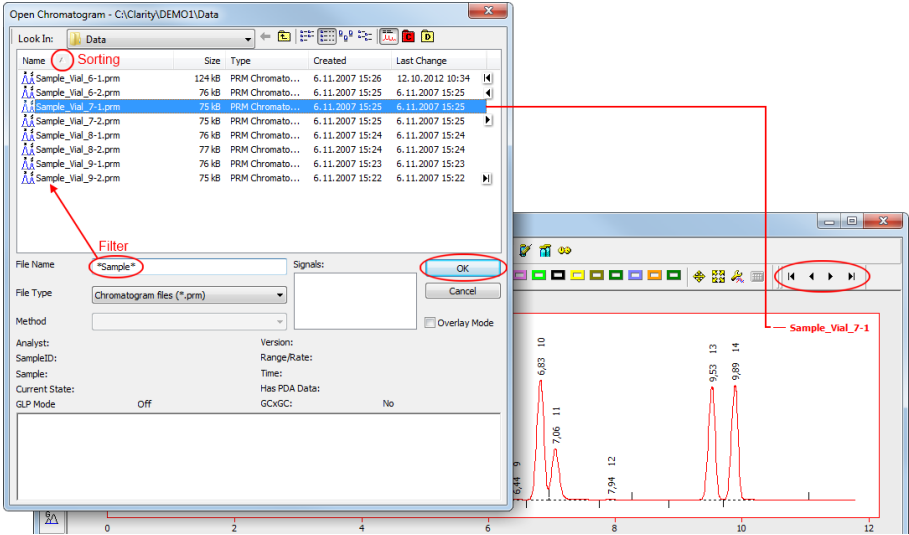


Fig 82: Open chromatogram sequentially

To apply sorting and filtering do the following:

- Check that the *Overlay* mode is **OFF**.
- Use the  *File - Open Chromatogram* command and navigate to the folder from which you want to open chromatograms.
- Specify the **filters** by writing partial name in the *File Name* field to limit the set to desired chromatograms only.

Note: It is possible to use the wildcards (*, ?) like in the example displayed in the image (*O91*).

- Click the desired column header to set ascending/descending **sorting** according to that column.
- Select a chromatogram from the defined set and open it using the *OK* button.
- In the **Chromatogram** window use the commands from *File - Open Chromatogram Sequentially* menu or icons     to browse through the chromatograms from the defined set.

The *Open Chromatogram Sequentially* function will always open the whole chromatogram (all of its signals).

3.1.4 Open Chromatogram from Sequence

Displays chromatograms measured according to sequence.

Caution: The *Open Chromatogram from Sequence* command will be active only if the *Overlay mode* is ON.

Using the command invokes the [Open Sequence](#) dialog. Once in the dialog, select the sequence file according to which the chromatograms you wish to display were measured. Only the chromatograms that have the checkbox in the *Run* column checked in the appropriate sequence will be opened. All checkboxes in the *Run* column are checked by default, but it might be useful to switch off some lines in the [Sequence](#) window prior to opening the chromatograms from sequence.

[Open Sequence](#) dialog (similar in appearance to the [Open Method](#) dialog described on pg 44.).

Method

Sets whether the opened chromatograms will be opened with the *Linked* calibration (might be changed from the time the chromatogram was opened for the last time) or with the *Stored* calibration.

Signals

Signal X number of opened signals

Only measured chromatograms will be displayed (i.e. chromatograms according to the rows of the sequence table where the *Sts.* field is colored blue).

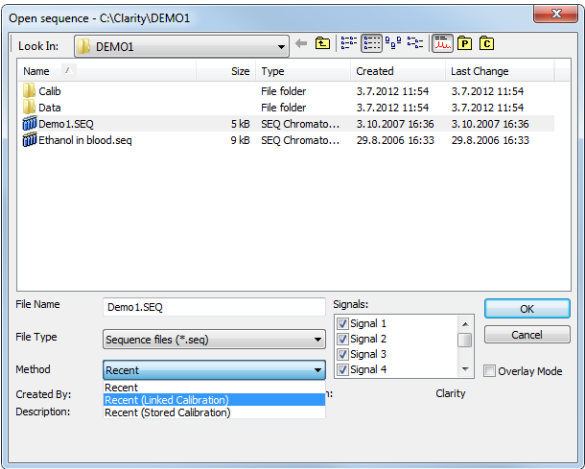


Fig 83: Open Sequence

If the same sequence was used several times, only the chromatograms from the last run of the sequence table will be remembered.

Note: This command is particularly suitable for a fast evaluation of a series of chromatograms in the summary table.

The sequence file must be saved after completing the measurement! Otherwise it will not be possible to use the *Open Chromatogram from Sequence* command on it, since the current state of the *Sts.* column has not been saved.

3.1.5 Save As...

Used to save the active chromatogram under a chosen name and/or to another directory (**Ctrl** + **Shift** + **S**). Saving always means that a copy of the active chromatogram is created.

The command opens the [Save As](#) dialog, similar to the [Open Method](#) dialog, where the new name can be entered or a new directory selected.

Note: If a file of that name already exists and has been opened on other Instrument or window, the name cannot be used. Moreover, in the protected mode (full version  only), a user may not overwrite a file he/she is not entitled to modify (see the *User Account* command).

Saving chromatogram in format of CSW station

In the *File Type* listbox select the version **CSW32** or **CSW17**.

3.1.6 Import Chromatogram...

Imports a chromatogram in **AIA**, **text** or **EZChrom** ASCII format. The Text files are supported with *.TXT, *.CHR and *.CSV suffixes. Select the [Open Files to Import](#) dialog using the *File - Import Chromatogram...* command. Select the desired directory and one or more files to import. The selection of files displayed may be influenced by using the filter in the *Files of Type* listbox.

Caution: Imported chromatograms are processed using the current method from the [Instrument](#) window and not according to the method of the potentially displayed chromatogram!

Import of several files simultaneously

You may indicate a number of files simultaneously in the [Open Files to Import](#) dialog. The [Import Text File](#) or [Import AIA File](#) dialogs will then open progressively for each imported file according to the type of imported file. If all the selected files share a same format, it is possible to import them all with the same settings.

Import of multi-detector chromatograms

The [Import Chromatogram...](#) dialog can import multi-detector chromatograms from *.TXT format.

3.1.6.1 AIA

The **AIA** file format is relatively strictly fixed in regards to the form, but not in the content. Thanks to this an **AIA** file may contain various custom items. The dialog for importing files in **AIA** format is designed as an unpacking tree, making it easier to view which items are contained in the imported file.

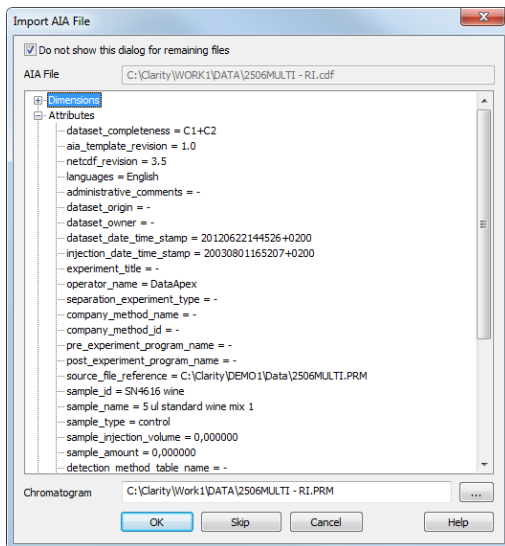


Fig 84: Import AIA File

Do not show this dialog for remaining files

With this checkbox checked, this dialog will not appear for other files in importing sequel, thus multiple files can be easily imported at once.

AIA File

Name of imported file including directory path.

Window with parameters

Items found in the current **AIA** file and their parameters. Expand collapsed items by clicking on the "+" sign with the left mouse button (the same as working with the **Explorer** application in **Windows**).

Clarity will correctly import only files with following parameters:

uniform-sampling-flag=Y

actual-sampling-interval value set

detector-minimum -value > -10000

detector-maximum-value < 10000

Note: Items and their values cannot be adjusted in the dialog.

Chromatogram

Name and location of the resultant chromatogram created after importing an **AIA** file. By design, the program will store the original name of the imported file along with the *.PRM suffix and the path to the directory of the current project.

To select another name and location of the new file, open the [Save As](#) dialog using the  button.

3.1.6.2 Text and EZChrom ASCII

Imports a chromatogram saved in a text file.

A chromatogram in text format may (but need not) contain a heading with a description of the conditions of measurement and must contain data of the chromatographic curve in one or more columns.

The station automatically interprets the headings of text files exported from the **Clarity** station and from the **EZChrom** program.

Note: The **EZChrom ASCII** text format uses the file suffix (*.ASC). It differs from the **text** format exported from **Clarity** station in the naming of individual items in the heading.

After selecting the chromatograms in the [Open Files to Import](#) dialog and pressing the *Open* button, the [Import Text File](#) dialog opens.

The conversion program determines the most suitable import setting according to the content of individual imported text files. Individual settings can subsequently be adjusted in the [Import Text File](#) dialog and saved for further use. Import is performed by the *OK* button.

Note: If a number of files was selected in the [Open Files to Import](#) dialog, a further [Import Text File](#) dialog or [Import AIA File](#) dialog will open (according to the type of the next of the indicated chromatograms) after clicking on the *OK* button. When *Apply Settings to* radio button is switched into the *All n Files* position, subsequent files will be proceed with the same settings. By checking *Prefer Detected Settings* checkbox you can force Clarity to use setting detected from file whenever it is possible.

Preset

When the same type of files are imported frequently, it can be useful and time-saving to save import settings for further use. In drop-down box on the left user can select from previously saved presets, by clicking  button right to list the presets management menu will open. The menu contain commands for saving current setting (*Create from the Current Settings*), *Rename* or *Remove* presets. Presets are stored as part of **Clarity** Desktop settings.

Apply Settings to

This radio button group indicates whether all the settings below will be applied to current file only, or to all files proceeded at once. In case multiple files are imported at a time, and this radio buttons are set at Current File, the Import Text File dialog will be invoked for every file in the sequel.

Current Text File

The name of imported file including directory path.

Preview

This button will open the current file in simple text viewer. This can help user to identify exact meaning of particular columns when it's not clear from file suffix.

Import From Line

Sets the line of the file on which the data start.

Import To Line

Sets the last line of the file that should be imported.

Separator

Selects the sign by which the individual items are separated in the text file. <TAB>, <SPACE>, <SEMICOLON> or <COMMA> are available to choose from.

Decimal Separator

Separator of decimal points - <DOT>, <COMMA> or <Windows Locale>. Separator indicated in Windows (as <Windows Locale>) is used as the default.

The 'Import Text File' dialog box contains the following fields and controls:

- Preset:** A dropdown menu set to '<None>'.
- Apply Settings to:** Radio buttons for 'Current File' (selected) and 'All 1 Files'. A checkbox for 'Prefer Detected Settings' is checked.
- Current Text File:** A text field showing 'C:\Clarity\WORK1\DATA\2506MULTI - RI.txt' with a 'Preview' button.
- Import From Line:** An empty text field.
- Import To Line:** An empty text field.
- Separator:** A dropdown menu set to '<TAB>'.
- Decimal Separator:** A dropdown menu set to '<DOT>'.
- Analyst:** A text field with 'DataApex'.
- Sample ID:** A text field with 'SN4616'.
- Sample:** A text field with '5'.
- Sample Date:** A text field with '1. 8. 2003' and a calendar icon.
- Amount:** A text field with '0'.
- Dilution:** A text field with '1'.
- ISTD Amount:** A text field with '0'.
- Injection Volume:** A text field with '0'.
- Imported Data:** A dropdown menu set to 'Time + Signal + Signal...'.
- Number of Detectors:** A dropdown menu set to '4'.
- Table:** A table with 7 columns: Sample Rate [Hz], Signal Units, Autoscale, Signal Name, Time Units, Y Multiplier, and Data Size. It contains 4 rows of data.
- Save as Chromatogram:** A text field with 'C:\Clarity\Work1\DATA\2506MULTI - RI.PRM' and a 'Load Defaults' button.
- Buttons:** 'OK', 'Skip', 'Cancel', and 'Help'.

Sample Rate [Hz]	Signal Units	Autoscale	Signal Name	Time Units	Y Multiplier	Data Size
10	V	☒		min	0,001	10142
10	V	☒		min	0,001	10142
10	V	☒		min	0,001	10141
10	V	☒		min	0,001	10142

Fig 85: Import Text File

Note: Parameters entered to *Import From Line*, *Import To Line*, *Separator* and *Decimal Separator* fields will be applied on all simultaneously imported chromatograms.

Analyst, Sample ID, Sample, Sample Date, Amount, Dilution, ISTD Amount, Injection Volume

These fields composing the chromatogram heading have the same meaning as described in the chapter "**Single**" on pg 59.

Imported Data

Time + Signal + Time + Signal - Every chromatographic curve is saved in a pair of columns, in which the first one represents a time and the second one a signal.

Signal + Time + Signal + Time - Every chromatographic curve is saved in a pair of columns, in which the first one represents a signal and the second one a time.

Signal only - Every chromatographic curve is saved in one column, where the values represent the signal only. The time basis is obtained from the *Sample Rate* field.

Time + Signal + Signal - Time base is contained in first column, every next column represents a signal of one detector.

Number of detectors

Specifies the number of detectors for the import of multi-detector chromatograms.

Sample Rate

Selects the sampling speed in the number of samples per second (Hz). Select one of the speeds from the list corresponding to the speeds of your A/D converter or enter your own value.

Signal Units

Determines the units in which the signal values are stated.

Note: Imported Data will not be parsed to recognize scaled units.

Autoscale

Sets the possibility to automatically scale the signal units in graphs.

Signal Name

Sets the signal name for further display.

Time Units

Determines the units in which time values are stated.

Possible values are: *ms* - milliseconds, *s* - seconds, *min* - minutes.

Note: Minutes are entered in decimal format (e.g. 1.5 min. = 1 min. 30 seconds).

Y Multiplier

Sets the multiplier for scaling the data on the signal axis.

Data Size

Number of points of a curve.

Note: This field cannot be adjusted. It serves only for referential checking as to whether the import takes place correctly.

Save as Chromatogram

Sets the name and location of the resultant chromatogram created after importing a text file. By default, the program will offer to use the original name of the imported file using the suffix *.PRM with a path to the directory of the current project.

To select another name and location for the new file, open the [Save As](#) dialog using the  button.

Load Defaults

Selects implicit values of all fields (i.e. returns the dialog into the same state it was in at the original moment the opening).

OK

Confirmation of import.

Cancel

Cancellation of import.

3.1.6.3 Import RAW...

Transforms raw data (from **CSW17** or the external **PAD** converter) to a chromatogram. Invoking the command opens the [Open RAW Data](#) dialog (similar in appearance to the [Open Method](#) dialog described on pg 44.). Here you will be invited to enter the name of the chromatogram to be created.

Caution: Raw data is processed using the current method from the [Instrument](#) window (not by the method of the currently displayed chromatogram, if any!).

3.1.7 Export Chromatogram

Exports chromatographic curve into **AIA** (*.CDF), **text** (*.TXT), **Multidetector text** (*.CHR) or **EZChrom ASCII** (*.ASC) format. Using the command invokes the **Export Chromatogram** dialog.

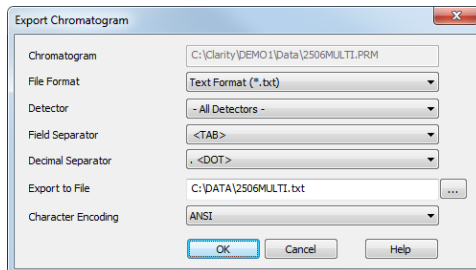


Fig 86: Export Chromatogram

Chromatogram

Name of the exported chromatogram including its full path. This field serves for informational purposes only and cannot be modified.

File Format

Selects one of the available formats for export:

AIA Format (*.CDF) - is one of the standard formats used for exchanging data among various chromatography systems. Multidetector chromatograms will be exported to multiple files, each signal into a separate file distinguished by the name of the detector.

Note: Further information about AIA format can be found for example in following article: Terrance A. Rooney - "STANDARDS FOR CHROMATOGRAPHY DATA SYSTEMS", Today's Chemist at Work, 1998, 7(5), 15-17.

Clarity only exports raw data and basic chromatogram description, no results, method data, calibration data or GLP data are exported into the *AIA Format*.

Following text formats (*Multidetector Text Format* (*.CHR), *Text Format* (*.TXT), *EZ Chrom ASCII Format* (*.ASC)) will contain a header and exported data in two columns separated by an arbitrary separator (TAB, COMMA, SEMICOLON).

Multidetector Text Format (*.CHR) - will always contain all available signals in a single exported file.

Text Format (*.TXT) - Multi-detector chromatograms will be exported to multiple files, each signal into a separate file distinguished by the name of the detector.

EZ Chrom ASCII Format (*.ASC) - is a text format used by EZ Chrom software. Comparing to the *Text Format* the data are exported in one

column only (voltage), separator can be <TAB> only and the structure of the header is slightly different.

Detector

Selects the signals for export. An individual detector, or all detectors at once can be selected.

Note: Only for *AIA* and *EZChromASCII Format*.

Field Separator

Selects the character to be used as data separator.

Note: Only for *Multidetector Text Format* and *Text Format*.

Decimal Separator

Selects the character to be used as decimal separator.

Note: Only for *Multidetector Text Format* and *Text Format*.

Export File

Invokes the [Save As](#) dialog in which the name and location of the file into which you wish to export the chromatogram is entered.

Note: If no name is entered, the name of the exported chromatogram is automatically used with the relevant suffix.

Character Encoding

You can chose to export Chromatogram using *ANSI* or *Unicode (UTF 8)* encoding.

Note: Only for *Multidetector Text Format* and *Text Format*.

Caution: Chromatogram containing unsaved modifications will be saved before exporting. Export is logged in the **Audit Trail**.

3.1.8 Sign ✓ Full version

Command for electronic signature of current version of chromatogram. After selection, the [Sign](#) dialog will be displayed.

Caution: Chromatograms cannot be signed unless you have set up a user account in the [User Accounts](#) dialog. It is also impossible to sign a chromatogram which has been opened using one of the methods from the history (*Method* field of the [Open Chromatogram](#) dialog).

Choose User

Selects the user who is to sign the current version. A list of users from the [User Accounts](#) dialog is displayed.

User Description

Content of the *Description* field of selected user from the [User Accounts](#) dialog.

Sign as Current User

The chromatogram will only be signed using the name and password entered within the framework of the **Clarity** station.

Sign Using Certificate

The chromatogram will be signed using the certificate installed in your computer. Select the certificate by using the *Select Certificate* command.

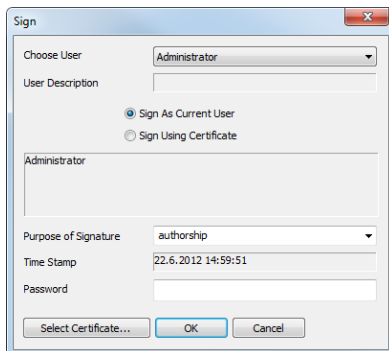


Fig 87: Sign

Purpose of Signature

Allows to enter the purpose of signature. Select from the preset options or supplement your own.

Time Stamp

Date and time of signature.

Password

Mandatory password of selected user.

Select Certificate

Selects certificate when selecting a signature using certificate (*Sign Using Certificate*).

3.1.9 Verify

Displays a list of the current chromatogram's electronic signatures. After selection, the [Verify Signatures](#) dialog will be displayed.

A list of all the signatures of all versions of the chromatogram is displayed in the window. The signatures are arranged according to all the signed versions. Individual versions are displayed in the list by the date that the chromatogram was saved.

Caution: Take note that this is not the date of signature but the date of saving of the signed version of the chromatogram.

Each version may be signed several times.

Each signature is represented in the list by the name of the current user and the purpose of signature. Double-clicking on the signature or clicking on the *Show Details* button will display all details about the currently selected signature:

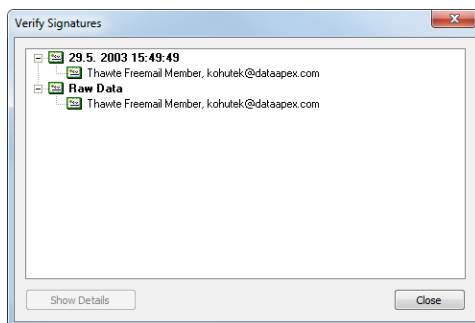


Fig 88: Verify Signatures

Signature is Valid

Confirms validity of signature (no attempt to forge the signature).

Content was not changed

Confirms validity of content (no attempt to change the signed content).

Purpose of signature

Displays the purpose of signature.

Time Stamp

Displays the date and time of signature.

3.1.10 PostRun Setting (From Chromatogram Window)

Post-processing of an already measured chromatogram. This dialog can be invoked using the *Perform Postrun...* command from the *File* menu of the [Chromatogram](#) window. When prompted, the [PostRun Setting \(From Chromatogram Window\)](#) opens, where selected actions can be manually performed on the active chromatogram.

The same function is done by the  icon (hidden by default), as long as the [PostRun Setting \(From Chromatogram Window\)](#) dialog is unset. When some actions are selected in the dialog, the  icon only performs the actions set, without opening of the dialog.

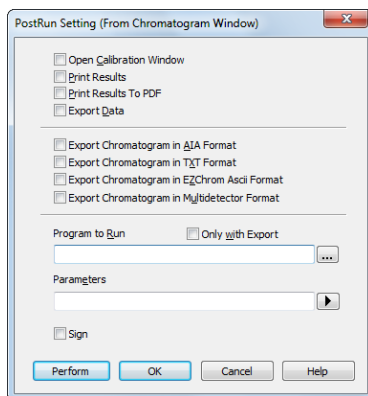


Fig 89: PostRun Setting (From Chromatogram Window)

Open Calibration Window

The created chromatogram will be automatically loaded into the [Chromatogram](#) window.

Print Results

A report will be printed according to the print style set in the [Instrument](#) window by using the *File - Report Setup* command. The default style is CHROMATOGRAM.STY.

Print Results To PDF

A report will be printed to *.PDF file according to the print style set in the [Instrument](#) window by using the *File - Report Setup* command. The default style is CHROMATOGRAM.STY.

Export Data

The created chromatogram will be exported according to settings made in the [Export Data](#) dialog (for more details see also chapter **Export Data** on pg 96).

Export Chromatogram in AIA Format

Exports a chromatogram in AIA format. This corresponds to the *File - Export Chromatogram* command in the [Chromatogram](#) window. The exported chromatogram will be saved under the same name as the source chromatogram (with the *.CDF suffix) and located in the same directory (for more details see also chapter **Export Chromatogram** on pg 140).

Export Chromatogram in TXT Format

Exports a chromatogram in the TXT format. Corresponds to the *File - Export Chromatogram* command in the [Chromatogram](#) window (for more details see also chapter **Export Chromatogram** on pg 140).

The exported chromatogram will be saved under the same name as the source chromatogram (with the *.TXT suffix) and located in the same directory.

Export Chromatogram in EZChrom ASCII Format

Exports a chromatogram in the ASC text format, which is used by the EZ CHROM program. This corresponds to *File - Export Chromatogram* command in the [Chromatogram](#) window (for more details see also chapter **Export Chromatogram** on pg 140).

The exported chromatogram will be saved under the same name as the source chromatogram (with the *.ASC suffix) and located in the same directory.

Export Chromatogram in Multidetector Format

Exports a chromatogram in the TXT format. Multiple signals will be stored in separate columns of the resulting file. Corresponds to the *File - Export Chromatogram* command in the [Chromatogram](#) window (for more details see also chapter **Export Chromatogram** on pg 140).

The exported chromatogram will be saved under the same name as the source chromatogram (with the *.TXT suffix) and located in the same directory.

Program to Run

Sets the program which will be run after the completion of the measurement.



Searches for the program which should be entered in the *Program to Run* field.

Only with Export

This item has two functions:

- The program specified in the *Program to Run* field will only be run if the *Automatic Export* checkbox has been selected.
- While manually exporting from the [Chromatogram](#) window using **Ctrl + E** shortcut this item enables the simultaneous running of the program specified in the *Program to Run* field.

Parameters

Parameters of the run program, which are normally entered in the command line behind the program name (e.g. name of loaded file, macro, etc.).

The following variables can be used in the *Parameters* field:

- Variable %f will be substituted by the name of the chromatogram at the moment of processing.
- Variable %e will be substituted by the name of the file to be exported at the moment of processing. (This does not apply when exporting multidetector chromatograms in TXT or AIA formats.)

Click the  button next to the *Parameters* field to enter the variables automatically.

For facilitating of exported chromatogram files copying you can use COPYFILES.EXE utility, located in UTILS subfolder. It allows you to copy all the exports from one chromatogram (.TXT, .PDF) at once. An utility is thoroughly described in COPYFILES.TXT file in the same folder.

Sign

Sign the chromatogram in which the changes are performed during the postrun according to the settings done in the [Sign](#) dialog. This function is only available in the Full version.

Caution: Chromatograms cannot be signed unless you have set up a user account in the [User Accounts](#) dialog. It is also impossible to sign a chromatogram which has been opened using one of the methods from the history (*Method* field of the [Open Chromatogram](#) dialog).

Use the *Perform* button to perform the actions set in this dialog, or the *OK* button to save the changes made in the dialog and leave it without performing the actions.

3.2 Edit

Sub-menu of commands for working with clipboard and tables.



Undo

Cancels adjustments in the integration table. This may be used repeatedly as the station remembers all adjustments made to tables from its initial start-up. The history of commands only applies within the framework of the table and each table maintains its own independent list of operations. The command can also be invoked using the **Ctrl** + **Z** shortcut.

Note:

Changes performed interactively in the chromatogram graph are recorded into the **Integration Table** on the [Integration](#) tab. The history of operations for the *Undo* function ceases to be accessible after switching to another table. If you wish to return to the given operation, first place the cursor into the area of the graph or switch to the [Integration](#) tab and then apply the *Undo* function.



Redo

Countermands the *Undo* command which can be invoked repeatedly. The station remembers all modifications made in tables during the last session. The history of commands applies within the framework of the table and table of each signal maintains its own independent list of operations. Commands can also be invoked using the **Ctrl** + **Shift** + **Z** shortcut.



Cut

Removes selected text and stores it in clipboard. The command can also be selected using the **Ctrl** + **X** shortcut.



Copy

Copies selected text into clipboard. The command can also be selected using the **Ctrl** + **C** shortcut.



Paste

Copies the clipboard content to the desired location (**Ctrl** + **V**). The *Paste* function automatically inputs the content of the clipboard into the desired fields of tables. If the area is larger than the content, the content will be tiled. You may make use of this characteristic if you wish to insert a larger quantity of identical fields into a table or several periodically repeating fields.

Select All

Selects all fields of table. This command can also be selected using the **Ctrl** + **A** shortcut.

Delete All	Deletes all fields of the table.
Delete	Deletes selected fields of the table. To access the edition of user columns the user must be authorized to edit methods. See the <i>Edit Method</i> checkbox in the User Accounts dialog.
Add User Column... ✓ Full version	Adds column as defined by the user.
Delete User Column... ✓ Full version	Deletes all user columns. To delete only some of the user columns use the <i>Setup Columns...</i> command. In the Setup Columns dialog then delete the selected user columns using the <i>Delete</i> button.
<i>Note:</i>	If you wish to remove a user column only temporarily, use the <i>Hide Column(s)</i> command (available from the local menu of the table or from the Setup Columns dialog).
Restore Default Columns...	Restores implicit columns, their visibility, order, and width. All user columns are hidden.
<i>Note:</i>	By contrast to the <i>Delete User Columns</i> command, user columns are only hidden and not deleted.
Setup Columns...	Sets the appearance of columns. Opens the Setup Columns dialog.
Hide Column(s)	Hides selected columns. Select the columns by left clicking the mouse on the column header. Multiple columns can be selected by pressing and holding the Ctrl key.
<i>Note:</i>	Hidden columns will not be exported from the Result Table .
Show Hidden Column(s)	Displays columns that have been hidden using the <i>Hide column(s)</i> command or Setup Columns dialog.
Caution:	When the <i>Show Hidden Columns</i> command is not active and a column is still not visible, it means that a zero width has been set for the given column. To restore the column, use the <i>Restore Default Columns</i> command.

Filter

Displays/hides duplicate rows in the **Integration Table**. For example, after a series of adjustments on the beginning of a peak has been made, it is not necessary to display its entire history, since only the last adjustment is valid. The filter attempts to combine all overlapping intervals when repeating interval operations (e.g. *Integration Interval*).

3.2.1 Add User Column ✓ Full version

Appends a new column with user-defined calculation and opens the [Add User Column](#) dialog:

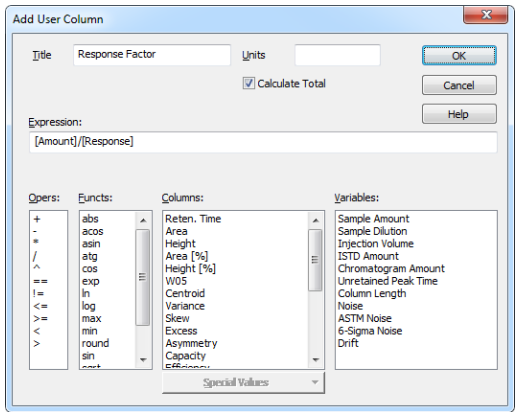


Fig 90: Add User Column

Note: To access the edition of user columns the user must be authorized to edit methods. See the *Edit Method* checkbox in the [User Accounts](#) dialog.

Title

Enter the name of the new column here. Only one line of text is allowed.

Units

Enter the units for the user column.


Calculate Total

Determines whether the *Total* row in the corresponding *User Column* should be displayed.


Expression

A calculation formula that can be entered directly or created progressively by double-clicking items in the provided *Columns*, *Ops*, *Funcs*, *Special Values* and *Variables* fields.

You will be notified if you have composed an expression that causes a recursion, after clicking the *OK* button.

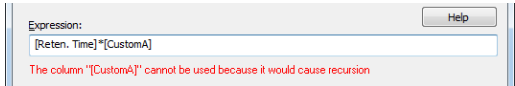

The *Details* button will appear below the *Expression* field. Click it to display a message box which will identify the columns that have caused the recursion.

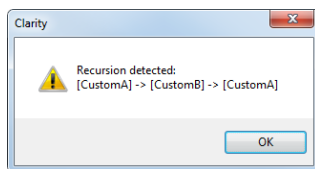

Opers

Contains a list of all mathematical operators available. Double-click an operator to insert it in the formula being compiled.

Functs

List of available mathematical functions. Insert a function into an formula by double-clicking the mouse.

Tab 6: List of available functions

Function	Explanation
abs	- Absolute value
acos	- Arcuscosinus
asin	- Arcussinus
atg	- Arcustangent
cos	- Cosinus
exp	- Exponential function
ln	- Natural logarithm
log	- Logarithm
max	- Maximum value of parameter of a function. E.g. MAX (1; 5; 3) returns value 5.
min	- Minimum value of parameter of a function. E.g. MIN (1; 5; 3) returns value 1.
round	- Rounding off
sin	- Sinus
sqrt	- Square root
switch	- SWITCH(n; H1; H2; H3;...) returns the value of the nth item Hn
tg	- Tangent

Columns

Lists all available columns as variables. Double-click a column to insert it in the formula (by default as a name of the column in square brackets).

Spec. Values

Calculated values related to the current (selected) column.

Tab 7: List of special values

Value	Description
Sum	- Sum of all values in the column
Average	- Average of the values in the column
Std. Deviation	- Standard Deviation of values in the column
Minimum	- Minimal value in the column
Maximum	- Maximal value in the column
First	- Value from the first line of the column.
Last	- Value from the last line of the column.
Previous	- Value from the previous column.
Next	- Value from the next column.
Compound	- Opens a table, from which is possible to pickup the name of compound (coined in Calibration or Integration table) and use it as a variable. (e.g. its retention time or amount)

The values of the *First*, *Last*, *Previous* and *Next* variables are not dependent on the current order of the **Result Table** and are derived from the retention time order.

Only peaks displayed in the **Result Table** are calculated. Due to this, it is possible that the values of these variables may change depending on the setting of the *Report in Result Table* item from the right side of the [Chromatogram - Results](#) tab.

Note: These functions enable the creation of inter-line calculations, e.g. calculation **RRT** - relative retention time. For calculating an individual RRT, one can input expression [Reten. Time]/[REFER\$Reten. Time], where REFER will be the name of selected reference compound.

Variables

Lists all available internal variables of the **Clarity** station. Double-click a variable to insert it in the formula.

3.2.2 Setup Columns...

Opens the [Setup Columns](#) dialog which is used to customize the table columns:

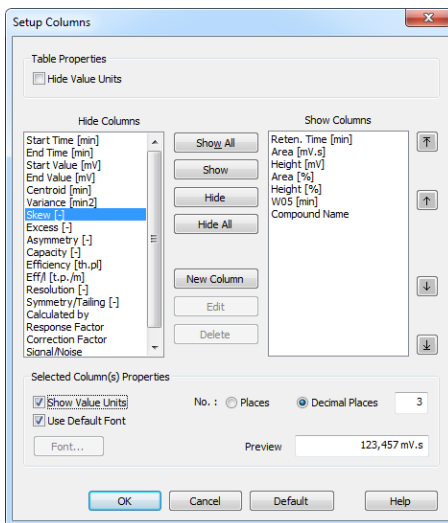


Fig 91: Setup Columns

Note: To access the edition of user columns the user must be authorized to edit methods. See the *Edit Method* checkbox in the [User Accounts](#) dialog.

Hide Value Units

Hides all value units in the whole table.

Hide Columns

Shows the list of all hidden columns.

Show Columns

Displays a list and specifications of currently displayed columns.

Show All

Displays all existing columns by transferring hidden columns from the left *Hidden Columns* list to the end of the right *Displayed Columns* list.

Show

Transfers the selected hidden column(s) from the left list to the list of displayed columns. The displayed columns are added to the *Show Columns* field in the fixed order they had in the *Hide Column* list, but may be moved forward/backwards via the arrow icons to the right of the *Show Columns* field.

Hide

Hides the selected displayed column(s) by transferring them from the right to the left list.

Hide All

Hides all columns by transferring them from the right to the left list.

New Column ✓ Full version

Appends a new column with user-defined calculation and opens the [Add User Column](#) dialog (for more details see the chapter **Add User Column** on pg 151).

Edit ✓ Full version

Modifies the selected column containing a user-defined calculation.

Delete ✓ Full version

Deletes the selected column containing a user-defined calculation.

Show Value Units

Shows the values with units for the selected column.

Use Default Font

Applies the default font, specified in the *Table Fonts - Default Fonts* item of the [User Options - General](#) dialog (available from the [Instrument](#) window by using the *View - Options* command), to the columns marked.

Font

Prescribe the font for the column selected. Invoking the command opens the [Font](#) dialog.

No. of Places

Specifies the number of valid digits to be displayed in the column selected.

No. of Decimal Places

Specifies the number of decimal places to be displayed in the column selected.

Preview

Displays a preview of the results in the selected column.

Default

Restores the default columns and their values and hides the user columns (for more details see the chapter **Edit** on pg 149).

3.3 Display

This is an entry point to a submenu of commands for displaying chromatograms.

Preview Graph	Displays/hides the orientation graph.
Zoom Out	Zooms out the current part of chromatogram. Also invoked by pressing  key on numeric keyboard or by turning mouse wheel clockwise with Ctrl pressed.
Zoom In	Zooms in the current part of chromatogram. Also invoked by pressing  key on numeric keyboard or by turning mouse wheel counter-clockwise with Ctrl pressed.
 Previous Zoom	Using the  icon or Ctrl +  shortcut on numeric keyboard restores the previously displayed cut. This function can be used repeatedly to go backwards through previous cuts. The station remembers all cuts effected during current session.
 Next Zoom	Restores the original cut abandoned by the <i>Previous Zoom</i> command. (Ctrl +  on numeric keyboard) It can be used repeatedly to go forwards through existing cuts. The station remembers all cuts effected during current session.
 Unzoom	Using the  icon or Ctrl +  shortcut on numeric keyboard cancels all cut-outs and displays the chromatogram in its original size. The original size may not always represent the whole chromatogram. If you have a fixed display range (<i>Range - Fixed</i>) set by the <i>Display - Properties...</i> command on the Time Axis or Signal Axis tabs, the chromatogram will be displayed in this range. Equally, the chromatogram may not be displayed whole if using the <i>Scale Y Mode</i> function from the Signals tab.
Tables Up	Mutually exchanges the locations of the chromatogram and the tabs with tables so that the tabs with tables are at the top and the graphs are below.
 Properties	Controls the displayed graph and chromatogram(s).

3.3.1 Preview Graph

Displays/hides an auxiliary graph. The auxiliary graph is designated for displaying the cut-out performed in the main graph. The command can also be found in the local menu invoked by right clicking above the area of the graph. The auxiliary graph is always displayed in its original size.

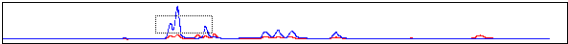


Fig 92: Auxiliary Graph

Caution: If the size of the graph pane part falls below certain dimensions, the auxiliary graph is no longer displayed, regardless of the activated *Preview Graph* command. In such a case the graph can again be displayed by shifting the splitter separating the graph from the tables or by enlarging the entire *Chromatogram* window.

Note: The original size means either the size where the chromatogram fills the entire graph or a fixed size preset by the *Range - Fixed* commands from the *Graph Properties - Signal Axis* and *Graph Properties - Time Axis* dialog.

Work with cut-outs in auxiliary graph

A cut-out is illustrated in the auxiliary graph by a dotted rectangle. A cut-out can be easily created and shifted and its dimensions changed here.

Tab 8: Manipulating preview graph using the mouse cursor

Action	Effect
Creation	Place the cursor in one of the corners of the envisaged cut-out. Click and hold the left mouse button, then draw the cursor diagonally to the opposite corner and release the button.
Shifting	Place the cursor inside the illustrated dotted rectangle. The cursor changes to  . Click and hold the left mouse button. Shift the cut-out to the required place and release the button.
Change of size	Place the cursor on the edge of the cut-out. The cursor changes to  ,  or  . Click and hold the left mouse button. Change the size of the cut-out and release the button.

Tab 9: Manipulating preview graph using the mouse wheel

Action	Effect
Wheel alone	Shifts cut-out of chromatogram up/down
Shift + wheel	Shifts cut-out of chromatogram left/right
Ctrl + wheel	Enlarges/reduces cut-out of chromatogram
Ctrl + Shift + wheel	Enlarges/reduces the cut-out of the chromatogram in the dimension of the Y-axis only.

Note: Navigation using the mouse wheel will function even when the auxiliary graph is hidden.

3.3.2 Properties

Sets the manner of the graph and chromatogram display. After selection, the [Graph Properties](#) dialog will be opened. The tabs of this dialog are very similar to the tabs of the [User Options](#) dialog described in the chapter "[Options...](#)" on pg 103. The difference is found primarily in the validity of set values. Here the set parameters are valid only for the main graph of the [Chromatogram](#) window. By contrast, parameters set in the [User Options](#) dialog are valid for all graphs of the current Instrument where the *Use User Options* checkbox has been checked.

3.3.2.1 Graph Properties - Graph tab

Except for the *Use User Options* and *Show Data Points* parameters, the tab is identical to the **Graph** tab of the **User Options** dialog described in the chapter "**Graph tab**" on pg 107.

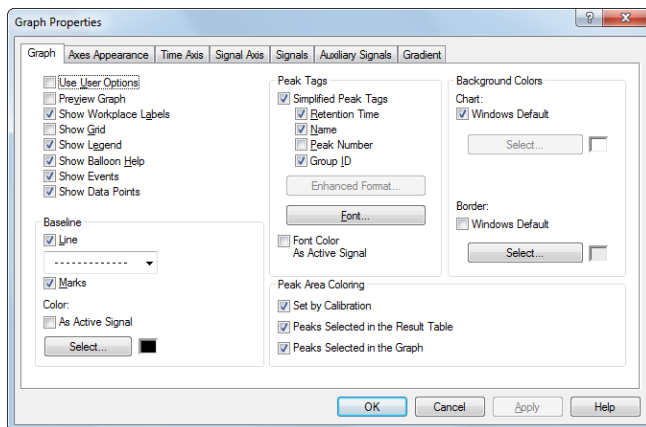


Fig 93: Graph Properties - Graph

Use User Options

All parameter settings will be taken from the global setting in the **User Options - Graph** dialog (accessible from the **Instrument** window using the *View - Options...* command).

Show Data Points

When checked, shows the individual data points in the graph. These data points may be inspected using the magnification of the area of the graph.

Note: This parameter setting is not saved in the desktop file and its selection is only valid until the **Instrument** window is closed.

3.3.2.2 Graph Properties - Axes Appearance tab

Displays graph axes. Except for the *Use User Options* parameter, the tab is identical to the [Axes Appearance](#) tab of the [User Options](#) dialog described in the chapter "[Axes Appearance tab](#)" on pg 110.

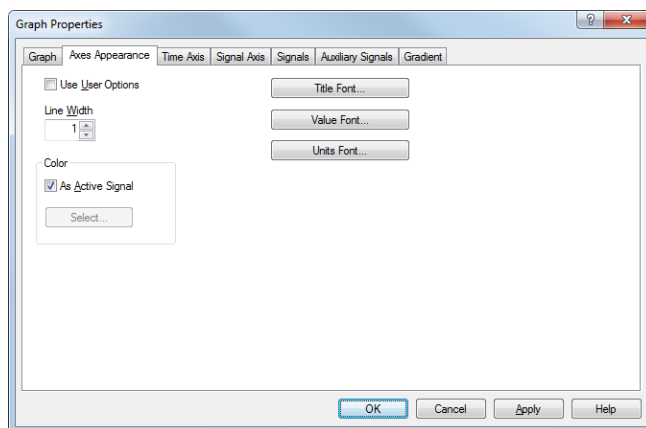


Fig 94: Graph Properties - Axes Appearance

Use User Options

All parameter settings will be taken from the global setting in the [User Options - Axes Appearance](#) dialog (accessible from the [Instrument](#) window using the *View - Options...* command).

3.3.2.3 Graph Properties - Time Axis tab

Sets the scale of the display and the units for time on the horizontal axis of graph.

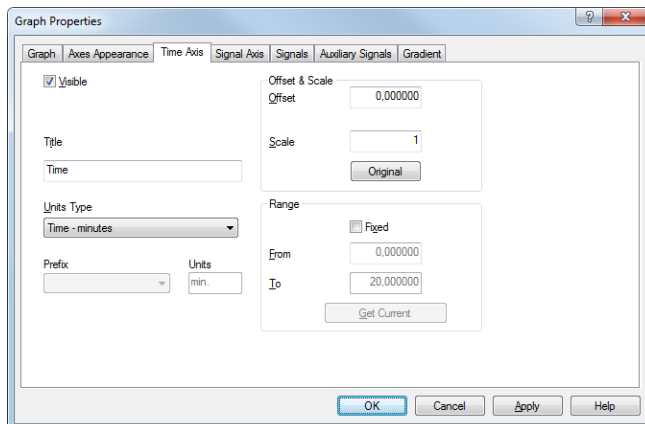


Fig 95: Graph Properties - Time Axis

Visible

Enables the display of time axis.

Title

Name of the axis.

Units Type

Type of units of the axis, displayed in square brackets. Following unit types are available:

Auto - Numbers will be displayed on the axis by the normalized basic units stated in the *Units* field. The axis is thus automatically normalized. This means that, when necessary, suffixes expressed in multiples of 1000 or fractions of the basic units are added to the basic units.

Fixed - Numbers will always be displayed on the axis in the units stated in the *Units* field without regard to size. Whenever a change of the range of the axis occurs, the numbers are not normalized and thus no prefix is added to the units.

Time (User Options) - Units will be taken from the [User Options](#) dialog.

Time - Seconds - The axis will be displayed in seconds.

Time - Minutes - The axis will be displayed in minutes and their hundredths.

Time H:M:S - The axis will be displayed in hours, minutes and seconds.

Units

By selecting *Auto* or *Fixed* in the *Units Type* field, you can enter the units in which you wish to display the horizontal axis. The conversion between the implicit units (minutes) and your own units is entered with the help of

the *Offset* and *Scale* fields. When the Units Type is switched to *Auto*, only basic units without prefixes can be entered.

Offset & Scale

Offset

Displayed data will be shifted by the value entered here.

Scale

Displayed data will be multiplied by the value entered here.

Original

Sets original values (i.e. *Offset* to 0 and *Scale* to 1).

Note: This settings applies to the graph axis only and not to the results tables. Consequently, only simplified peak tags will be shifted, while enhanced peak tags will keep the original values, as they are derived from result table.

Range

Fixed

Fixes range of horizontal axis of graph. This means that the range is determined by the *From* and *To* parameters and not by the actual size of the chromatogram.

From, To

Values for fixed range on horizontal axis.

Get Current

The *From* and *To* fields are entered by the minimum and maximum time of the active chromatogram (signal) according to the current selection.

3.3.2.4 Graph Properties - Signal Axis tab

Sets scale of display and units for signal (vertical) axis of graph. Both the signal name and signal units are exclusive to the signal displayed and are taken (by default) from the [Device Setup](#) dialog of the detector.

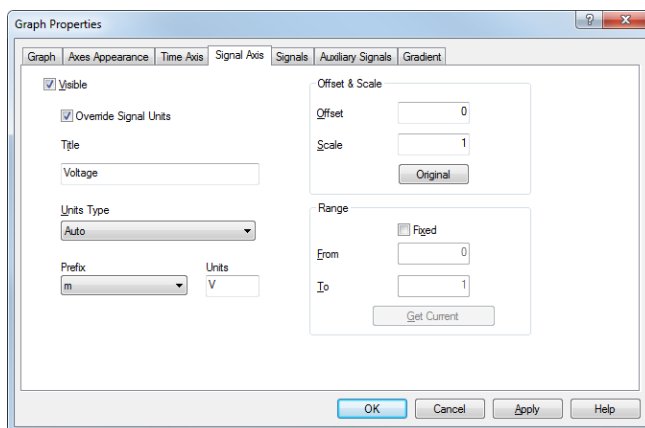


Fig 96: Graph Properties - Signal Axis

Visible

Enables the display of the vertical axis.

Override Signal Units

Allows to manually enter units which will be displayed on the y-axis of the graph. When unchecked (default), signal units set in the [System Configuration](#) dialog will be used.

Title

Name of the axis.

Units Type

Type of the units of the axis displayed in square brackets.

Fixed

Numbers will always be displayed on the vertical axis in the units stated in the *Units* field without regard to size. When changing the range of the axis the numbers are not normalized and thus no prefix is added to the units.

Auto

Numbers will be displayed on the vertical axis in the normalized basic units stated in the *Units* field. The vertical axis is thus automatically normalized. This means that in the case of necessity, prefixes expressing multiples of 1000 or fractions of the basic units are added to the basic units. For example, instead of the value *0.001 V* the value *1 mV* will be displayed.

Prefix

Sets the prefix for the *Units* if the *Auto* option for *Units Type* is selected.

Units

The type of units entered (e.g. V (Volts), RI (Refractive Index),...). The conversion between the implicit units (volts) and your own units is entered with the help of the fields *Offset* and *Scale* while the *Fixed* option for *Units Type* is selected. When *Auto* option is selected, only the basic units (without prefixes) should be entered.

Offset & Scale**Offset**

Displayed data will be shifted by the value entered here. This applies to the graph signal axis only and not to the result tables.

Scale

Displayed data will be multiplied by the value entered here. This applies to the graph signal axis only and not to the results tables.

Original

Sets original values (i.e. *Offset* to 0 and *Scale* to 1).

Range**Fixed**

Fixed range of the vertical axis of the graph. This means that the range is determined by the *From* and *To* parameters and not by the actual size of the chromatogram.

From and To

Values for fixed range on vertical axis.

Get Current

The *From* and *To* fields are filled in by the minimum and maximum voltage of the active chromatogram (signal) according to the current selection.

3.3.2.5 Graph Properties - Signals tab

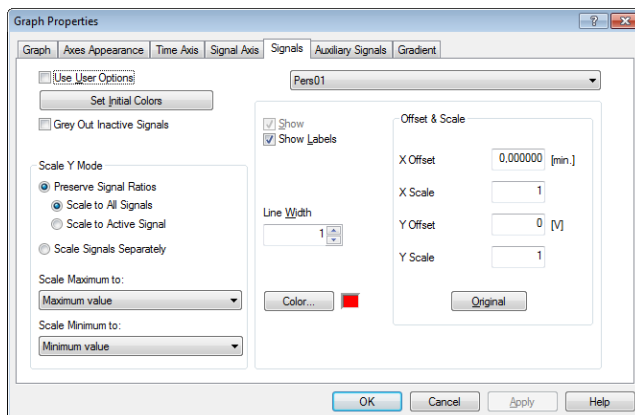


Fig 97: Graph Properties - Signals tab

Use User Options

All parameter settings will be taken from the global setting in the [User Options - Signals & Curves](#) dialog (accessible from the [Instrument](#) window using the *View - Options...* command).

Set Initial Colors

Sets colors back to those fixed at the initial installation of the station.

Grey Out Inactive Signals

When checked, all signals available in the [Chromatogram](#) window save the active one will be greyed out.

Scale Y Mode:

Automatically sets a scale on the vertical axis.

Preserve Signal Ratios

Preserves common axis for all displayed chromatograms (signals). Two alternatives may be selected:

Scale to All Signals - The scale will be set so that all the displayed signals fit into the window according to the minimum and maximum value of all displayed signals.

Scale to Active Signal - The scale will be set according to the active signal.

Scale Signals Separately

Enlarges or reduces each signal so that they fit, full height, into the graph. Using this method will eliminate the mutual proportions between the various signals and each signal will have its own scale on the vertical axis.

Scale Maximum to

The upper limit for scale for displaying the chromatogram (signal) will be derived from:

Maximum value - maximum value of signal.

Largest peak - highest integrated peak.

2nd largest peak - second highest integrated peak.

3rd largest peak - third highest integrated peak.

Maximum baseline value - maximum value not belonging to any identified peak (useful for negative chromatograms with positive spikes).

Scale Minimum to

The lower limit for scale for displaying the chromatogram (signal) will be derived from:

Minimum value - minimum value of signal.

Largest negative peak - highest negative integrated peak.

2nd largest negative peak - second highest negative integrated peak.

3rd largest negative peak - third highest negative integrated peak.

Minimum baseline value - minimum value not belonging to any identified peak (useful for chromatograms with positive peaks and negative spikes).

List of open chromatograms

Provides a list of all chromatograms (or signals when multi-detector chromatograms are in use) displayed in the right-hand section of the tab. If less than sixteen chromatograms are open, the list is supplemented by the **#X unassigned** fields, enabling the color of further chromatograms to be defined without having to open them.

Caution: All fields in the right-hand section are valid only for the chromatogram (signal) selected in this field.

Show

Displays/hides selected chromatogram. This does not close a chromatogram meaning that all the data remain valid (i.e. the results table). It is possible to hide an active chromatogram.

Show Labels

Displays/hides descriptions and lines joined to the chromatogram. Descriptions and lines joined to the working area (Workplace) remain unchanged (to hide these use the *Show Workplace Labels* command from the [Graph](#) tab).

Line Width

Width of chromatogram curve in points.

Color

Selects the color of a chromatogram

Default - A preset color will be used.

Select... - Selects user desired color.

Offset & Scale

X Offset (Y Offset)

Shifts chromatogram in currently set units.

X Scale (Y Scale)

Enlarges or reduces chromatogram in time or voltage axis. The permitted values range is between 0.01-100.

Note: The *X Offset* and *Y Offset* fields correspond to the *Move* command, and the *X Scale* and *Y Scale* fields to the *Scale* command. These are performed by right-clicking the mouse within the graph and then using the *Overlay - Move* or *Overlay - Scale* command from the context menu (for more details see the chapter **Overlay** on pg 195).

Note: The *Offset* and *Scale* operation are not stored in the current chromatogram, after opening the chromatogram again they will be reset to the default values. Only in case the chromatogram window is closed with some chromatogram opened, its offset and scale values will be restored after reopening the window.

Original

Sets original values (i.e. *Offset* to 0 and *Scale* to 1).

3.3.2.6 Graph Properties - Auxiliary Signals tab ✓ Full version

This window lets you to adjust the appearance of Auxiliary signals recorded together with the chromatogram. The tab copies part of its functionality from the [Auxiliary Signals](#) tab of the [User Options](#) dialog described in the chapter "[Auxiliary Signals tab](#)" on pg 112.

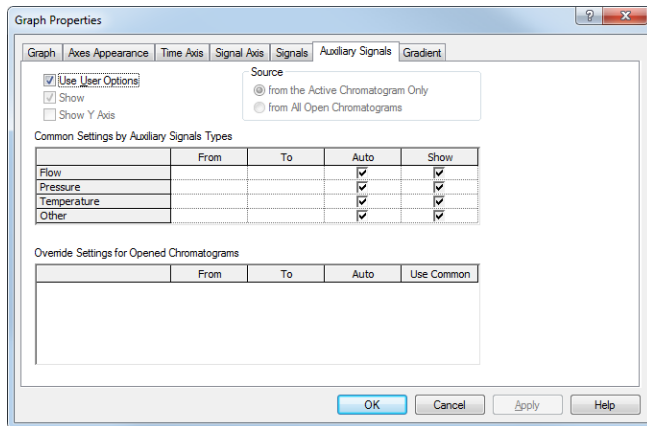


Fig 98: Graph Properties - Auxiliary Signals

Use User Options

All parameter settings will be taken from the global setting in the [User Options - Auxiliary Signals](#) dialog (accessible from the [Instrument](#) window using the *View - Options...* command).

Show

Auxiliary signals will be displayed.

Show Y Axis

Y axis will be displayed.

Source

Select for which chromatograms the auxiliary signals will be displayed.

Common Settings by Auxiliary Signals Types

Sets the auxiliary signal axis range for auxiliary detectors. When the *Auto* checkbox is checked, the auxiliary signal axis range starts at a minimal value and enlarges according to the auxiliary signal course. While unchecked, the axis range is fixed with the values entered.

Override Settings for Opened Chromatograms

Override the settings for any particular auxiliary signal from currently opened chromatogram.

3.3.2.7 Graph Properties - Gradient tab

Sets the display of gradient graph on directly controlled HPLC configurations. Except for the *Use User Options* parameter, the tab is identical to the [Gradient](#) tab of the [User Options](#) dialog described in the chapter "[Gradient tab](#)" on pg 113.

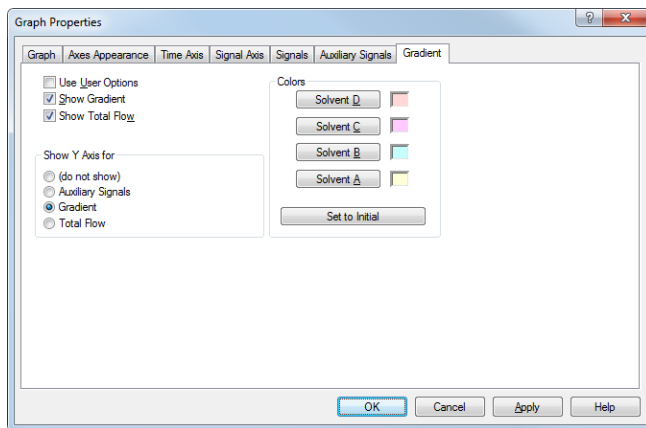


Fig 99: Graph Properties - Gradient

Use User Options

All parameter settings in the left column will be taken from the global setting in the [User Options - Gradient](#) dialog, accessible from the [Instrument](#) window using the *View - Options...* command.

3.4 Chromatogram

A submenu of commands for working with chromatograms.

	Global Peak Width	Use of this command or <i>Global Peak Width</i>  icon calculates the <i>Peak Width</i> parameter for the whole chromatogram from interactively selected interval. The <i>Global Peak Width</i> can also be set manually in the first row of the Integration Table found on the Chromatogram - Integration tab). Detailed description of the <i>Peak Width</i> parameter can be found in the chapter " Integration " on pg 183.
	Global Threshold	Use of this command or <i>Global Threshold</i>  icon calculates the <i>Threshold</i> parameter for the whole chromatogram from interactively selected interval. The <i>Global Threshold</i> can also be set manually in the second row of the Integration Table found on the Chromatogram - Integration tab). Detailed description of the <i>Threshold</i> parameter is in the chapter " Integration " on pg 1.
	Baseline	Commands for baseline correction.
	Peak	Commands for operations with peaks.
	Integration	Commands for adjustment of integration parameters.
	Separation	Commands for adjustment of separation parameters.
	Noise & Drift	Commands for evaluation of noise and drift.
	Overlay	Commands for simultaneous work with several chromatograms.
	Chromatogram list	Contains a chronological list of all currently displayed chromatograms. The active chromatogram is denoted by a point before its name. To set a chromatogram to active status, click on its name. Chromatograms can be displayed using the <i>Chromatograms...</i> command.
	Chromatograms...	Displays the Chromatogram List dialog with a list of all chromatograms/signals that are opened in the Chromatogram window.

Measure Distance	Shows the distance to the relative point in the graph. After invoking the command or clicking on the <i>Measure Distance</i> icon  (hidden by default, can be displayed from the <i>Overlay</i> toolbar icons group) the cursor changes to  and will prevent you from leaving the Graph pane. Click and drag to see the distance from the location you clicked on and the current cursor position. This distance is shown just out of the upper right corner of the Graph pane. Normally, absolute position of the cursor in the graph is shown.
Create Label	Adds labels and lines.
Remove Label(s)	Deletes labels and lines. Selected Deletes the selected label or line. From Workplace Deletes all labels and lines attached to the workplace. From Chromatogram Deletes all labels and lines attached to the active chromatogram. Remove All Deletes all labels and lines regardless of whether they are attached to the workplace or to the active chromatogram. Labels of other opened but inactive chromatograms will be retained.

Merge...

Creates multi-detector chromatograms. The command is commonly used when a multi-detector analysis was measured on one of the **CSW** stations which do not allow for the creation of multi-detector chromatograms. The procedure is following:

1. Open the chromatogram to which you would like to add further signals.
2. Prompt the *Merge* command and the [Open Chromatogram](#) dialog will open where you can select a chromatogram.
3. Select the required chromatogram and click the *OK* button.
4. Repeat the procedure to add more signals.

It is possible to create a chromatogram containing a maximum of 16 signals. Only the whole chromatograms will be merged, regardless of the number of signals selected in [Open Chromatogram](#) dialog. Chromatograms containing PDA data cannot be merged.

Set Signal Names... Edits names of signals.

3.4.1 Baseline

Processes the baseline. Individual commands can also be invoked with the icons in the *Baseline* toolbar.

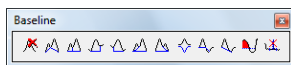


Fig 100: Baseline Toolbar

All operations involving the use of peaks are performed on an interval, which must include peak beginnings and ends of all participating peaks. Any incomplete peaks are ignored. Another essential condition of the success of an operation is that a new baseline can not intersect the signal.

Note: If done improperly, the process will either be completed without any changes to the actual baseline, effect some peaks only or only some of the beginnings and ends of peaks might be modified.

	Lock	Eliminates peaks within a selected interval from integration.
	Valley	The baseline will pass through all valleys that separate individual peaks.
	Together	An inverse operation to <i>Valley</i> - the baseline will not pass valleys, creating the perpendicular line from the valley.
	Forward Horizontal	The baseline is extrapolated horizontally forwards.
	Backward Horizontal	The baseline is extrapolated horizontally backwards.
	Front Tangent	Rider peaks on the front of the following larger peak are separated by the "tangent skimming method".
	Tail Tangent	Rider peaks on the tail of the preceding larger peak are separated by the "tangent skimming method".
	Clamp Negative	Changes the interpretation of the signal inside given area.
	Cut Negative	Discards negative peaks inside a given interval from integration without baseline modification.
	Reject Negative	Discards negative peaks inside a given interval from integration with baseline modification.

	Allow Crossing	Allows the baseline to cross the signal line. Only available in Clarity version 4.0 or higher.
	Spike Removal	Removes the spike from the chromatogram. Only available in Clarity version 4.0 or higher.

3.4.1.1 Lock

Eliminates peaks within the selected interval from integration ( icon). Only peaks lying entirely within the specified interval are eliminated this way.

3.4.1.2 Valley

The baseline will pass through all valleys separating peaks within the selected interval ( icon).

3.4.1.3 Together

Peaks inside the selected interval will be separated by a vertical line from the valley to the baseline ( icon). This operation is directly inverse to the *Valley* operation.

3.4.1.4 Forward Horizontal

The baseline is extrapolated horizontally forwards ( icon). In the selected interval, the baseline will proceed horizontally from the first identified peak beginning and will end in a vertical line at the peak end. Should the baseline intersect the signal prior to the peak end, the forward horizontal method may not be applied. The end of the previous peak is then identified before the point of intersection and the station will again attempt to project a horizontal baseline. In this instance extrapolation is understood to proceed from left to right.

3.4.1.5 Backward Horizontal

The baseline is extrapolated horizontally backwards (icon ). The baseline will proceed in the selected interval horizontally from the last identified peak end and will finish in a vertical line at the peak beginning. Should the baseline intersect the signal prior to the peak beginning, the backward horizontal method may not be applied. The end of the previous peak is then identified after the point of intersection and the station will again attempt to project a horizontal baseline. In this instance extrapolation is understood to proceed from right to left.

3.4.1.6 Front Tangent

This method (icon ) separates small peaks on the ascending part of the following larger peak by what is known as the tangent skimming method. Place the left interval borderline before the peak you still wish to separate and place the right interval borderline inside the fronting peak from which you wish to separate the rider.

3.4.1.7 Tail Tangent

This method (icon ) separates small peaks on the tail of the preceding larger peak by what is known as the tangent skimming method. Place the left interval borderline inside the peak with the riders you wish to separate. Place the right interval borderline behind the end of last rider to be separated.

3.4.1.8 Clamp Negative

Modifies interpretation (icon ) of the signal inside the specified interval by creating positive peaks from a group of negative ones inside the region (the original apexes then become valleys and *vice versa*). This works only for at least partially unresolved peaks, totally resolved negative peaks in the interval will just be eliminated (no positive peaks will be created).

3.4.1.9 Cut Negative

Negative peaks (icon ) inside the specified interval will be eliminated from integration. If an eliminated negative peak overlaps the following positive one, neither its beginning nor its end will be shifted. Use this option to eliminate negative baseline disturbances.

3.4.1.10 Reject Negative

Negative peaks (icon ) inside the specified interval will be eliminated from integration. If an eliminated negative peak is immediately followed by a positive one, the beginning and end of the latter are shifted to the original apex of the eliminated negative peak. Eliminated negative peaks are then considered to be proper part of baseline.

3.4.1.11 Allow Crossing

This function (icon ) causes the check for the baseline crossing the signal line inside the specified interval to be disabled. The peak starts and peak ends that were moved because of this check are returned to their original positions and further manual moves of peak starts and peak ends will ignore the crossing. Only available in **Clarity** version 4.0 and higher.

3.4.1.12 Spike Removal

This function (icon ) removes the spikes in the selected interval and replaces the missing area with a calculated signal line. Usage of this function needs careful placing of the start and end point of the selected interval, as **Clarity** will replace the "missing" data with a polynomial using the same slope at the start point and the end point as the original signal had. The usage of this function only available in **Clarity** version 4.0 and higher may be disabled by the use of the *Disallow Spike Filter and Spike Removal* parameter from the [GLP Options](#) dialog (for more details see the chapter **GLP Options...** on pg 23).

3.4.2 Peak

Set of commands for operations with peaks. All commands can also be invoked using icons from the *Peak* toolbar.

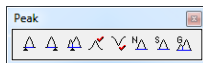


Fig 101: Peak Toolbar

	Start	Shifts the peak beginning.
	End	Shifts the peak end.
	Both	Shifts the common point of two peaks.
	Add Positive	Creates a new positive peak.
	Add Negative	Creates a new negative peak.
	Force Peak Name	Forces the peak a new name from calibration.
	Solvent Peak	Marks the peak as solvent peak.
	Peak Groups	Invokes a dialog for working with groups of peaks.

3.4.2.1 Start

Invoking this command or the *Start* icon shifts the beginning of a peak. If this shift would move the peak start ahead of the end of the preceding peak, that end will be shifted to the selected point and both peaks will be separated by a vertical line. If the beginning of the peak before shifting has the same retention time as the end point of the preceding peak and will be shifted closer to the peak apex, the previous peak's end will stay at the original point and both peaks will be separated. The peak beginning may be shifted only within the interval delimited by the apexes of the previous peak and peak being modified.

3.4.2.2 End

Invoking this command or the *End* icon  shifts the end of a peak. If this shift would move the peak end behind the beginning of the following peak, that beginning will be shifted to the selected point and both peaks will be separated by a vertical line. If the end of the peak before shifting has the same retention time as the start of the following peak and will be shifted closer to the peak apex, the following peak's beginning will remain at the original point and the peaks will be separated. The peak end may be shifted only within the interval delimited by the apexes of the following peak and the peak being modified.

3.4.2.3 Both

Invoking this command or using the *Both* icon  shifts the common point of two peaks (a valley or the separating vertical line). The point may be shifted only within the interval delimited by the apexes of the both separated peaks.

3.4.2.4 Add Positive

Invoking this command or using the *Add Positive* icon  creates a new, positive peak. Its beginning and end are defined by the interval borderlines, the apex is identified automatically. If the peak overlaps the neighboring peaks, their respective start and end points will be shifted and the peaks will be separated by vertical lines. New peaks cannot be created inside the interval of tangent skimming, nor close to the apex of an existing peak.

Note: A two-step process must be used to create a peak in the area of a monotonous signal. First, the beginning (end) must be placed at the point where the apex ultimately should be placed and the end (or beginning) placed in its final position. Only then can the peak beginning or end be shifted by the *Start* (or *End*, as the case may be) command to the required position.

3.4.2.5 Add Negative

Invoking this command or using the *Add Negative* icon  creates a new negative peak. Its beginning and end are defined by the interval borderlines, the apex is identified automatically. If the peak overlaps neighboring peaks, their respective start and end points will be shifted and the peaks will be separated by vertical lines. New peaks cannot be created inside the interval of tangent skimming, nor close to the apex of an existing peak.

Note: A two-step process must be used to create a peak in the area of a monotonous signal. First, the beginning (end) must be placed at the point where the apex should ultimately be placed and the end (or beginning) placed at its final position. Only then can the peak beginning or end be shifted by the *Start* (or *End*, as the case may be) command to the required position.

3.4.2.6 Force Peak Name

Invoking this command or using the *Force Peak Name* icon  lets the user to select a peak whose name should be changed and then opens the **Force Peak Name** dialog with a list of all peaks included in the calibration file. Alternatively, user can define its own name in the field on the top of the dialog.

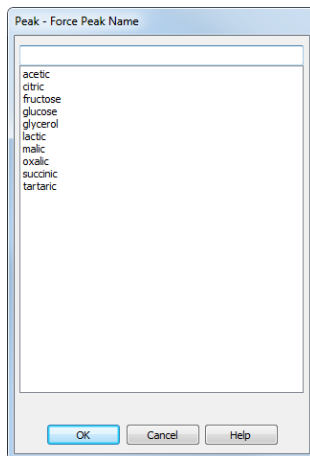


Fig 102: Force Peak Name

Selected peak name will be set to the given peak.

Note: A peak with the same name cannot appear twice in the **Result Table** - only the peak named so latest will hold its name, other peaks with the same name will become uncalibrated.

3.4.2.7 Solvent Peak

Invoking this command or the *Solvent Peak* icon  marks the selected peak as solvent peak; several peaks may be marked this way by repeatedly using the command. Solvent peak will be labeled by the letter S before the serial number (if displayed) in the **Result Table** and is excluded from the integration results.

3.4.2.8 Groups

Invoking this command or the *Group* icon  opens the *Groups* dialog used for the operations with peak groups.

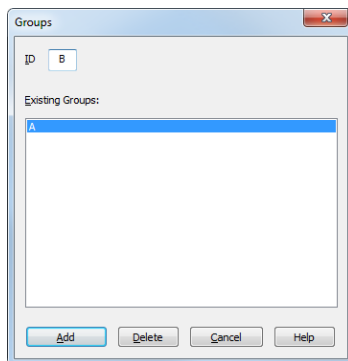


Fig 103: Groups

ID

A group is marked by a single capital letter symbol (26 letters of the English alphabet can be used). If the identifier of a group already containing at least one peak is entered here, the group will be highlighted in the *Existing Groups* list.

Existing Groups

All defined groups containing at least one peak are listed here. When a calibrated calculation is used, the name taken from the calibration file is appended after the group identifier (provided the calibration file contains the relevant group). To transfer the group identified to the *ID* field, click the corresponding line.

Add

Adds peaks to the group selected in the *ID* field or creates a new group, if the one set in the *ID* field does not exist. All peaks completely bracketed by the selected interval will be added to the group. A peak can only be included in a single group - only the last designation will thus apply.

Delete

Deletes peaks from the group selected in the *ID* field. Only peaks assigned to that group and completely bracketed by the selected interval will be deleted.

Note: If the *Group ID* field on the *Graph Properties* dialog is checked, the capital letter of a group that a peak belongs to will be indicated in the graph at the peak's description.

3.4.3 Integration

Menu for adjustment of integration parameters. All commands can also be invoked using icons from the *Integration* toolbar.

All integration parameters are executed as a part of the **Integration Table** (for more details see the chapter **Integration** on pg 212). All of these parameters are interval parameters, which means that it is possible to define the time frame of their validity and in particular to select various values for various parts of the chromatogram.

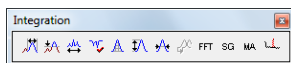


Fig 104: Integration commands in Baseline bar

	Local Peak Width	Defines minimum width of peak.
	Local Threshold	Describes the threshold of noise.
	Integration Interval	Sets the interval for the integration algorithm.
	Detect Negative	Detects negative peaks.
	Minimal Area	Sets the minimum peak area.
	Minimal Height	Sets the minimum peak height.
	Minimal Half Width	Sets the minimum peak width in half-height of the peak.
	Detector Delay	Shifts the signal on the time axis.
FFT	FFT Filter	Applies/suspends the FFT Filter (Fast Fourier Transformation) for noise reduction.
SG	Savitzky-Golay Filter	Applies/suspends the Savitzky-Golay Filter for noise reduction.
MA	Moving Average Filter	Applies/suspends the Moving Average Filter for noise reduction.
	Spike Filter	Applies a filter that removes spikes to the selected region of the chromatogram. Only available in Clarity version 4.0 or higher.

3.4.3.1 Local Peak Width

Invoking this command or using the *Local Peak Width* icon  opens the *Peak Width* dialog which defines the method of peak integration on a given time interval. **Peak Width** parameter should be set according to the narrowest peak of interest and influences the number of points used in the evaluation. The *Value* must be in the range of 0.001 - 10 minutes.

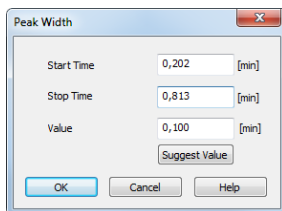


Fig 105: Peak Width

Caution: Setting an excessively high value will distort the results.

The Local Peak Width command procedure:

- You will be prompted to place the cursor lines to the beginning and the end of the interval and mark each by left-click of the mouse. The *Local Peak Width* parameter will be applied within this interval only.
- Then the *Peak Width* dialog will open in which you may adjust the graphically set interval in the *Start Time* and *Stop Time* fields and enter the required peak width into the *Value* field.
- Click the *Suggest Value* button to user is advised to interactively select the narrowest peak that still has to be integrated (again using the cursor lines). Suggested value will be transferred to the *Value* field.
- Click *OK* to apply the selected value to the specified interval.

Note: The value as well as the interval can be later edited in the **Integration Table**.

Start Time

Enables to edit the interactively set *Start Time* of the interval. The value set in the *Value* field will be applied within this interval only.

Stop Time

Enables to edit the interactively set *Stop Time* of the interval. The value set in the *Value* field will be applied within this interval only.

Value

Enables to edit the suggested value of the *Local Peak Width* parameter. The value can be suggested by **Clarity** using the *Suggest Value* button.

Suggest Value

Clicking the button will display the cursor lines again and the user will be prompted to select the interval containing narrowest peak that still has to be identified. The suggested **Peak Width** value will be transferred to the *Value* field where it can be confirmed or edited.

Note: The *Peak Width* value is not critical, so narrower peaks than the set value are also detected, but this may lead to an imprecise definition of the beginning and end of the peak and incorrect interlining of the baseline.

To adjust the **Peak Width** value for the entire chromatogram, use the *Global Peak Width* command (for more details see the chapter **Chromatogram** on pg 171).

3.4.3.2 Local Threshold

Invoke this command or press the *Local Threshold* icon  to open the **Threshold** dialog which determines the noise threshold value for the specified interval. The **Threshold** parameter determines the sensitivity of the integration algorithm while detecting the beginning and end of a peak. The height of all peaks (measured from the baseline to the apex) which you wish to detect should be at least twice this parameter. The *Value* must be in the range of $1\ \mu\text{V} - 10\ \text{V}$.

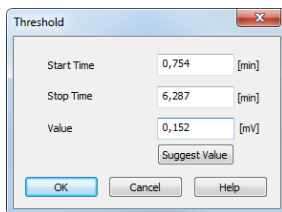


Fig 106: Threshold

The Local Threshold command procedure:

- You will be prompted to place the cursor lines to the beginning and the end of the interval and mark each by left-click of the mouse. The *Local Threshold* parameter will be applied within this interval only.
- Then the **Threshold** dialog will open in which you may adjust the graphically set interval in the *Start Time* and *Stop Time* fields and enter the required threshold value into the *Value* field.
- Click the *Suggest Value* button to interactively select the area with the "highest peaks" that still should not be integrated (using the cursor lines once more). Suggested value will be transferred to the *Value* field.

Note: Keep in mind that the suggested value will be half of these "highest peaks" because only peaks twice the size of the noise threshold will be detected.

- Click *OK* to apply the selected value to the specified interval.

Note: The **Threshold** value as well as the *Local Threshold* interval can be later edited in the **Integration Table**.

Start Time

Enables to edit the interactively set *Start Time* of the interval. The value set in the *Value* field will be applied within this interval only.

Stop Time

Enables to edit the interactively set *Stop Time* of the interval. The value set in the *Value* field will be applied within this interval only.

Value

Enables to edit the suggested value of the *Local Threshold* parameter. The value can be suggested by **Clarity** using the *Suggest Value* button.

Suggest Value

Clicking the button will display the cursor lines again and the user will be prompted to place an interval containing "highest peaks" that still should not be integrated. The suggested **Threshold** value will be transferred to the *Value* field where it can be confirmed or edited.

To adjust the **Threshold** value for the entire chromatogram use the *Global Threshold* command (for more details see the chapter **Chromatogram** on pg 171).

3.4.3.3 Integration Interval

Invoking this command or using the *Integration Interval* icon  allows to define the interval in which the integration algorithm is run. The course of the chromatogram outside of this interval will be ignored.

After using the function you are prompted to select the beginning and end point of the interval in the graph using the mouse. The starting time or ending time of the interval may be adjusted or directly set in the **Integration Table** on the *Integration* tab. When the *Integration Interval* command is not used, the integration algorithm is run on the whole chromatogram.

Caution: It is not true that the *Integration Interval* command is an opposite to the *Baseline - Lock* command. Although their logic is opposite (the first command will permit the integration in the interval and in the second will prohibit it) there is a significant difference between using the *Integration Interval* command (which directly defines where the integration algorithm will be used) and using the *Baseline - Lock* command (which only prohibits already integrated peaks). In other words: the *Baseline - Lock* command does not radically change the course of the baseline before and beyond the area with suppressed peaks.

3.4.3.4 Detect Negative

Invoking this command or using the *Detect Negative* icon  enables the detection of negative peaks in the defined interval. When the function is not active on the given part of the chromatogram, all negative peaks present are considered to be the correct course of the baseline (similar to the part of the chromatogram which uses the *Reject Negative* function).

3.4.3.5 Minimal Area

Invoking this command or using the *Minimal Area* icon  allows you to set minimum allowed area of the peak in millivolt-seconds. Peaks with areas smaller than or equal to this value are excluded from the integration.

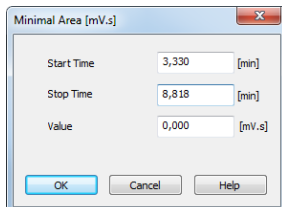


Fig 107: Minimal Area

Start Time

Enables to edit the interactively set *Start Time* of the interval. The value set in the *Value* field will be applied within this interval only.

Stop Time

Enables to edit the interactively set *Stop Time* of the interval. The value set in the *Value* field will be applied within this interval only.

Value

Enables to edit the suggested value of the *Minimal Area* parameter.

3.4.3.6 Minimal Height

Invoking this command or using the *Minimal Height* icon  allows to set the minimum height of the peak in millivolt-seconds. Peaks with heights smaller than or equal to this value are excluded from the integration. Parameters and fields in the *Minimal Height* dialog are the same as in the *Minimal Area* dialog.

3.4.3.7 Minimal Half Width

Invoking this command or using the *Minimal Half Width* icon  allows to set the minimum half-width of the peak in minutes. Peaks with *W05* parameter smaller than or equal to this value are excluded from the integration. Parameters and fields in the *Minimal Half Width* dialog are the same as in the *Minimal Area* dialog.

3.4.3.8 Detector Delay

Invoking this command or using the *Detector Delay* icon  allows to correct the time delay of the two or more detectors connected in a series. This command cannot be used on the signal of the first detector (it is inactive), but only on the second or any further signal. You are advised to point and click the cursor on the desired peak of the active signal and then on a corresponding peak in the first signal.

3.4.3.9 FFT Filter

Invoking this command or using the *FFT Filter* icon  allows to apply the **FFT Filter** (*Fast Fourier Transformation*) for noise reduction in the selected interval. If more than one type of the filter is set on the same time interval, only the last used filter on that interval is actually used.

3.4.3.10 Savitzky-Golay Filter

Invoking this command or using the *Savitzky-Golay Filter* icon  allows to apply the **Savitzky- Golay Filter** for noise reduction in the selected interval. If more than one type of the filter is set on the same time interval, only the last used filter on that interval is actually used.

3.4.3.11 Moving Average Filter

Invoking this command or using the *Moving Average Filter* icon  allows to apply the **Moving Average Filter** for noise reduction in the selected interval. If more than one type of the filter is set on the same time interval, only the last used filter on that interval is actually used.

3.4.3.12 Spike Filter

Invoking this command or using the *Spike Filter* icon  allows to remove the spikes on the chromatogram signal. The parameter of the command (that can be edited from the Integration table) describes how wide spikes will be removed (their missing points replaced by a calculated value). The usage of this function only available in **Clarity** version 4.0 and higher may be disabled by the use of the *Disallow Spike Filter and Spike Removal* parameter from the [GLP Options](#) dialog (for more details see the chapter [GLP Options...](#) on pg 23)..

3.4.4 Separation

3.4.4.1 Valley To Valley Slope

Specifies the maximum slope of the baseline. If the baseline slope is smaller than the specified value, the separating perpendicular line is canceled and the baseline passes through the valley. The default value of 0 prevents the baseline from passing through the valley.

3.4.4.2 Tangent Area Ratio

Constitutes the first condition imposed on tangential separation. The condition is satisfied if the ratio of areas of the main and the separated peak exceeds the specified value.

The default value of 0 prevents tangential separation. If there is no *Tangent Slope Ratio* item defined in the **Integration table**, its value is considered to be 1.5.

3.4.4.3 Tangent Slope Ratio

This is the second condition imposed on tangential separation. The condition is satisfied if the ratio of slopes of the second and the first peak exceeds the specified value. The slope ratio is defined as the tangent to the first peak at its end point divided into the slope of the line connecting the beginning and end of the second peak.

The default value of 0 prevents tangential separation. If there is no *Tangent Area Ratio* item defined in the **Integration table**, its value is considered to be 0.1.

3.4.5 Noise & Drift

The *Noise* (all variants) and *Drift* parameters are set as intervals. All values can be utilized in custom calculations in **User Columns**.

Result Table (ESTD - Wine_Sample - UV detector)						
Noise (1,35-4,85 min.): 960,7550 [mAU], ASTM Noise (1,92-4,90 min.): 245,9011 [mAU], 6-Sigma Noise (1,76-5,90 min.): 857,1555 [mAU]						
	Reten. Time [min]	Response	Amount [g/l]	Amount [%]	Peak Type	Compound Name
2	1,063	58,456	0,028	0,9	Ordnr	oxalic
3	1,703	3808,667	0,246	8,1	Ordnr	citric

Fig 108: Noise & Drift above the Result table

The noise evaluation determines the amount of noise in the specified interval of the chromatogram. The interval for any type of noise evaluation is set in the **Integration Table**.

After invoking the particular *Noise Evaluation* command, select part of the chromatogram that contains noise only (ideally part of the baseline). In case several intervals for the same noise calculation method are set, only the last one is valid. Previous entries can be filtered from the **Integration Table**, so that only the last entry will be displayed. Set the interval to zero to display the noise over the whole length of the chromatogram.

If the interval exceeds the length of the actual chromatogram, only the valid part of the interval will be taken into account.

The noise value will be displayed above the **Result Table** along with appropriate signal units and the interval, from which it was calculated.

Note: It is possible to calculate the noise using all three methods - linear interlining (*Noise Evaluation*), *ASTM Noise Evaluation* and *6-Sigma Noise Evaluation* at the same time.

Current noise during acquisition can also be displayed in the upper left corner of the graph in the **Data Acquisition** window.

Note: The **Noise** in the **Data Acquisition** window is calculated using the cubic interlining to reduce the influence of the trend to its values.

3.4.5.1 Noise Evaluation

The **Noise** value for the *Noise Evaluation* command is calculated using the linear interlining on the interval defined in the **Integration Table**. If there is a peak in this interval, it will influence the noise values.

3.4.5.2 ASTM Noise Evaluation

The **Noise** value for the *ASTM Noise Evaluation* command is calculated as an average from values counted for smaller intervals. The procedure goes as follows:

- The selected time interval is divided into parts of the same length (30s by default, may be changed in the **Integration Table**, *Value* column). The remaining part creates its own interval.
- Noise is counted for each of these intervals using the linear interlining and the least squares method.
- The values gained from these intervals are then averaged.

3.4.5.3 6-Sigma Noise Evaluation

The **Noise** value for the *6-Sigma Noise Evaluation* command is calculated from the whole interval set in the **Integration Table**. The value gained is calculated as a six times the square root of least square method for each point in the whole interval divided by the number of points:

$$Noise = 6 * \sqrt{\left(\frac{\sum_{i=1}^n (E_i - f(E_i))^2}{n - 1} \right)}$$

where n is the number of acquisition points, E_i is the signal value of the i -th point and $f(E_i)$ is the value corresponding to the linear regression performed on the selected interval at the i -th point.

3.4.5.4 Drift Evaluation

The drift evaluation determines the drift between the beginning and the end point of the specified interval, which is set in the **Integration Table**. After invoking the *Drift Evaluation* command, select part of the chromatogram where the **Drift** should be calculated. If more then one interval has been set, the latest one will be applied. Previous entries can be filtered from the **Integration Table**, so that only the last entry will be displayed. Set the interval to zero to get the drift over the whole length of the chromatogram.

If the interval exceeds the length of the actual chromatogram, only the valid part of the interval will be taken into account.

3.4.6 Overlay

Includes commands for working with several chromatograms simultaneously. The maximum number of displayed chromatograms is practically unlimited, the actual limit can be set in the [User Options](#) dialog. Most of the commands identified below can be invoked by the icons found in the [Overlay](#) toolbar.



Fig 109: Overlay Toolbar

	Move	Moves the active chromatogram.
	Scale	Scales the active chromatogram up or down.
	Original	Cancels all shift and scaling operations.
	3D View	Displays chromatograms in a three-dimensional perspective.
	Clear 3D	Restores the original perspective.
	Properties...	Opens the Graph Properties dialog on the Signals tab.
	Mathematics...	Performs mathematical operations with chromatograms.

The *OVERLAY* mode can be activated using the *File - Overlay Mode* command, double-clicking on the **Overlay** inscription in the status bar or clicking on the icon. An active *OVERLAY* mode is indicated by the depressed icon in the *Overlay* toolbar, highlighted **Overlay** inscription in the status bar and by the checkmark by the *Overlay Mode* in the *File* menu.

Note: Too many chromatograms opened at once may (especially on slower computers) notably slow down the response of the computer.

3.4.6.1 Move

Invoking this command or clicking the *Move* icon  enables you to move the active chromatogram. To use the function, left-click the mouse and hold down while dragging the chromatogram to the new location, then release the button. Exact values of the shift or the possibility to return to the original location are found on the [Graph Properties - Signals](#) tab (accessed by using the *Chromatogram - Overlay - Properties...* command or the *Signal Properties* icon ). In the shifted chromatogram (signal) all commands from the *Chromatogram* menu other than those in the *Overlay* group are disabled and **Result Table** and peak labels are not shown.

Caution: The performed moving cannot be stored in the original chromatogram. The only possible way to save such chromatogram is to make a copy of the modified chromatogram by the *Overlay - Mathematics - Copy* command and save the copy. After finished *Move* operation, all the values in the **Result table** will disappear as they are no longer valid.

Note: Hold **Shift** while clicking the menu command or  icon to apply the move to all signals from the current chromatogram. Hold **Ctrl** while moving the chromatogram, to limit the moving direction to vertical or horizontal.

3.4.6.2 Scale

Invoking this command or using the *Scale* icon  allows you to change the scale of the active chromatogram. After invoking the command, left click the mouse and hold it while dragging the chromatogram to the new scale and then release it. The chromatogram will be drawn to the new scale when released. Exact values of the scale operation or the possibility to return to the original state are found on the [Graph Properties - Signals](#) tab (accessed by using the *Chromatogram - Overlay - Properties...* command or the *Signal Properties* icon ). In the scaled chromatogram (signal) all commands from the *Chromatogram* menu other than those in the *Overlay* group are disabled and **Result Table** and peak labels are not shown. The scaled chromatogram is marked by the *MODIFIED* inscription in the title bar.

Caution: The performed scaling operation cannot be stored in the original chromatogram. The only possible way to save such a chromatogram is to make a copy of the modified chromatogram by the *Overlay - Mathematics - Copy* command and save the copy. After finished *Scale* operation, all the values in the **Result table** will disappear as they are no longer valid.

Note: Hold **Shift** while clicking the menu command or  icon to apply the scaling to all signals from the current chromatogram. Hold **Ctrl** while moving the chromatogram, to limit the moving direction to vertical or horizontal.

3.4.6.3 Original

Invoking this command returns the chromatogram to its original location and scale. The initial state is the setting in which the scale is unchanged (**Scale** = 1) and the chromatogram is not shifted from the start (**Offset** = 0).

Note: Hold **Shift** while clicking the menu command to apply the change to all signals from the current chromatogram.

3.4.6.4 3D View

Invoking this command enables to display the chromatogram(s) in three-dimensional perspective. After prompting a pair of lines that serve to set the angle and depth of the 3D rendering is shown. Click the mouse button to uniformly space out all displayed chromatograms along the specified line segment. The actual spacing can be modified by changing the X (Y) *Offset* parameters on the [Graph Properties - Signals](#) tab.

3.4.6.5 Clear 3D

Invoking this command clears the 3D representation mode and restores the original display mode. All chromatograms are restored to their original positions (corresponding to the 0 values of the X (Y) *Offset* parameters). These values can be set for each chromatogram (signal) separately using the *Original* button on the [Graph Properties - Signals](#) tab.

3.4.6.6 Properties...

Invoking this command or using the *Signal Properties* icon  allows you to set the explicit shift and scale parameters for the active chromatogram. Invoking the command opens the [Signals](#) tab of the [Graph Properties](#) dialog (for more details see the chapter **Graph Properties - Signals tab** on pg 166). All set values refer only to the active chromatogram. If any parameter from the *Offset & Scale* group differs from the default value, the baseline, all commands from the *Chromatogram* menu other than those in the *Overlay* group are disabled and **Result Table** and peak labels are not shown.

3.4.6.7 Mathematics... (Mathematical Operations dialog)

Invoking this button or using the *Mathematics* icon  enables to perform mathematical operations with chromatograms. After selection, the *Mathematical Operations* dialog will be displayed. It is necessary to work in *Overlay* mode to have this option available. Mathematical operations are only performed with chromatograms (signals) that are currently open in the *Chromatogram* window.

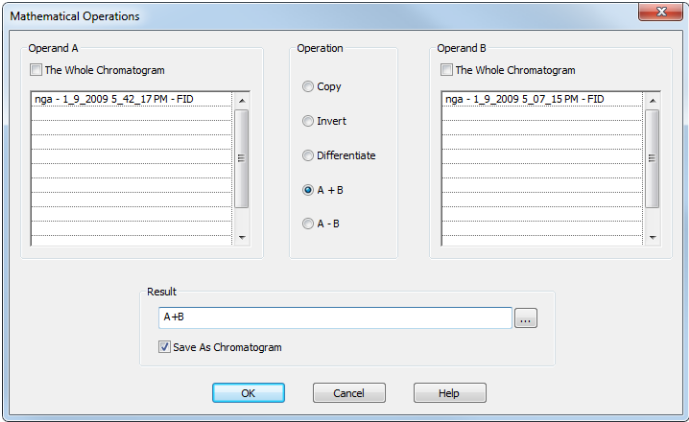


Fig 110: Mathematical Operations

Operand A (Operand B)

Select up to 16 signals in **Operand A** section in the listbox on the left. The whole chromatogram (all of its signals) will be included when the *The whole Chromatogram* checkbox has been checked. When unchecked, you may select individual signals (and change their order) from one chromatogram, or select signals from various chromatograms. This is done by simply clicking the row intended for editing and using the arrow icon  to the right to select the particular chromatogram or signal.

Operand B is selected analogically to the **Operand A**, if appropriate.

Operation

Type of required operation. The following operations are possible:

Tab 10: List of available mathematical operations

Operation	Effect
<i>Copy</i>	- Copies chromatogram (Selection of the Operand A only).
<i>Invert</i>	- Inverts signal or chromatogram (Selection of the Operand A only).
<i>Differentiate</i>	- Calculates first derivation (Selection of the Operand A only).
<i>A+B</i>	- Sum of signals or chromatograms.
<i>A-B</i>	- Difference of signals or chromatograms.

Result

This is the selection of the name under which the mathematical result will be displayed. The name offered is implicitly derived from the performed mathematical operation - e.g. $A+B$.



Searches for the file and directory that you would like to enter into the **Result** field.

Save As Chromatogram

If this checkbox is checked, the mathematical result will automatically be saved as a new chromatogram.

If this field is left unchecked, the mathematical result will only be displayed as a new curve, which you may save using the *Save* or *Save As* command from the *File* menu.

In the full version ☒ Full version, the resultant chromatogram contains information about the manner of its creation in [Chromatogram Audit Trail](#) described on pg 327.

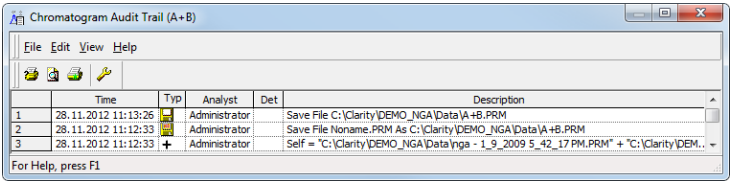


Fig 111: Chromatogram Audit Trail

3.4.7 Chromatograms

Displays the **Chromatogram List** dialog with a chronological list of all opened chromatograms/signals.

Note: Active signal is highlighted in bold letters, hidden signals are displayed in gray color.

The maximum number of signals that can be simultaneously displayed in the **Chromatogram** window ranges from 2 to 10 000 (the default setting is 20) is determined by the *Maximum Chromatograms in Overlay* item from the **User Options - General** dialog.

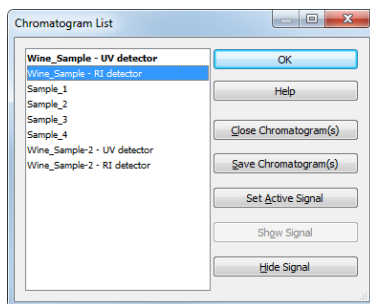


Fig 112: Chromatogram List

Close Chromatogram(s)

Closes the selected chromatogram(s) and all of its(theirs) signals.

Save Chromatogram(s)

Saves the entire selected chromatogram(s).

Set Active Signal

Sets the selected chromatogram (signal) as active.

Note: You can also double click a signal in the list to make it active.

Show Signal

Displays selected signals.

Note: The command will only be active if the selected signal has previously been hidden.

Hide Signal

Hides selected signals.

3.4.8 Create Label

Adds labels and lines.

Text

Appends a text label. Select the command and click on the desired location to open the **Text Label** dialog. The **Text Label** dialog may also be invoked by double-clicking on the actual text label to modify it.

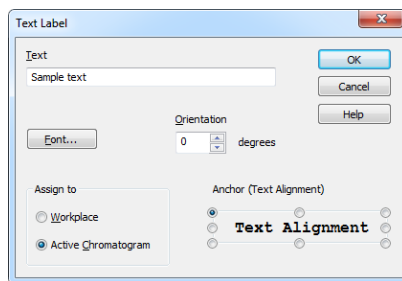


Fig 113: Text Label

Text

The text of the user's label is entered/edited on this line.

Font...

Selects the font, size and color of characters. Invoking the command opens the standard **Font** dialog.

Orientation

Select the orientation of the label text in degrees counted counterclockwise (e.g. enter 0 for horizontal or 90 for vertical orientation). Allowed values are 0 - 359.

Note: You can use *Orientation* field only when using **TrueType** fonts.

Delete

Deletes the selected label.

Assign to

Workplace

The label will lie, unattached to any specific chromatogram, in the background and will be saved in the current desktop.

Active Chromatogram

The label will be attached and become integral part of the chromatogram. When the chromatogram is scaled up or down, the label will be anchored to the nearest chromatogram point.

Anchor (Text alignment)

Defines the method of label anchoring. One of eight points on the label border or the label center may be selected as the anchor point. The label

is then firmly "bound" by that point to the nearest chromatogram point (or to the graph border when the label is attached to the **Workplace**). This can be of particular use when scaling or shifting a chromatogram where labels supplement peak identification etc.

A label can be shifted and/or scaled in the same way as in usual drawing applications. Left clicking on a label displays the controlling frame and handles for sizing it. The font size will change with the frame size. To move the frame containing the label, click inside the frame and drag.

Line

Adds connecting lines. To create a line label, invoke the command and then left click and hold the button on the desired starting point of the line, drag to the desired end point and release. The **Line Label** dialog will open to determine the properties of the line. The dialog can also be called by double-clicking on the actual line label to change its appearance.

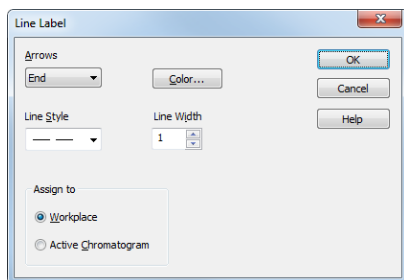


Fig 114: Line Label

Arrows

Selects optional line end:

None - No arrows will be placed on the line.

Beginning - An arrow will be added to the line beginning.

End - An arrow will be added to the line end.

Both - Both line ends will be provided with arrows.

Color...

Selects the line color.

Line Style

Defines the line style. If the *Line Width* parameter is greater than 1, the line will be solid. Possible variants for the *Line Style* are solid line, broken line, dotted line, dashed and dotted line or dashed and double dotted line.

Line Width

Defines the line width in pixels (1 to 6). Lines with the width greater than 1 have to be solid.

Assign to**Workplace**

The label will lie, unattached to any specific chromatogram, in the background and will be saved in the current desktop.

Active Chromatogram

The label will be attached and become integral part of the chromatogram. When the chromatogram is scaled up or down, the label will be sized along with the chromatogram (both ends are anchored to two nearest chromatogram points).

A line can be shifted or sized in the same way as in drawing applications. Left clicking on a line will display the controlling frame and handles the sizing of the line. To shift the line, left click and hold anywhere on the line except the end points, drag to the desired location and release.

3.4.9 Set Signal Names

The *Set Signal Names* command is used to invoke the [Set Signal Names](#) dialog where existing names of signals can be edited.

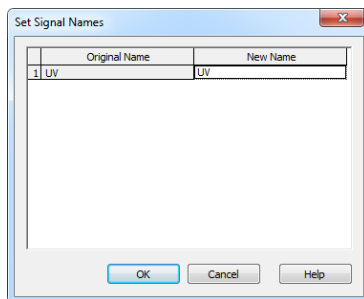


Fig 115: Set Signal Names

If the signal has no allocated name, the name **Signal n** (where n is the serial number of the signal in the chromatogram) will be displayed in the original name column.

Note: By default, individual signals are named in regards to the name of the detector used or upon import or connection by implicit names (1 .. 16). The default names of the signals can be changed in the [System Configuration](#) dialog by invoking the *Setup* button for the particular detector.

3.5 Method

An entry point to commands for modifying chromatograms and copying methods.

Calculation	Displays the parameters of chromatographic results calculation.
Integration	Displays integration parameters.
Measurement	Displays the conditions under which the chromatogram was created.
Save as Template...	The command saves the method used for the measurement of active chromatogram as template method, i.e., a method that can be used in the Instrument window for other analyses. Invoking the command displays the Save As dialog, where the name of the new template method should be specified. By default, template methods will be stored to the current project directory, but the path can also be changed. The method is a component of the the chromatogram file (files with the *.PRM suffix). The template method will be stored in a separate file with the *.MET suffix. It is not allowed to save the method into a file currently opened on the Instrument. Recommended workflow is to save template with different name, then open it on the Instrument.
Set Model 	Marks the method of the active chromatogram as model method. Any model method can be copied to another chromatogram by the <i>Copy from Model</i> command. The model method selection is valid until the Chromatogram window is closed or another model method is selected by the <i>Set Model</i> command.
Copy from Model 	Copy a Model Method to another chromatogram. Prior to invoking the command a method must have been marked as model method by the <i>Set Model Method</i> command. The model method can then be applied on any chromatogram opened in the Chromatogram window by using the <i>Copy from Model</i> command.

Copy from Template Method...

Copies Template Method into active chromatogram. Use this command to invoke the [Open Method](#) dialog (described in the chapter "**Open Method (Sequence, Chromatogram, Calibration)**" on pg 45.), which will display a list of all **Template Methods** from current project.

Copy from Chromatogram...

Copies the method from a selected chromatogram to active chromatogram. Use this command to open the [Open Chromatogram](#) dialog (described in the chapter "**Open Method (Sequence, Chromatogram, Calibration)**" on pg 45.), in which you can select the required chromatogram.

3.5.1 Calculation

Selects the method for calculating the results of the chromatogram. Once selected, the **Results** tab will be displayed in the lower part of the **Chromatogram** window. In its right-hand section it is possible to define the method of results calculation.

The parameters stated here are common for all signals when a multi-detector chromatogram is being used. The displayed values come from the template method; they are filled in on the **Calculation** tab of the **Method Setup** dialog as a part of the method according which the chromatogram is measured.

Calibration File (Peak Table)

Name of the calibration file used to obtain calibrated calculations.

Open with stored calibration

When checked, the chromatogram will be opened with calibration stored within the chromatogram file when it is opened next time. To open the chromatogram with the stored calibration this way it is necessary to open the chromatogram again using the **Open Chromatogram** dialog - it is not enough to just close and open the **Chromatogram** window. Similar behavior may be set directly when the chromatogram is created from **Sequence** window or **PostRun Setting** dialog.

Set...

Invokes the **Open Calibration** dialog where the calibration file may be selected.

None

Closes the opened calibration file and sets the calculation type in the **Calculation** field to *Uncal* (uncalibrated).

View

Opens the current calibration file in the **Calibration** window.

The screenshot displays the 'Results' tab of the 'Chromatogram - Results' window. On the left, a 'Result Table (ESTD - Wine_Sample - UV detector)' is shown with columns for Retention Time, Response, Amount (g/l), Amount (%), Peak Type, and Compound Name. The table lists 12 peaks and a total. On the right, the 'Calculation' settings are configured for 'Wine' with 'ESTD' as the calculation method. The 'Open with stored calibration' checkbox is checked. The 'Report in Result Table' section shows 'All Peaks in Calibration' selected. The 'Unidentified peaks' section shows 'Response Base' and 'Area' selected. The 'Scale' section shows 'Use Scale Factor' checked with a scale factor of 1. The 'Units' section shows 'ul' selected. The 'Amount (g/l)' and 'ISTD1 Amount' fields are set to 0. The 'Inj. Volume [mL]' is set to 0.1 and 'Dilution' is set to 1. The bottom of the window shows a tabbed interface with 'Results' selected.

	Reten. Time [min]	Response	Amount [g/l]	Amount [%]	Peak Type	Compound Name
2	1.063	58.456	0.028	0.9	Ordnr	oxalic
3	1.703	3808.662	0.246	8.1	Ordnr	citric
4	1.917	8486.434	0.440	14.5	Ordnr	tartaric
5	2.800	3331.770	1.146	37.8	Ordnr	malic
9	4.660	2730.333	0.243	8.0	Ordnr	succinic
10	5.050	2972.323	0.745	24.6	Ordnr	lactic
12	6.847	2304.521	0.185	6.1	Ordnr	acetic
	Total		3.033	100.0		

Fig 116: Chromatogram - Results

Calculation

Selects the type of calculation. Available options are shown **Tab 11** on pg **208**.

Tab 11: List of available calculation types

Calculation type	Way of calculation
<i>Uncal</i>	- Uncalibrated calculation without use of calibration file. The result of the integration is a percentage representation of individual areas and heights of peaks in the chromatogram. When using this option, all other sections on this tab will be inactive and dimmed.
<i>ESTD</i>	- External standard calculation. The result will be both a percentage and absolute representation of individual substances and their names.
<i>ISTD</i>	- Internal standard calculation. The result will be both a percentage and an absolute representation of individual substances and their names.
<i>NORM</i>	- External standard calculation with add-on checks. The result will be both a percentage and an absolute representation of individual substances and their names. In case any peak found in the Result Table is not present in a calibration and/or the <i>Amount</i> parameter differs from zero, the calculation of the <i>Peak Type</i> for all rows in the Result Table will be <i>Error</i> and the <i>Amount [%]</i> values will not be calculated.
<i>STDADD</i>	- Standard Addition calculation. The result will be both a percentage and an absolute representation of individual substances and their names.
<i>Note:</i> If only the <i>Peak Type</i> field for selected rows of the Result Table show the <i>Error</i> code, than the corresponding compounds do not have the <i>Amount</i> value filled in the calibration file.	

Report in Result Table

This field determines which peaks will be displayed in the **Result Table** and included in the sum for the **Total** row in the case of calibrated calculations.

All Peaks

All integrated peaks in the chromatogram will be displayed.

All Identified Peaks

Only peaks identified in the calibration file will be displayed.

All Peaks in Calibration

All peaks from the calibration file will be displayed without regard to how many peaks the chromatogram contains and how many were

identified. This provides what is known as a fixed table format given by the calculation of peaks in the calibration file.

Hide ISTD Peaks

The peaks identified as the internal standards will not be displayed in the **Result Table**.

Note: The command will be active only when the *ISTD* calculation type has been selected.

Unidentified Peaks

Response Base

Selects a base for calibrated calculation of unidentified substances.

Area - Calculations will be derived from peak areas.

Height - Calculations will be derived from peak heights.

Response Factor

Coefficient used for amount calculation for unidentified substances.

Amount is set as area (height) of peak multiplied by this coefficient.

This value serves in fact as an artificial single-point calibration for unidentified substances.

Scale

Area for entering the auxiliary calculation parameters.

Use Scale Factor

The whole area is activated by checking this checkbox.

Scale Factor

All values of the *Amount* column in the **Result Table** will be multiplied by the stated value.

Units

Units for application of scale. The original units taken from the calibration file are replaced by the units defined here in the results table.

Amount

Amount of the sample. This value is used as a basis for the calibrated calculation of the percentage representation of individual substances for *ESTD* and *ISTD* methods.

When the *ISTD* method is used, enter the amount of the sample without the internal standard. State the parameter in units used in the calibration file (*Units - Compound* parameter in the [Calibration Options](#) dialog). The implicit value is 0. In this case the percentage representation is calculated from the *Amount* sum in the *Total* row.

Note: If the *Amount* parameter is set to zero, the percentage representation of individual values will be calculated from the sum of all calibrated amounts (including substances calibrated using the *Uncal. Response* parameter). Otherwise the calculation is based on the *Amount* value entered.

ISTD Amount

Specifies the amount of internal standard (ISTD1) added when the results are calculated using the *ISTD* method. Similar settings are available for the ISTD2 - ISTD5 in the [Edit ISTD Amounts](#) dialog accessible through the  button.

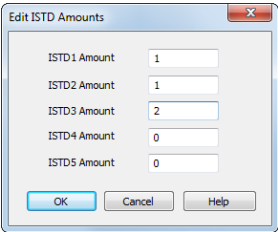


Fig 117: Edit ISTD Amounts

[Edit ISTD Amounts](#) dialog allows to set the ISTD Amounts for all internal standards used in particular chromatogram. It is also accessible from the [Single Analysis](#) dialog and amounts can be entered from the [Sequence](#) window (settings in the *ISTD1 Amount* - *ISTD5 Amount* columns, only *ISTD1 Amount* column is displayed by default). Enter the parameters in units used in the calibration file (*Units* - *Compound* parameter in the [Calibration Options](#) dialog). The implicit value is 0. If the same amount of the internal standard is used in all used calibration standards and unknown samples, this amount need not be entered and can be left at 0 in both calibration and chromatograms (sequence).

Caution: When the amount of the internal standard is entered either in the calibration file (**Amount** column) only or in the unknown sample (*ISTD Amount* parameter) only and zero will be left in the other place, the **Result Table** will not display any results and the header will show an inscription stating the cause of the problem. The compound will be also omitted from the **Summary Table**.

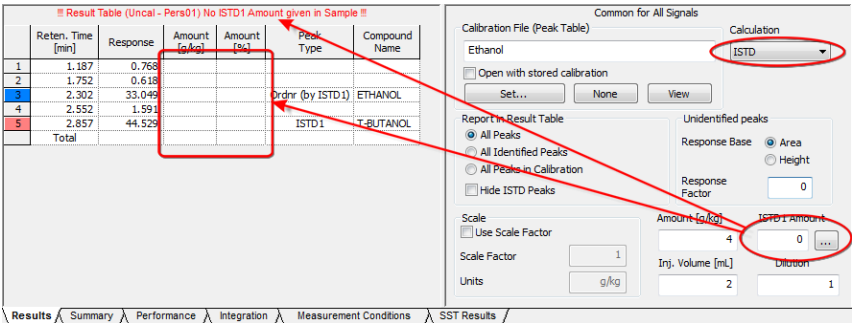


Fig 118: Chromatogram - Calculation - ISTD settings

Inj. Volume

Specifies the size of the injection. The response of all substances calculated using the calibrated calculation is corrected by the ratio of this value to the value of *Default Injected Volume* field set in the [Calibration Options](#) dialog. *Inj. Volume* parameter will influence the calculation only after the *Default Injected Volume* value is set. The implicit value is 0.

Note: The *Inj. Volume* parameter can be applied for injected amount correction while using autosamplers unable to determine precisely the quantity to be injected, but fully capable to determine (ex-post) the quantity which has been injected.

Dilution

Specifies the dilution of the sample. Each calibrated amount in the **Amount** column of the **Result Table** is multiplied by this value. The implicit value is 1.

3.5.2 Integration

Invoking this command switches the results panel of the **Chromatogram** window to the **Integration** tab, which contains the **Integration Table**. When using a multi-detector chromatogram, the tab always displays the **Integration Table** for the active signal. The displayed values are primarily taken from the **Integration** tab of the template method, according to which the chromatogram was measured, and will have integration adjustments added caused by operations performed in the **Chromatogram** window.

Integration table

The integration table records all operations performed on the baseline and peaks.

Chromatogram Operation

Selects the operations to be performed on the chromatogram. Clicking the right border of the field displays a list of all possible operations (identical to the list of commands from the *Chromatogram - Baseline, Peak, Integration, Separation* and other sub-menus in the **Chromatogram** window).

Integration Table (Sample_1, Integration Algorithm: 1.0)				
Chromatogram Operation	Grp.	Time A [min]	Time B [min]	Value
Global Peak Width				0.500 min
Global Threshold				2.0000 mV
Baseline - Valley		7.860	48.066	
Baseline - Lock		0.000	14.712	
Baseline - Lock		15.317	19.549	
Peak - End		14.900	0.570	
Peak - End		20.300	0.490	
Baseline - Together		19.807	20.339	
Baseline - Lock		24.675	25.697	
Peak - End		25.833	0.581	
Peak - End		25.833	0.600	
Baseline - Lock		29.359	31.642	
Peak - End		32.833	0.776	
Peak - Start		36.417	-0.495	
Peak - Start		37.983	-0.440	

ResultsSummaryIntegrationMeasurement ConditionsSST Results

Fig 119: Chromatogram - Integration

Grp

Group identifier for the *Add Group* and *Delete Group* commands.

Time A

- a) Indicates the starting time of the operation (for interval commands).
- b) Indicates the peak retention time (for commands that modify the peak start, peak end, etc.).

Time B

- a) Indicates the final time of the operation (for interval commands).
- b) Indicates the position of a new start, end, etc. relative to the peak apex (for commands that modify the start, end, etc.).

Note: Zero entered in the *Time B* will set the end of the interval to the end of the chromatogram.

Note: Indicating the relative position (i.e. time) is more suitable than indicating a mere absolute value, since in using the **Integration Table** the peak with the most similar retention time is then found and the peak start/end is shifted to the point calculated from the relative shift on the actual retention time. This substantially diminishes the differences in actual retention time of the same peak in different chromatograms.

Value

Value (if relevant for the operation) with appropriate units.

Description of fields of integration table

Global Peak Width / Local Peak Width [min]

This parameter defines the smallest peak width, which is to be detected. The value range is between 0,001 and 10 minutes.

The value is not critical, so even narrower peaks than the entered value will be detected, but if set incorrectly, the beginning and end of a peak may be inaccurately determined and the baseline could be incorrectly inset.

The *Global Peak Width* parameter is always displayed in the first row of the integration table and is applied to the whole chromatogram.

It is possible to set a different peak width for individual sections by adding the *Local Peak Width* parameters into further rows of the integration table which then override the *Global Peak Width* settings in the given area of chromatogram.

Global Threshold / Local Threshold [mV]

Determines the sensitivity the integration algorithm has to noise. The height of all peaks (measured from the baseline to the top) that should be detected has to be at least double this parameter. The value range must be 1 μV - 10 V.

The *Global Threshold* parameter is always displayed in the second row of the integration table and is applied to the whole chromatogram.

It is possible to set a different noise threshold for individual sections by adding the *Local Threshold* parameters into further rows of the integration table which then override the *Global Threshold* parameter in the given area of the chromatogram.

Integration Interval

Limits the time period over which the chromatogram will be integrated.

This means that the chromatogram profile will be ignored beyond the set time interval. This is the main difference compared to *Lock* command, which cancels already integrated peaks only in defined sections and does not reintegrate the whole chromatogram. If the parameter is not defined, integration is done on the whole chromatogram.

Detect Negative

Enables the detection of negative peaks. If this function has not been enabled, all negative peaks are considered to be correct course of the

baseline (in a similar way as removing the peak with the use of *Reject Negative* function).

Min. Area [mV.s]

This parameter sets the minimum peak area in mV.s. Peaks whose area is smaller than the specified value are not displayed.

Min. Height [mV]

This parameter gives the minimum peak height in mV. Peaks whose height is smaller than the specified value are not displayed.

Half Width [min]

This parameter gives the minimum peak half-width in minutes. Peaks whose *W05* is smaller or equal to the specified value are not integrated.

ValleyToValley Slope

Specifies the maximum slope of the baseline. If the baseline slope is smaller than the specified value, the separating perpendicular line is canceled and the baseline passes through the valley. The default value of 0 prevents the baseline from passing through the valley.

Tangent Area Ratio

Constitutes the first condition imposed on tangential separation. The condition is satisfied if the ratio of areas of the main and the separated peak exceeds the specified value.

The default value of 0 prevents tangential separation. If there is no *Tangent Slope Ratio* item defined in the **Integration table**, its value is considered to be 1.5.

Tangent Slope Ratio

This is the second condition imposed on tangential separation. The condition is satisfied if the ratio of slopes of the second and the first peak exceeds the specified value. The slope ratio is defined as the tangent to the first peak at its end point divided into the slope of the line connecting the beginning and end of the second peak.

The default value of 0 prevents tangential separation. If there is no *Tangent Area Ratio* item defined in the **Integration table**, its value is considered to be 0.1.

Filter

Commands from *Chromatogram - Integration* menu:

FFT, *Savitzky-Golay*, *Spike*, *Moving Average*. A detailed description of these functions is provided in the chapter "**Integration**" on pg 183.

Baseline

Commands from *Chromatogram - Baseline* menu:

Lock, *Valley*, *Together*, *Forward Horizontal*, *Backward Horizontal*, *Front tangent*, *Tail tangent*, *Clamp Negative*, *Cut Negative*, *Reject Negative*, *Allow Crossing*, *Spike Removal*. A detailed description of these functions is provided in the chapter "**Baseline**" on pg 174.

Peak

Commands from *Chromatogram - Peak* menu:

Start, End, Both, Add Positive, Add Negative, Force Peak Name, Solvent Peak. A detailed description of these functions is provided in the chapter "**Peak**" on pg 178.

Group

Adds (*Add group*) and removes (*Delete group*) a group of peaks. A detailed description of these functions is provided in the chapter "**Peak**" on pg 182.

Evaluation

Commands from *Chromatogram - Noise & Drift* menu:

Noise, Drift, ASTM Noise, 6-Sigma Noise. A detailed description of these functions is provided in the chapter "**Noise & Drift**" on pg 191.

Sequence of evaluation

The evaluation process does not proceed exactly in the same order as the parameters in the integration table are shown. The process is:

- Primarily initiated with the *Global (Local) Peak Width, Global (Local) Threshold, Integration Interval* and *Detect Negative* parameters.
- Secondly the *ValleyToValley Slope* parameter will be applied. After all groups of non-resolved peaks have been tested by this parameter the remaining two parameters *Tangent Area Ratio* and *Tangent Slope Ratio* are applied to all remaining non-resolved groups. Tangential separation is used whenever the relevant criteria are met.
- Only then is the **Integration Table** used, followed by the *Rejections* parameters.

It thus follows that the separation parameters are not applied (e.g. to manually added peaks) and can be suppressed by the *Together* and *Valley* commands included in the **Integration table**.

3.5.3 Measurement

Displays the conditions under which the chromatogram was measured. Once selected, the **Measurement Conditions** tab will be displayed in the lower part of the **Chromatogram** window, containing **Instrument**, **Event Table** and possible **Detector** sub-tabs; when using a chromatogram that has been measured on an instrument with directly controlled chromatograph, autosampler or any other device having its own tab in the **Method Setup** dialog, the **Sampler**, **GC**, **LC Gradient**, **LC** or other sub-tabs are also present. The displayed values are taken from the relevant tabs of the template method describing the way how the chromatogram was measured.

Instrument

This sub-tab is common to all signals in given chromatogram.

The following fields in the **Measurement Conditions** section are designated for the information purposes only and can be filled in or left blank according to your choice.

Description - brief description of method.

Column - description of the column used.

Mobile Phase - description of mobile phase.

Flow Rate - description of flow rate.

Pressure - description of pressure.

Detection - detection method.

Temperature - description of temperature.

Note - any comment primarily connecting to the method.

Common for All Signals

Measurement Conditions Description: Polymer IEX - wine analysis Column: WATREX 250x8 mm Polymer IEX H form Mobile Phase: 1,7 mM H2SO4 Flow Rate: 0,8 ml/min Pressure: 10 MPa Detection: Refractometer R40 I, 8x, 100mV/UV2 Temperature: 50 C Note: SDS 030 Guard PEEK 10x3 mm PS-DVB 10 um		Sample Identification Sample ID: Vial 101 Sample: Acquired: 16.7.2012 13:52 Analyst: Acquisition Parameters AutoStop: 12 min., enabled Ext. Start: Down Acq. method: DEMO_NGA Range: <varies> Sampling Rate: 12,5 samples/sec Subtraction Chromatogram: Matching: No Change		Chromatogram Comments: Edit... Audit Trail: Acquisition Messages Show... GLP Mode: On
---	--	---	--	--

Instrument | Event Table | GLP Info | Sampler 1 | GCs | U-PAD2 - 1 | U-PAD2 - 2 |
 Perf | Summary | Performance | Integration | **Measurement Conditions** | SST Results

Fig 120: Chromatogram - Measurement Conditions

Sample Identification

Fields from the header of the analysis - **Single Analysis** dialog or the relevant row of the **Sequence Table** in the **Sequence** window.

Sample ID

Sample identifier. This corresponds to the *Sample ID* field from the **Single Analysis** dialog or the *Sample ID* column from the **Sequence** window.

Sample

Description of the sample. This corresponds to the *Sample* field from the [Single Analysis](#) dialog or the *Sample* column from the [Sequence](#) window.

Acquired

Date and time of the end of the measurement. If the date is not accessible (e.g. in some imports), N/A will be displayed.

Analyst

Name of the user who measured the analysis.

Acquisition Parameters

Parameters of measurement common for all signals. The remaining parameters can be found on the [Method Setup - Measurement](#) tab.

AutoStop

Notifies whether the *Enable Autostop* checkbox was used and what was the *Run Time* value set.

Ext. Start

Notifies whether and by what method the external control of the measurement was defined. This corresponds to the *External Start/Stop* section of the [Method Setup - Measurement](#) tab (see chapter 2.5.2 on page 59 for more details).

Acq. Method

Name of the template method used for measurement.

Range

Specifies the range of the detector signal used. If more than one detector was used to gather data with different *Ranges* set (including signal units), the *<varies>* inscription will be displayed here.

Sampling Rate

Specifies the sampling rate of the detector used. If more than one detector was used to gather data with different *Sampling Rate* set, the *<varies>* inscription will be displayed here.

Subtraction ✓ Full version

Sets the automatic subtraction of selected chromatogram. Subtraction of the chromatogram (e.g. solvent) is indicated by the note *Mathematically reprocessed* in the printed report. The use of automatic subtraction is also logged into the chromatogram audit trail.

There are two places in **Clarity** where the subtraction chromatogram can be set - [Method Setup - Advanced](#) dialog accessible from the [Instrument](#) window and [Measurement Conditions - Instrument](#) tab of the [Chromatogram](#) window.

If you wish to automatically subtract a chromatogram from the newly measured chromatograms, specify it in the template method in the [Method Setup - Advanced](#) dialog.

If you wish to subtract a chromatogram from an already measured chromatogram, specify which on the [Chromatogram - Measurement Conditions](#) tab (described further in this chapter).

The subtracted chromatogram is saved into the measured chromatogram, it is not simply linked to it. The changes occurring with the subtraction chromatogram thus do not influence the measured chromatogram from which the background is subtracted. The subtraction, however, does not hold its history - only last subtracted chromatogram's data is saved when it is changed, all older subtraction chromatograms are discarded.

Chromatogram

Field specifying the file with subtraction chromatogram.

Set...

Displays the [Open Chromatogram](#) dialog used for selecting the subtracted chromatogram.

None

Cancels subtraction of chromatogram.

Note: If you only wish to display the chromatogram on the background of the acquisition window without automatic subtraction, use the *File - Set Background Chromatogram...* command in the [Data Acquisition](#) window.

Matching

Selects the subtraction method.

No change - The chromatogram is subtracted from the measured chromatogram without any additional adjustments.

Offset Subtraction Chromatogram - Before subtraction, the **Clarity** station attempts to “move” the measured chromatogram with the subtracted chromatogram by shifting along the time axis.

Scale Subtraction Chromatogram - Before subtraction, the **Clarity** station attempts to “align” the measured chromatogram with the subtracted chromatogram by interlining it or shortening the time.

Note: Alignment of chromatograms is achieved using the biggest peaks in the subtraction chromatogram, i.e. either to the highest peak or to the peak with the largest area.

Chromatogram

Section that contains the additional information on the sample or chromatogram.

Edit...

This button serves for adding some information about the chromatogram or additional info on the sample in the [Rich Text Edit](#) dialog. The button also allows to modify the comment entered prior to the analysis from the [Single Analysis](#) dialog. This field can then be added to the report setup on the [Report Setup - Results](#) tab described on pg 348.

Audit Trail

Using the *Show* button allows to display the [Chromatogram Audit Trail](#) dialog containing the log of events happening with the given chromatogram.

Event Table

This tab shows the setting of the **Event Table** defined in the method on the [Method Setup - Event Table](#) tab. It is common for all signals in case of the multi-detector chromatogram. More information on the **Event Table** is in the chapter "**Event Table**" on pg 73.

GLP Info

This tab shows the list of control modules (hardware) configured on the instrument and used for the acquisition of the chromatogram, along with the serial numbers of this hardware. The contents of this tab may be added to the printed reports by checking the *GLP Info* checkbox on the [Report Setup - Method](#) tab.

Detector

The tab displaying detailed measurement parameters. In the case of a multi-detector chromatogram there is a tab for every detector in the configuration of the Instrument.

Sampler

Tab displaying the injection method of the autosampler.

GC/LC/LC Gradient

Tab displaying instrument method of gas or liquid chromatograph.

3.6 Results

This section contains commands for displaying the result tables.

Result Table	Switches to the Result Table on the Results tab of the Chromatogram window.
Summary Table	Switches to the Summary Table of all displayed chromatograms on the Summary tab.
Performance Table 	Switches to the Performance Table characterizing the chromatographic system quality on the Performance tab of the Chromatogram window.
Summary Options...	Opens the Summary Table Options dialog which sets the layout options for the Summary Table .

All tables are very flexible; the order of particular items and the visibility and size of columns can be adapted and customized, columns appended, fonts defined separately for each column and the number of digits displayed specified for selected columns.

Tables can be modified by commands from the local menu accessible by right clicking the mouse over the table.

3.6.1 Local Table Menu

Local menus contain frequently used commands relevant for the given table. Local menus can thus vary, according to the table you are currently editing. In the following overview commands common for most tables are described:

Select All

Selects all cells in the table. Also available by using the **Ctrl** + **A** shortcut.

Copy

Copies selected cells to the clipboard. Also available by using the **Ctrl** + **C** shortcut.

Paste

Pastes the content of the clipboard into the selected cells of the table. Also available by using the **Ctrl** + **V** shortcut.

Restore Default Columns

Restores the default columns and fonts for a given table. A detailed description of this function is in the chapter "**Edit**" on pg 149.

Setup Columns...

Invoking the *Setup Columns...* command opens the [Setup Columns](#) dialog, where it is possible to change the appearance and order of table columns. Detailed description of this dialog can be found in the chapter "**Setup Columns...**" on pg 154.

Note: To access the setting of user columns the user must be authorized to edit methods. See the *Edit Method* checkbox in the [User Accounts](#) dialog.

User Columns ✓ Full version

Submenu of commands enabling to handle user columns in the full version. The available commands are:

Add...

Use this command to add a customized user column to the table. Invoking this command will open the [Add User Column](#) dialog, whose detailed description can be found in the chapter "[Add User Column](#)" on pg **151**.

Edit Selected...

Invoking this command will open the [Add User Column](#) dialog for the active user column. It is used for modifying of the content of the user columns.

Delete Selected

Invoking this command deletes the active user column or columns.

Delete All

Deletes all previously added user columns. A detailed description is in the chapter "[Edit](#)" on pg **148**.

Hide Column(s)

Hides the selected columns.

Show Hidden Column(s)

Shows all hidden columns.

3.6.2 Result Table

Invoking this command switches the lower part of the **Chromatogram** window to the **Results** tab, displaying the **Result Table** in the left half. When using a multi-detector chromatogram the **Result Table** is specific only to the active signal.

Note: If you use *Peak(s) Selected in Result Table* option in **Graph Properties** dialog, the active row(s) in **Result Table** will color the area of corresponding peak(s) in the **Main Graph**. The peak color is the same as the signal color and corresponds to the text color of the table heading.

Result Table (ESTD - C:\Clarity\DEMO1\Data\2506MULTI - UV)						
	Reten. Time [min]	Response	Amount [g/l]	Amount [%]	Peak Type	Compound Name
1	4,563	3,613	0,027	0,3	Ordnr	oxalic
2	5,203	253,325	1,609	16,8	Ordnr	otric
3	9,417	561,767	2,372	24,7	Ordnr	tartrac
4	6,300	217,400	1,181	12,3	Ordnr	malic
8	8,160	177,150	1,126	11,7	Ordnr	succinic
9	8,550	193,934	1,970	20,5	Ordnr	lactic
11	10,347	152,602	1,312	13,7	Ordnr	acetic
C11	12,710					methanol
	Total		9,596	100,0		

Fig 121: Chromatogram - Result Table

The **Result Table** consists of the **Heading**, **Upper Header**, **Left Header** and **columns**, which may be default or user defined. The appearance of the **Result Table** of chromatogram with linked calibration differs from the one without it. For chromatograms without calibration file linked some of the **columns** are not available.

Note: The user columns defined in the **Result Table** and in the **Summary Table** are independent.

Note: The Result table is visible only when chromatographic data are unchanged. After modifying primary data (Scale or Move operations) the Result table will disappear as the results are no longer valid.

Heading

Shows the name of the active chromatogram, its active signal in the case of multi-detector chromatograms and calibration calculation type. If it is not possible to use the selected type of calculation for whatever reason, the entire heading is marked with exclamation marks and supplemented with a description of the cause of the problem.
 E.g., the following report **!! Result Table (2506multi - UV) <-- Uncal No ISTD peak detected !!** means that the calculation cannot be performed using the internal standard, because the calibration file does not contain any peak which has been indicated as internal standard (ISTD). Uncalibrated calculation will be used instead. The heading is always in the color of the active chromatogram (signal).

Upper header

Shows the names of columns. A click on the field in the column header selects the column, a double-click arranges the table in ascending order according to the values in the column. A second double-click arranges the table in descending order. This column can be moved by clicking on the selected column header and dragging the cursor to the desired position. The column can be widened or narrowed by pressing the left mouse button above the right edge (when the cursor is changed to ) and dragging it. A double-click above the right edge sets the lowest possible width of the column for the data and column header to fit in.

Left header

Shows the serial number of particular peak and possibly also the indication of appurtenance to group of peaks. When using the *All peaks in Calibration* option for the *Report in Result Table* selection, the substances present in the calibration but not found in the active chromatogram are indicated with a red serial number taken from the calibration beginning with the letter **C**.

Reten. Time

Retention time of peaks in minutes. The peaks are by default ordered according to the retention time, so double-clicking *Reten. Time* column header returns the **Result Table** to the default arrangement.

Start Time/End Time [min]

These columns display the retention time of particular Peak Start/Peak End. They are hidden in the default layout.

Start Value/End Value [mV]

These columns display the height of the baseline in the Peak Start/Peak End. They are hidden in the default layout.

Area [mV.s]

Shows the peak area in millivolt-seconds (or units defined in the detector setup*seconds). This column is hidden by default when the calibrated results are used. Instead the *Response* column is shown.

Height [mV]

Shows the height of the particular peak in mV (or units defined in the detector setup). This column is hidden by default when the calibrated results are used. Instead the *Response* column is shown.

Response

Area or height of the peak used in the calibrated calculation. The type of response is selected according to the calibration file and could be shown for each peak in the *RB* column.

RB

Describes the response base for the particular peak when the calibrated calculation is used (*A* stands for area, *H* for height). Unidentified peaks

use the response base according to the *Response Base* parameter set in the *Unidentified Peaks* section.

Area% [%], Height% [%]

Percentage representation of peak area or height.

$$Area\%_i = \frac{A_i}{\sum_i A_i} \cdot 100 \quad Height\%_i = \frac{H_i}{\sum_i H_i} \cdot 100$$

W05 [min]

Shows the width of the peak in half of its height.

Amount [x]

Shows calculated amount of particular substance in units x stated in the calibration file by the *Units - Compound* parameter of the [Calibration Options](#) dialog. These units may be overridden by units entered due to scaling in the *Units After Scaling* field of the [Method Setup - Calculation](#) dialog or *Units* field of the *Scale* section on the [Chromatogram - Results](#) tab. See the **Appendix 1** below.

Amount% [%]

Shows the percentage amount of the corresponding substance. The total amount is given by the sum of the corrected amounts of particular compounds or by the *Amount* parameter from the heading of the chromatogram if this is other than zero. The *Amount* parameter can also be set additionally in the *Scale* section on the [Chromatogram - Results](#) tab. See the **Appendix 2** below.

Peak Type

Shows the type of identified substance. The type may be one of the following:

Tab 12: List of possible peak types

Peak Type	Description
<i>Ordnr</i>	- Peak which is neither a reference peak nor internal standard.
<i>Refer</i>	- Reference peak.
<i>ISTD</i>	- Peak of internal standard.
<i>RISTD</i>	- Reference peak which is also an internal standard.
<i>Grp</i>	- Indication of calibrated group.
<i>Free</i>	- Freely calibrated substance (Free Calibration). The calibration curve is not used for the calibration calculation but rather the response coefficient of the substance (<i>Response Factor</i>) from the main Calibration Table .

The information may be complemented by the failed result of the linearity check (as defined by the *Curve Check* item in the [Calibration Options](#) dialog), in which case the *Error (Curve Check)* inscription is shown prior to the *Peak Type* value. Successful checks are not shown.

The *Peak Type* column may also indicate the fact that the amount is less than limit of detection (**LOD**) or quantification (**LOQ**). In such case the column will also display "< LOD" or "< LOQ" inscription. The *LOD* and *LOQ* limits can be set in the *LOD* and *LOQ* columns of the global calibration table (for more details see the chapter **Calibration** on pg 239).

The last type of the information given in the *Peak Type* column is the way how the response for the peak was calculated. This type of information is shown in brackets after the peak type definition only in case the identification or calculation was not performed in the ordinary way. If the peak was identified using the *Force Peak Name* function from the [Chromatogram](#) window, the *Forced* label is shown in the brackets. In case the response of the peak was calculated using the response factor of another peak (as defined in the [Calibration](#) window), the *by [compound]* label is shown, where [compound] is substituted by the name of the peak whose response was used for the calculation. In case a *Correction Factor* for the peak differs from 1 (as defined in the [Calibration](#) window), the *CF* label is displayed. These three labels can be freely combined, according to the actual settings in the [Chromatogram](#) and [Calibration](#) windows.

Compound Name

Displays the name of the substance from calibration file corresponding to the relevant peak on the basis of the basic identification of retention times. If a certain peak is not identified, the *Peak Type* and *Compound Name* columns for that peak are empty and the *Amount* field value is zero (unless the *Response Factor* parameter in the *Unidentified peaks* section is not equal to zero).

Centroid [min], Variance [min²], Skew [-], Excess [-], Asymmetry [-], Capacity [-], Efficiency [th.pl], Eff/I [t.p./m], Resolution [-]

These columns belong to functions used in the **Performance Table** on the [Chromatogram - Performance](#) tab. More information can be found in the chapter "**Performance Table**" on pg 232.

Response Factor ✓ full version

This is a predefined **User Column** displaying ratio of **Response** and **Amount**. The column is hidden in the default desktop.

Response Factor = Amount / Response

It is possible to calculate response factor in inverse form (RF = Response / Amount) by checking *Response Factor as Response / Amount* checkbox in [Calibration Options](#).

Signal/Noise ✓ full version

This is a predefined user column displaying ratio of **Signal** and **Noise**. The column is hidden in the default desktop.

Signal/Noise = Signal / Noise

Group-X row

This row is visible only in the case that peak groups are used, once for each group. It shows the sum of values of individual peaks belonging to

the group in the given column (only for columns when such total makes sense).

Total

This row shows the total value of a given column (only for columns when such total makes sense).

Appendix 1:

The calculation is determined by the type of calibration:

Tab 13: Calculation formulas

Calculation	Formula
ESTD, NORM	$Amount_i = CF_i[R_i] \cdot ScaleFactor$
ISTD	If the calibration does not have ISTD amount filled on any level, this equation is used. Same amount of the internal standard is considered to be present both in samples and in calibration standards, ISTD Amount values everywhere in Clarity are set to 0. $Amount_i = CF_i \left[\frac{R_i}{R_{ISTD}} \right] \cdot ScaleFactor$
ISTD	If the calibration contains ISTD Amount filled on at least one calibration level, following equation is used. ISTD Amounts then must be filled in everywhere in Clarity, both in unknown samples and in the calibration file, and these amounts may differ. $Amount_i = RCF_i \left[\frac{R_i}{R_{ISTD}} \right] \cdot Amount_{ISTD} \cdot ScaleFactor$

where:

Tab 14: Explanation of used symbols

Symbol	Description
$Amount_i$	- Amount of component "i" in sample.
R_i	- Response of the "i" component (height or peak area).
R_{ISTD}	- Response of internal standard (height or peak area).
$CF_i(R_i)$	- Amount of "i"th component read out from corresponding calibration curve.
$CF_i \left[\frac{R_i}{R_{ISTD}} \right]$	- Amount of "i"th component read out from corresponding calibration curve displaying dependence of response ratio (R_i/R_{ISTD}) of calibration standard on "i"th component amount in that calibration standard for given substance.

Symbol	Description
$RCF_i \left[\frac{R_i}{R_{ISTD}} \right]$	- Value read out from corresponding relative calibration curve, displaying dependence of response ratio (R_i/R_{ISTD}) of calibration standard on "i"th component amount ratio ($Amount_i/Amount_{ISTD}$) in that calibration standard for given substance.
$Amount_{ISTD}$	- Amount of internal standard added to unknown sample. This value must be entered into both the ISTD Amount field in the chromatogram heading and the Amount column of the calibration, or to none of these places in case all ISTD Amount values are the same (optional, the calculation works the same if the same amounts are included on both places).
ScaleFactor	- Multiplication factor.

In the case that the corresponding peak is not found in the calibration file, the *Uncal. Response* parameter from the [Method Setup - Calculation](#) dialog is used for the calculation instead of deriving the value from the calibration curve, according to the formulas:

Tab 15: Calculation formulas for non-calibrated peaks

Calculation	Formula
ESTD	$CF_i[R_i] = Uncal. response \cdot R_i$
ISTD	$RCF_i \left[\frac{R_i}{R_{ISTD}} \right] = Uncal. response \cdot \frac{R_i}{R_{ISTD}}$
NORM	The $CF_i(R_i)$ value is calculated exactly like in the ESTD calculation, but with peaks not found in the calibration, no <i>Amount[%]</i> value may be given in this case and the <i>Peak Type</i> column shows the <i>Error</i> inscription.

Caution: When using a different type of calibration than the type you selected, the used type of calibration is indicated above the table, stating the reason for the change.

Appendix 2:

Calculation of percentage representation:

$$Amount\%_i = \frac{\frac{Amount_i}{SampleAmount} \cdot ScaleFactor}{Dilution} \cdot 100$$

3.6.3 Summary Table

Invoking this command switches the lower part of the [Chromatogram](#) window to the [Summary](#) tab, displaying the **Summary Table** on it. The **Summary Table** serves for easy display and printing of selected parameters from all displayed chromatograms in one table. The table only displays the results of calibrated peaks.

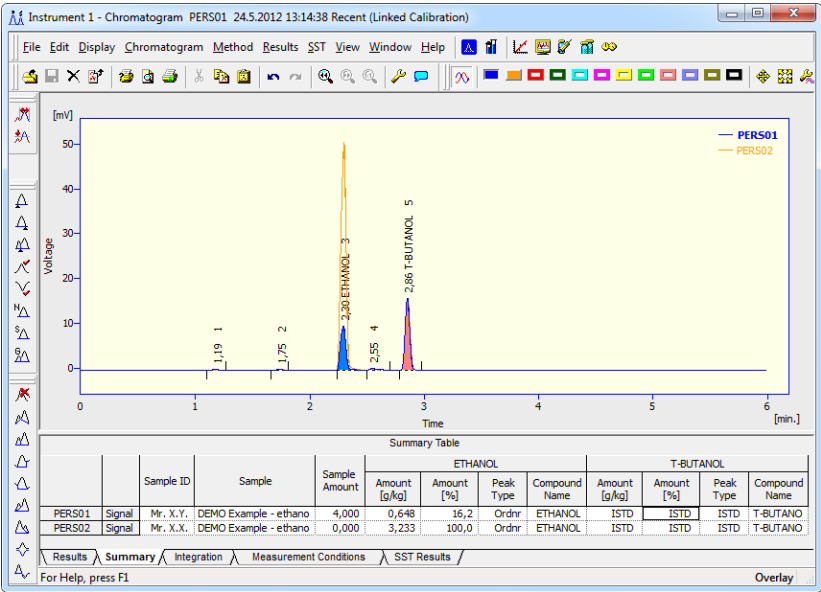


Fig 122: Chromatogram - Summary

Note: The user columns in **Result Table** and in **Summary Table** are independent.

Presentation of results

The results table can be added into the print configuration by checking the *Summary Table* checkbox in the [Report Setup - Results](#) dialog described on pg 348. The results table can also be exported using the *File - Export - Export Summary Table* command described in the chapter "**File**" on pg 123.

Selection of displayed parameters

Columns display the same parameters for each identified substance as in the [Results](#) tab for all calibrated compounds. The *Amount [x]* and *Amount [%]* parameters are implicitly offered. The displayed parameters are set using the *Edit - Setup Columns...* command from the main menu or *Setup Columns...* command from the local menu.

Selection of table layout

In the implicit setting, the rows correspond to individual chromatograms and their signals. The layout best corresponding to your requirements can easily be set using the *Table Inverted* and *Parameter Header* commands from the [Summary Options](#) dialog (see following chapter).

3.6.3.1 Summary Options

By right-clicking on the **Summary** tab and selecting the *Summary Options...* command, the **Summary Table Options** dialog with the settings of the **Summary Table** display appears.

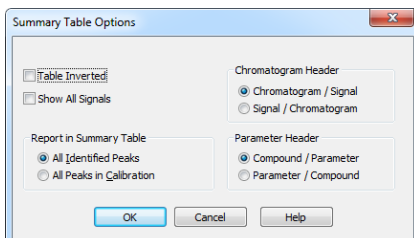


Fig 123: Summary Table Options

Table Inverted

When checked, chromatograms will be displayed in columns and individual compounds and their parameters will be placed in rows, inverting the table order. This display may be suitable for a smaller number of chromatograms and a large amount of compounds and/or parameters.

Show All Signals

The **Summary Table** implicitly hides signals that do not contain any identified peaks. When checked, all signals from opened chromatograms will be displayed. This can be used to identify problematic signals easily.

Chromatogram Header

Selects the grouping order of rows in the **Summary Table**. This serves for easy comparison of corresponding signals in a set of several multi-detector chromatograms.

Report in Summary Table

This option selects which compounds will be displayed in the **Summary Table**.

All Identified Peaks

Only compounds identified in the corresponding calibration files will be displayed.

All Peaks in Calibration

Summary table will contain all compounds from all used calibration files without regard to whether they were all identified.

Parameter Header

This section determines whether there will be main columns organized by compounds or parameters.

Compound / Parameter(s)

Main columns are organized by compounds. First compound column with all parameters of the first compound is followed by second compound column with all parameters of the second compound and so on.

Parameter / Compound(s)

The columns are organized by parameters. The values of the first parameter will be displayed for all compounds first, followed by the values of second parameter for all compounds and so no.

Available layouts of Summary Table:

Summary Table							
		Benzen		Toluen		Chlorbenzen	
		Amount [ug/l]	Amount [%]	Amount [ug/l]	Amount [%]	Amount [ug/l]	Amount [%]
SR 1_1710	Signal 1	0,082	12,2	0,085	12,6	0,110	16,2
	Signal 2	0,087	101,0	0,086	100,0	0,111	128,0
SR 1_1910	Signal 1	0,074	12,2	0,077	12,7	0,098	16,1
	Signal 2	0,087	101,0	0,086	100,0	0,111	128,0
SR 1_2510	Signal 1	0,074	11,9			0,099	15,9
	Signal 2	0,087	49,3	0,086	48,8		
Results Summary Performance Integration Measurement Conditions SST Results /							

Compound / Parameter

Summary Table							
		SR 1_1710		SR 1_1910		SR 1_2510	
		Signal 1	Signal 2	Signal 1	Signal 2	Signal 1	Signal 2
Benzen	Amount [ug/l]	0,082	0,087	0,074	0,087	0,074	0,087
	Amount [%]	12,2	101,0	12,2	101,0	11,9	49,3
Toluen	Amount [ug/l]	0,085	0,086	0,077	0,086		0,086
	Amount [%]	12,6	100,0	12,7	100,0		48,8
Chlorbenzen	Amount [ug/l]	0,110	0,111	0,098	0,111	0,099	
	Amount [%]	16,2	128,0	16,1	128,0	15,9	
Results Summary Performance Integration Measurement Conditions SST Results /							

Inverted Compound / Parameter

Summary Table							
		Amount [ug/l]			Amount [%]		
		Benzen	Toluen	Chlorbenzen	Benzen	Toluen	Chlorbenzen
SR 1_1710	Signal 1	0,082	0,085	0,110	12,2	12,6	16,2
	Signal 2	0,087	0,086	0,111	101,0	100,0	128,0
SR 1_1910	Signal 1	0,074	0,077	0,098	12,2	12,7	16,1
	Signal 2	0,087	0,086	0,111	101,0	100,0	128,0
SR 1_2510	Signal 1	0,074		0,099	11,9		15,9
	Signal 2	0,087	0,086		49,3	48,8	
Results Summary Performance Integration Measurement Conditions SST Results /							

Parameter / Compound

Summary Table							
		SR 1_1710		SR 1_1910		SR 1_2510	
		Signal 1	Signal 2	Signal 1	Signal 2	Signal 1	Signal 2
Amount [ug/l]	Benzen	0,082	0,087	0,074	0,087	0,074	0,087
	Toluen	0,085	0,086	0,077	0,086		0,086
	Chlorbenzen	0,110	0,111	0,098	0,111	0,099	
Amount [%]	Benzen	12,2	101,0	12,2	101,0	11,9	49,3
	Toluen	12,6	100,0	12,7	100,0		48,8
	Chlorbenzen	16,2	128,0	16,1	128,0	15,9	
Results Summary Performance Integration Measurement Conditions SST Results /							

Inverted Parameter / Compound

3.6.4 Performance Table ✓Full version

A table containing peak parameters characterizing the quality of the chromatographic system. Invoking the command opens the [Performance](#) tab.

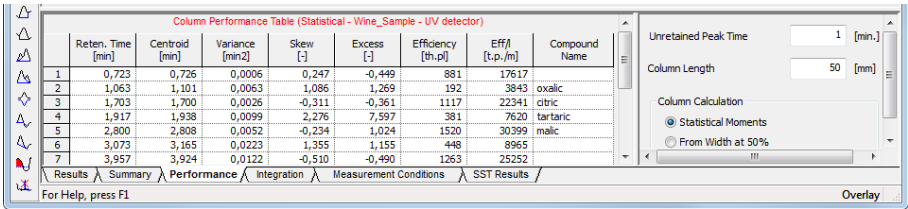


Fig 124: Chromatogram - Performance

The left-hand part contains the table, while on the right part the calculation type is selected.

Tab 16: Elements of the Performance Table

Element	Description
Upper Caption	- Shows column names. To display the values of the table in an ascending order, double-click the upper caption; double-click again to change the order to descending. Move the column to some other position by left clicking the mouse twice slowly and dragging the cursor above the desired column. To broaden or narrow a column, click, hold down and then drag the right border. Restore the default column width by double-clicking the right border.
Left caption	- Displays the peak number.
Reten. Time	- Retention time in minutes.
W05 [min]	- Peak half-width (peak width at half-height) in minutes. This column displays only if the Performance table is created using <i>Width at 50%</i> .
Compound Name	- Names of identified compounds.

The following columns vary depending on the type of calculation used:

Statistical Moments

Tab 17: List of columns in the Performance Table - Statistical Moments

Column	Equation
Centroid [min]	- Peak centroid: $M_1 = \int_{-\infty}^{+\infty} \frac{th(t)dt}{M_0}, \quad \text{where } M_0 = \int_{-\infty}^{+\infty} h(t)dt$
Variance [min ²]	- Peak variance: $\overline{M}_2 = \int_{-\infty}^{+\infty} \frac{(t-M_1)^2 h(t)dt}{M_0}$
Skew [-]	- A measure of peak skewness: $Skew = \frac{\overline{M}_3}{\overline{M}_2^{3/2}}$
Excess [-]	- A measure of Kurtosis: $Excess = \frac{\overline{M}_4}{\overline{M}_2^2} - 3$ where: $\overline{M}_n = \int_{-\infty}^{+\infty} \frac{(t-M_1)^n h(t)dt}{M_0}$
Efficiency	- Column efficiency measured by the plate number: $Efficiency = \frac{M_1^2}{M_2}$
Eff./l	- Relative efficiency: $Eff./l = Efficiency_i \cdot \frac{1000}{ColumnLength}$
Symmetry/Tailing [-]	- Symmetry/Tailing: $Symmetry = \frac{A05_i + B05_i}{2 \times A05_i}$

From Width at 50%:

Tab 18: List of columns in the Performance Table - Width at 50%

Column	Equation
Asymmetry[-]	- Peak asymmetry: $Asymmetry_i = \frac{b_i}{a_i}$
Capacity[-]	- The k' factor: $Capacity_i = \frac{RT_i - T0}{T0}$
Efficiency	- Column efficiency measured by the plate number: $Efficiency_i = 5,545 \cdot \left(\frac{RT_i}{W05_i} \right)^2$
Eff./l	- Relative column efficiency in plate numbers per column length: $Eff./l = Efficiency_i \cdot \frac{1000}{ColumnLength}$
Resolution[-]	- Represents resolution with regard to the preceding peak: $Resolution_i = 1,18 \frac{RT_i - RT_{i-1}}{(W05_{i-1} + W05_i)}$

Tab 19: Description of used symbols

Symbol	Description
RT_i	- is the retention time of the component.
$W05_i$	- is the peak width at half height.
a_i	- is the left half-width at 10 % height.
b_i	- is the right half-width at 10 % height.
$T0$	- is the unretained (dead) time of the column.
Column Length	- in mm.
$A05_i$	- is the left half-width at 5 % height.
$B05_i$	- is the right half-width at 5 % height.

Parameters influencing the calculation are stated in the right-hand part of the [Performance](#) tab:

Unretained Peak Time

The parameter defines the column dead time in minutes. The value is used to determine the capacity factor.

Column Length

The length of the column in millimeters. The value is used to calculate the theoretical column plate height.

Column Calculations

The method used to calculate some parameters in the table is specified here.

Statistical Moments - Calculations are based on statistical moments.

From Width at 50% - The peak width at 50% of the peak height and the corresponding half widths are used in the calculations.

3.7 SST

SST stands for the **System Suitability Test** characterizing the measurement quality. Invoking the command opens the [SST Results](#) tab. A detailed description of this **Clarity Extension** can be found in a separate **SST Extension** manual.

3.8 View

This section contains commands used to set up the appearance of the [Chromatogram](#) window.

Toolbars	Used to display/hide all toolbars(for more details see the chapter View on pg 30).
Customize	A command used to customize the toolbars. The command is similar to the <i>Customize...</i> command in the main Clarity window.
Reset All	Returns all menus and toolbars in the Chromatogram window into their initial state.
Maximize Graph	The command maximizes the graph to cover the entire window (hides the bottom tabs). Double click near the splitter (the cursor will change to ) to achieve the same effect.
Show Both	Used to display both parts of the window (the chromatogram and the tabs).
Maximize Tables	The command maximizes the bottom tabs to cover the entire window (hides the chromatogram). Double click near the splitter (the cursor will change to ) to achieve the same effect.
Show Tables	The submenu for displaying/hiding of the tabs of Chromatogram window. Checked tabs are displayed. The Results tab cannot be hidden this way.

3.9 Window

This is an entry point to a submenu of commands for displaying other main windows.

Main 	Displays the Clarity main station window described on pg 1 .
Instrument	Displays the Instrument window described on pg 40 .
Calibration	Displays the Calibration window described on pg 238 .
Single Analysis	Displays the Single Analysis window described on pg 59 .
Sequence	Displays the Sequence window described on pg 294 .
Data Acquisition	Displays the Data Acquisition window described on pg 279 .
Device Monitor	Displays the Device Monitor window described on pg 55 .
Station Audit Trail 	Displays the Audit Trail window described on pg 318 .
Chromatogram Audit Trail 	Displays the Chromatogram Audit Trail window described on pg 327 .

4 Calibration

The name of the current calibration file is displayed in the window header of the **Calibration** window. When displaying calibration curves saved in the chromatogram, the header contains the name of the chromatogram and the date the displayed calibration curves were saved.

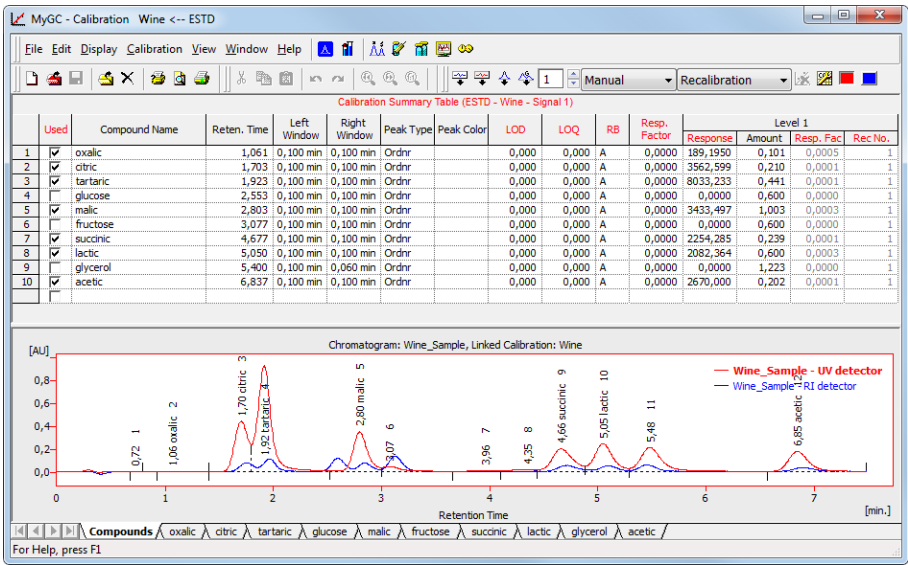


Fig 125: Calibration - Compounds

The window contains the main **Compounds** tab as well as tabs of every compound which is defined in the displayed calibration file. The **Compounds** tab contains the main calibration table and, eventually, the current calibration standard. The **Individual Compound** tabs contain a table for the displayed compound, several parameters and its calibration curve.

Compounds tab

Main calibration table

When using a multi-detector measurement, the main calibration table in the header contains the name of the current signal (detector) for which the data is being displayed.

Certain columns (e.g. compound name, retention time, amount) are **common for all signals** and are displayed in **black**, other columns which are **dependent on the corresponding signal** are displayed in the **color of the active signal**.

Serial number of compound according to retention time. This column is common for all signals.

Used Possibility to prohibit the use of a selected compound. This column is individual for each signal.

Note: This column is particularly useful in multidetector chromatograms, where each compound is assessed on one signal and can be removed from the results table on the remaining signals using this selector.

Compound Name Name of the compound. This column is common for all signals.

Reten. Time Retention time of the compound. This column is common for all signals.

Search Window Defines whether the *Left Window* and *Right Window* content will be inserted in *Abs* (absolute - set in minutes) or *Rel* (relative - set in percents of the *Reten. Time*) values.

Left (Right) Window Defines the width of the left (right) window for identification of compounds (the value is set in minutes). Maximum width is 50 minutes. This column is common for all signals.

Peak Selection Defines the way the correct peak is matched with the Retention time when several peaks are present in the same identification window. The options available are *Nearest* (closest possible to the selected retention time - default for standard peaks), *First*, *Last* and *Biggest* (default for Reference peaks).

Note: The *Peak Selection* choice is overridden whenever a peak is forced-linked to a compound in the calibration by the *Set Peak Name* function in the [Chromatogram](#) window. For more details see the chapter **Peak** on pg 178.

Peak Type	Type of the calibrated compound. This column is common for all signals. <i>Ordnr</i> - Peak which is not a reference peak. <i>Refer</i> - Reference peak. <i>Grp_</i> - Group of peaks.
Is ISTD	Defines whether the given compound is ISTD peak for the purposes of ISTD calculation and assigns the ISTD peak number to ISTD peaks. Up to 5 compounds in the calibration file may be indicated as internal standards from Clarity version 5.0 - before that only one ISTD peak was possible.
Use ISTD	Sets which ISTD peak will be used for the calculations of the given compound. When first ISTD peak is selected in the <i>Is ISTD</i> column of the calibration file, the <i>Use ISTD</i> column of all other compounds is automatically pre-filled with that peak name. ISTD peaks themselves have the <i>Use ISTD</i> column empty.
<i>Note:</i>	A group of peaks can be neither an internal standard nor reference peak.
Peak Color	Sets the coloration of the peak area to the desired color. Whenever the given peak will be identified in the chromatogram using this calibration, it will be colored accordingly.
Reten. Index	Retention (Kovats) indexes.
<i>Note:</i>	This column is hidden in the default layout. For more details see the chapter Setup Columns... on pg 154 .
LOD	Limit of Detection (LOD) - the exceeding of the limits can be displayed in the result table. This column is individual for each signal.
LOQ	Limit of Quantification (LOQ) - the exceeding of the limits can be displayed in the result table. This column is individual for each signal.
RB	Response Base - Calculations are displayed for height - <i>H (Height)</i> or area - <i>A (Area)</i> . This column is individual for each signal.
Resp. Factor	Global response coefficient for free calibration. The value must be entered manually (not automatically calculated) and is used upon a calibrated calculation for compounds which have the flag filed <i>Curve Fit Type</i> set to <i>Free Calibration</i> . This column is individual for each signal.

Correction Factor	Allows for the correction of amount due to the lower recovery etc. The <i>Amount</i> value is multiplied by the <i>Correction Factor</i> to calculate the final <i>Amount</i> .
Calculate By	Clicking the field of this column on a given row opens the Calculate By dialog, which allows to set a compound whose <i>Response</i> will be used for <i>Amount</i> calculation for a given compound.
Curve Fit Type	Determines the type of the curve used to fit the data - the available options are described in the section Curve Fit Type on pg 246
Origin	Specifies the method for including origin zero point into the calibration curve. This value is individual for each signal and is described in the section Origin on pg 247
Ending Point	Specifies the maximum interpolation of <i>Sigmoid</i> and <i>Half Sigmoid Curve Fit Type</i> . The column is only editable when either of these <i>Curve Fit Type</i> options is selected and is described in the section Ending Point on pg 248
Weighting Method	Determines the value of the weight of individual points when calculating the interposing curve. The function and its possible options are described in the section Weighting Method on pg 248 described on pg 248.
Linearization X	Sets the X-axis linearization for the calibration curve calculation. The function and its possible options are described in the section Linearization X on pg 248 described on pg 248.
Linearization Y	Sets the Y-axis linearization for the calibration curve calculation. The function and its possible options are described in the section Linearization Y on pg 249 described on pg 249.
Level X	Number of the calibration level to which the following columns relate. This column is common for all signals.
<i>Note:</i>	Calibration table can contain up to 21 calibration levels including the blank.
Response	Area or height of peak as a response. Type is determined by the <i>RB</i> column. This column is individual for each signal.
Amount	Amount of compound on the level displayed. Units are determined by the <i>Units - Compound</i> field from the Calibration Options dialog and are common for the entire calibration file. This column is common for all signals.

Note: On any level other than 21 (blank), zero in the response or amount columns excludes the level from the calculation. Amount for blank level is *Blank* (0 for all other purposes).

Resp. Fact. Response coefficient for each level. A more detailed description is provided below with the description of the compound table. This column is individual for each signal.

Rec. No. Number of recalibrations on the level displayed. After the recalibration of *Replace* type or after the manual adjustment of the response, the number is set to one. This column is individual for each signal.

Note: A double-click on the row header with the number of compound (gray field to the left next to the name of the compound) displays the tab of the corresponding compound.

Local menu of the Calibration Summary Table

In addition to the common commands for the majority of tables, the local menu contains the following items:

Delete Compound

Removes the actual substance from the calibration.

Delete All Compounds

Removes all compounds from the calibration (from all calibration levels).

Graph of the calibration standard

The graph refers to the calibration standard loaded by the *File - Open Standard* command. Its name is shown at the bottom right of the **Status Bar** and in the upper right corner of the graph. The graph facilitates the transfer of retention times and responses from the calibration standard to the calibration table, and is the same as the chromatogram graph in the [Chromatogram](#) window, including the characteristics set up by the *Properties...* command from the *Display* menu or the local menu. The only difference is the **Peak Area Coloring** area referring to the **Calibration Table**.

Note: To see the colored peaks in the [Calibration](#) window, the calibration standard must have the current calibration linked to it (this can be done in the [Chromatogram - Results](#) tab).

Tabs of individual compounds:

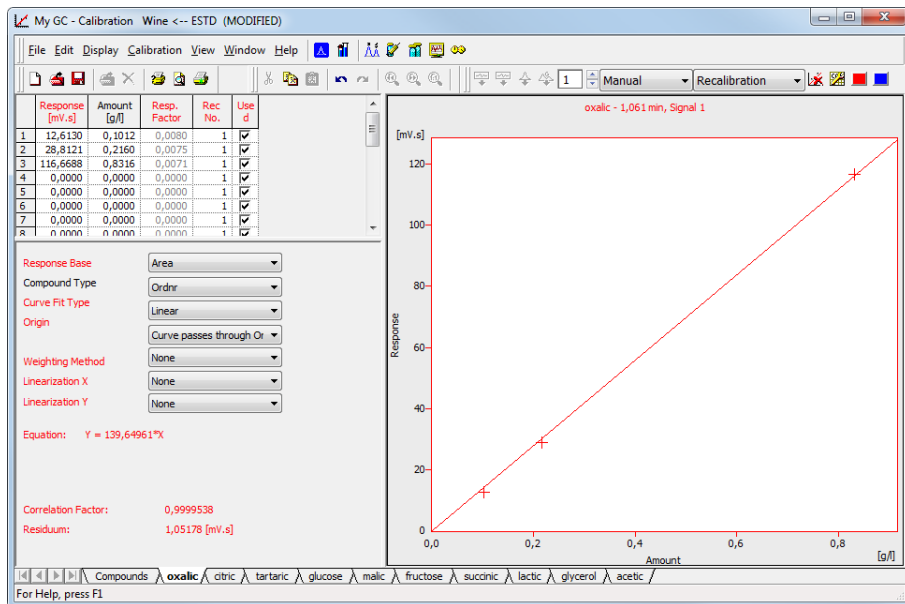


Fig 126: Calibration - Individual Compound (oxalic)

Compound Table

The table can be found in the top left corner of the relevant compound tab.

First column

Contains the level number of each displayed line.

Note: Calibration table can contain up to 21 calibration levels, including blank.

Response

Represents the response at a given level. Whether it will be area or height is defined in the *Response Base* field.

Amount

The amount of the compound at a given level.

Response Factor

The response factor at a given level is equal to the reciprocal value of the slope of a straight line passing through the origin and the calibration point at the given level. It is calculated as the amount of the compound divided into the response at the relevant level. The value serves for information only.

$$LRF_i = \frac{A_i}{R_i}$$

where:

A_i - is the amount of the compound at level i.

R_i - is the response (height or area) at level i .

It is possible to calculate response factor in inverse form ($RF = \text{Response} / \text{Amount}$) by checking *Response Factor as Response / Amount* checkbox in [Calibration Options](#)

Rec No.

Shows the number of times the given calibration point was recalibrated.

Used

Sets validity/invalidity of a point of the calibration curve. The point for which you cancel the indication will not be used in the calibration curve. It will be illustrated by a circle in the graph.

Deviation [%]

Deviation of the *Amount* of the calibration point from the calculated *Amount* given by the calibration curve.

$$\text{Deviation}[\%] = ((\text{Amount} - \text{Amount}_C) / \text{Amount}_C) * 100$$

where:

Amount - The actual Amount on the concentration level.

Amount_C - The calculated Amount corresponding to the concentration level (calculated from the calibration curve).

RSTD [%]

Displays the relative standard deviation of averaged points used during the recalibration of the respective concentration level.

Local menu of the Compound Table

In addition to common commands available for the majority of tables, the local menu contains the

Clear All Levels	Erases all entered values on all levels of the given substance.
Clear Selected Level	Erases all values on the given level.
Clear Selected Value	Erases the value in the current field of the table.
Delete Compound	Removes the current substance from the calibration.
Show History	Invokes the History of Calibration Point dialog with a history log of the selected calibration point. Stored information includes: the name of the calibration standard, retention time, area, height and date of a (re) calibration.

Resettable parameters

The section with resettable parameters is positioned in the bottom left part of each compound tab, below the **Compound Table**.

Response Base

Sets whether the peak height or area will be used as the base for calculation of the calibration curve, calibration dependency equation and two response factors and specifies the base to be displayed in the *Response* column. This setting is individual for each signal. Please note that the height is calculated as the distance from the peak maximum vertically to the peak baseline, so for the small peaks on very steep baselines, the peak height may be imprecise.

Area - Peak area constitutes the base.

Height - Peak height is the base.

Note: Other fields from the tabs of individual compounds are dependant on the *Response Base* selected - switching the *Response Base* will thus switch the content of all editable fields to the set of saved parameters connected to the other *Response Base* option, the graph display will change accordingly and the content of the *Response* column in the **Compound Table** will change too.

Compound Type

This setting is individual for each compound, but common for all signals.

Ordnr - Designates a peak or group that is neither a reference peak nor an internal standard.

Refer - Designates a reference peak.

ISTD - The internal standard peak.

RISTD - The internal standard peak that is at the same time a reference peak.

Note: Only one peak in the calibration file may be marked as the internal standard. A group of peaks can neither be the internal standard nor a reference peak.

Curve Fit Type

Determines the type of the curve used to fit the data:

Free Calibration - No calibration curve is construed and no calibration equation is calculated. Instead of a value read off the calibration curve the calculations use the corresponding global *Response Factor* from the global calibration table.

Point to Point - The calibration points are connected by a broken line. No common calibration equation is calculated, the equation of the relevant line segment is used instead in the calibration calculations.

Linear - A straight line is passed through the points.

Quadratic - The calibration curve will be interlined by a quadratic polynomial.

Cubic - The calibration curve will be interlined by a cubic polynomial.

Sigmoid - The calibration curve will be interlined by a sigmoid.

$$z = 2 \cdot b \sqrt{\frac{\text{Response}}{d}} - 1$$

$$\text{Amount} = c \cdot \sqrt[3]{\frac{1}{2} - \frac{1}{2k} \cdot \text{sign}(z) \cdot \ln(1 - \text{abs}(z) \cdot (1 - e^{-k}))}$$

c - Maximum Amount

d - Maximum Response

The **c** and **d** parameters can be increased up to 125% using the *Ending Point* field (both parameters are changed at once).

b = 1

a, **k** - are coefficients that are calculated using iteration from the interval $<0, 1 - 100>$ so that the curve will best fit the given calibration points.

ln - The calibration curve will be interlined by a logarithm function.

log10 - The calibration curve will be interlined by a decimal logarithm function.

exp - The calibration curve will be interlined by an exponent line.

pow10 - The calibration curve will be interlined by a decimal exponent line.

Hyperbola - The calibration curve will be interlined by a hyperbola.

Half Sigmoid - The calibration curve will be interlined by a half-sigmoid.

If the number of available points does not suffice for the required curve type, the curve will automatically be changed to the next lower type. The *Curve Fit Type* selection remains unchanged.

Note: The *Curve Fit Type* selection is disabled in the Standard Addition display mode - only *Linear* curve fit is available in this mode.

Origin

Specifies the method for including origin zero point into the calibration curve. This value is individual for each signal.

Ignore Origin - Origin will not be included.

Compute with Origin - Considers the Origin to be one of the calibration points.

Curve passes through Origin - Always passes the Curve through the origin.

Note: The *Ignore Origin* option is recommended for the *Origin* field when the blank response is included in the calibration curve calculation.

When interposing using the sigmoid, the *Origin* field will be replaced by the *Ending Point* field. In the Standard Addition display mode, the Origin Field is replaced with the Unknown field.

Ending Point

Enables the calculation of sigmoid interposing outside of the end points. The end calibration points are taken as limit values, increased by the relevant percentage according to the *Ending Point* field. Best used when the possibility to measure the calibration points within the sufficient range is limited.

Unknown

Specifies the method for including the unknown sample into the calibration curve calculation in the Standard Addition mode. This value is individual for each signal.

Without Unknown Sample - Unknown sample will not be included in the calculation.

Compute with Unknown - Considers the Unknown sample to be one of the calibration points.

Curve passes through Unknown - Always passes the calibration curve through the origin.

Weighting Method

Determines the value of the weight of individual points when calculating the interposing curve. When weighting is not applied, points of higher concentrations (larger quantity and response) have a greater influence over the calibration curve than points with a lower concentration.

Weighting enables to apply the same relative influence to each point ($1/x$ weighting method) or assign greater influence to points of lower levels ($1/x^2$ weighting method).

Weighting may be related to either the amount of the compound in the given point of the calibration curve ($1/Amount$ or $1/Amount^2$) or to the response of the compound in the given point of the calibration curve ($1/Response$ or $1/Response^2$).

Linearization X

Linearizes the values of the X-axis using the selected formula, and uses these linearized values to calculate the calibration curve equation. The X-axis and all calibration points are still displayed for the non-linearized values, so the interleaved curve will usually not match the calibration points.

Note: To display the curve as a linear line, it is also necessary to check the *Show Linearized Values* checkbox in the [Graph Properties](#) dialog. For more details see the chapter **Graph Properties** on pg 258.

Possible linearization formulas are *None* (no linearization), $1/Amount$, $\ln(Amount)$, $\sqrt{Amount}$, $1/\sqrt{Amount}$, $Amount^2$, $1/(Amount^2)$ and $\log(Amount)$.

Linearization Y

Linearizes the values of the Y-axis using the selected formula, and uses these linearized values to calculate the calibration curve equation. The Y-axis and all calibration points are still displayed for the non-linearized values, so the interleaved curve will usually not match the calibration points.

Note: To display the curve as a linear line, it is also necessary to check the *Show Linearized Values* checkbox in the [Graph Properties](#) dialog. For more details see the chapter **Graph Properties** on pg 258.

Possible linearization formulas are *None* (no linearization), *1/Response*, *ln(Response)*, *sqrt(Response)*, *1/sqrt(Response)*, *Response^2*, *1/(Response^2)* and *log(Response)*.

Equation

Contains the equation of the calibration curve, which determines the relationship between the response and the actual amount of the compound according to the curve type fit selected.

Correlation Factor

A number from the interval <0, 1> that characterizes the closeness-of-fit of the calibration curve. The closer is the number to 1, the better is fit of particular points to the calibration curve. The calculated value is individual for each signal.

$$Cor.Coeff = \sqrt{\frac{(\sum (x_i - \bar{x}) * (y_i - \bar{y}))^2}{\sum (x_i - \bar{x})^2 * \sum (y_i - \bar{y})^2}}$$

where:

- x_i - the actual response at level i
- y_i - the calculated response at level i
- $\bar{x}$ - the mean actual response
- $\bar{y}$ - the mean calculated response

Residuum

The calculated value displays the square root counted from the sum of squares of the difference between the actual responses and the calculated responses.

Curve Check

The calculated value displays the result of the linearity check for the calibration curve. The item is only displayed when the *Curve Check* option is set in the [Calibration Options](#) dialog. The result may be either *OK* (if all performed *Curve Check* tests are passed) or *ERROR* (if any of the performed tests is failed).

Calibration curve

The calibration curve of any particular compound is located in the right-hand section of each compound tab.

Name of compound and retention time.

Shows the data which identify the displayed compound. These data come from the main calibration table and include *Compound Name*, *Reten. Time* and the name of the active signal.

Graph

Displays the calibration points of the current compound along with the calibration curve. The graph has identical properties to the other graphs, including cut-outs, auxiliary lines and optional setting using the *Properties* command from the local menu.

A calibrated point is indicated with a cross, a recalibrated point indicated with a star. In the *STDADD* calibration type, *Blank* sample is displayed as a vertical line along with multiplication cross sign to signal that the amount of the sample is not calculated from the intersection of the calibration curve with the X-axis on zero response, but from the intersection of the calibration curve with the blank horizontal line.

When using the *ISTD* method, a relative calibration curve will be displayed, depending on whether the same or a different amount of the internal standard has been used:

Same amount of internal standard

Relative calibration curve as dependence of response ratio of compound and internal standard on amount of injection of compound.

$$\frac{R_i}{R_{ISTD}} = f(A_i)$$

R_i - response of i-th compound

R_{ISTD} - response of internal standard

A_i - amount of i-th compound

Different amount of internal standard

Relative calibration curve as dependence of response ratio of compound and internal standard on ratio of amount of compound and internal standard.

$$\frac{R_i}{R_{ISTD}} = f\left(\frac{A_i}{A_{ISTD}}\right)$$

R_i - response of i-th compound

R_{ISTD} - response of internal standard

A_i - amount of i-th compound

A_{ISTD} - amount of internal standard

4.1 Calculate By

The **Calculate By** dialog allows to select the compound whose *Response Factor* will be used for the calculation of the *Amount* value instead of own response. Only peaks already added to the calibration are valid for selection.

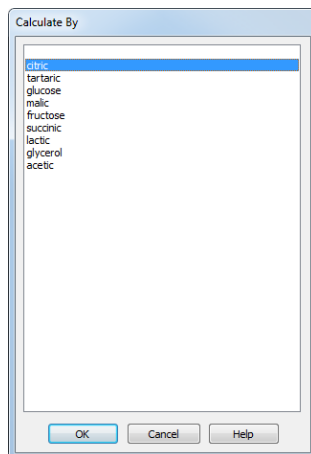


Fig 127: Calculate By

In case of ISTD calculation type, The ISTD compound may be used for calculation of the amounts for different peak, in which case the ISTD Amounts must be filled in both in the calibration file and in the unknown sample. It is not possible to use the *Calculate By* option for an ISTD peak (using the response of another compound for such peak).

4.2 File

This is an entry point to a submenu of commands for working with calibration files:

	New	Creates a new calibration file and at the same time closes the currently opened file. May also be invoked by pressing the Ctrl + N shortcut. New sequence will be based on the template calibration file (TEMPLATE.CAL) from the common directory (C:\CLARITY\COMMON\ by default). If the template calibration file does not exist in the project, the new calibration will be based on values internally preset in Clarity.
	Open	Selects and opens a calibration file. Invoking the command or using the Ctrl + O shortcut opens the Open Calibration dialog, which is identical to the Open Method dialog described on pg 44. The selected calibration file is loaded and its name, along with the calculation method, is displayed in the dialog title bar.
	Save	This command or the Ctrl + S shortcut saves the current calibration file. If a new file is to be saved, the Save As dialog will open where the name of the new file has to be filled in. The default name NONAME.CAL is reserved only for files that have not been saved yet.
Save As		Saves the calibration file under a chosen name (Ctrl + Shift + S).
Send Calibration by email		Sends the displayed calibration file as an email attachment. It is analogous to the <i>Send Method by Email</i> command (for more details see the chapter Send Method by Email on pg 47).

Export Calibration Curve as Picture to Clipboard	This command will copy the calibration curve picture into the Windows Clipboard, from which it can be instantly copied by using standard <i>Paste</i> command (Ctrl+V). This allows user to quickly utilize picture in another software, e.g. spreadsheet or word processor. This command is also available from the context menu invoked by right mouse-click in the calibration curve area.
Export Calibration Curve as Picture...	This command will open Save As dialog, which will save the picture of current calibration curve into Enhanced Metafile Format (.EMF). This format is extensively supported by Microsoft Windows and Microsoft Office. This command is also available from the context menu invoked by right mouse- click in the calibration curve area.
 Open Standard	Selects and opens a chromatogram that will be used as a calibration standard. Invoking the command opens the Open Chromatogram dialog similar to the Open Method dialog described on pg 44., which offers (by default) the content of the calibration subdirectory. The selected chromatogram is subsequently loaded and its name shown in the right-hand side of the status bar.
 Close Standard	Closes the displayed calibration standard.
Open Standard as Chromatogram	Opens the Standard in the Chromatogram window. Available only if the Standard is opened.
Link Calibration To Standard	Links currently opened chromatogram of standard to calibration file in which this standard is used. This will enable showing the names of peaks, as they were coined in the calibration file.
 Report Setup	Selects a report style (Ctrl + Alt + P). The command opens the Report Setup dialog described on pg 333.

 Print Preview	Displays a preview of the printout prior to actual printing (Ctrl + Shift + P). The command opens the Print Preview dialog described on pg 352 .
 Print to PDF	Prints the report to *.PDF file.
 Send Printed PDF by Email	Opens new email message in default email client, prints the report to *.PDF and adds created *.PDF file as an email attachment.
 Print	Selects a printer and starts actual printing (Ctrl + P). Invoking the command opens the Print dialog described on pg 354 .
-----	List of recently opened files so that you can easily get to the method files you are currently working with. The number of files in the list is determined by the <i>Recent Files</i> parameter in the User Options - General tab available from the Instrument window by <i>View - Options...</i> command.
Exit	Ends the work with the calibration standard and closes the Calibration window. A message will be displayed if the current calibration file has been changed and not saved yet. Modifications that are not saved are always marked by the inscription MODIFIED after the filename.

4.2.1 Save As...

This command (or the **Ctrl** + **Shift** + **S** shortcut) saves the opened calibration file under a newly desired name. The command opens the **Save As** dialog, similar to the **Open Method** dialog described on pg 44., where the new name can be entered or a new directory selected.

If a file of that name already exists and has been opened in some other Instrument or window, the name cannot be used. Moreover, in the protected mode (in full version ☒ Full version only) a user may not overwrite a file he/she is not entitled to modify (for more details see the chapter **User Accounts** on pg 3).

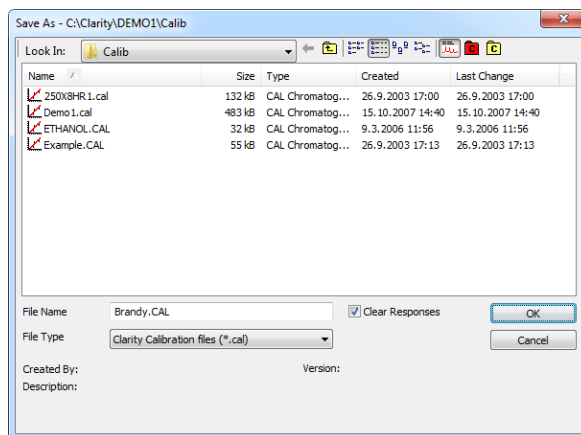


Fig 128: Save As (Calibration)

Clear Responses

The *Clear Responses* checkbox enables to clear all responses in the new calibration file.

4.3 Edit

Commands for working with the clipboard and individual tables.

 Undo	Cancels the modifications effected in tables (Ctrl + Z).
 Redo	Countermands the <i>Undo</i> command (Ctrl + Shift + Z).
 Cut	Moves the selected text to the clipboard (Ctrl + X).
 Copy	Copies the selected text to the clipboard (Ctrl + C).
 Paste	Copies the clipboard content to the indicated place (Ctrl + V).
Fill Down	Fills down all fields in the actual column with the value set in the active cell (Ctrl + F).
Fill Series...	Displays the Fill Series... dialog for filling up incremental values to cells.
Select All	Selects all cells of the table (Ctrl + A).
Restore Default columns	Restores the default view of columns.
Setup Columns...	Using this command invokes the Setup Columns dialog, where you can set the display of columns.
Hide Column(s)	Hides selected columns. You can select columns by left-clicking the mouse over the column header, more columns can be selected while holding the Ctrl key.
Show Hidden Column(s)	Shows columns hidden by <i>Hide Column(s)</i> command or Setup Columns dialog.

Detailed description of the commands from the *Edit* menu can be found in the chapter "**Edit**" on pg 148.

4.4 Display

This is an entry point to a submenu of commands for displaying chromatograms.

	Preview Graph	Displays/hides the orientation graph.
	Zoom Out	Zooms out the current part of chromatogram. Also invoked by pressing  key.
	Zoom In	Zooms in the current part of chromatogram. Also invoked by pressing  key.
	Previous Zoom	Restores the previous chromatogram cut. ( + )
	Next Zoom	Restores the original cut abandoned by the <i>Previous Zoom</i> command. ( + )
	Unzoom	Cancels all cuts. ( + )
	Tables Up	Swaps the placement of the graph with tables.
	Properties	Controls the displayed graph and chromatogram(s). If you are on the Compounds tab, this command will invoke the Graph Properties - Graph dialog described on pg 160. If you are on the Individual Compound tab, the Graph Properties dialog will be opened. This command is also available from the context menu invoked by right mouse-click in the calibration curve area.

Detailed description of the *Preview Graph*, *Previous Zoom*, *Next Zoom*, *Unzoom* and *Tables Up* commands is in the chapter "**Display**" on pg **156**.

4.4.1 Graph Properties

Sets the characteristics for displaying the graph of the calibration curve on the tabs of individual compounds.

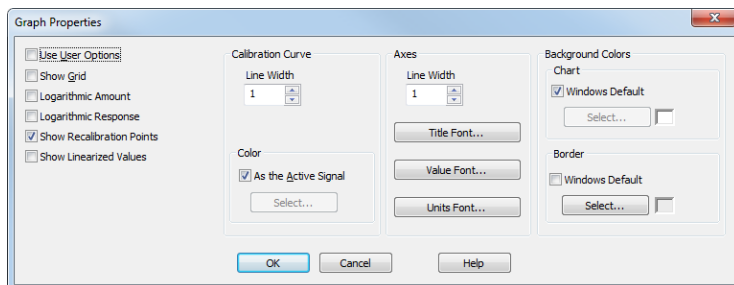


Fig 129: Graph Properties

Use User Options

All parameter settings will be taken from the global setting in the [User Options](#) dialog - [Graph](#), [Axes Appearance](#) and [Signals & Curves](#) tabs. These tabs are accessible from the [Instrument](#) window by using the [View - Options...](#) command.

This option is also available directly from the context menu invoked by right mouse-click in the calibration curve area.

Show Grid

Displays/conceals an auxiliary grid to help reading the values from a graph.

Logarithmic Amount

Sets the logarithmic axis for Amount.

Logarithmic Response

Sets the logarithmic axis for Response.

Show Recalibration Points

Displays calibration points along with all of their older versions.

Show Linearized Values

Displays the axes (both X and Y axis) in a linearized form of particular values of calibration points instead in the non-linearized form. Only effective when any linearization (*Linearization X* or *Linearization Y*) is set on the tab of a particular compound.

Calibration Curve

Line Width - Defines the width of calibration curve in points.

Color - Selects the color of the calibration curve.

As the Active Signal - Uses the color of the current signal.

Select... - Selects any desired color.

Axes

Line Width - Sets the width of a line in points.

Title Font... - Selects the font for the graph axes title.

Value Font... - Selects the font for the numerical figures on the axes.

Units Font... - Selects the font for the displayed units of the axes.

Background Colors

Chart - Sets the graph background color (the surface between axes).

Windows Default - Sets the Windows system initial color.

Select... - Selects the desired color.

Border - Sets the window background color (the surface of the outside axes).

Windows Default - Sets the Windows system initial color.


Select... - Selects the desired color.

4.5 Calibration

Commands for creation of calibration and setting parameters.

Set Level	Sets the level to be displayed in the Calibration Summary Table. This serves the identical purpose to the one for level setting in the <i>Calibration</i> toolbar. The selected level is indicated by the heading above the last four columns of the Calibration Summary Table. While the global Calibration Summary Table displays all substances on one of the selected levels, the tabs of individual substances will display all levels of that one particular substance.
Set Signal	Selects the active signal (when using multi-detector calibration). The sub-menu contains a list of available signals in the form of names taken from the signal names used for calibration in the first calibration standard. You can also select the active signal by clicking on the square with the color of the required signal in the <i>Calibration</i> toolbar. The name of the active signal is displayed in the header of the calibration table and in the header of the calibration curve next to the name and retention time of the active compound.
 Add All	Adds or updates all peaks and groups in the calibration standard into the calibration.
 Add Existing	Updates all peaks and groups of peaks that are already in the calibration with the data in the calibration standard.
 Add Peak	Adds or updates the peak selected in the calibration standard into the calibration.
 Add Groups	Adds or updates a group of peaks selected in the calibration standard into the calibration.

	Delete Compound	Deletes a selected compound from the calibration file. After invoking the <i>Calibration - Delete Compound</i> command or the  icon, use the cursor to select the peak of the compound to be deleted. The command can be aborted anytime by using the Esc key. Compounds may also be deleted from a calibration by deleting the corresponding line from the global table of compounds by using the Delete key or the <i>Delete Compound</i> command from the local menu.
Delete All Compounds		Deletes all compounds from the calibration file. Use of the commands in the Compounds tab cannot be undone using the <i>Undo</i> command. Before using the command a warning message will appear informing you of this condition.
Clear All Responses		Clears responses on all levels. Use of the commands in the Compounds tab cannot be undone using the <i>Undo</i> command. Before using the command a warning message will appear informing you of this condition.
Clear All Levels		Clears all responses and amounts on all levels. Use of the commands in the Compounds tab cannot be undone using the <i>Undo</i> command. Before using the command a warning message will appear informing you of this condition.
Clear Selected Level		Clears all responses and amounts on the selected level. Use of the <i>Clear All Responses</i>, <i>Clear All Levels</i> or <i>Clear Selected Level</i> command on the Compounds tab cannot be undone using the <i>Undo</i> command. Before using the command a warning message will appear informing you of this condition.
Show History		Invokes the History of Calibration Point dialog with a history log of the selected calibration point. Stored information includes: the name of the calibration standard, peak retention time, area, height and date of a (re)calibration.

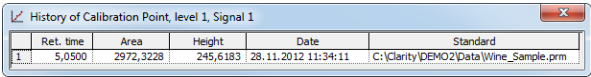

	Ret. time	Area	Height	Date	Standard
1	5,0500	2972,3228	245,6183	28.11.2012 11:34:11	C:\Clarity\DEMO2\Data\Wine_Sample.prm

Fig 130: History of Calibration Point

Note: Changing the response of a compound manually on any level that includes at least one recalibration completely clears the history of a given calibration point. You will be warned by a warning message before doing so.



Options...

Sets the calibration or recalibration method and implicit values.

4.5.1 Add All

Adds or updates all peaks and groups from the calibration standard into the calibration. Updating refers to recalibration or calibration on further levels. Unlike the *Add Existing* command, the *Add All* command will also add new compounds to the calibration during an update.

When using a multi-detector configuration, the command is always valid only for the active signal.

In the **Calibration Summary Table**, peaks are searched for according to retention time and identification windows. Groups are identified by a letter.

Dependence on type and method of calibration:

- When using automatic calibration, all compounds will be brought into the first available calibration file without regard to the current level.
- When using automatic recalibration, compounds are recalibrated on the current level. Further adjustments and supplements to the data can be performed later at any time.
- When using manual calibration or recalibration, the **(Re)Calibrate Peak** or **(Re)Calibrate Group** dialog will be displayed for direct checking and amendment of data for each compound.

Note: If you hide the part of the **Calibration** window with the calibration standard, for example by reducing the **Calibration** window or enlarging the **Compound Summary Table**, the *Add All*  command will not be active.

4.5.2 Add Existing

Adds all (or updates existing) peaks and groups from the calibration standard into the calibration. The behavior of the command depends on whether or not the calibration file already contains compounds:

- When creating a new calibration file, data on all peaks will be transferred from the calibration standard.
- When using calibration of a further level or recalibration, only data from already calibrated peaks and groups will be transferred. If you wish to add new compounds to the calibration file, you must use the *Add All*, *Add Peak* or *Add Group* commands.

Caution: When using a multi-detector configuration, only the active signal will be calibrated.

In the **Calibration Summary Table**, peaks will be searched for according to retention time and identification windows. Groups are identified by a letter.

Dependence on type and method of calibration:

- When using automatic calibration, all compounds are brought into the first completely free level in the calibration file without regard to the current level.
- When using automatic recalibration, compounds are recalibrated at the current level. Further adjustments and supplements to the data can later be performed at any time.
- When using manual calibration or recalibration the **(Re)Calibrate Peak** or **(Re)Calibrate Group** dialog for the direct inspection and amendment of data will be displayed for each compound.

Note: If you hide the part of the **Calibration** window with the calibration standard, for example by reducing the **Calibration** window or enlarging the **Compound Summary Table**, the *Add Existing*  command will not be active.

4.5.3 Add Peak

Use the *Calibration - Add Peak* command (or  icon) to add the selected peak to the calibration file. Invoking the command or icon moves the cursor to the graph of the calibration standard, where the desired peak should be selected. Use the **Esc** key to exit the command at anytime.

When using automatic calibration, the first available level of the calibration file will be calibrated regardless of the current level. When using automatic recalibration, the current level is always recalibrated. Data can be modified and supplemented at any later time. In manual (re)calibration, the **(Re)Calibrate Peak** dialog will open once a peak has been selected to enable immediate control and editing of the data.

In the **Calibration Summary Table** peaks are identified by their retention times and identification windows. If a peak is not identified in automatic (re)calibration, it is inserted as a new compound at the first level.

Note: If you hide the part of the **Calibration** window with the calibration standard, for example by reducing the **Calibration** window or enlarging the **Compound Summary Table**, the *Add Peak*  command will not be active.

4.5.3.1 Calibrate Peak

This dialog is opened by using the *Calibration - Add All* () or *Add Peak* () commands in manual calibration. It serves for editing/supplementing data transferred from the chromatogram to the calibration. The retention time of the calibrated peak will be displayed in the title bar.

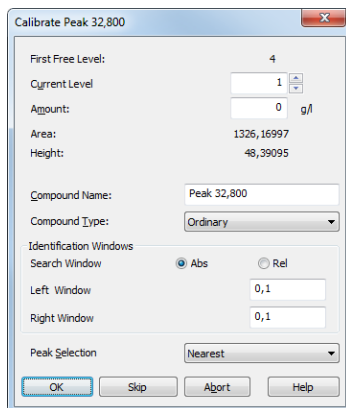


Fig 131: Calibrate Peak

First Free Level

Indicates the first free level for the displayed compound.

Current Level

Displays the level as it has been currently set. Click the arrows to select another level. The default value will differ depending on which command was used to invoke this window:

- **Add Peak** ( icon): By default, the *Current Level* command will be set to the value of the **First Free Level** parameter.
- **Add All** ( icon): By default, the *Current Level* command will be set to the first entirely free level. This new level will be also remembered for all remaining compounds.

Amount

Indicates the amount of the calibration compound. The units are taken over from the *Quantity Units - Compound* field in the [Calibration Options](#) dialog and are common to the entire calibration file.

Caution: In multi-detector calibration, the *Amount* column is common to all detectors. The amount can be modified on any level, it will always apply to all signals.

Area

Peak area in mV.s as the response.

Height

Peak height in mV as the response.

Compound Name

Refers to the calibrated peak and is common to all levels of the calibration file.

Compound Type

Defines the function of the peak in the calibration. The compound type may be:

Ordinary - Designates a peak that is neither a reference peak nor an internal standard.

Reference - Designates a reference peak.

Internal Standard - The given peak is an internal standard.

Reference ISTD - Designates a reference peak that is at the same time the internal standard.

Note: Only one compound in a calibration file may be marked as the internal standard. A group of peaks can neither be the internal standard nor a reference peak.

Identification Windows**Search Window**

Set *Abs* if the search window will be set absolutely or *Rel* if it will be set according to *Calculated By* field.

Left (Right) Window

Refers to the width of the left/right identification window and is common to all levels of the given peak. For an identified peak the value is taken from the global calibration table, otherwise the default value set by the *Identification Window* fields of the [Calibration Options - Defaults](#) dialog will be displayed. The value set there is used for identification purposes in automated or manual (re)calibrations and in all calibration calculations.

Peak Selection

Defines the way the correct peak is matched with the Retention time when several peaks are present in the same identification window. The options available are *Nearest* (closest possible to the selected retention time - default for standard peaks), *First*, *Last* and *Biggest* (default for Reference peaks).

Skip

Skips the current substance when calibrating using the *Calibration - Add All* (📈 icon) command.

Abort

When a calibration has been effected by the *Calibration - Add All* (📈 icon), this command aborts calibration of the remaining compounds.

4.5.3.2 Recalibrate Peak

This dialog is opened by using the *Calibration - Add All* (📈 icon) or *Add Peak* (📈 icon) commands in manual recalibration. It serves for editing/adding data transferred from a chromatogram to the calibration file. The retention time of the recalibrated peak is displayed in the title bar. The recalibration type is governed by the items in the *Recalibration* section of the [Calibration Options](#) dialog.

Recalibrate Peak 1,837

Current Level: 1 Recalibration: 1

Compound Name: malic

	Old Value	New Value	Units	Departure	
				Units	%
Area	4004,496	4030,496	[mV.s]	0,000	0,0
Height	908,457	915,457	[mV]	0,000	0,0
Amount	0,000	0,000	ng/mL		

OK Skip Abort Help

Fig 132: Recalibrate Peak

Current Level

Contains the current level that is set on the toolbar above the **Calibration Summary Table** in the [Compounds](#) tab of the [Calibration](#) window by

default. You can edit the number or use the up and down arrows near it to select any other level; in case of the *Add All* command that level is then retained for all compounds.

Recalibration

Indicates the number of recalibrations of the calibrated compound at the current level. In a recalibration that has been effected according to the *Replace* option set in the [Calibration Options](#) dialog or when the manual modification of the response value has been performed, the number of recalibrations will be set to 1.

Compound Name

Specifies the component name corresponding to the calibrated peak. The name is common to all levels of the calibrated compound.

Recalibration Table

Displays and nominally modifies the recalibration data items. All values are displayed for height, area, and amount.

Old Value

The last value prior to recalibration.

New value

The current value which will be used for the recalibration. The resulting value for a given recalibration point might be different if the recalibration has *Average* or *Weight* option switched in the [Calibration Options](#) dialog.

Units

Units of all displayed values.

Departure

Indicates the absolute and relative deviation of the value gained from the calibration standard and the last value in the calibration file.

Skip

Skips the current substance when a recalibration has been made using the *Calibration - Add All* command.

Abort

When a recalibration has been effected by the *Calibration - Add All* command, the recalibration of the remaining compounds will not be performed.

4.5.4 Add Group

Use the *Calibration - Add Group* command (or  icon) to add the selected peak group to the calibration file. Invoking the command or icon moves the cursor to the graph of the calibration standard, where the desired group should be selected. Use the **[Esc]** key to exit the command at anytime. When using a multi-detector configuration, the command will only effect the current signal.

When using automatic calibration, all data items will be assigned to the first free level of the group regardless of the currently set level. In automatic recalibration, the current level is always recalibrated. Data can be modified and supplemented at any later time. When using manual (re) calibration, the **(Re)calibrate Group** dialog will open once the group has been selected and enable immediate control and modification of data. Groups are identified by the corresponding letter.

Note: If you hide the part of the **Calibration** window with the calibration standard, for example by reducing the **Calibration** window or enlarging the **Compound Summary Table**, the *Add Group*  command will not be active.

4.5.4.1 Calibrate Group

Adds data from the calibration standard to the calibration file and possibly modifies them. This dialog is either opened by the *Calibration - Add All* ( icon) or *Add Group* ( icon) commands in manual calibration mode. The identifier of the selected group will be displayed in the title bar. The dialog contains items analogous to those shown in the **Calibrate Peak** dialog described on pg 265.

4.5.4.2 Recalibrate Group

Adds data from the calibration standard to the calibration file and possibly modifies them. This dialog is either opened by the *Calibration - Add All* ( icon) or *Add Group* ( icon) commands in manual recalibration mode. The identifier of the selected group will be displayed in the title bar. The dialog contains the (meaningful) items analogous to those shown in the **Recalibrate Peak** dialog described on pg 267.

4.5.5 Options

This section contains commands for selecting the (re)calibration mode. Invoking the command or using the  icon opens the [Calibration Options](#) dialog. Settings performed here are then stored in the calibration file. When a new calibration file is created, default values for all of its options are loaded from here.

Calibration Options - Calibration Options

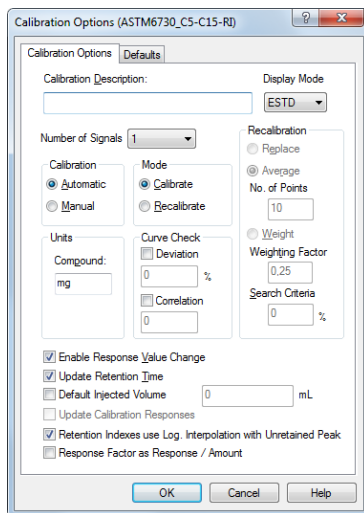


Fig 133: Calibration Options - Calibration Options

Calibration Description

Contains a brief description of the calibration file. When a calibration file is loaded, this text is displayed in the *Description* field of the [Open Method](#) dialog; it is also shown in the *Description* field of the [Method Setup - Calculation](#) dialog when a calibration is loaded there.

Display Mode

Sets the method of the calibration curve display (suitable for different types of calibration curve use):

ESTD

Displays the dependence of a compound's response to its quantity.

ISTD

Relative calibration curve:

- If the amount of the internal standard has not been entered, displays the dependence of compound/internal standard response ratio to amount of compound.

- If the amount of the internal standard has been entered, displays the dependence of compound/internal standard response ratio to the compound/internal standard amount ratio.

NORM

Displays the dependence of a compound's response to its quantity, just like for the *ESTD* type.

STDADD

Displays the dependence of a compound's response to its quantity, but the graph is shifted so that the calibration curve crosses the x-axis.

Caution: This parameter only indicates the display method. The actual type of calculation upon evaluation of unknown compounds can be selected in the [Chromatogram](#) window in the [Result](#) tab using the *Calculation* parameter.

Number of Signals

Selects the number of displayed signals when a multi-detector calibration is used. The implicit value is 1. The change is manifested in the number of commands in the *Calibration - Set Signal* menu and in the number of colored squares on the [Overlay](#) toolbar.

Note: The number of signals will also be adjusted without regard to the setting of *Number of Signals* field if the currently set calibration standard contains more signals than is set here.

Note: You do not risk deleting the calibration values when you reduce the number of signals, since the file always remembers the values for all 16 possible signals.

Calibration

Sets the calibration or recalibration method for the *Add All*, *Add Peak* and *Add Group* commands. Possible options are:

Automatic

Automatically transfers all data related to compounds from a chromatogram to the calibration file. The data can be modified and supplemented later on.

Manual

Enables to enter the amount of the component or to modify or add the remaining data for each compound by opening one of the [\(Re\) Calibrate Peak](#) or [\(Re\)Calibrate Group](#) dialogs.

Mode

Sets the mode of the calibration process, unless in the Standard Addition calibrations, where it is disabled.

Calibrate

Calibration mode. Data from the calibration standard will be transferred into the first empty calibration level. After all 20 calibration levels have been filled in, Clarity will rewrite the responses in the last (20th) calibration level.

Recalibrate

Recalibration mode. Unlocks the *Recalibration* section for setting up recalibration conditions.

Caution: The *Calibrate* mode applies only in the Single Analysis. In the Sequence the calibration always recalibrates applying the setting from *Recalibration* options.

Recalibration

Specifies the recalibration mode conditions:

Replace

The original response is replaced by the new one and the number of recalibrations is set to one.

Average

The resulting value is calculated as the arithmetic mean of the new value and all previous values (their number is taken over from the *Rec. No.* column of the *Compound Summary Table*).

No. of Points

Sets the number of recalibrations used for the Averaging. When the number of recalibrations exceeds this number, averaging calculates floating average.

Available range is 0 - 100, default value is 10.

Weight

The resulting value is calculated as a weighted average of the original and the new value. The statistical weight assigned to the new value is given by the content of the *Weight* field.

Weighting Factor

The statistical weight used to calculate the weighted average when the recalibration mode is *Weighted*.

Search Criteria

The maximum percent deviation (departure) of the new values for the recalibration to be effected. In the event it has been exceeded in manual recalibration the inscription *Over* replacing the numerical value of the deviation will notify you. When using automatic recalibration components whose deviation in height or area exceeds the pre-set value are not recalibrated. The allowed range is 1 to 50 %, enter the 0 % to disable the condition.

Units

Specifies the units for compound quantities.

Compound

Contains the units of component amounts. Units are stated for all displayed amounts in calibration tables and graphs as well as in the table of integration results in the [Results](#) tab, unless other units are specified by *Scale Factor* and *Units after Scaling* fields.

Curve Check

Enables to check the linearity of the calibration curve based on the two following parameters. The result of the linearity check is displayed on the tabs of each individual compound in the [Calibration](#) window and in case the *Curve Check* test is failed also in the *Peak Type* column of the **Result Table** in the [Chromatogram](#) window. Fail in any *Curve Check* test part that is switched on fails the whole *Curve Check*.

Deviation

Sets the value for the maximal allowed standard deviation of each point in the calibration curve of the given compound.

Correlation

Sets the value for the minimal allowed correlation factor as calculated on the tab of the individual compounds of the [Calibration](#) window.

Enable Response Value Change

Enables the manual modification of responses (in the *Response* or *Area* and *Height* fields).

Update Retention Time

Enables the automatic update of retention time after each calibration or recalibration. The retention time is calculated as the arithmetic mean of retention times established in all (re)calibrations of the given component at all levels.

Default Injected Volume

The value is used to perform automatic correction for differences in injected volume. In calibration, recalibration and calibration calculations the value of *Inj. Volume* from the chromatogram header is compared with *Default Injected Volume* and a new response is calculated from their ratio. This correction method is only applicable to small deviations since it assumes that the amount is proportional to the response. If the checkbox is left unchecked, the last recalculated values remain in effect regardless of potential changes effected in *Injected Volume* and *Default Injected Volume* fields. In this instance the last two parameters serve only for information.

Update Calibration Response

If the *Default Injected Volume* is changed from non-zero value to another non-zero value, then, provided the *Update Calibration Response* checkbox has been checked, all responses will be recalculated in the proportion of the new and old value of *Default Injected Volume*. These changes occur when closing the [Calibration Options](#) dialog.

Retention Indexes use Log. Interpolation with Unretained Peak

Use *Unretained Time* value (set on the [Method Setup - Advanced](#) tab) for retention index calculation (for more details see the chapter **Calculation** on pg 88).

Caution: When this option is unchecked, then the retention indexes are calculated using the **linear interpolation** and when it is checked, then the **logarithmic interpolation with unretained peak time** is used.

Equation used for the linear calculation of the Retention Indexes:

$$I = 100 \times \left[n + (N - n) \frac{t'_{r(\text{unknown})} - t'_{r(n)}}{t'_{r(N)} - t'_{r(n)}} \right]$$

Where:

- I = Kovats retention index,
- n = the number of carbon atoms in the smaller n-alkane,
- N = the number of carbon atoms in the larger n-alkane,
- z = the difference of the number of carbon atoms in the smaller and larger alkane,
- t'_r = the retention time.

Equation used for the logarithmic calculation of the Retention Indexes:

$$I = 100 \times \left[n + (N - n) \frac{\log(t'_{r(\text{unknown})}) - \log(t'_{r(n)})}{\log(t'_{r(N)}) - \log(t'_{r(n)})} \right]$$

Where:

- I = Kovats retention index,
- n = the number of carbon atoms in the smaller n-alkane,
- N = the number of carbon atoms in the larger n-alkane,
- t'_r = the adjusted retention time ($t'_r = t_r - t_0$).

Response Factor as Response / Amount

Enables calculating Response Factor as RF = Response / Amount (on default : RF = Amount / Response). This setting changes the format of RF in [Calibration](#) window as well as on [Results](#) tab of the [Chromatogram](#) window.

4.5.5.1 Options

Calibration Options - Defaults

Defines parameters that will be used for new compounds.

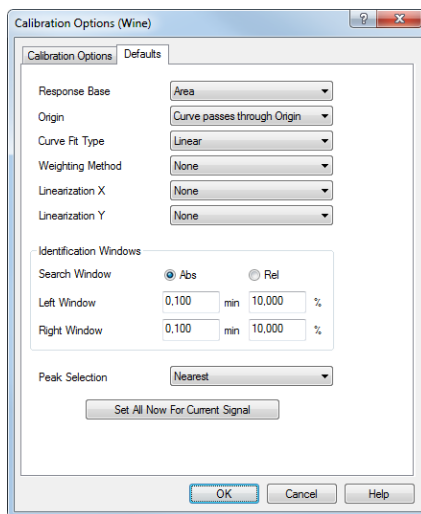


Fig 134: Calibration Options - Defaults

Response Base

Indicates the type of response used: *Area*, *Height*.

Origin (Ending Point)

Defines the treatment of the origin point in the calibration graph (for more details see the chapter **Calibration** on pg 247).

Curve Fit Type

Indicates the type of curve fitted to the data (for more details see the chapter **Calibration** on pg 246).

Weighting Method

Defines the weighting method of calibration points upon interlining of the curve (for more details see the chapter **Calibration** on pg 248).

Linearization X

Defines the default X-axis display on the tabs of individual compounds. Possible values are *None*, *1/Amount*, *ln(Amount)*, *sqrt(Amount)*, *1/sqrt(Amount)*, *Amount^2*, *1/Amount^2* and *log(Amount)*.

Linearization Y

Defines the default Y-axis display on the tabs of individual compounds. Possible values are *None*, *1/Response*, *ln(Response)*, *sqrt(Response)*, *1/sqrt(Response)*, *Response^2*, *1/Response^2* and *log(Response)*.

Identification Windows

This section handles the defaults for the identification windows under the *Abs* and *Rel Search Window* conditions. These conditions are more closely described in the chapter "**Calibration**" on pg 239.

Search Window

Defines the default calculation type of both Left and Right identification window boundaries. Defines whether the Left Window and Right Window content will be inserted in *Abs* (absolute - set in minutes) or *Rel* (relative - set in percents of the Reten. Time) values.

Left (Right) Identification Window

Defines the width of the left (right) identification window (in minutes or percents).

Peak Selection

Defines how the target peak will be selected in a given identification window. The available options are the peak *Nearest* to the exact retention time set in the calibration, *First* peak in the identification window, *Last* peak in the identification window and the *Biggest* peak in the identification window.

Note: The peaks that are identified must still follow the trend of rising retention times, so when two identification windows are selected so that they overlap each other, the first compound must define a peak with lower retention time than the second compound. Moreover, each peak in the calibration must be defined using a unique retention time.

Set All Now For Current Signal

Sets the selected parameters for all compounds in current calibration and current signal.

Caution: Pressing the *Set All Now For Current Signal* button is different from pressing the *OK* button, which will cause the use of set parameters for future signals and compounds.

4.6 View

This section contains commands that govern the appearance of the **Calibration** window.

Toolbar	Displays/hides all toolbars. The command is similar to the <i>Toolbar</i> command in the main Clarity window (for more details see the chapter View on pg 30).
Customize...	Used to customize the toolbars. The command is similar to the <i>Customize...</i> command in the main Clarity window (for more details see the chapter Customize... on pg 31).
Reset All	Sets all menus and toolbars in the Calibration window into the initial state.
Maximize Graph	Enlarges the chromatogram over the whole window or conceals the Compound Summary Table underneath. This may also be performed by double-clicking on the edge of the graph close to the splitter where the cursor arrow changes to  . This command is only valid for the Compounds tab.
Show Both	Displays both parts of the window (chromatogram and tabs). This command is only valid for the Compounds tab.
Maximize Table	Enlarges the part of the window with the Compound Summary Table over the whole window or conceals chromatogram. This action may also be performed by double-clicking on the edge of the table section close to the splitter where the cursor arrow changes to  . This command is only valid for the Compounds tab.

4.7 Window

This is an entry point to a submenu of commands for displaying other main windows.

Main 	Displays the Clarity main station window described on pg 1 .
Instrument	Displays the Instrument window described on pg 40 .
Chromatogram	Displays the Chromatogram window described on pg 118 .
Single Analysis	Displays the Single Analysis window described on pg 59 .
Sequence	Displays the Sequence window described on pg 294 .
Data Acquisition	Displays the Data Acquisition window described on pg 279 .
Device Monitor	Displays the Device Monitor window described on pg 55 .
Station Audit Trail 	Displays the Audit Trail window described on pg 318 .
Calibration Audit Trail 	Displays the Calibration Audit Trail window described on pg 327 .

5 Data Acquisition

Displays the signal from detectors configured on the given Instrument. The **Data Acquisition** window can be opened using the *Monitor - Data Acquisition* command or using the  icon in the **Instrument** window. During the measurement the name of the measured chromatogram is displayed in the window header.

Note: It is possible to switch between signals and hide/display them in the same way as in the **Chromatogram** window.

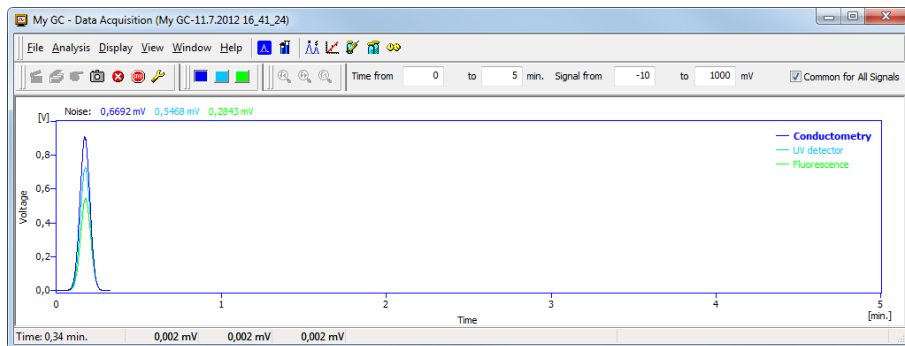


Fig 135: Data Acquisition

Note: When the analysis is started, it takes some time to initialize the detector (200 ms for **DataApex** Converters). In this time, the data is not valid.

The window contains two fundamental parameters, **Time** and **Signal**. In the default setting (if you have the *Context Axes* item set in the *View* menu) the function of these two parameters will be dependent on whether an analysis is currently taking place:

a) Analysis not running (state *STOP*, *WAITING* or *IDLE*):

The axes are set as *Floating Axes*.

Time Range

Sets the width of display window in minutes (0.5 - 20 min). Should the signal exceed the set time, the window will be reset and the signal will once again be displayed from the start.

Signal Range

Sets the height of display window in units set for the given active signal. (in range 0.01 - 1000000000 mV or equivalent units). The window dynamically monitors the course of the signal and is automatically redrawn after reaching the upper or lower edge for the given signal.

b) Analysis running (state *RUNNING*):

The axes are set as *Fixed Axes*. This means that if the current time or the level of the signal exceeds any of values set, the window is not redrawn and the signal is not displayed.

Time: from... to...

Sets the fixed range of displayed time in minutes.

Signal: from... to...

Sets the fixed range of the signal in units selected for the given active signal.

Note: Automatic switching between the fixed and floating range of axes, depending on whether or not analysis is taking place, may be canceled by switching the axes display in the *View* menu from *Context Axes* to *Fixed Axes* or *Floating Axes* (for more details see the chapter **View** on pg 289).

Regardless of the axes display type selection, there is the *Common for All Signals* checkbox on the toolbars. When checked, the axes range settings will be common for all signals displayed in the **Data Acquisition** window. Otherwise, the change will be performed for the active signal only.

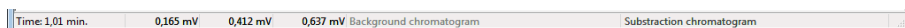
Status Bar

Fig 136: Data Acquisition - Status Bar

The **Data Acquisition** window contains the following items in the status bar:

Time [min.]

Displays the current time of an analysis in minutes (only during the measurement).

Current Signals

Displays the immediate value of all signals connected to the Instrument, along with appropriate signal units.

Background Chromatogram ✓ Full version

If it has been defined, using the *File - Set Background Chromatogram* command, the name of the background chromatogram will appear in the status bar. If the workstation is in the *WAITING* state, the chromatogram will not be displayed and its title in the status bar will be shaded gray.

Subtraction Chromatogram ✓ Full version

If it has been defined in the **Method Setup - Advanced** dialog and the display has been set using the *File - Show Subtraction Chromatogram* command, the name of the subtraction chromatogram will appear in the status bar.

5.1 File

A submenu of commands used for displaying chromatograms in the background.

Set Background Chromatogram 	Displays the background chromatogram on the background of the Data Acquisition window. Invokes the Open Chromatogram dialog, in which you may select which of the measured chromatograms you would like to display on the background of the running analysis. The background chromatogram is colored gray in order to prevent it from being mistaken with the measured signal. It will only be displayed after the measurement has been started.
Close Background Chromatogram 	Switches the display of the background chromatogram off.
Show Subtraction Chromatogram 	Determines whether the subtracted chromatogram defined in the <i>Subtraction</i> section of the method file (Method Setup - Advanced) will be displayed in the background of the Data Acquisition window. The checkmark will appear before the command if the display is activated.
Exit	Closes the Data Acquisition window.

5.2 Analysis

Submenu with commands for controlling the running of the analysis.



Send Method

It sends the method to all connected control modules/hardware. This button is not available during data acquisition.



Run Single

Starts individual measurement. A change of color in the background of the graph in the [Data Acquisition](#) window and the commencement of time counting in the status bar will indicate a running measurement. This command can be selected during the measurement only when *External Start* setting in the [Method Setup - Measurement](#) dialog is set to *Start-Restart*. Invoking the command terminates measurement and at the same time starts another one. This feature is often called "Continuous Measurement". Invoking the command during analysis is possible only when there are analog detectors (INT7, Net-PAD, ...) on instrument. If there is one or more digital detectors on instrument, invoking the command during analysis is not possible. This command is inactive during sequence measurement.



Run Sequence

Starts the sequence measurement. Measurement will be performed according to the current sequence file in the [Sequence](#) window. A change of color in the background of the graph in the [Data Acquisition](#) window and the commencement of time counting in the status bar will indicate a running measurement. This command is inactive during analysis.



Resume Sequence

Resumes the operation of a suspended sequence. This function may also be invoked by pressing the **Ctrl** + **Space bar** shortcut (for more details see the chapter [Sequence](#) on pg 310).

**Snapshot**

Saves the data from the running analysis to a chromatogram without terminating the measurement. The snapshot chromatogram will be automatically replaced by the final chromatogram after the analysis is finished.

If the modifications of a snapshot chromatogram have to be preserved, it is necessary to save the snapshot chromatogram under a different name (using the *Save As* command) before the end of the analysis.

As of **Clarity** version 5.0, *Snapshot* is also available outside of the analysis to save the data measured from the point when the last analysis run ended. In case more than 300 minutes has passed from the end of the last analysis, only the last 300 minutes of the data will be stored.

**Stop**

Immediately terminates an individual measurement and processes it. The Instrument will return to the *STOP* or *WAITING* state. If the control program for HPLC pumps or other devices is still running, the *CONTROL* state will be displayed.

When running a sequence measurement, it will be terminated after the length of analysis set in the *Enable Autostop - Run Time* field of the [Method Setup - Measurement](#) dialog has elapsed.

**Abort**

Terminates the measurement without storing and processing any data. The Instrument will return to the *STOP* or *WAITING* state. Any control program for HPLC pumps will be stopped. Running gradient will be also stopped immediately and the device should be properly conditioned by the user before next analysis.

The sequence will be terminated without processing the current analysis. The sequence run can be continued from the actual point using the *Resume* command.

Set Autostop

Opens the [Method Setup - Measurement](#) dialog where the *Enable Autostop* and *Run Time* parameters can be set.



Set Zero

Sets the active signal value as offset. That means that the current value of the signal will be subtracted from the signal from now on. The offset value can be reset using *Reset Zero* command. The *Set Zero* function is only active when the measurement is not running.



Reset Zero

Resets the subtracted offset value set using the *Set Zero* command to 0.
The offset value is automatically reset when an Instrument is closed.

Note: The above commands are similar to the commands in the [Single Analysis](#) dialog described on pg 59.

5.3 Display

A submenu of commands for setting up the graph display.

Zoom Out	Zooms out the current part of chromatogram. Also invoked by pressing  key on numeric keyboard.
Zoom In	Zooms in the current part of chromatogram. Also invoked by pressing  key on numeric keyboard.
 Previous Zoom	Invoking the command, using the  icon or the Ctrl +  shortcut on numeric keyboard displays the previous cut-out. Use the command repeatedly to browse backwards through the history of the cut-outs - the workstation remembers cut-outs which have been performed since the Data Acquisition window was opened.
 Next Zoom	Invoking the command, using the  icon or the Ctrl +  shortcut on numeric keyboard progressively returns to original cut-outs, which had been abandoned using the <i>Previous Zoom</i> command. Use the command repeatedly to browse forwards through the history of the cut-outs. The workstation remembers cut-outs which have been performed since the Data Acquisition window was opened.
 Unzoom	Invoking this command, using the  icon or Ctrl +  shortcut on numeric keyboard cancels all cut-outs and displays chromatogram in its original size.
 Properties...	Opens the Graph Properties dialog for displaying the graph display settings.

5.3.1 Properties...

Sets the characteristics of the [Data Acquisition](#) window display. Invoking the command opens the [Graph Properties](#) dialog.

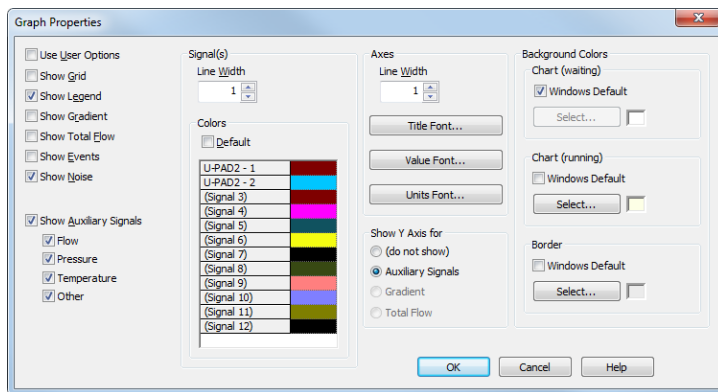


Fig 137: Graph Properties

Use User Options

All parameter settings will be taken from the global settings of the [Graph](#), [Axes Appearance](#) and [Signals & Curves](#) tabs of the [User Options](#) dialog, accessible from the [Instrument](#) window by using the *View - Options...* command.

Show Grid

Displays/conceals auxiliary grid for easier reading of values in graph.

Show Legend

Displays/conceals names of the individual signals including the eventual subtraction chromatogram and background chromatogram.

Show Gradient

Permits the display of gradient when using direct control of LC pumps.

Show Total Flow

Permits the display of total flow when using direct control of LC pumps.

Show Events

Permits the display of events from the **Event Table** in the [Data Acquisition](#) window.

Show Noise

Permits the display of **Noise** (along with appropriate units) in the upper left corner of the graph. Noise in the [Data Acquisition](#) window is calculated using the cubic interlining to reduce the influence of the trend to its values. In the [Chromatogram](#) window and [Results](#) tab the **Noise** values are calculated using the linear interlining on the interval defined in the **Integration** table. If there is a peak in this interval, it will influence the

noise values. If there are multiple intervals defined in the **Integration** table the last one will be applied.

Show Auxiliary Signals

This section determines which auxiliary signals will be displayed in the graph in the [Data Acquisition](#) window. Each auxiliary signal is included in one of following categories: *Flow*, *Pressure*, *Temperature*, *Other*. The display of all signals in each category is allowed together.

Signal(s)

Line Width

Selects the curve width in points.

Colors

Selection of color of the signal

Default - When checked, sets the display colors to the default state.

Select... - Click the particular color to change it.

Axes

Line Width

Selects the axes width in points.

Title Font...

Selects the font for graph axes title.

Value Font...

Selects the font for numerical figures on axes.

Units Font...

Selects the font for displayed units of axes.

Show Y Axis for

Determines whether the second (right) Y axis will be displayed and for which signal. The axis may be displayed for the **Gradient** (the percentage rate of individual compounds), **Total Flow** or **Auxiliary Signals**. The *Show Gradient* and *Show Total Flow* options must first be checked in order to display the Y axis for this particular signal. Applies only to the directly controlled devices with auxiliary signals.

Background Colors

Chart (waiting)

Sets the color of the graph background (surface of the graph between axes).

Windows Default - Sets the initial color of Windows system.

Select... - Serves for selecting of the desired color.

Chart (running)

Sets the graph background color (surface between axes) when the analysis is running. Setting a different colored background to the graph enables you to quickly differentiate as to whether an analysis is

actually running or if the detector signal is merely being monitored in the [Data Acquisition](#) window.

Windows Default - Sets the initial color of Windows system.

Select... - Serves for selecting of the desired color.

Border

Sets the color of the window background (surface outside axes).

Windows Default - Sets the initial color of Windows system.

Select... - Serves for selecting of the desired color.

5.4 View

Toolbars	Displays/conceals the toolbars in the Data Acquisition window.
Customize...	User setting of toolbars. The command is analogous to the <i>Customize...</i> command located in the Clarity window (for more details see the chapter Customize... on pg 31).
Reset All	Resets the menu and toolbars back to their initial state as they were from the initial installation of the Clarity .
Grid Lines	Displays/conceals auxiliary grid for easier reading of values in the graph.
Fixed Axes	Sets the fixed range axes display. The range of signal axis is determined by entering the lower and upper limit (e.g. <i>15 - 30mV</i>) both in the analysis and outside of it, the range of the time axis is determined by entering the time range (outside of the analysis) or lower and upper limit (during the analysis - e.g. <i>0 - 20min</i>).
Floating Axes	The range is defined by its size (e.g. <i>± 20 mV; 5 min</i>) While being displayed, the display will "float" so that the signal always stays in the displayed part of the graph. If the signal exceeds the y-axis range, the graph is moved vertically so the actual signal value will be in the center of the left axis. If the signal reaches end of the time range, the graph is redrawn from beginning.
Context Axes	Automatically switches between <i>Fixed Axes</i> and <i>Floating Axes</i> setting according to whether the workstation is measuring the analysis (in the <i>RUNNING</i> state) or is merely monitoring the signal (<i>WAITING</i> state). When measuring, the <i>Fixed Axes</i> display will automatically be used, while when in the <i>WAITING</i> state, the <i>Floating Axes</i> will be used.

Auto Axes	For the time axis, automatically switches between <i>Fixed Axes</i> and <i>Floating Axes</i> setting according to whether the workstation is measuring the analysis (in the <i>RUNNING</i> state) or is merely monitoring the signal (<i>WAITING</i> state). When measuring, the <i>Fixed Axes</i> display will automatically be used, while when in the <i>WAITING</i> state, the <i>Floating Axes</i> will be used. The signal axis behaves the same as when set to <i>Floating Axes</i> while the analysis is not <i>RUNNING</i>. While it is <i>RUNNING</i>, the range of the signal axis starts at a minimal value and enlarges according to the signal course; thus, the signal is always displayed in the graph and the view is not overly large at any moment of the analysis - the signal always fits in it.
Reset Axes Ranges	Sets the default values into all fields of the Axes Ranges dialog.
Set Axes Ranges...	Opens the Axes Ranges dialog used for defining the ranges of all signals in the Data Acquisition window, both in the run and outside of the analysis run, including ranges used for auxiliary signals.

5.4.1 Axes Ranges

The [Axes Ranges](#) dialog, serving for setting the axes ranges both during the acquisition run and outside of it, can be invoked using the *Set Axes Ranges...* command from the *View* menu of the [Data Acquisition](#) window.

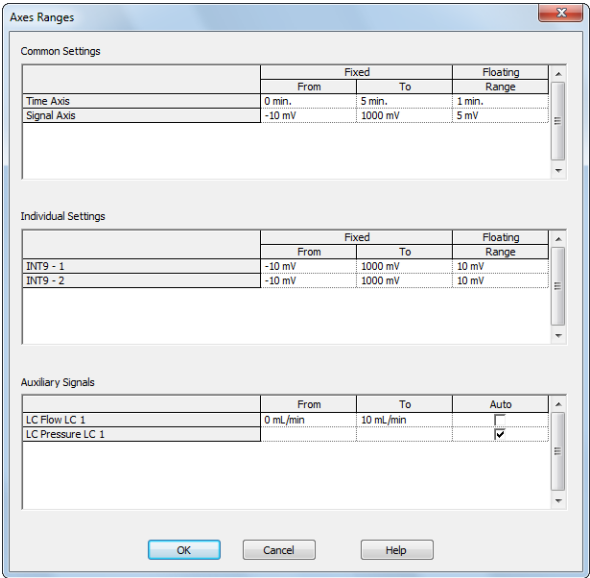


Fig 138: Axes Ranges

The settings performed in the [Axes Ranges](#) dialog are applied to all of the display possibilities in the [Data Acquisition](#) window - *Fixed Axes*, *Floating Axes*, *Context Axes* and *Auto Axes*. Only some of the settings are usually valid for each axes display type. Default values in the [Axes Ranges](#) dialog may be set by using the *Reset Axes Ranges* command from the *View* menu of the [Data Acquisition](#) window.

Common Settings

Sets the ranges for time axis and for the signal axes in case when *Common for All Signals* checkbox in the [Data Acquisition](#) window is selected. The values are set in appropriate value units for the given detector.

Time Axis

Sets the range of the time axis both in the mode where start display time (*From*) and end display time (*To*) are set and where the total range (*Range*) of the time axis is set. The time values are set in minutes.

Signal Axis

Sets the range for the signal axes, both for the regular and auxiliary signals.

Individual Settings

Sets the signal axis ranges for individual detectors.

Auxiliary Signals

Sets the auxiliary signal axis range for auxiliary detectors. When the *Auto* checkbox is checked, the auxiliary signal axis range starts at a minimal value and enlarges according to the auxiliary signal course. While unchecked, the axis range is fixed with the values entered.

5.5 Window

This is an entry point to a submenu of commands for displaying other main windows.

Main 	Displays the Clarity main window described on pg 1 .
Instrument	Displays the Instrument window described on pg 40 .
Chromatogram	Displays the Chromatogram window described on pg 118 .
Calibration	Displays the Calibration window described on pg 238 .
Single Analysis	Displays the Single Analysis window described on pg 59 .
Sequence	Displays the Sequence window described on pg 294 .
Device Monitor	Displays the Device Monitor window described on pg 55 .
Station Audit Trail 	Displays the Audit Trail window described on pg 318 .

6 Sequence

This Chapter describes the window which is used for consecutive measurement of multiple samples. Open the window using the  icon or the *Analysis - Sequence* command from the *Instrument* window.

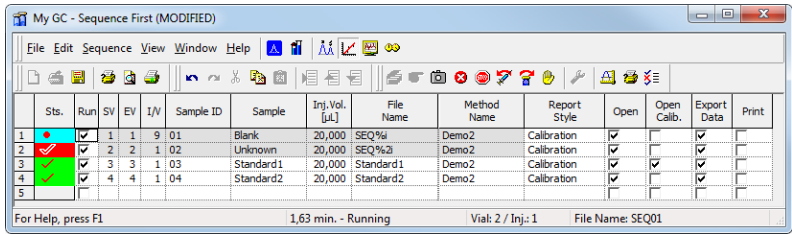


Fig 139: Sequence

The window includes a toolbar, sequence table and status bar.

Table description

Each line defines the setting for processing one or more injections. To process several injections according to a single line enter a different value for the initial = start (SV) and last = end (EV) vial. The I/V parameter indicates that there will be multiple injections taken from a single vial.

Sts.

Status of the row in the sequence. Each line can be in one of the following four states:

Tab 20: List of possible states of sequence rows

Status	Description
Green field	No analysis of this line has been (or is being) carried out.
Red field	Analysis from this line is being measured.
Dark blue field	All analyses on the line have been measured out. Already measured lines will be blocked against changes (dimmed).
Light blue field	Lines that have been measured but the <i>Skip Vial</i> command was used so the line is incomplete.

Moreover, each row is marked with a symbol indicating its correctness:

Tab 21: List of possible line correctness check results

	Status	Description
✓	Check mark	Indicates a line that doesn't contain any errors which could prevent a sequence from being run.
•	Bullet	Indicates the line where the validity check wasn't performed (e.g. line that has already been measured, omitted from a sequence etc.)
✗	Cross mark	Indicates the row with one or more errors on it. Such row is disabled from running in a sequence.

Note: Analysis will be marked as finished after the analysis time elapsed. Note that the instrument method may be still running.

The *Sts.* fields will return to their initial status (all fields green) in the following cases:

- Running of the sequence (by *Run* command, the  icon or **F5** or **Ctrl** + **Q** shortcut).
- Using a *Sequence - Reset Status* command (**Ctrl** + **E**).

Note: A text description will be printed in the printouts instead of colored tags in the *Sts.* column.

Run

Selects whether the particular line in the **Sequence Table** will be measured. The column corresponds to the *Run Lines* field in [Sequence Options](#) dialog. Working with *Run* column is more suitable for discontinuous selection of lines or for exclusion of individual lines from measurement.

SV

Number of the first vial used in the processing of the given row.

EV

Number of the last vial used in the processing of the given row. Its number must be the same or higher than the number entered in the SV field. The difference between the end and start vial defines how many vials will be processed by selected line.

If *Options - Format - Automatically* setting has been applied, then for more than one vial or injection, the *File Name* pattern will automatically be verified and corrected to ensure that it is in the correct format. In other words, %v and %i variables will be added to the pattern.

If automatic formatting has been applied, the variables in the filename will be supplemented with zeros from the left in order to correspond with defined number of injected samples (e.g. names will be supplemented by numbers ranging from 001 to 100 for SV 1 and EV 100).

Note: Some autosamplers may support alphanumeric or graphical setting of the *SV* or *EV* parameter. In such case, an arrow will appear in the right side of the given field, and clicking it will open particular [Select Vial](#) dialog.

IV

Indicates how many injections will be processed from each vial based on this line in the **Sequence Table**.

If the *Format - Automatically* item from [Sequence Options](#) dialog has been activated, then, for multiple vials or injections, the *File Name* pattern will automatically be verified and corrected to ensure that it is in the correct format. In other words, the %v and %i variables will be added to the template.

Sample ID

The sample identification field will be transferred to the headers of all created chromatograms. Maximum number of characters in this field is limited to 64. The parameter corresponds to the *Sample ID* field from the [Single Analysis](#) dialog.

This field serves for information only. After clicking on the cell in this column the  icon will appear. Using it invokes the list where you can pick from the available variables for composing the name. Variables that can be used are described in the **Chromatogram File Name** section of the [Single Analysis](#) dialog described on pg 59.

Sample

Sample description - will be transferred to the headers of all created chromatograms. Maximum number of characters in this field is limited to 64. The parameter corresponds to *Sample* field from the [Single Analysis](#) dialog.

This field serves for information only. After clicking on the cell in this column the  icon will appear. Using it invokes the list where you can pick from the available variables for composing the name. Variables that can be used are described in the **Chromatogram File Name** section of the [Single Analysis](#) dialog.

Sample Amount

The value is used in calibration calculations of the *ESTD* and *ISTD* methods. When using the *ISTD* calculation, enter the sample amount without the internal standard. Enter values in units used in the calibration file (as defined in the *Units - Compound* field of the [Calibration Options](#) dialog). The default value is 0. The parameter corresponds to the *Amount* field of the [Single Analysis](#) dialog and the value will be transferred to the chromatogram header.

ISTD1 Amount

The amounts of the internal standards for the *ISTD* method. Other columns for entering amounts of ISTD2 - ISTD5 are hidden by default. This value has to be entered in the same units as those used in the calibration file (as

defined in the *Units - Compound* field of the [Calibration Options](#) dialog). The default value is 0. The parameter corresponds to the homonymous field from the [Single Analysis](#) dialog and the value will be transferred to the chromatogram header.

Sample Dilut.

This field holds the sample dilution factor - its value is used to multiply each calibrated amount in the *Amount* column of the **Result Table**. The default value is 1. The parameter corresponds to *Dilution* field in the [Single Analysis](#) dialog and the value will be transferred to the chromatogram header.

Inj. Vol.

Defines the volume of injection. The response of all compounds in a calibration or when using calibrated calculations is modified by the ratio of this value and *Default Injected Volume* value from the calibration file. This parameter corresponds to the analogous field from [Single Analysis](#) dialog and its value will be transferred to the chromatogram header. The default volume is 0 and default units are μl .

Caution: When using a directly controlled autosampler this value will define the injection volume! The station then automatically verifies whether the entered volume corresponds to the current autosampler setting.

When the value set in this column is 0, no correction to the *Default Injection Volume* value will be made. However, when *Inj.Vol.* of 0 is entered into the sequence when using a directly controlled autosampler, a warning will be given.

File Name

Defines the chromatogram name pattern. The pattern may use a combination of fixed text and variables (usually entered in the form of % and a character). The same variables may be used as in the [Single Analysis](#) dialog, plus %v and %i variables for automatic calculations of vials and injections when measuring multiple analysis from one line. After clicking on the cell in this column the  icon will appear. Using it invokes the list where you can pick from the available variables for composing the file name. You can also enter a subfolder in the chromatogram name. Variables and how to enter the subfolder name are described in the **Chromatogram File Name** section of the [Single Analysis](#) dialog described on pg 59.

It is possible to include also the name of a subfolder in the name to store chromatograms into subfolders of the [Data Subdirectory](#). E.g. the *MySubProjectMyFileName* in the DEMO1 project will result in the MYFILENAME.PRM stored in the C:\CLARITY\DEMO1\DATA\MYSUBPROJECT folder. If the subfolder doesn't exist, then it will be created when the chromatogram will be stored.

Note: If variable %n is being used, it will be independent from the variable using the same name in the [Single Analysis](#) dialog.

Std

Sets the type of the sample being measured. The possible options are:

Unknown - denotes an unknown sample.

Standard - specifies the calibration standard. Samples marked as *Standard* must have the calibration level filled in the *Lvl* column. Chromatograms created on the sequence rows marked as calibration standards will be stored in the CALIB subdirectory instead of the DATA subdirectory.

Bypass - this *Std* type allows for performing an analysis in the *ACTIVE* sequence with controlled autosampler without actually injecting a sample. This may be useful for system clean-up etc..

Blank - in fact a calibration standard with no amount of added sample. The *Blank* is thus one point of a calibration curve (with the *Amount* of 0).

Lvl

Indicates the number of the calibration level when the resulting chromatogram is to be used for automatic recalibration when the *Std* column of the sample is set to *Standard*. The file has to be set as the calibration standard in the *Std* column and a proper calibration file has to be set in the current method as a prerequisite for automatic recalibration.

Method Name

Indicates the name of the method governing all analyses set on a given line. The active template method name from the [Instrument](#) window will automatically be inserted in the first line of the new sequence table.

Note: If you point the mouse cursor to the title of the method, the yellow tooltip will open, in which the title of the method including the entire path will be displayed.

Report Style

Sets the report style to be used for printing. If no name has been entered, no printing on a given row of the **Sequence Table** will be performed. After clicking on the cell in this column the  icon will appear. Using it opens the [Insert File Name](#) dialog for specifying the print method. Another necessary condition for printing the report is checking the *Print Results* field in the [PostRun Setting](#) dialog and/or setting the direction mark next to the  icon in the [Instrument](#) window to .


Caution: Print style cannot be selected while a sequence is running.

If you point the mouse cursor to the title of the print style, the yellow tooltip will open, in which the title of the print style including the entire path will be displayed.

Postrun functions

The commands *Open*, *Open Calib.*, *Print*, *Print to PDF*, *Export Data*, *Export AIA*, *Export TXT*, *Export EZChrom*, *Export Multidetector*, *Run Program*, *Program to Run* and *Parameters* generally correspond to analogous items in the **PostRun Setting** dialog (for more details see the chapter **Postrun...** on pg **100**). All of these columns contain checkboxes, with the exception of the *Program to Run* column (corresponding to the *Program to Run* field of the **PostRun Setting** dialog) and the *Parameters* column.

Note: When there is a printing set in the *Postrun functions* which includes the chromatogram or results, the *Open* chromatogram *Postrun function* has to be selected too. Only chromatograms opened in the **Chromatogram** window can be reported.

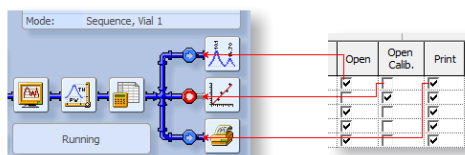

When the sequence is running, the postrun switches in the **Instrument** window will be locked to the state that has been set on the currently running row. Also, when any change in *Postrun functions* on still-not-measured row is performed, even when the sequence itself is saved, no changes will actually take place until the **Sequence** window is closed and reopened.

Note: In the implicit layout only the *Open*, *Open Calib.* and *Print* columns are visible, the rest is hidden. Use the *Edit - Setup Columns...* command to display the desired columns.

Include in SST

This checkbox allows user to set whether the particular chromatogram will be included in the SST calculations or not.

Stored Calib.


This checkbox saves the chromatogram with the stored version of the calibration file, as opposed to the linked calibration when not checked. The checkbox corresponds to the *Open with stored calibration* checkbox in the right part of the Chromatogram - Results tab.

Close All

This checkbox allows user to close all currently opened chromatograms. Opening of the chromatogram measured in the same line of sequence is not influenced by this setting, it follows setting of *Open* checkbox only.

Status Bar

The status bar provides the following information:

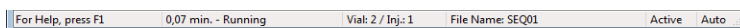

State

If the sequence is running, the runtime of the actual sample is displayed in the first half of the sequence status section. The status itself is shown in its second part. Possible sequence states are described in the **Appendix 1** of the **User Guide**.

Vial No.

The vial number / injection number from a given vial.

File Name

The name of the chromatogram currently being measured.

Active (Passive)

Active/passive sequence flag.

Auto (Manual)

Chromatogram name control flag showing the settings of the *Format* section in the [Sequence Options](#) dialog described on pg **313**.

6.1 File

This is an entry point to a submenu of commands for working with sequence files.



New

Creates a new sequence and closes the sequence previously opened. An empty sequence table will be created with default parameter values based on the template sequence file (TEMPLATE.SEQ) saved in the common directory (C:\CLARITY\COMMON\ in the default projects). New sequence is *ACTIVE* by default (or set according to the setting in TEMPLATE.SEQ). If the template sequence file does not exist in the project, the new sequence will be based on internal preset values in Clarity. The first row in a new sequence takes its settings from the Method and Report setup currently loaded to the Instrument, the **PostRun** information is loaded from the actual setting of the [PostRun Setting](#) dialog. For further rows, the settings are copied from previous line.



Open

Selects and opens a sequence. The command opens the [Open Sequence](#) dialog which is identical to the [Open Method](#) dialog described on pg 44.

In the full version , the same particular sequence cannot be opened on multiple Instruments at once.



Save

Saves the currently displayed sequence.

If you forget to save the sequence file after measuring, the measured chromatograms may not be automatically opened using the *Open Chromatogram from Sequence* command in [Chromatogram](#) window, nor be batch processed using *Batch* command from the [Instrument](#) window, because the current status of *Sts* column will not have been saved.

Furthermore, file names of measured files will not be remembered and their listing will not be available.

Save As

Saves the current sequence under a selected name and/or to another directory. The command invokes the [Save As](#) dialog similar to [Open Method](#) dialog where you can set a new name or select a different directory.

When another file of the same name already exists, you will not be able to use this name. Also, in the protected mode (full version  only), you are not allowed to change a file without the appropriate access rights (see *System - User Accounts...* command).

Send Sequence by Email	Sends the displayed sequence as an email attachment. It is analogous to the <i>Send Method by Email</i> command (for more details see the chapter Send Method by Email on pg 47). The sequence will automatically be saved prior to sending.
Import...	Imports the sequence table from a text file (for more details see the chapter Import... on pg 303).
Export...	Exports a Sequence Table to a text file. Values are exported in tab delimited format, first row contains column headers. Hidden columns are not exported.
 Report Setup	Selects and configures the report style. The command opens the Report Setup dialog described on pg 333.
 Print Preview	Displays a preview of the printout prior to actual printing. The command opens the Print Preview dialog described on pg 352.
 Print to PDF	Prints the Sequence Table to *.PDF file.
 Send Printed PDF by Email	Opens a new email message in default email client, prints the Sequence Table to *.PDF and adds created file as an email attachment.
 Print...	Selects a printer and starts actual printing. Invoking the command opens the Print dialog described on pg 354.
-----	Displays the list of recently opened sequence files so it is easily possible to get to the files you are currently working with. The number of files in the list is determined by the <i>Recent Files</i> parameter in the User Options - General dialog available from the Instrument window by <i>View - Options</i> command.
Exit	Closes the Sequence window.

6.1.1 Import...

Imports a **Sequence Table** that has been stored in a text file. Values have to be in delimited format and separated by an arbitrary delimiter.

Import of the sequence is one of the functions that enables an interconnection between the **Clarity** and **LIMS** systems. It can be performed in three steps:

Select text file and set import settings

Invoking the command will open the [Import Sequence Step 1](#) dialog, that is used to select which text file will be imported, set the delimiters and preview the values.

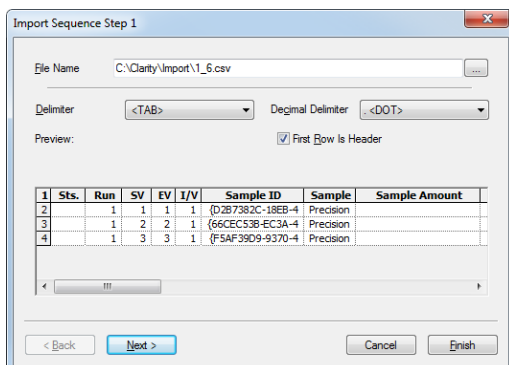


Fig 140: Import Sequence Step 1

File Name

Serves for setting the filename and directory path of the imported file. Use the  button to invoke the [Open Files To Import](#) dialog and locate the desired file. Press OK to confirm the selection of the file.

Delimiter

Selects the character used as field delimiter. Possible options are **<TAB>**, **<SPACE>**, **<COMMA>** or **<SEMICOLON>**.

Decimal Delimiter

Defines the character used as decimal delimiter. Options for the decimal delimiter are **<Window's Locale>** (the station will use the setting specified in the local settings of MS Windows), **<COMMA>**, **<DOT>** or **<SEMICOLON>**.

First Row is Header

If the text file to be imported contains column descriptions in the first row, use this checkbox to immediately transfer it to the column headers of the *Preview* section and also use this row for mapping in the [Import Sequence Step 2](#) dialog.

Preview

Preview of the first five rows of the imported sequence table.

Column mapping

After clicking the *Next>* button the dialog changes to the [Import Sequence Step 2](#). This section is used to map the columns of the imported file with the corresponding columns in the **Sequence Table**.

Import can only be finished after the required columns (SV, File Name, Method Name - all marked in bold) have been mapped.

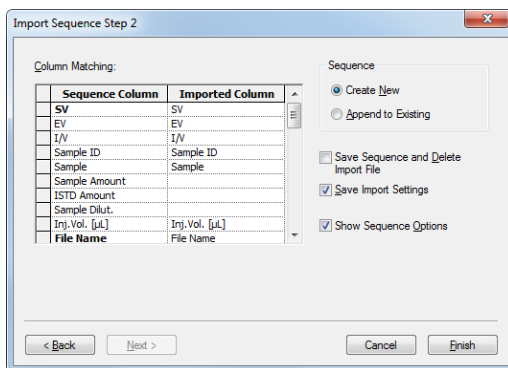


Fig 141: Import Sequence Step 2

Column Matching

Set the mapping so that the columns used in **Clarity** (*Sequence Column*) will correspond to the columns from the imported file (*Imported Column*). Columns that are highlighted in bold letters are required.

When the imported table contains column headers (the *First Row is Header* checkbox from the [Import Sequence Step 1](#) dialog has been checked), **Clarity** will attempt to automatically map the columns according to their names. Manual mapping will override the default name mapping.

Clarity will store these names for future imports.

Sequence

Determines whether the imported table will be appended to the end of the currently opened sequence table in the [Sequence](#) window (*Append to Existing*) or whether a new sequence table will be created (*Create New*).

Save Sequence and Delete Import File

When checked, automatically saves the imported sequence file and deletes the source text file. When the *Create New* checkbox has been checked, the [Save As](#) dialog for selecting the filename and path of the new **Sequence Table** will open automatically after import.

Save Import Settings

Saves the current import settings for future imports. The setting will be stored in the station configuration and is thus shared by all users and

Instruments.

Show Sequence Options

Automatically opens the [Sequence Options](#) dialog after the import.

Note: When importing into a running sequence, this checkbox will be disabled.

Adjust the Sequence Options

If you checked the *Show Sequence Options* checkbox, the [Sequence Options](#) dialog will automatically open after the import. This is where the sequence settings can be adjusted. A detailed description of the [Sequence Options](#) dialog can be found in the chapter "**Sequence Options**" on pg 313.

6.2 Edit

Submenu containing commands for work with clipboard and tables.

Note: Commands may also be found in the local tables menu, accessed by right clicking the mouse on the table.

Note: The procedure for selecting lines is detailed in the **User Manual**.

	Insert New Line	Inserts a new line ahead of a currently selected one. The inserted line will be filled with the same values as the selected line.
	Line Up	Moves selected rows one position up. Moving can be also easily done using the mouse - left click and hold the mouse on the first column of the Sequence Table (with line numbers), drag the line up to the desired position and release. Selected lines do not have to form a continuous block.
	Line Down	Moves selected rows one position down. Moving can be also easily done using the mouse - left click and hold the mouse on the first column of the Sequence Table (with line numbers), drag the line down to the desired position and release.
	Undo	Cancels the modifications effected in tables (Ctrl + Z).
	Redo	Countermands the <i>Undo</i> command (Ctrl + Shift + Z).
	Cut	Moves the selected text to the clipboard (Ctrl + X).
	Copy	Copies the selected text to the clipboard (Ctrl + C).
	Paste	Copies the clipboard content to the indicated place (Ctrl + V). If you are pasting to an area that is larger than the clipboard content, the content will be tiled (<i>Fill</i> function). This is useful if you want to insert a large number of the same fields or several periodically repeating fields into the table.

Paste Insert	Pastes clipboard content to automatically inserted new lines (Ctrl + I). Unlike the <i>Paste</i> command, the original table content is not replaced, but new lines are instead inserted. New lines will be inserted above (ahead of) the currently selected line.
Fill Down	Copies the content of current cell to all cells below it or to selected range (Ctrl + F). The values in the <i>Start Vial</i> column of the Sequence Table will automatically be incremented.
Fill Series...	Displays the Fill Series... dialog for filling up incremental values to cells.
Delete	Deletes the selected table fields. Deleted fields will be reset to their default values. If you have selected entire lines, the lines will be removed.
Mark for Run	Checks the checkbox in the <i>Run</i> column for selected lines. These rows will be processed when the sequence will be run.
Clear Run Marks	Unchecks the checkbox in the <i>Run</i> column for selected lines. These rows will not be processed when the sequence will be run.
Invert Run Marks	Inverts the marks in the <i>Run</i> column of the Sequence Table . All selected rows with the <i>Run</i> column not checked will be checked (marked for run) and all rows currently checked will be excluded from the run (unchecked).
Reset Status (Sts.)	Invoking this command (or using the Ctrl + E shortcut) resets the measured line flags on the selected lines. The command will change the color of <i>Sts.</i> column fields of all selected rows to green (not measured yet). The status of all lines will be automatically reset while running a sequence using the <i>Run</i> command.
Restore Default Columns	Restores the default columns setting (for more details see the chapter Edit on pg 149).
Setup Columns...	Opens the Setup Columns dialog for the Sequence Table .
Select All	Invoking the command (or using the Ctrl + A shortcut) selects the whole Sequence Table .

Remove All	Deletes the entire Sequence Table , only the empty first row will remain.
Hide Column(s)	Hides all selected columns.
Show Hidden Column(s)	Shows all hidden columns.

6.2.1 Fill Series

Dialog for filling up a series in selected cells of the sequence or other table. When selected one cell or multiple cells in one row, the changes will apply from the first top selected line to the last. To apply changes in selected cells only, select multiple cells (rows and columns) as you are used to do in your favorite spreadsheet.

This dialog is also available using the popup menu in the [Gradient Table](#).

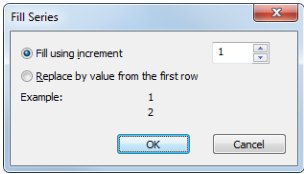


Fig 142: Fill Series dialog

Fill using increment

Sets this option if you want to fill up the series incrementally using the entered step in the box on the right. Starting value will be taken from the first top cell of the each selected column. The number is also detected at the end of the string in the cell.

Replace by value from the first row

Replaces all selected cells with the value from the top selected cells.

Example

Contains preview how the first and second cell in the first selected column will look like.

6.3 Sequence

Summarizes commands for the control of the running sequence. The majority of these commands is also placed on the toolbar by default.

Note: If icons representing these commands are disabled, the tooltip may help in discovering the reason of the disablement.



Run

Invoking this command or using the  icon or **Ctrl** + **Q** shortcut runs the sequence. The command is only active when the sequence is not in the *RUNNING* state. Before the sequence is run, the validity check is performed. All methods used in *ACTIVE* sequence must:

- have *External Start/Stop* option in the [Method Setup - Measurement](#) dialog checked.
- have *Enable Autostop* checkbox in the [Method Setup - Measurement](#) dialog selected and *Run Time* field filled.

The *Run* command automatically resets the status of all lines in the **Sequence Table** (all fields in the *Sts.* column will turn green).

Running a partially measured sequence using the *Resume* command starts the measurement from the first non-measured line, while running it using the *Run* command the entire sequence will be measured again. The station automatically checks the entire sequence upon its start. During the measurement only the currently used line is checked in this way, which means that the sequence can be modified during the run.



Pause

Using the *Pause* command,  icon or **Ctrl** + **Space bar** shortcut suspends the active sequence. A suspended sequence does not allow injections and does not react to an external start signal issued.

The status of the sequence after the use of the *Pause* command depends on the status of the sequence before invoking the command:

WAITING, *IDLE* sequence is turned to *PAUSED* status. *RUNNING* sequence is turned to the *PAUSED* status only after the data acquisition of the current analysis is completed (the intermediate status is *RUNNING/PAUSE*).

	Resume	Using this command, icon or the Ctrl + Space bar shortcut resumes the operation of a suspended sequence. The sequence status becomes <i>IDLE</i> or <i>WAITING</i> and, accordingly, the signal that allows the injection defined on the next row will either be sent immediately or after the Idle time has passed. When returning to the <i>IDLE</i> state, the sequence counts the time from the moment it was set to the <i>PAUSED</i> state. When the <i>Resume</i> command is used while in the <i>RUNNING/PAUSE</i> state, the planned pause of the sequence will be canceled. If you run a partially measured sequence using the <i>Resume</i> command, measurement continues from the first line that wasn't measured yet. If you run it using <i>Run</i>, the entire sequence will be measured again.
	Abort Sequence or Acquisition	Invoking this command or using the icon or Ctrl + B shortcut immediately aborts the running sequence. Any data acquisition in progress will be interrupted, the results will not be processed, and the sequence will be terminated.
	Stop Sequence or Acquisition	Using the command, icon or Ctrl + T shortcut halts the running sequence. Currently running analysis will be completed (the actual state will change to <i>RUNNING/STOP</i>). If the <i>Stop</i> command is used for the second time, it will stop the running analysis.
	Repeat Injection	Invoking this command or using the icon or Ctrl + shortcut interrupts the actual measurement and repeats the injection (runs the analysis according to the row in the Sequence Table). If the sequence is in the <i>PAUSED</i> state before using the <i>Repeat Injection</i>, triggering the command will repeat the measurement of last completed injection.

**Skip Vial**

Using this command,  icon or **Ctrl** + **J** shortcut terminates the analysis of active vial and also omits all remaining injections using the current vial defined on the same row of the **Sequence Table**. This command works principally in the same way as the *Abort* command does, but only for the currently measured sample. The sequence continues with the measurement of the next vial.

The line containing skipped injections will be indicated in the *Sts.* column by light blue color.

Potential re-measurement of skipped vials has to be done manually (e.g. by creating a new line in the **Sequence Table**).

**Check Sequence**

This command performs a check of every row in the **Sequence Table** for potential errors. These may be attempt to use a non-existent method or print style, incorrectly set control parameters, prohibited chromatogram name entered etc. The check same as the one invoked by the *Check Sequence* command is automatically performed on the start of the sequence measurement.

During a measurement, only the line currently being measured is checked. Manual use of the *Check Sequence* command allows the user to test the whole sequence, for example after modifying the sequence while it is running to verify that all changes are correct.

**Snapshot**

Invoking the command,  icon or **Ctrl** + **H** shortcut processes and saves the data from currently measured sample without terminating the measurement. Sample counter *%n*, if used, will not be incremented in the chromatogram name. This command is identical to the *Snapshot* command from the [Single Analysis](#) dialog described on pg 59.

**Options...**

When invoked, opens the [Sequence Options](#) dialog used for defining the sequence parameters (for more details see also chapter **Sequence Options** on pg 313).

**Edit Current Method...**

Displays and modifies the method defined on the row of the **Sequence Table** with currently active field. The command opens the [Method Setup - Measurement](#) dialog described on pg 80. The operation may also be performed by the *Edit Method...* command from the local menu for the given row of the **Sequence Table**.

**Edit Current Report...**

Displays and modifies the report style set on the row of the **Sequence Table** with currently active field. The command opens the [Report Setup](#) dialog described on pg 333. The operation may also be performed by the *Edit Report...* command from the local menu for the given row of the **Sequence Table**.

6.3.1 Sequence Options

Sequence Options dialog serves for defining the sequence parameters. It is opened by using the  icon or the *Sequence - Sequence* command from the **Sequence** window.

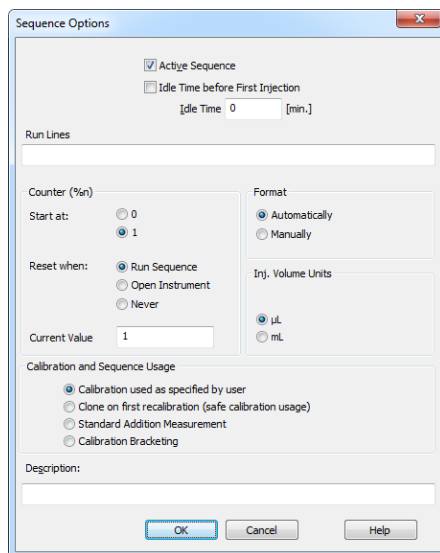


Fig 143: Sequence Options

Active Sequence

This checkbox serves for setting the operation mode of the sequence. *ACTIVE* sequence differs from the *PASSIVE* sequence in the fact that the operation is governed by **Clarity** (instead of the autosampler). After the active sequence is started, the station sends the permission signal to the autosampler and waits (in the *WAITING FOR INJECTION* state) until the sampler acknowledges the injection. Data acquisition is started after the confirmation signal is sent back and the permission to inject is disabled.

Caution: When the *ACTIVE* sequence is used, *Enable Autostop*, *External Start/Stop* must be set in the **Method Setup - Measurement** dialog and the *Start-Restart* option of the *External Start/Stop* behave as *Start Only*.

Caution: When the *PASSIVE* sequence is used, *External Start/Stop* must be set in the **Method Setup - Measurement** dialog.

Once the analysis time of the method in question has elapsed (as set in the *Run Time* field of the **Method Setup - Measurement** dialog with *Enable Autostop* field enabled), the analysis will be completed and injection will be permitted once again.

Note: Only the **Start In** and **Ready Out** synchronization cables are needed for the *ACTIVE* sequence to function. The *ACTIVE* sequence can thus be used even with an autosampler not governed by a specialized control module.

Note: The confirmation signal is represented by a change against the initial value of the selected digital output. The default assignment of outputs is described in the corresponding hardware manual.

The assignment of digital inputs and outputs used for the synchronization in the full version  can be changed on the tab of appropriate Instrument in the [System Configuration](#) dialog, where *Ext. Start Dig. Input* and *Ready Dig. Output* fields are placed.

Idle Time before First Injection

Determines whether the *Idle Time* parameter will be used even before the first injection in the *ACTIVE* sequence. This field can only be edited when the sequence is set to *ACTIVE*.

Idle Time

Delays the sending of the permission to inject. The next injection after the end of an analysis will always be permitted only after the defined time has elapsed. Possible values are in the 0.01-100 min range. This field can only be edited when the sequence is set to *ACTIVE*.

Run Lines

Sets the line numbers of the **Sequence Table** to be used in the measurement. Particular line numbers should be separated with a comma (e.g. 1,3,5), ranges should be entered using a dash (e.g. 1-8).

A list of lines that should be run defined interactively by the checkboxes in the *Run* column of the **Sequence Table** will be automatically filled to this field.

Counter (%n)

Sets the value and properties of the %n variable (analysis counter) used in a **Sequence Table**.

Start at:

Defines if the %n variable will be reset to 0 or 1 at the moment defined in the following **Reset when** field.

Reset when:

Sets the condition when the counter (%n variable) is reset. Available options are:

Run Sequence - The counter is reset every time the sequence is run.

Open Instrument - The counter is reset every time the [Instrument](#) window is opened.

Never - The counter is never reset.

Current Value

This field shows (and allows to modify) the current value of the sequence counter (%*n* variable).

Caution: The station keeps two separate counters: one for the sequence and another one for the single analysis. This means that two %*n* variables with possibly different values exist simultaneously.

Format

Sets the verification of the chromatogram names pattern (set in the *File Name* field) when measuring more than one chromatogram from particular line.

Automatically

The pattern will be automatically modified to correspond to the number of measured chromatograms and to avoid possible loss (overwriting) of chromatograms.

Note: When using the automatic format, variables will be complemented with zeros from the left in order to correspond with the defined number of samples (e.g. in a sequence where SV = 1 and EV = 100, filenames will automatically be numbered from 001 to 100).

Manually

The pattern will not be modified. Once the sequence has been run, you will be notified of a possible overwrite of chromatograms within one line.

Inj. Volume Units

Sets the units for the *Inj. Vol.* field. Select the units your autosampler is working with (μ l, ml).

Caution: It is important to select the correct units when the autosampler is controlled by the direct control module. In such a case, values in the *Inj. Vol.* field are checked against the syringe volume used by autosampler and incorrectly set values would actually cause wrong injections.

The standard units used throughout the whole **Clarity** station is *ml*. The **Sequence** window is the only place where other units than *ml* can be set and used.

Calibration and Sequence Usage

Selects the mode of the calibration and sequence usage.

Calibration used as specified by user

Sets the sequence and calibration to the standard **Clarity** use (the only way available since version 3.0). The calibration set in the **Method Setup - Calculation** dialog is used for calibration and recalibration, no calibration cloning is performed.

Clone on first recalibration (safe calibration usage)

Sets the sequence and calibration to the Safe Calibration Usage mode. The calibration will be cloned at each start of the sequence and at each start of a new sequence. The option is not available on EA Instruments.

Note: This is because of the possibility to devalue the already measured samples because the calibration with the same name, but different responses, is linked to them.

Standard Addition Measurement

Sets the sequence and calibration to the Standard Addition mode use. As each unknown sample will use one calibration file, calibration will be cloned from the original calibration file whenever an unknown sample or blank follows the calibration standard. All responses are cleared from a newly cloned calibration. The option is not available on EA Instruments.

Calibration Bracketing

Sets the sequence and calibration to the Calibration Bracketing mode use. As every unknown sample series is demarcated by calibration standards and as such uses a single calibration file, calibration will be cloned from the previous calibration clone whenever an unknown sample or blank follows the calibration standard. In a newly cloned calibration file, all responses other than responses of the last series of calibration standards (only those standards that follow the unknown samples) are cleared.

In the EA mode, the name of the cloned calibration file is hard-set to the original name incremented by the sequence line number (%L parameter).

Description

Describes a displayed sequence.

6.4 View

Commands for setting up the [Sequence](#) window.

Toolbar	Displays/hides all toolbars. The command is similar to the <i>Toolbar</i> command in the main Clarity window (for more details see the chapter View on pg 30).
Customize...	Customizes the toolbars. The command is similar to the <i>Customize...</i> command in the main Clarity window (for more details see the chapter Customize... on pg 31).
Reset All	Restores all menus and toolbars in the Sequence window to their default state.

6.5 Window

This section contains commands allowing to open other base windows.

Main 	Displays the Clarity main station window described on pg 1 .
Instrument	Displays the Instrument window described on pg 40 .
Chromatogram	Displays the Chromatogram window described on pg 118 .
Calibration	Displays the Calibration window described on pg 238 .
Single Analysis	Displays the Single Analysis window described on pg 59 .
Data Acquisition	Displays the Data Acquisition window described on pg 279 .
Device Monitor	Displays the Device Monitor window described on pg 55 .
Station Audit Trail 	Displays the Audit Trail window described on pg 318 .
Sequence Audit Trail... 	Displays the Sequence Audit Trail window described on pg 327 .

7 Audit Trail ✓full version

Audit Trail window is used for logging of individual operation of the station. By allowing backward control of all performed operation including information of who and when performed it, the logging of operations supports higher reliability and reproducibility of results in the sense of Good Laboratory Practice (GLP) rules.

Audit Trail										
File Edit Filter View Window Help										
OK	Time	Type	Analyst	Instrument	Area	Description	User	Computer		
1	ok	9.7.2012 11:53:32		System	System	System Options: Generate Chromatogram After Abort has been c	lenka	LENKA	Ver: 4.0.1.696,	
2	ok	9.7.2012 11:53:32		System	System	System Options: Chromatogram Files Overwriting has been chang	lenka	LENKA	Ver: 4.0.1.696,	
3	ok	9.7.2012 11:53:32		System	System	System Options: Chromatogram GLP Mode has been changed fro	lenka	LENKA	Ver: 4.0.1.696,	
4	ok	9.7.2012 11:53:15		System	Configuration	Control module 542 added.	lenka	LENKA	Ver: 4.0.1.696,	
5	ok	9.7.2012 11:53:03		MyGC	Instrument	Close Instrument	lenka	LENKA	Ver: 4.0.1.696,	
6	ok	9.7.2012 11:53:02		MyGC	Acquisition and	Waiting: Instrument was aborted by user	lenka	LENKA	Ver: 4.0.1.696,	
7	ok	9.7.2012 11:53:02		MyGC	Files	Save File C:\Clarity\Projects\MyMethod.PRJ	lenka	LENKA	Ver: 4.0.1.696,	
8	ok	9.7.2012 11:51:50		MyGC	Detector	INT9 - 2 (SN DemoSN): Method sent to sub-device.	lenka	LENKA	Ver: 4.0.1.696,	
9	ok	9.7.2012 11:51:50		MyGC	Detector	INT9 - 1 (SN DemoSN): Method sent to sub-device.	lenka	LENKA	Ver: 4.0.1.696,	
10	ok	9.7.2012 11:51:50		MyGC	Acquisition	Method 'C:\Clarity\MyMethod\MyMethod.MET' has been sent	lenka	LENKA	Ver: 4.0.1.696,	
11	ok	9.7.2012 11:51:49		MyGC	Files	Open File C:\Clarity\Common\Instrument.sty	lenka	LENKA	Ver: 4.0.1.696,	
12	ok	9.7.2012 11:51:49		MyGC	Files	Open File C:\Clarity\MyMethod\MyMethod.MET	lenka	LENKA	Ver: 4.0.1.696,	
13	ok	9.7.2012 11:51:49		MyGC	Files	Open File C:\Clarity\Projects\MyMethod.PRJ	lenka	LENKA	Ver: 4.0.1.696,	
14	ok	9.7.2012 11:51:49		MyGC	Instrument	Open Instrument	lenka	LENKA	Ver: 4.0.1.696,	
15	ok	9.7.2012 11:51:47		Administ	My LC	Instrument	Close Instrument	lenka	LENKA	Ver: 4.0.1.696,
16	ok	9.7.2012 11:51:47		Administ	My LC	Acquisition and	Not ready to run acquisition: Instrument was aborted by user	lenka	LENKA	Ver: 4.0.1.696,
17	ok	9.7.2012 11:51:47		Administ	My LC	Files	Save File C:\Clarity\Projects\MyMethod.PRJ	lenka	LENKA	Ver: 4.0.1.696,
18	ok	9.7.2012 11:51:43		Administ	My LC	Files	Save File C:\Clarity\MyMethod\Data\My LC - 9.7.2012 11_51_18.	lenka	LENKA	Ver: 4.0.1.696,
19	ok	9.7.2012 11:51:43		Administ	My LC	Chromatogram	C:\Clarity\MyMethod\Data\My LC - 9.7.2012 11_51_18.PRM - Det	lenka	LENKA	Ver: 4.0.1.696,
20	ok	9.7.2012 11:51:43		Administ	My LC	Chromatogram	C:\Clarity\MyMethod\Data\My LC - 9.7.2012 11_51_18.PRM - Det	lenka	LENKA	Ver: 4.0.1.696,
21	ok	9.7.2012 11:51:43		System	Messages	My LC - 9.7.2012 11_51_18.PRM has been changed. Save before	lenka	LENKA	Ver: 4.0.1.696,	
22	ok	9.7.2012 11:51:38		Administ	My LC	Files	Open File C:\Clarity\DEMO1\Calib\250038R1.lcal	lenka	LENKA	Ver: 4.0.1.696,
23	ok	9.7.2012 11:51:38		Administ	My LC	Files	Open File C:\Clarity\DEMO1\Calib\250038R1.lcal	lenka	LENKA	Ver: 4.0.1.696,
24	ok	9.7.2012 11:51:21		Administ	My LC	Chromatogram	C:\Clarity\MyMethod\Data\My LC - 9.7.2012 11_51_18.PRM - Op	lenka	LENKA	Ver: 4.0.1.696,
25	ok	9.7.2012 11:51:21		Administ	My LC	Chromatogram	C:\Clarity\MyMethod\Data\00000000.PRM - Opened method 4.7.	lenka	LENKA	Ver: 4.0.1.696,
26	ok	9.7.2012 11:51:21		Administ	My LC	Files	Open File C:\Clarity\MyMethod\Data\00000000.PRM	lenka	LENKA	Ver: 4.0.1.696,
27	ok	9.7.2012 11:51:21		Administ	My LC	Files	Open File C:\Clarity\Common\Chromatogram.sty	lenka	LENKA	Ver: 4.0.1.696,

Session

Daily Audit Trail: 2012_07_09

For Help, press F1

Fig 144: Audit Trail - Daily

Audit Trail window is accessible from the main **Clarity** window via **System - Audit Trail** command or from all other station windows using **Window - Station Audit Trail** command. Other windows, such as **Calibration** or **Chromatogram**, have their own **Local Audit Trails**.

The **Clarity** station keeps two global tables with logs of operations, each of them corresponds with one tab in **Audit Trail** window - these are **Session** audit trail and **Daily audit trail**.

Audit Trail files are stored in YYYY-MM-DD.AUDIT file in the C:\CLARITY\LOG directory.

7.1 Station Audit Trail

Session

When **Clarity** is running, logs of all operations performed since the last startup can be found in the Station Audit Trail. Once the **Clarity** station is closed, the **Session** table will be reset and upon next run it will start to record from the beginning.

Note: The Session audit trail will also be reset when the date changes - sessions longer than one day will be divided into several files, each holding its 1-day part of the session.

Daily Audit Trail

Logs all operations performed during a single day (24 hours). It will form something of a daily sum of all the **Session** logs. Daily logs can be found in the LOG subdirectory of the main station directory (C:\CLARITY\LOG by default).

7.1.1 File

Open Audit Trail (Append)	Opens audit trail file. An opened file will be appended to the end of current listing.
Export	Exports data from the audit trail to the text or database file.
Print Preview...	Prints a preview of the audit trail (Ctrl + Shift + P).
Print...	Prints the audit trail (Ctrl + P) from a specified printer.
Exit	Exit from Audit Trail window.

Note: The print format of the audit trail is in the WYSIWYG (What You See Is What You Get) format - all columns displayed on the screen will be displayed accordingly in the report.

7.1.2 Open Audit Trail (Append)

When invoked, displays the selection of all previously saved **Daily Audit Trail** files. Newly opened files will be displayed together within the current **Daily Audit Trail** table, appended to its upper part. Opening previous day's tables will not threaten the process of logging data into the current audit trail file.

Other than current **Daily Audit Trail**s, that have been opened, will be indicated on the **Daily Audit Trail** tab by their date separated by the "+" sign.



Fig 145: Daily Audit Trail tab - several audit trails open

7.1.3 Export

Exports a record of logged operations of the station into the text or database file. The submenu contains two fields:

Text...

Exports data into a text file with the *.TXT or *.CVS extension. Individual items in the export will be separated by semicolons. When invoked, the [Save As](#) dialog will appear where you must enter the name of the new file or select some existing file.

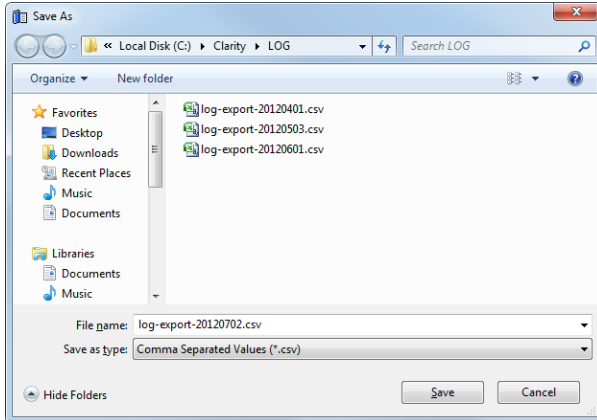


Fig 146: Audit Trail - Export - Save As dialog

When trying to use an existing file a dialog will appear where you need to select whether exported data will be appended to the end of file (*Append*), or if the data will overwrite any existing content of the file (*Overwrite*).

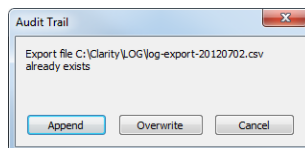


Fig 147: Audit Trail System Message

Note: You need to select the extension to be used to filter the list of available files or the extension of a newly saved file in the *Save as type* field.

DBF...

Exports the logged data in database format. The procedure is similar to the export into text file (it is necessary to select whether to append the exported data to the end of the existing file or overwrite the file).

Note: MS Access supports import of DBF files with **DOS** filename (eight characters max) format only.

7.1.4 Edit

Submenu of commands for working with the clipboard and tables.

Copy	Invoking the command or using the Ctrl + C shortcut copies the selected text into clipboard.
Select All	Invoking the command or using the Ctrl + A shortcut selects all fields of the table.
Restore Default Columns	Invoking the command returns the display of the given table to default settings (for more details see the chapter Edit on pg 149).
Setup Columns...	Using this command opens the Setup Columns dialog which allows to set the visibility and order of appearance of columns in the given table
Hide Column(s)	Invoking the command hides all selected column(s) from display (for more details see the chapter Edit on pg 149).
Show Hidden Column(s)	Using this command displays columns hidden by the <i>Hide Column(s)</i> command (for more details see the chapter Edit on pg 149).

7.1.5 Filter

This menu helps to define which operations will be displayed in the table on the [Session](#) and [Daily Audit Trail](#) tabs. Using this setting will not impact the logging of items into file, just their display. To set the items to be logged, use the [Audit Trail Settings](#) dialog (accessible by using the *View - Properties* command).

Individual menu items will work like checkboxes. If the item is checked, the corresponding logs will be displayed.

Show All	Displays all items of the <i>Filter</i> menu.
Show Instrument X	Displays items from individual instruments.
Show System	Displays system logs and items from the main Clarity window.

7.1.6 View

Displays the setting for the [Audit Trail](#) window.

Toolbar	Shows/hides all toolbars (for more details see the chapter View on pg 30).
Customize...	Customizes toolbars according to a user specification of toolbars (for more details see the chapter Customize... on pg 31).
Reset All	Returns all menus and toolbars in the Audit Trail window into their initial state.
Properties...	Invokes the Audit Trail Settings dialog where the user can specify the operations to be logged.

7.1.7 Properties...

The [Audit Trail Settings](#) dialog is used to specify which operations will be automatically logged by **Clarity** station. The operations are classified into five categories: *System Messages*, *Chromatogram*, *Calibration*, *Sequence* and *Method* operations and *Other*. This corresponds to six tabs.

Note: Optional extensions, like the GPC extension, may use and display additional tabs. For its description see the manual of the particular extension.

You can always select whether the events should be logged to *Daily Audit Trail* or to current audit trail (*Session*) table and, in case of the events concerning the [Chromatogram](#), [Calibration](#) or [Sequence](#) windows, to their local audit trails: *Chromatogram*, *Calibration* and *Sequence*.

Local audit trails are displayed from the respective window through the *Window - Audit Trail* command and their description is in the chapter **"Local Audit Trails"** on pg 327.

The default setting will log all operations into all relevant files.

You can easily disable the access to the [Audit Trail Settings](#) dialog for selected users in the *Open Audit Trail Setting* field of the [User Accounts](#) dialog.

Messages

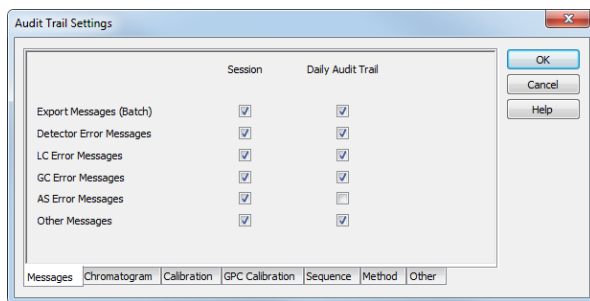


Fig 148: Audit Trail Settings - Messages

Export Messages (Batch)

Sets whether to log messages concerning export of files and batch operations into the *Session* audit trail and/or *Daily Audit Trail*.

Detector Error Messages

Defines whether to log messages concerning problems with acquisition into the *Session* audit trail and/or *Daily Audit Trail*.

LC Error Messages

Defines the logging of messages concerning communication problems with a directly controlled LC chromatograph into the *Session* audit trail and/or *Daily Audit Trail*.

GC Error Messages

Defines the logging of messages concerning communication problems with a directly controlled GC chromatograph into the *Session* audit trail and/or *Daily Audit Trail*.

AS Error Messages

Defines the logging of messages concerning communication problems with a directly controlled autosampler into the *Session* audit trail and/or *Daily Audit Trail*.

Other Messages

Sets whether to log any other system messages into the *Session* audit trail and/or *Daily Audit Trail*.

Chromatogram

Selects which chromatogram operations will be logged (for more details see the chapter **Properties...** on pg 330).

Calibration

Selects which calibration file operations will be logged (for more details see the chapter **Properties...** on pg 331).

Sequence

Selects which sequence operations will be logged (for more details see the chapter **Properties...** on pg 331).

Method

Selects which method operations will be logged (for more details see the chapter **Properties...** on pg 332).

Other

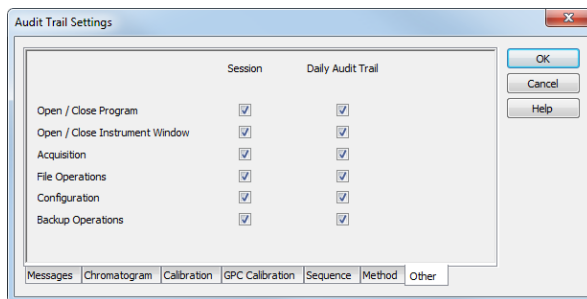


Fig 149: Audit Trail Settings - Other

Open / Close program

Defines whether the data concerning the running and shutdown of **Clarity** station will be logged into the *Session* audit trail and/or *Daily Audit Trail*.

Open / Close Instrument Window

Defines whether the data concerning the running and shutdown of the [Instrument](#) window will be logged into the *Session* audit trail and/or *Daily Audit Trail*.

Acquisition

Defines whether the data concerning the running and shutdown of data acquisition will be logged into the *Session* audit trail and/or *Daily Audit Trail*.

File Operations

Defines whether the operations with files (opening, closing, deleting) will be logged into the *Session* audit trail and/or *Daily Audit Trail*.

Configuration

Defines whether the changes of the station's configuration will be logged into the *Session* audit trail and/or *Daily Audit Trail*.

Backup Operations

Defines whether the use of the *Backup* command from the [Instrument](#) window will be logged into the *Session* audit trail and/or *Daily Audit Trail*.

7.1.8 Window

A submenu of commands for working with the station windows.

Main	Displays the main Clarity window of the station described on pg 1.
Instrument 1 .. 4	Displays the relevant Instrument window described on pg 40.

7.2 Local Audit Trails

Besides the events logged in the global audit trail, the **Clarity** station may also log all operations with chromatograms, calibrations and sequences. The audit trail always forms the integral part of a given file. A list of logged operations is selectable using the *View - Properties* command either from the global [Audit Trail](#) window on the relevant tab or by using the same command directly in window of corresponding local audit trail.

Chromatogram Audit Trail

The audit trail of the active chromatogram can be displayed using the *Window - Chromatogram Audit Trail* command from [Chromatogram](#) window.

Chromatogram Audit Trail also keeps log of errors that occurred during its acquisition (Pump Errors, missing spectra, etc.)

Calibration Audit Trail

An opened calibration log can be displayed using *Window - Calibration Audit Trail* command from [Calibration](#) window.

Sequence Audit Trail

An opened sequence log can be displayed using the *Window - Sequence Audit Trail* command from [Sequence](#) window. The sequence audit trail is reset when the sequence itself is reset.

Method Audit Trail

An opened method log can be displayed using the *Audit Trail* button from the [Method Setup](#) dialog.

7.2.1 File

Export	Invoking this command exports logged operations into a text or database file (for more details see the chapter Export on pg 321).
Report Setup...	Invoking the command opens the Report Setup dialog which allows you to set the print style and the content of the printed report.
Print Preview...	Displays a preview of the print output. The printing format of the audit trail table is WYSIWYG (What You See Is What You Get) and thus depends on the current display of the table.
Print...	Opens the Print dialog used for specifying the printer and starting the actual printing.
Exit	Exits the Audit Trail window.

7.2.2 Edit

Submenu of commands for working with the clipboard and tables.

Copy	Invoking the command or using the Ctrl + C shortcut copies the selected text into clipboard.
Select All	Invoking the command or using the Ctrl + A shortcut selects all fields of the table.
Restore Default Columns	Invoking the command returns the display of the given table to default settings (for more details see the chapter Edit on pg 149).
Setup Columns...	Using this command opens the Setup Columns dialog which allows to set the visibility and order of appearance of columns in the given table (for more details see also chapter Setup Columns... on pg 154)
Hide Column(s)	Invoking the command hides all selected column(s) from display (for more details see the chapter Edit on pg 149).
Show Hidden Column(s)	Using this command displays columns hidden by the <i>Hide Column(s)</i> command (for more details see the chapter Edit on pg 149).

7.2.3 View

Submenu of commands for displaying and modifying of the [Local Audit Trail](#) window settings.

Toolbar	Shows/hides all toolbars (for more details see the chapter View on pg 30).
Customize...	Invoking this command allows to customize the toolbars (for more details see the chapter Customize... on pg 31).
Reset All	Invoking this command returns all menus and toolbars in the Local Audit Trail window into their initial state.
Properties...	Invoking this command or using the  icon opens the Audit Trail Settings dialog where you can specify operations to be logged.

7.2.4 Properties...

The [Audit Trail Settings](#) dialog specifies the operations that will be logged automatically by the **Clarity** station. When invoked from any [Local Audit Trail](#) window, only the tab for changing the settings for the given local audit trail is displayed.

Note: Default setting will log all operations.

Note: The option to change the settings for specified users can easily be disabled using the [User Accounts](#) dialog in the *Open Audit Trail Setting* field.

According to the window from which the local audit trail was opened, one of the tabs described below will appear:

Chromatogram

Sets the logging of chromatogram operations into particular audit trails.

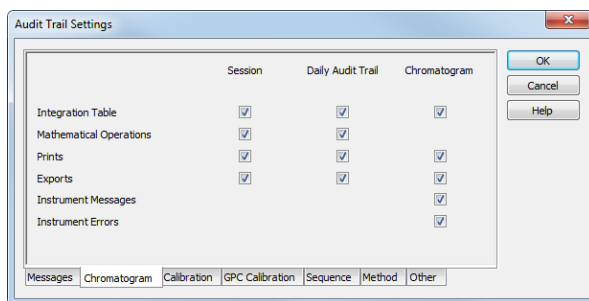


Fig 150: Audit Trail Settings - Chromatogram

Integration Table

Defines whether the changes in the integration table will be logged into the current Global audit trail (*Session*), Daily audit trail (*Daily Audit Trail*) or Chromatogram audit trail (*Chromatogram*).

Mathematical Operations

Defines whether a performed mathematical operations will be logged into the current Global audit trail (*Session*) or Daily audit trail (*Daily Audit Trail*).

Note: Logging of mathematical operations into the chromatogram audit trail is not performed because they create a new chromatogram instead of modifying the current one.

Prints

Defines whether print operations will be logged into the current Global audit trail (*Session*), Daily audit trail (*Daily Audit Trail*) or Chromatogram audit trail (*Chromatogram*).

Exports

Defines whether the exporting operations will be logged into the current Global audit trail (*Session*), Daily audit trail (*Daily Audit Trail*) or Chromatogram audit trail (*Chromatogram*).

Instrument Messages

Defines whether the communication messages sent by and to the controlled instrument will be logged into the current Chromatogram audit trail (*Chromatogram*).

Instrument Errors

Defines whether the errors sent from the controlled instruments will be logged into the current Chromatogram audit trail (*Chromatogram*).

Calibration

Sets the logging of calibration operations into particular audit trails.

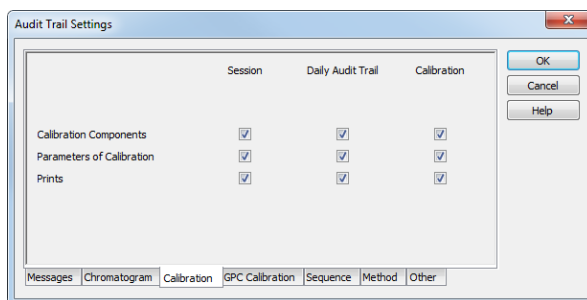


Fig 151: Audit Trail Settings - Calibration

Calibration Components

Defines whether the changes in calibrated compounds will be logged into current Global audit trail (*Session*), Daily audit trail (*Daily Audit Trail*) or Calibration audit trail (*Calibration*).

Parameters of Calibration

Defines whether the changes to calibration parameters will be logged into the current Global audit trail (*Session*), Daily audit trail (*Daily Audit Trail*) or Calibration audit trail (*Calibration*).

Prints

Defines whether print operations will be logged into the current Global audit trail (*Session*), Daily audit trail (*Daily Audit Trail*) or Calibration audit trail (*Calibration*).

Sequence

Sets the logging of sequence states into particular audit trails.

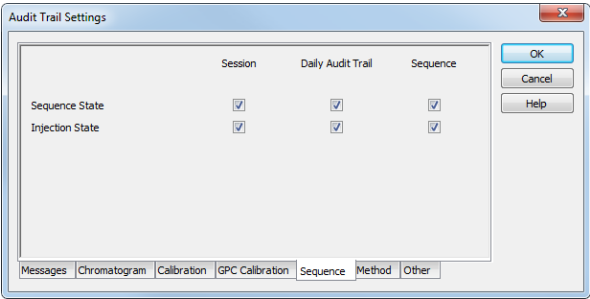


Fig 152: Audit Trail Settings - Sequence

Sequence State

Defines whether the changes of the sequence status will be logged into the current Global audit trail (*Session*), Daily audit trail (*Daily Audit Trail*) or Sequence audit trail (*Sequence*).

Injection State

Defines whether the information about all injections performed during the run of a sequence will be logged into the current Global audit trail (*Session*), Daily audit trail (*Daily Audit Trail*) or Sequence audit trail (*Sequence*).

Method

Sets the logging of method changes into particular audit trails.

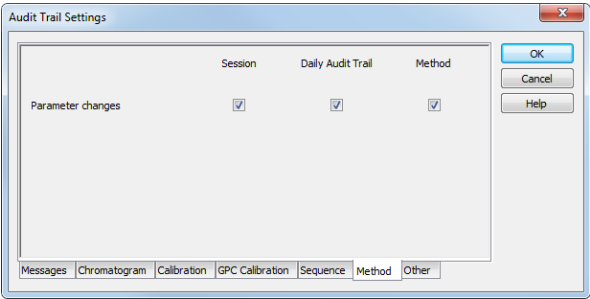


Fig 153: Audit Trail Settings - Method

Parameter Changes

Defines whether the changes in the method will be logged into the current Global audit trail (*Session*), Daily audit trail (*Daily Audit Trail*) or Method audit trail (*Method*).

8 Report Setup

This dialog serves for selecting and adjusting the report style, setting of a printer and initiating of the print itself. The dialog opens either by using the *Report Setup* command or  icon from any major window or by using the  button in the workflow diagram in the *Instrument* window. The displayed report style depends on the window from which the *Report Setup* window was opened - report styles set may differ between different windows.

Note:

When the mouse cursor is placed over the  icon (*Report Setup*) and left still for a moment, the tooltip will appear with the name of the current report style. When placed over the  icon (*Print*) the tooltip will display the current report style and currently selected printer.

The station is supplied with pre-set report styles that are set to be used from all windows that contain printing report commands. These styles may be modified, copied, relocated or replaced by user specified styles.

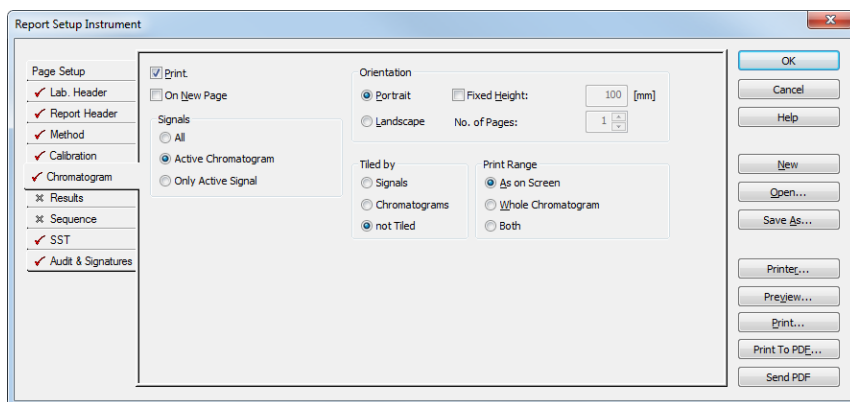


Fig 154: Report Setup

The report is printed in order respective to tabs order on the left. By right click on the tab you can move the tab up or down. The order of *Page Setup* and *Audit & Signatures* can't be changed since they don't represent the print section.

All tabs of the *Report Setup* dialog use the same set of buttons aligned to the right side. These buttons perform following actions:

OK

Saves any changes made in the *Report Setup* dialog and exits it.

Cancel

Exits the [Report Setup](#) dialog. If there were any changes made to some parameter, the user is asked if he wants to save the new values.

New

Creates a new, empty report style. The values of all variable items of the report style will be erased or reset to default values.

Open...

Selects and opens a report style. Invoking the command will open the [Open Report Style](#) dialog similar to the [Open Method](#) dialog described on pg 44.

Save As...

Saves the print style under a newly specified name. When invoked, the [Save As](#) dialog will be displayed with the name of the original file. Change the file name to save the file as a new report style, or leave the name as it is to save the changes to the same report style. When the file name selected is already used, you will be notified.

Printer...

Selects the printer to be used. When invoked, the standard [Print Setup](#) dialog for selecting and setting up the printer will open.

Preview

Displays a preview of the printout (for more details see the chapter [Print Preview](#) on pg 352).

Print

Initiates the printing of the report. When invoked, the command will open the standard [Print](#) dialog. The date and time of the report will be printed on each page. For tabs where the *OVERLAY* mode is possible only information about the active chromatogram will be printed.

Print To PDF

Prints the report to *.PDF file. Using the button invokes the [Print to PDF](#) dialog where the user can select filename and location of the resulting *.PDF file.

Note: *.PDF files can be read by freely available Adobe Acrobat Reader.

Send PDF


Sends the *.PDF file as an email attachment. The command will automatically print the report to *.PDF file, create new message in your default email client and enclose the created *.PDF file as an attachment. Then you can fill in the remaining information (recipient, text of the message, etc.) into the email message and send it.

8.1 List of tabs:

- [Page Setup](#) - Sets up the printed page.
- [Lab. Header](#) - Sets up the page header.
- [Report Header](#) - Reported data about the analysis.
- [Method](#) - Sets up the method printout.
- [Calibration](#) - Sets up the calibration printout.
- [Chromatogram](#) - Sets up the chromatogram printout.
- [Results](#) - Sets up the results printout.
- [Sequence](#) - Sets up the sequence printout.
- [SST](#) ✓ full version - Sets up the system suitability report.
- [PDA](#) ✓ full version - Sets the printout for the **PDA** method.
- [NGA](#) ✓ full version - Sets the printout for the **NGA** method.
- [Audit & Signatures](#) ✓ full version - Sets up the audit trail and electronic signatures report.

The report itself is ordered in the order of the tabs shown - to change the order of the item groups in the resulting reports, re-order the tabs in the [Report Setup](#) dialog by using the *Move to the Top/Move Up/Move Down/Move to the Bottom* commands from the local menu on each tab. The [Page Setup](#) tab cannot be moved down, same as the [Audit & Signatures](#) tab cannot be moved up.

Most of the tabs contain the following two checkboxes:

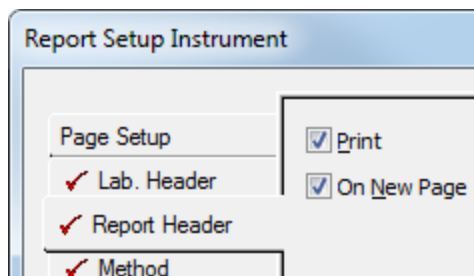

Print

Allows/disallows the printing of the relevant part. If checked, the symbol ✓ will be shown before the tab name and other fields on the tab will become editable.

On New Page

If checked, the relevant part will be printed on a new page.

8.2 Page Setup

Specifies the setup of the printed page.

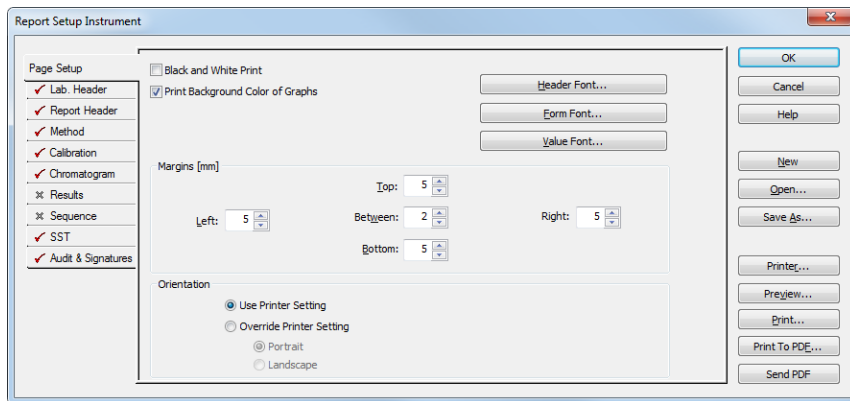


Fig 155: Report Setup - Page Setup

Black and White Print

Prints the report in black and white color (even on color printers).

Note: Some black&white printers in MS Windows are identified as color printers. In such case it is necessary to use the *Black and White Print* checkbox to ensure that the B&W optimization will be applied. These printers would otherwise, for example, not print background colors of the graph.

Print Background Color of Graphs

Prints the background color of the graphs. If not checked, the graph background will be white.

Header Font...

Selects the font for the report header. Invoking the command will open the standard [Font](#) dialog.

Note: There is always the method name, date and time of the printing and the page number displayed in the header. To hide this header, use white color for the *Header Font*.

Form Font...

Selects the font for item names and table headers. Invoking the command will open the standard [Font](#) dialog.

Value Font...

Selects the font for values. Invoking the command will open the standard [Font](#) dialog.

Caution:

In case you want to print PDF Reports in Chinese:

- You have to install True Type font supporting Chinese signs on your system.
- You have to set this font to use in **Clarity** reports.

Margins [mm]

Sets the margins. The allowed range is 0 to 50 mm.

Left, Right, Top, Bottom

Specifies the corresponding margin.

Between

Indicates the space left blank between individual parts of the printed report.

Orientation

Sets the orientation of the printed page. This may be either left on the selected printer setting or overridden by custom page orientation, which is advantageous when using two report styles with different page orientation on the same printer.

Use Printer Setting

Uses the printer page orientation selection for the given report style.

Override Printer Setting

Overrides the printer page orientation selection by one of the available options - *Portrait* or *Landscape*. During the selection and preparation of the report style, you might be prompted to select the printer that will be used for the print.

The size and print quality of the printed page are set in the [Print Setup](#) dialog of the selected printer (opened by the *Printer* command in the [Report Setup](#) dialog).

8.3 Lab. Header

Specifies the page header.

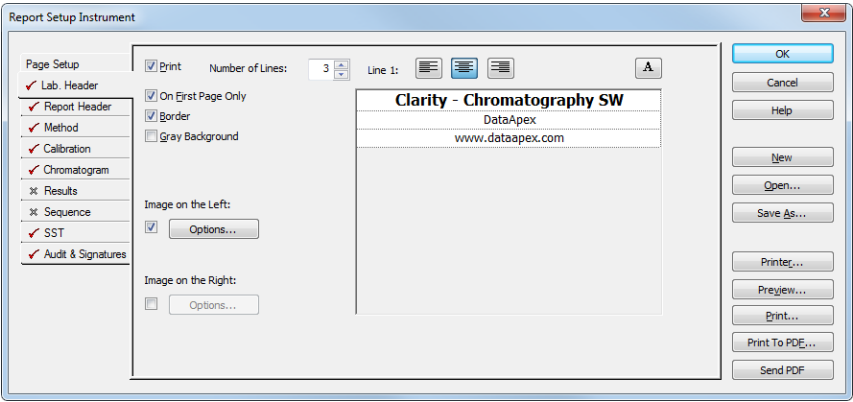


Fig 156: Report Setup - Lab. Header

On First Page Only

Sets the print of the header on the first page only.

Border

When checked, places the header in a frame made of thin lines.

Gray Background

While checked, prints the header with a gray background (blue when using a color printer).

Number of Lines

Defines the number of lines in the header. New lines will be appended/removed to/from the end of the current list. The maximum number is 9.

Line 1(..9)

Indicates the current line for the purpose of formatting. Use arrows or click a line to make it current, then use one of the icons to change the alignment or format of the text:

Tab 22: Description of icons on the Report Setup - Lab. Header tab

Icon	Description
	Left-aligns the text on the current line.
	Centers the text on the current line.
	Right-aligns the text on the current line.
	Selects the font for the current header line. Invoking the command will open the standard Font dialog.

The changes in the header are then shown in the preview. Click the corresponding line and enter the required text, or click the existing text to edit it.

Image on the Left (Right)

Mark the checkbox to print an image on the left (right) side of the page header. Use the *Option...* button to define the picture.

Options...

Invoking this button opens the [Set Image Options](#) dialog where the desired image and its dimensions can be set.

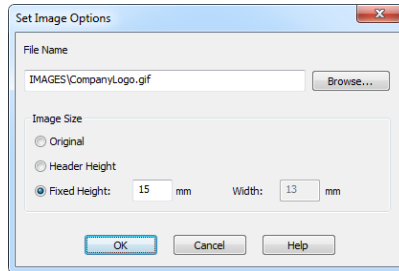


Fig 157: Set Image Options

File Name

Displays the filename of the image including its path (relative to the **Clarity** root directory). **Clarity** supports image files of following formats: *.BMP, *.GIF, *.JPG, *.ICO, *.TIF, *.TGA, *.PCX, *.WMF, *.PNG.

Image Size

The image size can either correspond to the specifications in the file (*Original*) or can be resized to the height of the header (*Header Height*) or to explicitly specified dimensions (*Fixed Height*).

If the image and text do not fit to the page width, **Clarity** will provide a warning when opening the [Print Preview](#) dialog, divide the header in thirds (halves when only one image is selected) and print its contents without respect to the actual dimensions.

Caution: Check the page header using the *Preview* button to see whether the image and text fits on the page.

Caution: The size of the page depends on the selected printer. Re- check the page when you switch to another printer.

8.4 Report Header

Selects the information that will be used in the report header. These information may include the origin of the chromatogram, date and time of analysis, user account used, calibration file name, name of the report style, data from the chromatogram header or directory where the chromatogram is stored.

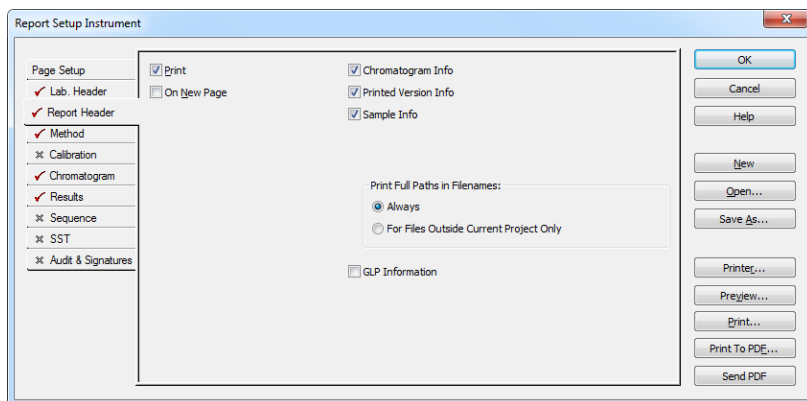


Fig 158: Report Setup - Report Header

Chromatogram Info

Adds the detailed info about the origin of the chromatogram to the printed report. Following parameters are added:

Filename - shows the filename of the chromatogram.

Origin - defines whether the chromatogram was *Acquired*, *Imported* or *Self* created by mathematical operations.

Project - shows the name of the project used to create the chromatogram.

File Created - shows the date and time when the chromatogram file was created (according to the Windows file system).

Acquired Date - shows the date and time when the chromatogram was acquired (the end of acquisition as stored in the chromatogram).

By - displays the name of the user account used to acquire the chromatogram.

Printed Version Info

Adds the detailed info about the print to the printed report:

Printed Version - shows the date and time in which the chromatogram version being printed was saved, thus enabling to trace the given chromatogram version.

Report Style - shows the report style used to print the report.

Calibration File - shows the name of the calibration file used.

Printed Date - shows the date and time when the report was printed.

By - displays the name of the user account used to print the report.

Sample Info

Adds the detailed info about the sample to the printed report. This info comes from the chromatogram header and is defined in the [Sequence](#) or [Single Analysis](#) dialog. The fields reported are **Sample ID**, **Sample**, **Inj. Volume**, **Amount**, **ISTD Amount** and **Dilution**.

Print Full Paths in Filenames

Determines the format of filenames by switching on and off the full path in the name of the files in the given project directory. For files outside of the project directory the full path is always printed.

Always

The full path will be printed for all files.

For Files Outside Current Project Only

- The files that are in the appropriate directory specified by the project will be printed without path.
- Files that are in any other directory or subdirectory will be printed with full path.

GLP Information

Adds the information about the computer, **Clarity** version and methods used for obtaining the printed information to the *Chromatogram Info* and *Printed Version Info* sections. If none of these sections is selected for display, the *GLP Information* checkbox will be disabled.

8.5 Method

Determines the elements of the method that will be printed.

Note: Parts of the method report dedicated to a particular device are printed only when this device is configured on the instrument in the moment of printing.

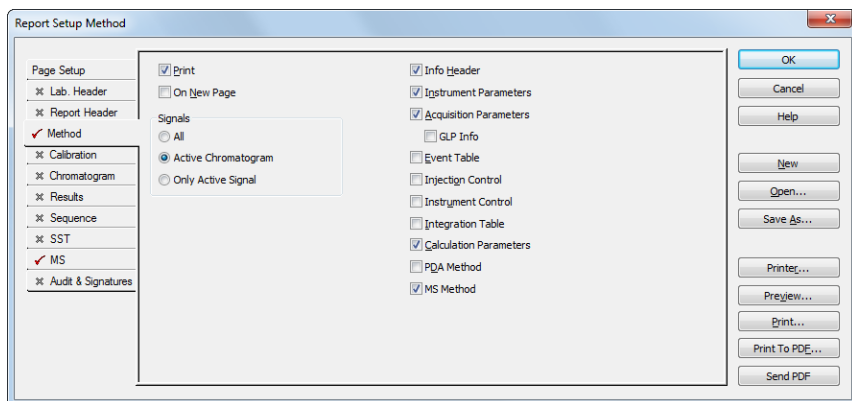


Fig 159: Report Setup - Method

Signals

Determines which signals (detector) information will be printed when using a multi-detector configuration. You can choose to include either method parameters from all signals of all chromatograms (*All*), all signals from active chromatogram (*Active Chromatogram*) or only the parameters from the active signal (*Only Active Signal*).

Info Header

Adds the section containing method **Name**, method **Description**, the name of the **User Account** used to create the method and the time of the **Creation** and **Last Modification** parameters to the report.

Instrument Parameters

Adds the content of the text fields (**Column**, **Mobile Phase**, **Flow Rate**, **Pressure**, **Detection**, **Temperature** and **Note**) from the [Method Setup - Measurement](#) dialog to the report.

Acquisition Parameters

Adds *Autostop* and *External Start/Stop* sections from the [Method Setup - Measurement](#) tab, measurement scale from the [Method Setup - Acquisition](#) tab and the *Subtraction* section from the [Method Setup - Advanced](#) dialog to the report.

GLP Info

Adds the content of the [GLP Info](#) sub-tab of the [Measurement Conditions](#) tab in the [Chromatogram](#) window.

Event Table

Adds the **Event Table** from the [Method Setup - Event Table](#) dialog to the report.

Injection Control

Adds the injection control parameters from the [Method Setup - AS](#) tab to the report. This checkbox adds no parameters to the report if there is no directly controlled autosampler used.

Instrument Control

Adds any parameters set on the tabs of individual directly-controlled chromatographs ([Method Setup - LC](#) tab, [Method Setup - GC](#) tab) and other directly controlled instruments to the report in case these instruments are used. This section also includes settings of the signals from A/D converters used on the given Instrument when analog acquisition is used.

Integration Table

Adds the **Integration Table** from the [Method Setup - Integration](#) dialog to the report.

Calculation parameters

Adds the parameters from the [Method Setup - Calculation](#) tab and from the **Column Calculations** section of the [Method Setup - Advanced](#) tab to the report.

PDA Method

Adds the parameters from the [Method Setup - PDA Method](#) tab to the report. This section is only active when the method has **PDA** part. Detailed description is provided in a separate **PDA** manual.

MS Method

Adds the parameters from the [Method Setup - MS Method](#) tab to the report. This section is only active when the method has **MS** part. Detailed description is provided in a separate **MS** manual.

8.6 Calibration

Determines the contents and layout of the calibration printout.

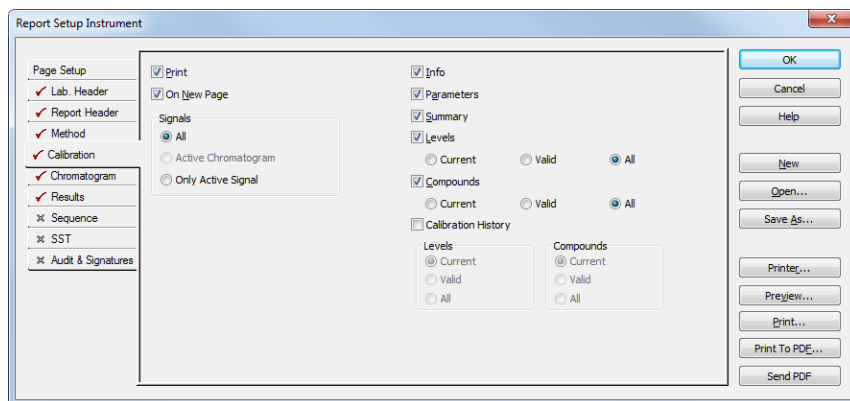


Fig 160: Report Setup - Calibration

Signals

Determines which signals (detectors) will be printed when using a multi-detector configuration. Include either calibration parameters from all signals (*All*) or only the parameters from the active signal (*Only Active Signal*).

Info

Prints the name and description of the calibration.

Parameters

Prints the parameters from the [Calibration Options](#) dialog.

Summary

Prints the **Calibration Summary Table** from the [Compounds](#) tab of the [Calibration](#) window.

Levels

Prints Calibration Summary Table(s) for specified levels including the section related to the active level and signal only.

Current

Prints only the table which refers to the current level.

Valid

Prints only tables which contain at least one valid compound. A compound will be valid when the *Response* and *Amount* columns differ from zero and the *Used* checkbox in the **Calibration Summary Table** has been checked.

All

Prints the table for all calibration levels (including blank) irrespective of the fact if the level has any valid data.

Compounds

Prints calibration tables from tabs of individual compounds including the calibration graphs. Information concerning individual compounds is printed on separate pages.

Current

Prints only the table and graph of the current compound.

Valid

Only tables and curves of compounds having at least one valid level will be printed. A level is valid when columns *Response* and *Amount* in the calibration table of compound differ from zero and on the same level the parameter *Used* has been checked.

All

Prints calibration tables and graphs for all compounds.

Calibration History

Allows to print the history of calibration points for the selected combination of calibration levels compounds.

Levels

Defines which levels will be used for the display of calibration point history. Options are *Current*, *Valid* and *All*.

Compounds

Defines which compounds will be used for the display of calibration point history. Options are *Current*, *Valid* and *All*.

8.7 Chromatogram

Sets the sections and layout of printed chromatograms.

All tags, descriptions and labels in the graphs are WYSIWYG (*What You See Is What You Get*). That means that the printed report will correspond to the display set on the screen.

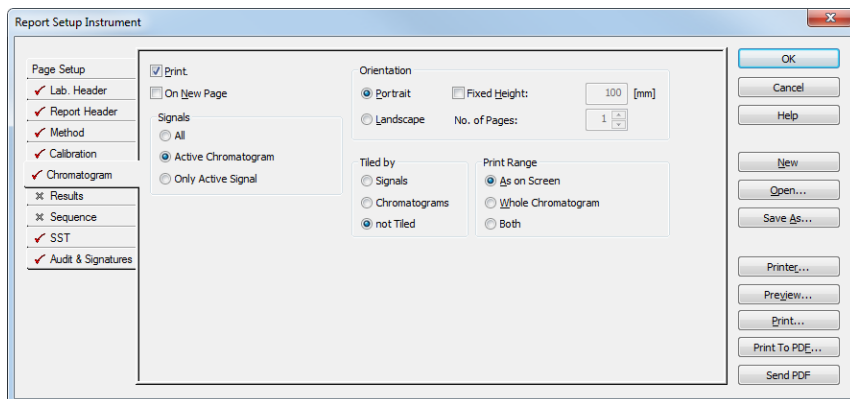


Fig 161: Report Setup - Chromatogram

Signals

Determines which signals (detector) information will be printed when using a multi-detector configuration. You can choose to include either method parameters from all signals of all chromatograms (*All*), all signals from active chromatogram (*Active Chromatogram*) or only the parameters from the active signal (*Only Active Signal*).

Orientation

Sets the type of the chromatogram orientation:

Portrait

Prints the chromatogram crosswise (the time axis is parallel to the printed lines). The width will always be given by page width and pre-set borders. Unless *Fixed Height* has been selected, the chromatogram will be printed with a fixed ratio of sides (approximately 2:3) on a single page.

Landscape

Prints the chromatogram on a separate page with the time axis perpendicular to the printed lines. The size will be given by page size and pre-set borders.

Fixed Height

If the graph orientation is set to *Portrait*, the chromatogram will be printed with a fixed height instead of the fixed ratio 2:3. The permitted

lower height range is *30 mm*, the upper height range is not set - the graph will be scaled to the page height if larger than the page itself.

No. of pages

If the graph orientation is set to *Landscape*, this field determines the chromatogram length display by number of pages. The permitted range is from 1 to 9.

Tiled by**Signals**

Prints multiple curves in separate graphs tiled by signals (each signal will have its own graph in the report).

Chromatograms

Prints multiple curves in separate graphs tiled by chromatograms (each chromatogram will have its own graph in the report, possibly consisting of more than one signal in case of multi-detector chromatograms).

not Tiled

Prints all curves overlaid in one graph.

Print Range**As on Screen**

Prints only the portion of the chromatogram which is visible on screen.

Whole Chromatogram

Prints the entire chromatogram(s) regardless of the size of the cut displayed in the [Chromatogram](#) window.

Both

Prints both the whole chromatogram and the active cut-out zoomed, in this order.

8.8 Results

Determines the content and layout of the printed analysis results.

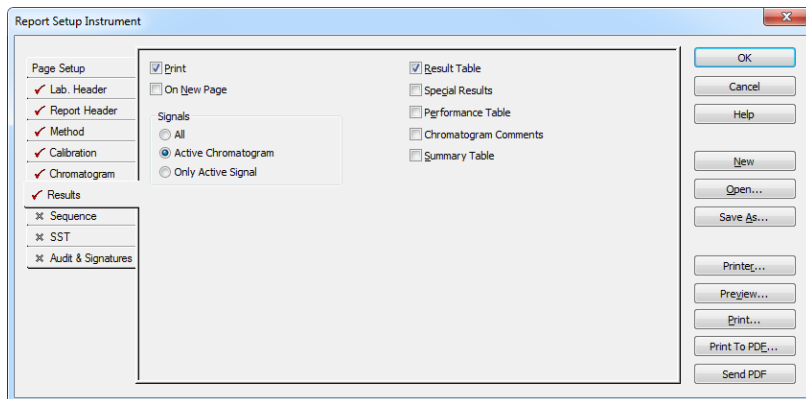


Fig 162: Report Setup - Results

Signals

Determines which signals (detector) information will be printed when using a multi-detector configuration. You can choose to include either method parameters from all signals of all chromatograms (*All*), all signals from active chromatogram (*Active Chromatogram*) or only the parameters from the active signal (*Only Active Signal*).

Result Table

Prints the **Result Table** from the [Chromatogram - Results](#) tab.

Special Results

Prints special **Result Tables** generated by some control modules such as Virtual FC module.

Performance Table

Prints the **Performance Table** from the [Chromatogram - Performance](#) tab.

Chromatogram Comments

Adds the chromatogram comments entered into the [Rich Text Edit](#) dialog from the [Single Analysis](#) dialog using the *Comments...* button or from the [Measurement Conditions](#) tab of the [Chromatogram](#) window using the *Edit...* button to the report.

Summary Table

Prints the **Summary Table** from the [Chromatogram - Summary](#) tab.

8.9 Sequence

Determines the parts of the sequence displayed in the report.

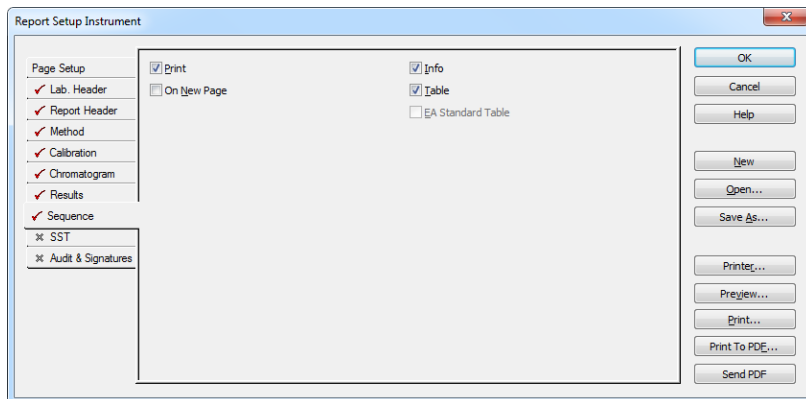


Fig 163: Report Setup - Sequence

Info

Prints the description of the sequence - namely the **Sequence Name** and **Description**, name of the **User Account** used to create the sequence, date and time when the sequence was **Created** and **Modified**, the **Type** of the sequence (*ACTIVE* or *PASSIVE*) and the **Idle Time** parameter set. All these parameters are described in the chapter describing the [Sequence](#) window (for more details see the chapter **Sequence** on pg 294).

Table

Prints the sequence table.

8.10 Audit & Signatures ✓ Full version

Governs the print of local audit trails and electronic signatures.

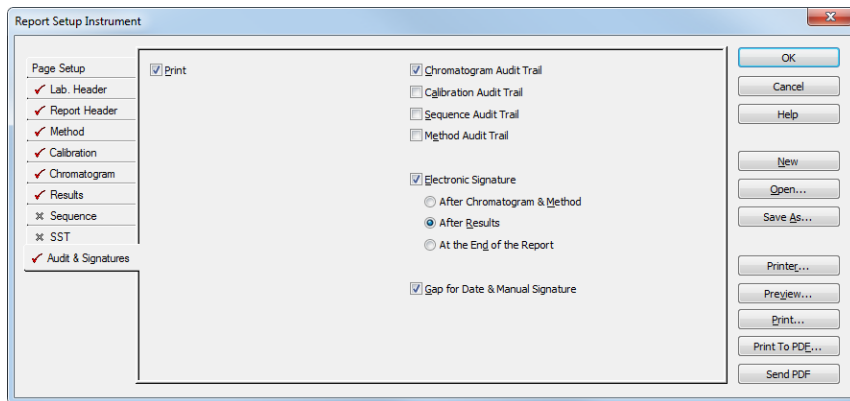


Fig 164: Report Setup - Audit & Signatures

Chromatogram Audit Trail

Prints the operations made with the **Active Chromatogram** logged in the [Chromatogram Audit Trail](#) window.

Calibration Audit Trail

Prints the operations made with the **Calibration** file logged in the [Calibration Audit Trail](#) window.

Sequence Audit Trail

Prints the operations made with the **Sequence** logged in the [Sequence Audit Trail](#) window.

Method Audit Trail

Prints the operations made with the **Method** file logged in the [Method Audit Trail](#) window.

Electronic Signature

Determines the placement of the electronic signature:

After Chromatogram & Method

Places the signature after the chromatogram and its method.

After Results

Places the signature after the result tables.

After the End of the Report

Places the signature at the end of the report.

Gap for Date and Manual Signature

Prints a pre-printed blank line at the end of the report so that the date and signature may be filled in by hand after the printout.

Note: Printing the entire station's logs (**Session Audit Trail** and **Daily Audit Trail**) can only be done directly and manually from the [Audit Trail](#) window using the *File - Print* command.

8.11 Print Preview

Invoking the *Preview* button in the [Report Setup](#) dialog or using the *Print Preview...* command opens the [Print Preview](#) dialog which displays a preview of the report prior to printing it.

Note: If there is no printer installed in MS Windows the *Print* and *Print Preview* commands will be disabled.

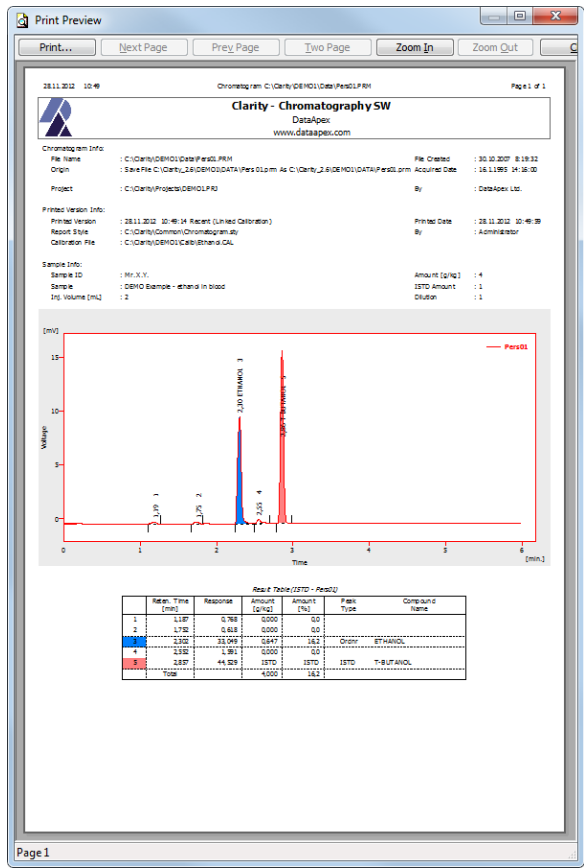


Fig 165: Print Preview

Print...

Initiates the print of the report. When invoked, the standard [Print](#) dialog is opened.

Next Page

Displays the next page of the report. Only active for reports consisting of more than one page and when the last page is not already displayed.

Prev Page

Display the previous page of the report. Only active for reports consisting of more than one page and when the first page is not already displayed.

Two Page/One Page

This button switches between the display of one page and the display of two pages of the preview on one screen. Only active for reports with more than one page.

Zoom In

Enlarges the size of the preview by one step. If the maximum enlargement has already been reached, switches to the least enlargement possible instead.

Zoom Out

Reduces the size of the preview by one step. If the maximum reduction of the size has already been reached, switches to the maximally enlarged view instead.

Close

Closes the preview.

Print to PDF

Prints the report to the *.PDF file. Using the button invokes the [Print to PDF](#) dialog (similar to [Open Method](#) dialog) where the user can select filename and location of the resulting *.PDF file.

Note: *.PDF files can be read by freely available Adobe Acrobat Reader.

Send PDF

Sends *.PDF file as an email attachment. The command will automatically print the report to *.PDF file, create new message in your default email client and enclose the created *.PDF file as an attachment. Then you can fill in the remaining information (recipient, text of the message, etc.) into the email message and send it.

8.12 Print

Invoking the *Print...* command, pressing Ctrl + P combination or using the *Print...* button from the [Report Setup](#) or [Print Preview](#) dialogs will opens the [Print](#) dialog, which is used for specifying the printer and starting the actual printing. The layout of this dialog will vary depending on which version of Windows operating system is installed and the local settings being used.

Pressing the  directly from toolbar will print out the report immediately, according to actual settings.

Note: If there is no printer installed in MS Windows the *Print* and *Print Preview* commands will be disabled.

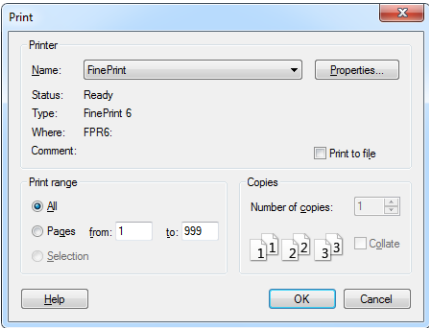


Fig 166: Print

9 Troubleshooting

This section briefly overviews possible error messages and the procedures which will lead to the identification of the problem and its solution.

9.1 Report error after Clarity crashed

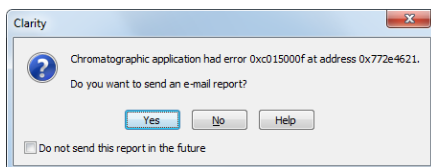


Fig 167: Send Report dialog

If **Clarity** encounters serious problems and crashes, it will invoke a dialog during the crash that offers the possibility to send report about the problem to **DataApex**. Such report would help us locate the conditions leading to crash and improve **Clarity** accordingly.

However, in some cases, the **Clarity** may crash without producing any message. In such case, similar error message is prompted at the next start of the **Clarity** station.

More details about sending an e-mail report can be found in the chapter "**Send Report by E-Mail**" on pg 38. If you check *Do not send this report in the future* checkbox, **Clarity** will not open this dialog in case of next crashes.

9.2 How to get help

Invoke the help in **Clarity** and try to use fulltext search.

Consult the **Troubleshooting** chapter of the **Getting Started** manual.

If you will not find your answers here, use the www.dataapex.com website where the Support menu will navigate you to frequently asked questions (FAQ), **Clarity** forum or contact to **DataApex** helpdesk.

9 Appendices

9.3 List of terms

To prevent any errors or misunderstandings, the following list of terms (often-used or differing) has been provided.

21 CFR Part 11

A directive which specifies the conditions under which a company may use electronic records and signatures. The directive is mandatory for companies that use computer system designed for the manufacturing or distribution of products which are subject to regulation by the FDA (known as, "regulated products").

Access rights

A system which governs the activities that specified users are allowed to perform. Among other things, a user may be restricted in their authorization to change fundamental station settings, access certain instruments, inspect or change the data of other users, etc.

Active chromatogram

The chromatogram name which is displayed in the header of the [Chromatogram](#) window. All data being displayed and operations performed will only relate to this active chromatogram. The name of the active chromatogram is highlighted in the key by bold letters.

Active chromatogram signal

When performing a multi-detector measurement, each chromatogram file may contain a record from up to 16 detectors (this is what we call "signals"). The active signal in the chromatogram is the signal whose name is being displayed behind the name of the chromatogram in the window header and header of the tables (divided by a hyphen).

The name of the active signal together with the name of the chromatogram are highlighted in the graph key by bold letters.

Active calibration signal

When performing a multi-detector measurement, each calibration file may contain calibration curves for up to 16 signals, the same as with chromatograms. Only one of these signals may be active at a time and this signal is modified.

The active calibration signal (i.e. its table or curve) is recognizable by its name highlighted in the header of the **Calibration Summary Table** or its curve. At the same time, all values that are signal specific will be displayed in the same color as the active signal.

Active detector

A detector whose measurement and calculation parameters are displayed in the [Acquisition](#) or [Calculation](#) tabs in the [Method Setup](#) dialog with the

help of the *Select Detector* field. The field is visible only when performing a multi-detector measurement.

Analogous *Select XXX* fields may be present for other devices than detectors - for example valves on the [Method Setup - Valve](#) tab if more than one valve is present.

AIA format

Also named NetCDF format, export format serving as a common format for chromatographic data. The data are saved in the vector format with the header containing the information on the data type and chromatogram information. **Clarity** only exports and imports raw data, not the results.

Calibrated group

A group of peaks calibrated as a single component.

Calibration file (Calibration)

Contains calibration curves and further data for calibrated calculations. When performing a multi-detector measurement, each substance will have its own calibration curves for all detectors.

Calibration standard

A chromatogram containing known amounts of some compounds that is used to calibrate/recalibrate at a specified calibration level.

Channel

Part of the converter which enables the independent collection of data from one single detector. **DataApex** produces A/D converters that has 1, 2 or 4 channels.

Clarity Offline (used to be Clarity EVAL)

A version of the **Clarity** station that, although otherwise fully functional, does not allow the data acquisition.

Clipboard

A **Windows** tool that uses the *Copy*, *Cut* and *Paste* commands for the simple transfer of data between items, including those in other applications.

Closed window

A closed window is invisible and its activity has been terminated. A window is usually closed using the *Close* command (see also the **Hidden window** term).

Configuration of the station

Sets the number of Instruments, the allocation of the A/D converters, directly controlled chromatographs, autosamplers, pumps and other devices.

Data processing string

A series of operations comprising the entire procedure used for everything from processing data from an analysis to printing the report. This

procedure is shown graphically in the Instrument window using icons representing the main commands.

Desktop file - extension DSK

A file used to store the size, location and appearance of all windows and parameters of the Instruments.

Direct control of chromatograph (autosampler, pump)

An integrated additional module for selected chromatographs (pumps, autosamplers etc.) enabling the direct control and monitoring of selected parameters. For example, directly controlled detectors (GC or LC chromatographs with detectors) might not need A/D converter as they gather digital data directly from the controlled instrument.

Electronic signature

An electronic signature is an substitution of the handwritten signature for electronic documents. These electronic signatures thus must fulfil the same function and prerequisites. Meaning that it is a similar process of marking a document for protection and that it must be unique, clear and ensure the immutability of the document.

File sharing

Clarity enables file sharing between multiple stations. Any changes (saving) of such files are only enabled in the station that first opened it. Modifications performed with the shared file will take effect in the other stations only after the file has been reloaded. **Clarity** does not enable file sharing between Instruments of a single station. On the other hand,

Clarity does enable projects to be shared between multiple Instrument of the same station.

Note: It is still not possible to use the same file on multiple Instruments even if they share the same project. The same project on two Instruments of the same station are useful when one instrument is *OFFLINE* and used to prepare methods for the other Instrument.

GLP

Good Laboratory Practice (GLP) is an internationally agreed upon system for ensuring and monitoring the quality of laboratory work. This is verified and its fulfillment confirmed by the issuance of a certificate.

Hidden window

A window which is not visible or has been “hidden”. Its activity has not been terminated but rather suspended or is proceeding in the background. Click the minimize button  to hide a window (see also the **Closed window** term).

Information table

A table in the [Instrument](#) window displaying the Instrument name, the measurement methods and the identification of the measured chromatogram.

Instrument

(a) A part of the program designated to measure and evaluate analyses that share a common time base (from one chromatograph). The Instrument is indicated by the [Instrument](#) window described on pg 40., from which further windows and dialogs can be opened, e.g.

[Chromatogram](#), [Calibration](#), etc. Detectors and control modules can be configured into individual Instruments. **Clarity Lite** has only one Instrument available. In this meaning, the word "Instrument" is written with the capital I.

(b) Sometimes also a chromatograph or generally any piece of equipment in the laboratory.

Instrument directory

The directory identified in the main [Clarity](#) window for each Instrument, where projects (i.e. project directories) are sought.

Integration Table

Constitutes a part of the method and contains a list of all modifications of the baseline or peaks that have been or will be effected in the chromatogram measured by that method.

Locked file (READ ONLY)

A file that can only be read and never modified.

Locked instrument

An Instrument in the protected mode with locked control. Lock mode can be activated using a command from the *Lock* submenu in the main [Clarity](#) window. An Instrument is unlocked with the same password as was used to log on to the Instrument.

Main station directory

The directory where the executable file of the station (CLARITY.EXE, CLARITYLITE.EXE in case of **Clarity Lite**) is stored (C:\CLARITY by default). The same directory contains all configuration files, the directories of individual projects, PROJECTS directory holding the project files, COMMON directory with print styles and other files.

Method file - extension MET

A template method file with a description of all parameters affecting the analysis.

Model Method

A chromatogram method marked as **Model Method** that can be easily used with other chromatograms by invoking the *Copy from Model* command.

Multi-detector configuration (measurement)

Setting of the station so that it measures, saves and evaluates one analysis by simultaneously using a signal from multiple detectors.

Password

A string of several characters which are necessary for working in the **Protected Mode** of the station. Each user has his own password. The minimal length of the password may be set centrally in the [User Accounts](#) dialog.

Printout

The layout of individual parts of analysis results and the setting of parameters controlling the resulting report. The layout is defined by the employed report style.

Project directory

Directory used to store all working files of an Instrument in which the homonymous project has been opened.

Project file - extension PRJ

The file that specifies the directories in which all user files will be stored.

Protected mode

A regime which enables only authorized users (users who have an account assigned in the [User Accounts](#) dialog) to work with the station and access data.

Raw data file - RAW

A file used in previous versions of the **CSW10 - 17** and for the external PAD unit to save non-analyzed and non-adjusted data from the analysis. The **CSW32** and **Clarity** stations save raw data directly into the chromatogram file, the *.RAW data file is used only for the gathering of the data during the acquisition (chromatogram file (*.PRM) is created only after the acquisition ends).

Report style

An outline determining which analytical results will be incorporated into a report and subsequently printed. Each report style will be stored in a separate report style file.

Report style file - extension STY

A file storing information which determines what to print and how to print it.

Sequence file - extension SEQ

A file that defines an automatic sequence of analyses.

Splitter

A resettable bar dividing certain windows (e.g., the [Chromatogram](#) or [Calibration](#) windows) into several panes. The bar can be moved to change the size of individual panes.

SST

The **System Suitability Test** is an extension module designated to validate the chromatographic system on the basis of evaluation of chromatograms obtained from the **Clarity** station.

Station configuration

Determines the number and arrangement of Instruments and their interconnection to possible A/D converter board(s) and directly controlled chromatographic equipment. Also allows to name the connected detectors and set signal units for them.

Status table

A table displayed in the [Instrument](#) window containing the names of system files, the status of the analysis currently in progress, potentially the retention time and identification number of a sample.

System files

The decisive files for working with the station and individual data files: **project**, **method**, **report style**, **sequence** and **calibration** files.

Tabs

Tab is the term for a section of the window or dialog where several layers of control elements are "stacked" on itself. The system is frequently used in dialogs and also in the [Chromatogram](#) and [Calibration](#) windows. Click the only visible part of the tab - that with the tab name - to switch to the corresponding tab.

Template method

A method file that serves as a template for new analyses. Its name is displayed in the **Information Table** of the [Instrument](#) window and its contents are always copied to the newly created chromatogram.

Toolbar

A block of icons symbolizing a set of commands. Each toolbar can be situated anywhere on the screen and be configured by accessing the *View - Customize...* command in the window or by right clicking the mouse button on any toolbar to access the local menu.

Unprotected mode

A regime of the station that enables any user to work with the station and access its data. No passwords are necessary for opening the Instruments.

Window title bar

Title bar is the upper band of each window. It is comprised of the button which activates the control menu of the window, the window name, the minimize button and the maximize button.

9.4 Shortcuts

MS Windows shortcuts:

Enter

Applies or implements the selected command. The command that has been selected will be highlighted in the menu; in a dialog the borderline of the button will be emphasized by a continuous or broken line).

From the menu the required command can be selected using the cursor keys, and from a window using the Tab key.

Esc

In a dialog, the Esc key substitutes for the Cancel key and will close a dialog without saving the changes.

Hides an expanded menu.

Tab

Gradual selection of commands, parameters, edit lines and buttons in the active window. A selection is usually completed using the **Enter** key.

Space bar

Rapidly switches between selected parameters by checking or unchecking them.

Alt

Selects the first menu item.

Alt + letter

Rapidly selects a command or parameter that has the selected letter underscored.

Tab 23: Clarity shortcuts:

Shortcut	Function	Validity
F1	Help	Everywhere
F2	Edit table mode	
Ctrl + N	New file	
Ctrl + A	Select all items in the list/table	
Ctrl + O	Open file (<i>Open</i>)	
Ctrl + S	Save file (<i>Save</i>)	
Ctrl + W	Close current file (<i>Close</i>)	
Ctrl + Shift + W	Close all opened files (<i>Close All</i>)	
Ctrl + P	Print (<i>Print</i>)	
Ctrl + X	Cut (<i>Cut</i>)	
Ctrl + C	Copy (<i>Copy</i>)	
Ctrl + V	Insert (<i>Paste</i>)	

Shortcut	Function	Validity
Ctrl + Z	Cancel the last command (<i>Undo</i>)	Tables
Ctrl + Shift + Z	Repeat the last command (<i>Redo</i>)	
Ctrl + I	Insert and create lines (<i>Paste Insert</i>)	
Delete	Erase the table field	
Ctrl + E	Automatic export of data from Chromatogram window according to settings in Export Data dialog.	Chromatogram
Ctrl + -	Display the previous cut-out (<i>Previous Zoom</i>)	Chromatogram Calibration Data Acquisition
Ctrl + +	Display the next cut-out (<i>Next Zoom</i>)	
Ctrl + *	Display in original size (<i>Unzoom</i>)	
Ctrl + R	Start an analysis (<i>Run Single</i>)	Data Acquisition
Ctrl + Q	Start a sequence (<i>Start, Run Sequence</i>)	Sequence Data Acquisition
Ctrl + Space bar	Pause a sequence (<i>Pause</i>) / Resume running of a set sequence (<i>Resume</i>)	
Ctrl + B	Immediately terminate a sequence (<i>Abort</i>)	
Ctrl + H	Process the sample currently being measured without terminating the analysis (<i>Snapshot</i>)	
Ctrl + T	Terminate a sequence after the completion of a sample measurement (<i>Stop</i>)	
Ctrl + ↓	Skip the current sample (<i>Skip</i>).	Sequence
Ctrl + ←	Repeat an injection from the current vial (<i>Repeat Injection</i>)	
Ctrl + E	Reset the flags in the Sts. column (<i>Reset Status (Sts.)</i>)	

10 Technical Specification

10.1 Declaration of Software Validation

You can find **Declaration of Software Validation** certificate on **Clarity** Installation DVD-ROM in \DOC PDF\DATASHEETS\D021- CLARITY-DECLARATION-OF-SW-VALIDATION.PDF.

Note: Actual version of this certificate is available for download as the D021 datasheet from www.dataapex.com.

10.2 Command line parameters

It is possible to specify command line parameters for automating certain functions. If not specified otherwise, CLARITY.EXE file can be run with the same or different set of parameters again and again. This creates new instance, which is temporary and just passes new parameters into **Clarity** station that is already running.

Some of the command line parameters may be also defined in a DATAAPEX.INI file. This file is not present at default, but when placed into the C:\WINDOWS directory, settings in it will override any of the same parameters used in the command line. DATAAPEX.INI file supports the use of following command line parameters: *bad_trace_cnt*, *create_emf_during_pdf*, *dump_memory*, *enable_asserts*, *mail_dump*, *report_cmdline_in_bad_trace*.

All command line parameters are processed in the order they are set on the command line. It is thus usually necessary to open the file first, before making any changes to it (for example, open calibration, open chromatogram, recalibrate the calibration on the given level, close chromatogram and close calibration, in this order).

Caution: The list contains complete set of parameters. In **Clarity Lite**, some of them cannot be used due to product limitations (e.g. **Clarity Lite** has only one Instrument etc.).

No wildcards (*, ?) are allowed in the command line.

All parameters (but not file paths etc.) are case-sensitive, unless stated otherwise. No spaces are allowed between the command and the argument and in the argument itself.

Example: *clarity.exe command=value* is correct while *clarity.exe command = value* is wrong.

If you want to use spaces or other white space in parameters, you may use the quotation marks to do so.

Example: *clarity.exe c="C:\Documents and Settings\Smith\Desktop\Ethanol in blood.met"* or *clarity.exe user="John Doe"*.

10.2.1 List of commands

abort

description:

Stops the single run/sequence measurement on the specified Instrument. This parameter is suitable for automatic testing. Does the same as **stop** command line parameter, but it does not store the chromatogram measured up to the time of abort.

syntax:

abort=VALUE

possible values:

1 (for Instrument 1), 2 (for Instrument 2), 4 (for Instrument 3), 8 (for Instrument 4), and their binary combinations.

implicit value:

None.

example:

clarity.exe abort=7 (will abort single analysis/sequence on Instruments 1, 2 and 3)

bad_trace_cnt

description:

Sets the number of lines in the BADTRACE.TXT file. This parameter may be also used (and is overwritten by settings) in DATAAPEX.INI file.

syntax:

bad_trace_cnt=VALUE

possible values:

0 - virtually any number

implicit value:

100

example:

clarity.exe bad_trace_cnt=1000

cal_apply

description:

Applies the specified calibration to the active chromatogram. It is recommended to use absolute paths because the calibration file will not be searched for in the working directory but in the calibration directory of the active project.

syntax:

cal_apply=PATH\FILENAME

possible values:

A valid path to the calibration (*.CAL) file.

implicit value:

None.

example:

clarity.exe cal_apply=c:\tmp\1.cal

cal_save*description:*

Saves the active calibration. Use together with the **instrument** parameter (which is ignored by **Clarity Lite**) to specify the Instrument whose **Calibration** window holds the calibration file you are interested in. Calibration file can be opened by specifying its path and filename. It is recommended to use absolute paths.
This command is ignored in GPC mode - use **gal_save** instead.

syntax:

cal_save

possible values:

None.

implicit value:

None.

example:

clarity.exe i=2 cal_save - saves the calibration opened in the second Instrument.

cal_save_as*description:*

Saves the active calibration under the specified name. Use together with the **instrument** parameter (which is ignored by **Clarity Lite**) to specify the Instrument whose **Calibration** window holds the calibration file you are interested in. Calibration file can be opened by specifying its path and filename. It is recommended to use absolute paths.
This command is ignored in GPC mode - use **gal_save_as** instead.

syntax:

cal_save_as=PATH\FILENAME

possible values:

A valid path to the calibration (*.CAL) file.

implicit value:

None.

example:

clarity.exe cal_save_as=c:\tmp\results.cal

cal_type*description:*

Changes the type of calibration calculation for the active chromatogram. The parameter allows to specify calibration type for each signal of the chromatogram -the first character is corresponding to the first signal, second character to the second signal and so on.

syntax:

cal_type=TYPE

any combination of _, U, I and E. For each signal in the chromatogram, a separate symbol is needed.

Characters used mean:

possible values:

- _ (underline) - do not change
- U - uncalibrated
- I - ISTD
- E - ESTD

implicit value: _ (underline) for each signal.

example: `clarity.exe cal_type=_ UIE` (the first signal is left unchanged, the second is set to UNCAL, the third to ISTD, the fourth to ESTD).

cfg or g

description: Opens the **Clarity** with the specified *.CFG file. Must be used when opening **Clarity** for the first time, otherwise it is ignored.

syntax: `g=PATH\FILENAME`

possible values: The path to *.CFG file.

Empty.

If the path is not specified, the system will look for the *.CFG file in the current directory. If not found, the system will look for the *.CFG file in the same directory where the CLARITY.EXE file is stored.

implicit value: When the path is specified absolutely, the **Clarity** will start with the *.CFG file specified.

When the path is specified relatively, the *.CFG file will be searched for in the directory relative to the current directory. Working directory will then be set to the new location.

example: `clarity.exe g=C:\CLARITY\user1.cfg`

clear_cal

description: Clears the responses in the opened calibration file. When used with **recal_level**, the calibration is cleared first and recalibrated after that.

syntax: `clear_cal`

possible values: None.

implicit value: None.

example: `clarity.exe test.cal clear_cal`

clear_gpc_cal*description:*

Clears the responses in the opened GPC calibration file. When used with **gpc_recal_stdno** and **gpc_recal_type**, the calibration is cleared first and recalibrated after that.

*syntax:**clear_gpc_cal**possible values:*

None.

implicit value:

None.

*example:**clarity.exe test.gal clear_gpc_cal***create_emf_during_pdf***description:*

Forces the creation of the *.EMF file from which the *.PDF files are converted. This parameter may be also used (and is overwritten by settings) in DATAAPEX.INI file.

*syntax:**create_emf_during_pdf**possible values:*

None.

implicit value:

None.

*example:**clarity.exe create_emf_during_pdf***delayed_exit***description:*

Exits **Clarity** after the period of time specified in seconds if it has no running acquisition. This parameter is suitable for automatic testing.

*syntax:**delayed_exit=VALUE**possible values:*

Numeric.

implicit value:

None.

*example:**clarity.exe delayed_exit=120***disable_message_boxes***description:*

Suppresses the showing of message boxes. Must be used when opening **Clarity** for the first time, otherwise it is ignored. This parameter is suitable for automatic testing.

*syntax:**disable_message_boxes**possible values:*

None.

implicit value:

None.

*example:**clarity.exe disable_message_boxes*

disable_send_report*description:*

Suppresses the sending of error reports. Must be used when opening **Clarity** for the first time, otherwise it is ignored. This parameter is suitable for automatic testing.

*syntax:**disable_send_report**possible values:*

None.

implicit value:

None.

*example:**clarity.exe disable_send_report***dsk or k***description:*

Opens the **Clarity** with the specified desktop (*.DSK) file. Must be used when opening **Clarity** for the first time, otherwise it is ignored.

*syntax:**k=PATH\FILENAME**possible values:*

The path and filename of the *.DSK file.

Empty.

implicit value:

If the path is not specified, the system will look for the *.DSK file in the current directory. If not found, the system will look for the *.DSK file in the same directory where the CLARITY.EXE file is stored.

*example:**clarity.exe k=C:\CLARITY\Peter.dsk***dump_memory***description:*

Causes to dump the memory into a dump file (*.DMP) in case **Clarity** crashes. If the memory is full, the dump file is about 200MB in size, while empty memory creates dump files about 200kB. This function can also be set via the DATAPEX.INI file that may be stored in WINDOWS directory. Must be used when opening **Clarity** for the first time, otherwise it is ignored. This parameter is suitable for troubleshooting. This parameter may be also used (and is overwritten by settings) in DATAPEX.INI file.

*syntax:**dump_memory**possible values:*

None.

implicit value:

None.

*example:**clarity.exe dump_memory*

enable_asserts*description:*

Specifies what to do if an ASSERT (software developer check point) emerges. Must be used when opening **Clarity** for the first time, otherwise it is ignored. This parameter is suitable for troubleshooting. This parameter may be also used (and is overwritten by settings) in DATAAPEX.INI file.

syntax:

`enable_asserts=VALUE`

0, 1, 2, ..., 7

The meaning of particular values when ASSERT emerges are:

0 - nothing happens

1 - note is written into BAD_TRACE.TXT

2 - note with stack unwind is written into BAD_TRACE.TXT

3 - note is written into BAD_TRACE.TXT and crash dump file is created - it can be used in

Debugging tools for Windows to solve the problem.

4 - note with stack unwind is written into BAD_TRACE.TXT and crash dump file is created

5 - ASSERT dialog box is shown, where the **Retry** button can be pressed. That causes the debugging of the application in **Visual Studio**.

6 - ASSERT dialog box is shown and note is written into BAD_TRACE.TXT

7 - ASSERT dialog box is shown, note is written into BAD_TRACE.TXT and crash dump file is created.

possible values:

0 when **enable_asserts** command is omitted in the command line, 6 when there is **enable_asserts** command without a value.

*implicit value:**example:*

`clarity.exe enable_asserts=7`

err*description:*

Stores the audit trail records also in the specified text file. File name is valid until it is replaced by a new one.

syntax:

`err=PATH\FILENAME`

possible values:

A valid path and filename.

implicit value:

None.

example:

`clarity.exe err=C:\tmp\error.txt`

exit*description:*

Immediately exits **Clarity**. Does not work if acquisition is running or if any modal window is opened. This parameter is suitable for automatic testing.

*syntax:**exit**possible values:*

None.

implicit value:

None.

*example:**clarity.exe exit***export_emf***description:*

Exports the currently opened chromatograms into enhanced Windows metafile.

*syntax:**export_emf=PATH\FILENAME**possible values:*

A valid path and filename.

implicit value:

None.

*example:**clarity.exe export_emf="c:\temp\test.emf"***export_results***description:*

Exports the results of active chromatogram into the specified *.TXT file (for more details see the chapter **Export Chromatogram** on pg 140). Setting of the export is taken from the active *.CFG or *.DSK file.

*syntax:**export_results=PATH\FILENAME**possible values:*

A valid path and filename.

implicit value:

None.

*example:**clarity.exe export_results=c:\tmp\results.txt***export_results_dbf***description:*

exports results of active chromatogram into specified database (*.DBF) file (for more details see the chapter **Export Chromatogram** on pg 140). Setting of the export is taken from the active *.CFG or *.DSK file.

*syntax:**export_results_dbf=PATH\FILENAME**possible values:*

A valid path and filename.

implicit value:

None.

*example:**clarity.exe export_results_dbf=c:\tmp\results.dbf*

export_run_time*description:*

Exports the *Run Time* value from the method currently opened on the Instrument as ANSI text into the specified file. Decimal separator is always ".". In case of any error the text "No current method." is exported instead.

syntax:

export_run_time=PATH\FILENAME

possible values:

Any valid file name, including the path.

implicit value:

None.

example:

clarity.exe export_run_time="c:\runtime.txt"

gal_apply*description:*

Applies the specified GPC calibration to the active chromatogram. It is recommended to use absolute paths because the calibration file will not be searched for in the working directory but in the calibration directory of the active project.

syntax:

gal_apply=PATH\FILENAME

possible values:

A valid path to the GPC calibration (*.GAL) file.

implicit value:

None.

example:

clarity.exe gal_apply=c:\tmp\1.gal

gal_save*description:*

Saves the active GPC calibration. Use together with the **instrument** parameter to specify the Instrument whose [Calibration](#) window holds the GPC calibration file you are interested in. GPC calibration file can be opened by specifying its path and filename. It is recommended to use absolute paths. This command is ignored when not in the GPC mode - use **cal_save** instead.

syntax:

gal_save

possible values:

None.

implicit value:

None.

example:

clarity.exe i=2 gal_save saves the calibration opened in the second Instrument.

gal_save_as*description:*

Saves the active GPC calibration under the specified name. Use together with the **instrument** parameter to specify the Instrument whose **Calibration** window holds the GPC calibration file you are interested in. GPC calibration file can be opened by specifying its path and filename. It is recommended to use absolute paths.

This command is ignored when not in the GPC mode - use **cal_save_as** instead.

syntax:

gal_save_as=PATH\FILENAME

possible values:

A valid path to the calibration (*.GAL) file.

implicit value:

None.

example:

clarity.exe gal_save_as=c:\tmp\results.gal

gpc*description:*

Switches the Instrument into the **GPC** or **Standard** mode.

syntax:

gpc=VALUE

possible values:

0 or 1.

implicit value:

None.

example:

clarity.exe gpc=0

gpc_recal_stdno*description:*

Together with the **gpc_recal_type** parameter enables to recalibrate an opened GPC calibration. The calibration is recalibrated by the active chromatogram at specified standard number and standard type. For successful recalibration it is necessary to specify the calibration file in the same command line as **gpc_recal_stdno** and **gpc_recal_type** parameters, which has to be both set to nonzero value.

syntax:

gpc_recal_stdno=VALUE

possible values:

0, 1, .. 20, 21 (Blank). Value 0 does nothing.

implicit value:

0

example:

clarity.exe test.gal gpc_recal_stdno=5 gpc_recal_type=1

gpc_recal_type*description:*

Together with the **gpc_recal_stdno** parameter enables to recalibrate an opened GPC calibration. The calibration is recalibrated by the active chromatogram at specified standard number and standard type. For successful recalibration it is necessary to specify the calibration file in the same command line as **gpc_recal_stdno** and **gpc_recal_type** parameters, which has to be both set to nonzero value.

syntax:

gpc_recal_type=VALUE

0, 1, 2.

possible values:

The meaning of values are as follows:

0 - does nothing (no recalibration).

1 - narrow calibration type.

2 - broad calibration type.

implicit value:

0

example:

clarity.exe test.gal gpc_recal_stdno=5 gpc_recal_type=1

hide_splash*description:*

Suppresses the splash screen that emerges when starting **Clarity**.

syntax:

hide_splash

possible values:

None.

implicit value:

None.

example:

clarity.exe hide_splash

instrument or i*description:*

Automatically opens an Instrument specified by the number. The command line parameter for user (and password) should follow, otherwise the opening is stopped on the [Login](#) dialog.

syntax:

i=VALUE

possible values:

1, 2, 3, 4. To open multiple Instruments, the CLARITY.EXE file has to be called separately for each Instrument.

implicit value:

1

clarity.exe i=3

example:

To work with two or more Instruments at once, you should use the MS Batch "START" command. See "Examples" on page 386 (for more details see the chapter **Command line parameters** on pg 386). section below the commands for more information.

mail_dump

description:

Disables the sending of memory dumps via e-mail. This function can also be set via the DATAPEX.INI file that may be stored in WINDOWS directory. Must be used when opening **Clarity** for the first time, otherwise it is ignored. This parameter is suitable for troubleshooting, when we certainly do not want to mail 200MB files. This parameter may be also used (and is overwritten by settings) in DATAPEX.INI file.

syntax:

mail_dump=VALUE

possible values:

0 or 1

implicit value:

1

example:

clarity.exe mail_dump=0

met_apply

description:

Applies the specified method to the active chromatogram (for more details see the chapter **Batch** on pg 67). It is recommended to use absolute paths because the method file will not be searched for in the working directory but in the default directory of the active project.

syntax:

met_apply=PATHFILENAME

possible values:

A valid path to the method (*.MET) file.

implicit value:

None.

example:

clarity.exe met_apply=c:\tmp\1.met

Open file *.PRM, *.MET, *.CAL, *.GAL, *.SEQ, *.PRJ, *.SST, *.NGA, *.AUDIT, *.CRG and *.LIB files

description: Opens the specified file(s) on the specified Instrument (using the **instrument** parameter).

syntax: **PATH\FILENAME**

possible values: The path and filename of the selected file.

implicit value: Empty.

example: **clarity.exe C:\CLARITY\PROJECTS\Work1.prj** or
clarity.exe "C:\CLARITY\DEMO1\Ethanol in blood.seq"

password or p

description: Tries to log the user on using the password specified.

syntax: **p=STRING**

possible values: String.

implicit value: Empty.

If empty, the system will deem that the user has an empty password.

example: **clarity.exe user=someuser p=mypassword**

prm_close

description: Closes the active *.PRM file opened. If any changes were performed on it, creates the dialog box asking whether these changes should be saved or discarded.

syntax: **prm_close**

possible values: None.

implicit value: None.

example: **clarity.exe prm_close**

prm_close_discard

description: Closes the active *.PRM file opened. If any changes were performed on it, discards the changes in the chromatogram without asking.

syntax: **prm_close_discard**

possible values: None.

implicit value: None.

example: **clarity.exe prm_close_discard**

prm_close_save*description:*

Closes the active *.PRM file opened. If any changes were performed on it, saves the changes to the chromatogram without asking.

*syntax:**prm_close_save**possible values:*

None.

implicit value:

None.

*example:**clarity.exe prm_close_save***prm_export***description:*

Exports the active chromatogram into the specified file. The type of the exported file is determined by its file name extension - supported file types are *.CDF, *.CHR, *.TXT, *.ASC. To learn more about file types See "Export Chromatogram" on page 140.

*syntax:**prm_export=PATH\FILENAME**possible values:*

A valid path and filename.

implicit value:

None.

*example:**clarity.exe prm_export=c:\tmp\1.chr***prm_import***description:*

Imports the specified file as a chromatogram. The type of a file imported is determined by its name extension - supported file types are *.CDF, *.CHR, *.TXT, *.ASC. To learn more about file types See "Export Chromatogram" on page 140. Imported chromatogram is stored in the TMP directory of the chromatographic software under the name IMPORT.PRM. If there is no such directory, it is stored in windows TMP directory under the same name. Existing *.PRM file is overwritten without warning.

A method must be used on and saved to the resulting chromatogram during import - we use the method currently opened on the Instrument, so its **Integration Table** parameters, linked calibration etc. are defined by it. Imported chromatogram is made active.

*syntax:**prm_import=PATH\FILENAME**possible values:*

A valid path and filename.

implicit value:

None.

*example:**clarity.exe prm_import=c:\tmp\1.chr*

prm_import_ex*description:*

Imports the specified file as a chromatogram. The type of a file imported is determined by its name extension - supported file types are *.CDF, *.CHR, *.TXT, *.ASC. To learn more about file types See "Export Chromatogram" on page 140. Imported chromatogram is stored in the DATA directory of the current project under the name FNAME.PRM, where FNAME is file name part of the imported file. Possible existing *.PRM file with the same name is overwritten without warning.

A method must be used on and saved to the resulting chromatogram during import - we use the method currently opened on the Instrument, so its **Integration Table** parameters, linked calibration etc. are defined by it. Imported chromatogram is made active.

syntax:

prm_import_ex=PATH\FILENAME

possible values:

A valid path and filename.

implicit value:

None.

example:

clarity.exe prm_import_ex=c:\tmp\1.chr

prm_save*description:*

Saves the active chromatogram. Use together with the **instrument** parameter (which is ignored by **Clarity Lite**) to specify the Instrument whose **Chromatogram** window holds the chromatogram file you are interested in. Chromatogram file can be opened by specifying its path and filename. It is recommended to use absolute paths.

syntax:

prm_save

possible values:

None.

implicit value:

None.

example:

clarity.exe prm_save

prm_save_as*description:*

Saves the active chromatogram under the specified name. Use together with the **instrument** parameter (which is ignored by **Clarity Lite**) to specify the Instrument whose **Chromatogram** window holds the chromatogram file you are interested in. Chromatogram file can be opened by specifying its path and filename. It is recommended to use absolute paths.

syntax:

prm_save_as=PATH\FILENAME

possible values:

A valid path to the calibration (*.PRM) file.

implicit value:

None.

example:

clarity.exe prm_save_as=C:\CLARITY\Work1\Test.PRM

recal_level*description:*

Recalibrates the opened calibration file at the specified level with an active chromatogram. The calibration file must be specified in the same command line as **recal_level**. Value 0 does nothing.

syntax:

recal_level=LEVEL

possible values:

0, 1, .. 20, 21 (Blank)

implicit value:

0

example:

clarity.exe test.cal recal_level=5 Recalibrates the TEST.CAL calibration on the 5th level.

register*description:*

Registers **Clarity** file extensions.

syntax:

register

possible values:

None.

implicit value:

None.

example:

clarity.exe register

reintegrate*description:*

Any chromatogram opened to a Clarity session which was run with the **reintegrate** command line parameter will be reintegrated. By default, the opened chromatograms are not reintegrated.

syntax:

reintegrate

example:

clarity.exe reintegrate

warning: In Clarity 4.0, the same effect can be achieved by using the **trace_ia** command line parameter with a variable z. When these two are used together in one command, the last written parameter takes effect.

report_cmdline_in_bad_trace

description: Reports the start of a session in the BADTRACE.TXT file and corresponding command line parameters. Must be used when opening **Clarity** for the first time, otherwise it is ignored. This parameter is suitable for troubleshooting. This parameter may be also used (and is overwritten by settings) in DATAAPEX.INI file.

syntax: `report_cmdline_in_bad_trace`
`report_cmdline_in_bad_trace=2`

possible values: 0 - no command line in BADTRACE.TXT. Default value (nothing is added to BADTRACE.TXT without this parameter)
 1 - only basic traces are written into BADTRACE.TXT. Has the same meaning as report_cmdline_in_bad_trace without parameters.
 2 - also internal traces are written into BADTRACE.TXT.

implicit value: 1

example: `clarity.exe report_cmdline_in_bad_trace`

report_last_only

description: Suppresses the BADTRACE.TXT file at the start of **Clarity** and sends only one dump file. This function can also be set via the DATAAPEX.INI file that may be stored in WINDOWS directory. Must be used when opening **Clarity** for the first time, otherwise it is ignored. This parameter is suitable for troubleshooting.

syntax: `report_last_only`

possible values: None.

implicit value: None.

example: `clarity.exe report_last_only`

resume_seq

description: Resumes the sequence on specified Instrument (the Instrument and sequence file must be opened separately). This parameter is suitable for automatic testing or for LIMS interfacing in combination with seq_import.

<i>syntax:</i>	<i>resume_seq=VALUE</i>
<i>possible values:</i>	1 (for Instrument 1), 2 (for Instrument 2), 4 (for Instrument 3), 8 (for Instrument 4), and their binary combinations.
<i>implicit value:</i>	None.
<i>example:</i>	<i>clarity.exe resume_seq=3</i> (will resume sequence measurement on Instruments 1 and 2)
run	
<i>description:</i>	Runs a single run analysis on specified Instrument (which must be opened separately). Single run has higher priority than sequence run. This parameter is suitable for automatic testing.
<i>syntax:</i>	<i>run=VALUE</i>
<i>possible values:</i>	1 (for Instrument 1), 2 (for Instrument 2), 4 (for Instrument 3), 8 (for Instrument 4), and their binary combinations.
<i>implicit value:</i>	None.
<i>example:</i>	<i>clarity.exe run=7</i> (will run single analysis on Instruments 1, 2 and 3)
run_seq	
<i>description:</i>	Runs the sequence run on specified Instrument (the Instrument and sequence file must be opened separately). This parameter is suitable for automatic testing.
<i>syntax:</i>	<i>run_seq=VALUE</i>
<i>possible values:</i>	1 (for Instrument 1), 2 (for Instrument 2), 4 (for Instrument 3), 8 (for Instrument 4), and their binary combinations.
<i>implicit value:</i>	None.
<i>example:</i>	<i>clarity.exe run_seq=7</i> (will run sequence measurement on Instruments 1, 2 and 3)
seq_export	
<i>description:</i>	Exports the active sequence into the specified *.TXT file.
<i>syntax:</i>	<i>seq_export=PATH\FILENAME</i>
<i>possible values:</i>	A valid path and filename.
<i>implicit value:</i>	None.
<i>example:</i>	<i>clarity.exe seq_export=c:\tmp\seq.txt</i>

seq_export_utf8

<i>description:</i>	Exports the active sequence into the specified *.TXT file using UTF8 encoding.
<i>syntax:</i>	<code>seq_export_utf8=PATH\FILENAME</code>
<i>possible values:</i>	A valid path and filename.
<i>implicit value:</i>	None.
<i>example:</i>	<code>clarity.exe seq_export_utf8=c:\tmp\seq.txt</code>

seq_import

<i>description:</i>	Imports the specified *.TXT file as a sequence. Current sequence is replaced by imported one, Sequence window is opened. The command is ignored when the sequence is running. Sequence Options dialog is never opened after command line import. Import settings are the same as the last saved ones during manual import. File name of imported sequence is NONAME, so it is necessary to use seq_save_as to store imported sequence before run.
<i>syntax:</i>	<code>seq_import=PATH\FILENAME</code>
<i>possible values:</i>	A valid path and filename.
<i>implicit value:</i>	None.
<i>example:</i>	<code>clarity.exe seq_import=c:\tmp\seq.txt</code> . See "Examples" on page 386 (for more details see the chapter Command line parameters on pg 386).section below the commands for more information.

seq_import_append

<i>description:</i>	Imports the specified *.TXT file as a sequence. Current sequence will be appended with lines from the imported file, Sequence window will be opened and the sequence file will be saved. Can be performed when the sequence is running. Sequence Options dialog is never opened after command line import. Import settings are the same as the last saved ones during manual import.
<i>syntax:</i>	<code>seq_import_append=PATH\FILENAME</code>
<i>possible values:</i>	A valid path and filename.
<i>implicit value:</i>	None.
<i>example:</i>	<code>clarity.exe seq_import_append=c:\tmp\seq.txt</code>

seq_save*description:*

Saves the active sequence. Use together with the **instrument** parameter (which is ignored by **Clarity Lite**) to specify the Instrument whose **Sequence** window holds the sequence file you are interested in. Sequence file can be opened by specifying its path and filename. It is recommended to use absolute paths.

syntax:

`seq_save`

possible values:

None.

implicit value:

None.

example:

`clarity.exe i=2 seq_save` - saves the sequence opened in the second Instrument.

seq_save_as*description:*

Saves the active sequence under the specified name. Use together with the **instrument** parameter (which is ignored by **Clarity Lite**) to specify the Instrument whose **Sequence** window holds the sequence file you are interested in. Sequence file can be opened by specifying its path and filename. It is recommended to use absolute paths. Command is ignored when sequence is running.

syntax:

`seq_save_as=PATH\FILENAME`

possible values:

A valid path to the sequence (*.SEQ) file.

implicit value:

None.

example:

`clarity.exe seq_save_as=c:\tmp\results.seq`

set_sample_name*description:*

Sets sample name parameter for the next single run. It is possible to set the Instrument by the `i=(number of Instrument)` parameter.

syntax:

`set_sample_name=text`

possible values:

Any non-empty text.

implicit value:

None.

example:

`clarity.exe set_sample_name="sample name"` or
`clarity.exe i=2 set_sample_name="sample name"`

show_help_id*description:*

After calling the context help (via the **F1** key or *Help* button) a message box with the help filename and topic ID will be displayed. Hardware and Control modules need to be configured to the Instrument, otherwise a classical help will be displayed for its dialogs.

*syntax:**show_help_id**possible values:*

None.

implicit value:

None.

*example:**clarity.exe show_help_id***sno or o***description:*

Uses the specified serial number (*.SNO) file. Must be used when opening **Clarity** for the first time, otherwise it is ignored.

*syntax:**o=PATH\FILENAME**possible values:*

The path and filename of the serial number (*.SNO) file.

Empty path.

implicit value:

If the path is not specified, the system will look for the serial number (*.SNO) file in the current directory. If not found, the system will then look for the file in the same directory where the *.CFG and CLARITY.EXE files are stored and in the current working directory.

*example:**clarity.exe o=C:\CLARITY\clarity.sno***stop***description:*

Stops the single run/sequence measurement on the specified Instrument. This parameter is suitable for automatic testing. In case of sequence it stops both current acquisition and sequence itself.

*syntax:**stop=VALUE**possible values:*

1 (for Instrument 1), 2 (for Instrument 2), 4 (for Instrument 3), 8 (for Instrument 4), and their binary combinations.

implicit value:

None.

example:

clarity.exe stop=7 (will stop single analysis/sequence on Instruments 1, 2 and 3)

unregister

description: Unregisters **Clarity** file extensions.

syntax: *unregister*

possible values: None.

implicit value: None.

example: *clarity.exe unregister*

user or u

description: Logs on the specified user.

syntax: *u=STRING*

possible values: String.

implicit value: Empty.

If empty, the system will show the [Login](#) dialog.

clarity.exe u="John Doe"

example: **Clarity** will automatically log in the user account called "John Doe".

10.2.2 Examples

10.2.2.1 Example 1:

This batch file will subsequently open **Instrument 1** and **Instrument 2**. On **Instrument 1** it will open the DEMO1 project, load the 2506MULTI.PRM chromatogram and ETHANOL IN BLOOD.SEQ sequence. On **Instrument 2** it will open PAH_EPA.CAL calibration file.

This batch file can be found in the UTILS subdirectory of the **Clarity** station main directory under the name CLARITY_COMMANDLINE_EXAMPLE.BAT.

```
start C:\CLARITY\clarity.exe i=1 u=Administrator
C:\CLARITY\PROJECTS\Demo1.prj
C:\CLARITY\DEMO1\DATA\2506multi.prm "C:\CLARITY\DEMO1\Ethanol
in blood.seq"
C:\CLARITY\UTILS\delay.exe 5
start C:\CLARITY\clarity.exe i=2 u=Administrator
C:\CLARITY\PROJECTS\Demo2.prj C:\CLARITY\DEMO2\CALIB\Pah_
epa.cal
```

10.2.2.2 Example 2:

This batch file will open the **Instrument 1**, import 1.CHR file as a chromatogram, apply a calibration on it, export the chromatogram and chromatogram results and close the chromatogram (single line command), then it closes **Clarity** itself. This can be useful when **Clarity** has to cooperate with other programs, such as LIMS.

```
start C:\CLARITY\clarity.exe i=1 u=Administrator
C:\CLARITY\UTILS\delay.exe 5
start C:\CLARITY\clarity.exe prm_import_ex=c:\tmp\1.chr cal_
apply=C:\CLARITY_Data\1\Calib\1.cal prm_export=C:\tmp\export.chr
export_results=C:\tmp\export.txt prm_close_discard
C:\CLARITY\UTILS\delay.exe 5
start C:\CLARITY\clarity.exe exit
```

10.2.2.3 Example 3:

This batch file will open **Instrument 1** as *Administrator*, import sequence file stored in C:\CLARITY\SEQUENCE.TXT and save it in the same directory in **Clarity** file format (*.SEQ) and finally start the sequence. Before trying this example create the SEQUENCE.TXT file in the given folder, e.g. using *File - Export* command in the [Sequence](#) window to save existing **Clarity** sequence to *.TXT.

```
start C:\CLARITY\clarity.exe i=1 u=Administrator
C:\CLARITY\UTILS\delay 15
C:\CLARITY\clarity.exe seq_import=C:\CLARITY\sequence.txt
C:\CLARITY\UTILS\delay 3
```

```
C:\CLARITY\clarity.exe seq_save_as=C:\CLARITY\sequence.seq  
C:\CLARITY\UTILS\delay 10  
C:\CLARITY\clarity.exe run_seq=1
```

10.2.2.4 Notes:

It is possible to call-up the CLARITY.EXE file when another instance of **Clarity** is already running. The new instance must quit (because only one running instance of **Clarity** is allowed at a time). Before quitting this new instance passes parameters to the instance already running. If CLARITY.EXE is already running, the **sdk**, **sno**, and **cfg** parameters will be ignored. If CLARITY.EXE is already running and the specified Instrument is opened, parameters **user**, **password** and **dsk** will be ignored. For all opening files some restrictions apply:

- A sequence must not be running.
- There must not be an opened modal dialog.
- Permission must be obtained to open a method.
- It must be allowed to change project.

Delay is not a regular batch command, it is a utility that waits a specified amount of time (set in seconds). DELAY.EXE file can be found in the UTILS directory.

Note: The necessary delay time that has to be set in the batch file is dependent on the configuration of the station and command line parameters used.

Several restrictions apply to the command line parameters used in the command line or a batch file:

- Only one project file (*.PRJ) can be accepted from the command line (actually the first one) for one Instrument.
- Only one method file (*.MET) can be accepted from one command line (actually the first one).
- Only one sequence file (*.SEQ) can be accepted from one command line (actually the first one).
- Multiple chromatograms (*.PRM) can be accepted for opening from one command line.

When opening an Instrument, the first project, method, sequence, calibration and then the chromatogram(s) are opened. The location of the working directory will be changed to where the *.CFG file is stored.

10.3 clarity.ini file

The CLARITY.INI file located in the root directory of MS Windows (C:\WINDOWS by default) contains keys for in-depth tuning of the **Clarity** Chromatography Station. The file in plain text format is structured into sections introduced in brackets, e.g. [General]. Each section can contain keys with values. Some sections are dependent on specific control modules and description of their keys can be found in appropriate manuals.

Note: Use e.g. Notepad program to edit this file. Using MS Word may result in destroying the proper format of the file.

Lines starting with semicolon (;) are comments.

[General]

The parameters common to all acquisition modules are set in this section.

PulseTime

Sets the length of the pulse in ms. Pulse on selected digital output of the A/D converter can be triggered by specified event from the **Event Table** found in the [Method Setup - Event Table](#) dialog (for more details see the chapter **Event Table** on pg 73).

Use60Hz

Set the *Use60Hz* parameter to 1 if the power supply frequency in your country is **60Hz** (countries in both Americas). Default setting 0 indicates the **50Hz** power supply frequency (rest of the world).

EventThreshold = 1

Sets the minimal necessary change of the signal (both the detector signal and auxiliary signals) as an input of the event enabling to trigger the event again. For the detector signal, the value is set in mV (or equivalent units as set in configuration of acquisition device). Events are usually defined in the **Event Table** found in the [Method Setup - Event Table](#) dialog (for more details see the chapter **Event Table** on pg 73).

TestLostPackets=0

Tests the completeness of data from the A/D converter. When set to 1 **Clarity** will report any lost packets to the BADTRACE.TXT file located in the **Clarity** root folder (C:\CLARITY by default).

FirstRun=dd.mm.yyyy

Shows the date of the first running of the station after installation.

[PDA]

VertexCountLimit

Specifies how many vertices will 3D view of PDA data use for display. The lower the value, the faster the PDA 3D view will render but the more degraded will be resulting image quality. Higher value will give better

quality but slow responsiveness. Warning: setting value too high can cause communication failures with hardware and/or "Virtual memory too low." error. Default value is 120000.

Index

1	
1/Amount	275
1/Amount^2	275
1/Response	275
1/Response^2	275
1/sqrt(Amount)	275
1/sqrt(Response)	275
2	
21 CFR Part 11	356
3	
3D View	196
6	
6-Sigma Noise Evaluation	193
A	
A-B	198
A/D converter	82
A+B	198
Abort	55, 62, 267-268, 283, 310
Abort Sequence or Acquisition	310
About	11, 39
About dialog	39
Clarity tab	39
Readme tab	39
System Files tab	39
What's New tab	39
Abs	276
Access Rights	356
Access To	5
Acq. Method	217
Acquired	217
Acquisition	326
Acquisition Parameters	217, 342
Acquisition Signals	111
Acquisition tab	82
Active	300
Active calibration signal	356
Active Chromatogram	342, 346, 348, 356
Active chromatogram signal	356
Active detector	356
Active Sequence	313
Add	12
Group	182
Add All	263, 266
Add button	12
Add Existing	264
Add Group	269
Add Negative	180
Add Peak	265-266
Add Positive	179
Add User Column dialog	151
Advanced date and time formatting	65
Advanced tab	92
After Chromatogram & Method	350
After Each Method Change	19
After Results	350
After the End of the Report	350
AIA	134
AIA File	134
AIA format	357
All	342, 345-346, 348
All As Instrument 1	21
All Data	98
All Identified Peaks	90, 208, 230
All Peaks	90, 208
All Peaks in Calibration	90, 208, 230
All Possible Instruments	27
Allow Automated Export of Audit Trail	24
Allow Chromatogram GLP Mode	23
Allow Crossing	176
Always	341
Always on Top	30
Amount	59, 209, 224, 241, 243, 266
Amount%	224
Amount^2	275
Analysis	59
menu	58, 282
Analysis Subdir	48
Analyst	217
Anchor (Text alignment)	201
Animated Arrows	106
Append	98, 321
Append to Existing	304
Apply Settings to	136

Calibration Audit Trail	327	tab	330
Calibration Bracketing	316	window	118
Calibration Cloning In Sequence	91	Chromatogram Audit Trail	219, 327
Calibration Components	331	Chromatogram Comments	218
Calibration curve	250	Chromatogram Cursors	106
Calibration Curve	111, 258	Chromatogram File Name	62, 64
Calibration Description	270	Chromatogram Header	97, 230
Calibration File	88, 207, 357	Chromatogram Info	340
Calibration History	345	Chromatogram list	171
Calibration Level	60	Chromatogram List dialog	200
Calibration options dialog		Chromatogram Operation	84, 212
Calibration options tab	270	Chromatogram Signals	111
Defaults tab	275	Chromatogram tab	346
Calibration Options dialog	270	Chromatograms	67, 200
Calibration options tab	270	Chromatograms in Overlay	105
Calibration point - history	261	Clamp Negative	176
Calibration standard		Clarity	1
Graph	242	Clarity Offline	357
Calibration Standard	60, 357	Clarity Registration dialog	33
Calibration Standards	51, 53, 67	Clarity shortcuts	362
Calibration Subdir	48	Clarity Switch	2
Calibration tab	331, 344	Clarity tab	39
calibration template	252	clarity.cfg	9, 21
Calibration used as specified by	315	clarity.exe	359
user		clarity.psw	3
Calibration window	238	clarity.sno	36
Calibration Point - history	262	Clear 3D	197
Cancel	13, 51, 53, 72	Clear All Levels	245, 261
Cancel button	72	Clear All Responses	261
Capacity	234	Clear Certificate	7
Caption	232	Clear Responses	89, 255
Centroid	233	Clear Run Marks	307
Certificate	6	Clear Selected Level	245, 261
Certificate for Exporting to PDF	7	Clear Selected Value	245
Change Password	6	clipboard	256
Channel	357	Clipboard	99, 357
Character Encoding	98, 141	Clone	89
Chart	259	Clone on first recalibration (safe	316
Chart (running)	287	calibration usage)	
Chart (waiting)	287	Close All	299
Check for Updates	35	Close Chromatogram	200
Check Sequence	311	Close Standard	253
Choose User	142	Closed Window	357
Chromatogram	92, 97-98, 140, 218	Color	108, 110, 167, 258
AIA import	135	Colors	113, 287
description	118	Column	80, 97, 216
menu	171	Column Calculations	93, 234

Column Length	93, 234	Current Signals	280
Column Matching	304	Current State	25
Columns	152, 154	Current Text File	136
Command line parameters	364	Current Value	315
Commands		Curve Check	249, 273
Customize	31	Correlation	273
Comment	13	Deviation	273
Comments...	60	Curve Fit Type	241, 246, 275
Common for All Signals	280	Curve passes through Origin	247
Common Settings		Curve passes through Unknown	248
Axes Ranges	291	Customize	236, 277, 289
Common Settings by Auxiliary	169	Customize dialog	31
Signals Types		Commands tab	31
Complete Processing	68	Toolbars tab	31
Compound	273	Cut	306
Compound / Parameter(s)	231	Cut Negative	176
Compound Name	225, 266, 268		
Compound Table	243		
Compound Type	246, 266		
Compounds	345	D	
Compute with Origin	247	Daily Audit Trail	319
Compute with Unknown	248	Data Acquisition	279
Configuration	9, 326, 357, 361	Data Acquisition window	279
Configuration of the station	357	Data Processing String	357
Content was not changed	144	Data Recovery	63
Context Axes	289	Data Size	138
Control	56, 61	DataApex A/D Converter Setup	15
Control Module Is Not Installed	14	dialog	
Copy	220, 306, 322, 328	dBase File	99
Mathematical Operation	198	Decimal Delimiter	303
Copy from Chromatogram	206	Decimal Separator	99, 137, 141
Copy from Model	205, 359	Default	155
Copy from Template Method	206	Default Font	106
Correction Factor	241	Default Injected Volume	60, 273
Correlation	273	Defaults tab	275
Correlation Factor	249	Delete	4, 155, 307
Counter	63-64, 314	Group	182
Counter (%n)	105	Delete All Compounds	261
Create Archive	50	Delete Compound	245, 261
Create Label	201	Delete Selected Files	51, 53
Create New Project	49	Delimited By	99
Created	90	Delimiter	303
Created by	4, 46	Departure	268
Cubic	247	Description	5, 46, 48, 90, 216, 316
Current	344-345	Descriptions	25
Current co-ordinates of cursor	120	Desktop file	46
Current Level	266-267	Desktop File	4, 358
		Detail Info	128

Detailed View	45	Edit Calibration	5
Detect Negative	187, 213	Edit Chromatogram	5
Detection	80, 216	Edit Current Method	312
Detector	141	Edit Current Report	312
Detector Delay	189	Edit ISTD Amounts	210
Detector Error Messages	324	Edit Method	5
Detector tab	219	Edit Report Style	6
Detector Units	15	Edit Sequence	6
Detectors	18	Eff./I	233
Deviation	244, 273	Efficiency	233-234
Device Monitor	55	Electronic Signature	350, 358
Different amount of internal standard	250	Enable Autostop	81
Differentiate	198	Enable File Overwrite	63
Digital Output Control	25	Enable Response Value Change ...	273
Digital Outputs	25	Enabled	82
Digital Outputs for external Start/Stop	19	End	179
Dilution	60, 211	Ending Point	241, 248, 275
Direct control of chromatograph ...	358	Enhanced Format... ..	108
Directories	21	Equation	249
Directories tab	115	Error Opening Project	49
Disallow Calibration Response Change	23	ESTD	89, 208, 270
Disallow Chromatogram Files Overwriting	23	EV	295
Disallow Empty Reason of File Change	24	Evaluation	215
Disallow Spike Filter and Spike Removal	24	Evaluation process	87
Disallow User List in Login Dialog ...	24	Event Table	343
Display		Event Type	74
menu	156, 257, 285	EventThreshold	388
Display Mode	270	Excel	99
Displayed Data	98	Excess	233
Do Not Send Method to Instrument ...	20	Existing Groups	182
Do not show this dialog for remaining files	134	Exit	43
Down	81	exp	247
Download	35	Expiration Warning	4
Drift	191	Export	99, 302, 321
Drift Evaluation	193	Export Calibration Curve as Picture	253
		to Clipboard	
		Export Calibration Curve as Picture... ..	253
		Export Chromatogram	140
		Export Chromatogram dialog	140
		Export Chromatogram in AIA .	101, 146
		Export Chromatogram in AIA Format	69
		Export Chromatogram in EZChrom ASCII Format	69, 101, 146
		Export Chromatogram in Multidetector Format	
E			
Edit	155, 218		
menu	148, 256, 306, 322, 328		

Export Chromatogram in TXT	101, 146	CLARITY.EXE	21
Export Chromatogram in TXT Format	69	clarity.psw	3
Export Content	97	clarity.sno	36
Export Data	69, 96, 101, 145	IQ.html	39
Export Data dialog	96, 123	SEND.COM.EXE	77
Export DBF	321	SYSTEMINFO.TXT	38
Export directory	115	TEMPLATE.CAL	252
Export File	141	TEMPLATE.SEQ	301
Export Messages (Batch)	324	WHATSNEW.TXT	39
Export Sequence	302	Fill Series	308
Export text	321	Fill using increment	308
Exports	331	Filter	12, 214
Expression	151	menu	323
Ext. Start	217	First	276
Ext. Start Dig. Input	19	First Free Level	266
External Start/Stop	81, 217	First Row is Header	303
EZ Chrom ASCII Format	140	FirstRun	388
		Fixed	163, 165
	F	Fixed Axes	289
FFT Filter	189	Fixed Height	339, 346
Field Separator	141	Fixed Width	98
File	56	Floating Axes	289
menu	42, 122, 281, 301, 328	Flow Rate	80, 216
NONAME	42	Font	155, 201
File Format	140	Font Color As Active Signal	108
File list	50, 52	Font...	108
File	46, 99, 295, 297, 300, 303, 339	For Files Outside Current Project	341
Na-		Only	
me		Force Peak Name	180
File Operations	326	Form Font	336
File Sharing	358	Format	315
File Type	50, 52, 67	Forward Horizontal	175
Files		Free	224
*.ASC	140	Free Calibration	224, 246
*.CHR	140	From	165
*.DGZ	50, 53	from All Open Chromatograms	112
*.DSK	46	From Common	51
*.PRJ	49	from the Active Chromatogram	112
*.TXT	140	Only	
*LAST.RAW	63	From Width at 50%	93, 234-235
*RUN.RAW	63	from... to...	280
BADTRACE.TXT	38	Front Tangent	176
BEEP.EXE	77	Full Format	98
clarity.cfg	21	Functs	152
CLARITY.CFG	9		
clarity.exe	359		

G		Height%	
Gap for Date and Manual Signature	350	formula	224
GC	18	Help	57, 72
GC Control	95	menu	32, 117
GC Error Messages	325	Help button	72
GC Gradient	95	Hidden window	358
GC tab	219	Hide	155
General tab	104	Hide All	116, 155
Generate Chromatogram After Abort	23	Hide Column	221
Get Current	163, 165	Hide Column(s)	322, 328
Global Peak Width	85, 171, 213	Hide Columns	154
Global Threshold	85, 171, 213	Hide ISTD Peaks	90, 209
GLP	358	Hide Signal	200
GLP Info	343	Hide Value Units	154
GLP Info tab	219	History of Calibration Point	261-262
GLP Information	341	dialog	
GLP Options	23	Hold	55
Gradient tab	113, 170	Hziperbola	247
Graph	250		
Graph legend	120	I	
Graph Properties	159	I/V	296
Auxiliary signals tab	169	ID	182
Gradient tab	170	Identification Windows	266, 276
Signals tab	166	IDLE	279
Graph Properties dialog	258, 286	Idle Time	314
Axes Appearance tab	161	Idle Time before First Injection	314
Graph tab	160	Ignore Origin	247
Signal Axis tab	164	Image on the Left	339
Time Axis tab	162	Image on the Right	339
Graph tab	107, 160	Image Size	339
Gray Background	338	Import AIA File dialog	134
Grey Out Inactive Signals	166	Import Chromatogram	6, 133
Grid Lines	289	Import directory	115
Group	215	Import EZChrom ASCII dialog	136
Group-X row	225	Import From Line	137
Groups	182	Import of multi-detector	133
Grp	212, 224	chromatograms	
		Import of several files	133
H		simultaneously	
Half Sigmoid	247	Import RAW	139
Half Width	86, 214	Import Sequence	303
Header Font	336	Import Sequence Step 1 dialog	303
Header Height	339	Import Sequence Step 2 dialog	304
Height	246, 266	Import Text File dialog	136
		Import To Line	137
		Imported Data	138

In Fixed Format	97		
Include Chromatogram in SST	101	L	
Include in SST	69, 299	Lab. Header tab	338
Including Common	51, 53	Landscape	337, 346
Individual Settings		Language Selection	37
Axes Ranges	292	Last	276
Info	344	Last Login	6
Info Header	342	LC	18
Information Table	40, 44, 358	LC Control	95
Initial password	28	LC Error Messages	324
Initial State	25	LC Gradient	95
Inj. No.	300	LC Gradient tab	219
Inj. Vol.	297	LC tab	219
Inj. Volume	60, 211	Left Button Doubleclick Means	106
Inj. Volume Units	315	Left Identification Window	276
Injection Control	343	Left Window	239, 267
Injection State	332	Level X	241
Injections in Sequence	105	Levels	344-345
Insert New Line	306	Life Time	4
Instrument	40, 359	Limit of Detection	240
Instrument Control	343	Limit of Quanification	240
Instrument Directories for Projects	21	Limits	105
dialog		Line	107, 202
Instrument Directory	359	Line Down	306
Instrument Errors	331	Line Label dialog	202
Instrument Images	18	Line Style	202
Instrument Messages	331	Line Up	306
Instrument Method Sending dialog	19	Line 110-111, 167, 202, 258-259, 287	
Instrument Parameters	342	W-	
Instrument tab	216	i-	
Instrument Type	17	d-	
Instrument window	40	t-	
Instrument X	5	h	
Integration	89, 183	Linear	246
Integration Interval	86, 187, 213	Linearization X	241, 248, 275
Integration tab	84, 212	Linearization Y	241, 249, 275
Integration Table	84, 330, 343, 359	Link Calibration To Standard	253
Integration toolbar	183	List of files	45, 67
Invert	198	List of terms	356
Invert Run Marks	307	List of the drivers	13
IQ.html	39	List of the Equipment Connected to	18
Is ISTD	240	the Instrument	
ISTD	89, 208, 224, 246, 270	In	247
ISTD Amount	60, 210	In(Amount)	275
ISTD1 Amount	296	In(Response)	275
		Load Defaults	139
		Local Audit Trails	327

Local Peak Width	85, 184, 213	Messages tab	324
Local Table Menu	220	Method	127-128
Local Threshold	85, 186, 213	menu	71
Lock	175	Method Audit Trail	327
menu	29	Method Description	80
Locked File	359	Method File	359
Locked Instrument	29, 359	Method Name	298
LOD	225, 240	Method setup dialog	
log(Amount)	275	Calculation tab	88
log(Response)	275	Method Setup dialog	
log10	247	Acquisition tab	82
Logarithmic Amount	258	Advanced tab	92
Logarithmic Response	258	AS tab	95
Login		GC Gradient	95
menu	27	Integration tab	84
Login Dialog	27	LC Gradient	95
Logout	43	Measurement tab	80
menu	28	Method tab	332, 342
Logout All	28	Method...	60
LOQ	225, 240	Min. Area	86, 214
Lvl	298	Min. Height	86, 214
		Min. Length	4
	M	Minimal Area	188
Main Clarity window	1	Minimal Half Width	188
Main graph	119	Minimal Height	188
Main station directory	359	Mobile Phase	80, 216
main window	1	Mode	271
Main window as Clarity Switch	30	Model Method	359
Manual	271, 300	Modified	90
Manual Flow...	55	MODIFIED	28, 42, 44, 122, 125
Manually	315	Module Info	13
Margins	337	Moments	97
Margins [mm]	337	Monitor	
Mark for Run	307	menu	54
Marks	107	Mouse Wheel Step	105
Matching	93, 218	Move	195
Mathematical Operations	330	Move Down	335
Mathematical Operations dialog	198	Move from Archive	53
Mathematics	198	Move to Archive	51
Maximize Graph	236, 277	Move to the Bottom	335
Maximize Table	277	Move to the Top	335
Maximize Tables	236	Move Up	335
Measure Distance	172	Moving Average Filter	189
Measurement	216	MS Method	343
Measurement Conditions tab	216	Multi-detector configuration	359
Measurement tab	80	Multidetector Text Format	140
Merge	173		

N		Open / Close Instrument Window ...	326
Name	13, 18	Open / Close program	325
Nearest	276	Open Audit Trail	320
NetCDF format	357	Open Audit Trail Settings	6
New	3, 49, 88, 301, 334	Open Calibration dialog	252
New Column	155	Open Calibration Window ..	68, 100, 145
New Password dialog	6, 28	Open Chromatogram	127
New value	268	Open Chromatogram dialog	127
Next Zoom	156, 257, 285	Open Chromatogram from	131
NGA Result Table	97	Sequence	
NGA Summary Table	97	Open Chromatogram Sequentially ..	129
No Access	5	Open Chromatogram Window ..	68, 100
No change	93, 218	Open Chromatogram with Stored ..	101
No. of Decimal Places	155	Calibration	
No. of pages	347	Open Configuration	5
No. of Places	155	Open Files to Import dialog	133
No. of Points	272	Open Method	
Noise	191	Calibration	45
Noise & Drift	191	Chromatogram	45
Noise Evaluation	192	Sequence	45
NONAME	42	Open Method dialog	44
None	93, 218	Open RAW Data dialog	139
NORM	89, 208, 271	Open Sequence dialog	131
Note	81, 216	Open Standard	253
Number of detectors	138	Open User Accounts	5
Number of Instruments	16	Open with calibration	69
Number of Lines	338	Open with stored calibration	207
Number of Signals	271	Open... ..	48
O		Opened Instrument Image dialog ..	18
oem.Clarity.exe	21	Operand	198
Offline	357	Operation	198
Offset	163, 165, 168, 196	Ops	152
Offset & Scale	163, 165, 168	Options	103, 270, 275, 313, 339
Offset Subtraction	93, 218	Ordnr	224, 246
Chromatogram		Orientation	201, 337, 346
OK	51, 53, 71, 147	Orientation graph	119
OK button	71	Origin	241, 247, 275
Old Value	268	Original	163, 165, 168, 196, 339
On First Page Only	338	Other Messages	325
On New Page	335	Other tab	325
Only Active Signal	342, 346, 348	Other Users Can	5
Only Vertical Line	106	Output no.	25
Only with Export	70, 102, 146	Overlay	194
Open	301, 334	Overlay Mode	128, 194
		Overlay toolbar	126, 194
		Override Printer Setting	337

Override Settings for Opened Chromatograms	169	Postrun Setting dialog	100
Override Signal Units	164	pow10	247
Overwrite	321	Prefix	165
		Preserve Integration	68
		Preserve Signal Ratios	166
		Preset	136
		Pressure	80, 216
	P	Preview	137, 155, 304, 334
Page Setup tab	336	Preview Graph	107, 119, 156-157, 257
Parameter / Compound(s)	231	Previous Zoom	156, 257, 285
Parameter Changes	332	Print	34, 254, 334, 354
Parameter Header	230	Print Background Color of Graphs	336
Parameters	102, 147, 344	Print dialog	354
Parameters of Calibration	331	Print Full Paths in Filenames	341
Parameters	70	Print Preview	254
	Sets	Print Preview dialog	352
Passive	300	Print Range	347
Password	6, 142, 360	Print Results	68, 100, 145
Password Changed	6	Print Results to PDF	69
Password Restrictions	4	Print Results To PDF	100, 145
Password Reuse	4	Print to PDF	353
Paste	220, 306	Print To PDF	334
Paste Insert	307	Print to PDF Directory	115
Pause	309	Printed Version Info	340
PDA Method	343	Printer	334
pdf	115	Printout	360
Peak	178, 215	Prints	330-331
Peak Area Coloring	108	Proceed	70
Peak Color	240	Program to Run	70, 102, 146
Peak Selection	239, 267, 276	Project	48
Peak Tags	108	Project directory	360
Peak Tags Format	109	Project file	360
Peak Type	224, 240	Project Name (Directory)	48
Peak Width	85, 213	Project Setup dialog	48
Peak Width dialog	184	Projects	6
Perform	147	Prompt for Confirmation	20
Performance tab	232	Properties	159, 197, 286
Performance Table	232	Propertis	324, 330
Plate number	233-234	Protected Mode	360
Play sounds assigned to selected events	104	PulseTime	388
Point to Point	246	Purge... ..	55
Portrait	337, 346	Purpose of signature	144
Post Run Settings	6	Purpose of Signature	142
Postrun	100		
Postrun functions	299		
PostRun Setting (From Chromatogram Window)	145	Q	
		Quadratic	246

R	
Range	82, 163, 165, 217
RAW	360
Raw data file	360
RB	223, 240
Read	5
Read & Write	5
Readme tab	39
Ready Dig. Output	19
Rec No.	244
Rec. No.	242
Recalibrate	272
Recalibrate Group	269
Recalibrate Peak dialog	267
Recalibration	244, 268, 272
Recalibration Table	268
Recent Files	105
Redo	306
Refer	224, 246
Reference	246
Register	32
Register Clarity	34
Reject Negative	176, 214
Rejection section	86
Rel	276
Remove	11
Remove All	308
Remove Label(s)	172
Repair Directories and Open This Project	49
Repeat Injection	310
Replace	272
Replace by value from the first row	308
Report	72
Report button	72
Report Header tab	340
Report in Result Table	90, 208
Report in Summary Table	230
Report Setup	253
Report Setup dialog	333
Audit & Signatures tab	350
Calibration tab	344
Chromatogram tab	346
Lab. Header tab	338
Method tab	342
Page Stup tab	336
Report Header tab	340
Results tab	348
Sequence tab	349
Report Style	298, 360
Report style file	360
Reprocess by Instrument Method	68
Request confirmation when opening old file formats	104
Rescue Data	63
Rescue Data dialog	63
Reset All	236, 277, 289, 317
Reset Axes Ranges	290
Reset Status	307
Reset when	314
Reset Zero	284
Residium	249
Resolution	234
Resp. Fact.	242
Resp. Factor	240
Response	223, 241, 243
Response Base	209, 240, 246, 275
Response Factor	209, 225, 243
Response Factor as Response / Amount	274
Response^2	275
Restore	52-53
Restore Archiv	52
Restore Default Columns	322, 328
Result	199
Result Table	97, 222
Results	
menu	220
Results panel	120
Results tab	207, 222, 348
Resume	310
Resume Idle	55
Resume Sequence	282
Reten. Index	240
Retention Indexes use Log.	274
Interpolation with Unretained Peak	
Retention time	232, 239
Retention Time	108, 223
Rich Text Edit	66
Right Identification Window	276
Right Window	239, 267
RISTD	224, 246

RSTD	245	Search Criteria	272
Run	61, 295, 309	Search Window	239, 266, 276
Run Lines	314	Select	109
Run Sequence	55, 282	Select All	68, 220, 322, 328
Run Single	55, 282	Select All Files	50, 52
Run Time	81	Select Another Directory	49
RUNNING	61, 280	Select Another Project	49
S			
S/N	36	Select Certificate	143
Same amount of internal standard	250	Select Certificate dialog	7
Sample	59, 65, 217, 296	Select Control	26
Sample Amount	296	Select Control dialog	25
Sample Dilution	297	Select Detector	82, 84
Sample ID	59, 65, 216, 296	Select Method	6
Sample Identification	216	Selected Items	109
Sample Info	341	Selected Size	50, 52
Sample Rate	83, 138	Send Calibration by email	252
Sampler tab	219	Send Calibration by Email	47
Sampling Rate	217	Send Chromatogram by Email	47
Save	42, 49, 301	Send Method	61, 72, 282
Save as	49	Send Method button	72
Save As	42, 124, 132, 255, 301, 334	Send Method by Email	47
Save as Chromatogram	139	Send Method Confirmation	20
Save As Chromatogram	199	Send Method to Instrument	20
Save as Template	205	Send PDF	334, 353
Save Chromatogram	200	Send Printed PDF by Email	254
Save Import Settings	304	Send Report by Email	38
Save Method	42	Send reports about unsuccessfully	104
Save Method As	42	finished sessions	
Save Sequence and Delete Import	304	Separation	86, 190
File		Separator	137
Saving chromatogram in format of	132	Sequence	66, 304
CSW station		Description	294
Savitzky-Golay Filter	189	menu	309
Scale	90, 163, 165, 168, 196, 209	Sequence Audit Trail	327
Scale Factor	90, 209	Sequence file	360
Scale Maximum to	167	Sequence Files	67
Scale Minimum to	167	Sequence Options	313
Scale Signals Separately	166	Sequence Options dialog	313
Scale Subtraction	93, 218	Sequence State	332
Chromatogram		Sequence tab	331, 349
Scale to Active Signal	166	Sequence window	294
Scale to All Signals	166	Serial Number	36
Scale Y Mode	166	Session	319
Scheme of Data Processing	40-41	Set	88, 93, 218
		Set Active Signal	200
		Set All Now For Current Signal	276
		Set Autostop	283

Set Axes Ranges...	290	Show Value Units	155
Set Background Chromatogram	280	Show windows on the taskbar	104
Set Image Options	339	Show Workplace	107
Set Initial Colors	111, 166	Show Y Axis	112, 169
Set Level	260	Show Y Axis for	113
Set Model	205	Shutdown	55
Set Signal	260	Sigmoid	247
Set Signal Names	204	Sign	70, 142, 147
Set Signal Names dialog	204	Sign as Current User	142
Set to Initial	114	Sign dialog	142
Set Zero	284	Sign Using Certificate	142
Setting		Signal	280
menu	96	Signal Axis	
Setting of digital inputs/outputs for	19	Axes Ranges	292
external Start/Stop		Signal Axis tab	164
Setup	11	Signal Name	138
Setup Columns	154, 322, 328	Signal Range	279
Setup Columns dialog	154	Signal Units	15, 138
Setup Control Modules	10	Signal(s)	287
Shortcuts	362	Signal/Noise	225
Show	154, 167, 169	Signals	127, 342, 344, 346, 348
Show All	116, 154, 323	Signals & Curves tab	111
Show All Signals	230	Signals tab	166
Show Auxiliary Signals	287	Signature is Valid	144
Show Balloon Help	107	Simplified Peaks Tags	108
Show Both	236, 277	Single	59
Show Columns	154	Single Analysis dialog	59
Show Data Points	160	Skew	233
Show Details	144	Skip	267
Show Events	107, 286	Skip Vial	311
Show Gradient	113, 286	Slope Ratio	190
Show Grid	107, 258, 286	Snapshot	62, 283, 311
Show Hidden Column	221	Solvent D-A	114
Show Hidden Column(s)	322, 328	Solvent Peak	181
Show History	245, 261	Sort - Backword	68
Show Instrument X	323	Sort by Name	68
Show Labels	167	Sort by Time	68
Show Legend	107, 286	Sort Normal	68
Show Linearized Values	258	Source	52-53, 169
Show Noise	286	Spec. Values	152
Show Recalibration Points	258	Special Results	97, 348
Show Sequence Options	305	Spike Filter	189
Show Signal	200	Spike Removal	177
Show Subtraction Chromatogram	280	Splitter	360
Show System	323	sqrt(Amount)	275
Show Tables	236	sqrt(Response)	275
Show Total Flow	113, 286	SST	235, 360

SST Result tab	235	System Files	361
Standard	298	System Files tab	39
Standard Addition Measurement	316	System Suitability Test	360
Start	178	SYSTEMINFO.TXT	38
Start - Restart	81		
Start - Stop	81		
Start at	314		
Start Only	81	T	
Start Time	184, 186, 188	Table Fonts	106
Station Audit Trail	319	Table Headers	9
Station configuration	361	Table Inverted	230
Statistical Moments	93, 234	Tables Up	156, 257
Status	13	Tabs	361
Status Bar	280, 300, 361	Tabs of individual instruments	17
Status Line	40	Tabs of Instruments 1 - 4	17
Status table	361	Tail Tangent	176
Std	298	Tangent Area Ratio	86, 190, 214
Blank	298	Tangent Slope Ratio	87, 190, 214
Bypass	298	Target	51
Standard	298	Temperature	81, 216
Unknown	298	template calibration	252
STDADD	89, 208, 271	Template method	361
Stop	55, 61, 283, 310	template sequenc	301
STOP	279	testLostPackets	388
Stop Flow	55	Text	201
Stop Sequence or Acquisition	310	Text File	99
Stop Time	184, 186, 188	Text Format	98, 140
Store	94	Text Label dialog	201
Stored Calib.	299	Threshold	186
Sts.	294	Threshold dialog	186
Subtraction	92, 217	Tiled by	347
Subtraction Chromatogram	280-281	Time	280
Successful registration	34	Time A	84, 212
Suggest Value	185, 187	Time Axis	
Summary	344	Axes Ranges	291
Summary Options	230	Time Axis tab	162
Summary tab	228	Time B	85, 212
Summary Table	97, 228	Time Format	106
Summary Table Options dialog	230	Time Range	279
SV	295	Time Stamp	142, 144
Switch	30	Time Step	98
Switch panel	30	Time Units	138
Symmetry/Tailing [-]	233	Title	151, 162, 164
System		Title Font	110, 259, 287
menu	2	To	165
System Configuration	9	To Common	53
System Configuration dialog	9	Together	87, 175, 215
		Toolbar	277, 361
		Baseline	174

Integration	183	Use	160-161, 166, 169-170, 258, 286
Overlay	126, 194	U-	
Toolbars	236, 289	s-	
Customize	31	e-	
Total row	226	r	
Threshold	85, 213	O-	
Threshold Value	77	p-	
Trial Mode	36	ti-	
TTL output	25	o-	
Type of Chromatograph	17	n-	
		s	
		Use60Hz	388
		Used	239, 244
		User	4
		User Access Rights	5
		user accounts	3
		User Accounts dialog	3
		User Code	36
		User Column	151
		User Description	142
		User Details dialog	8
		User Details for	4
		User Info	6
		User List	3
		User Name	4
		User Options dialog	103
		Auxiliary Signals tab	112
		Axes Appearance tab	110
		Directories tab	115
		General tab	104
		Gradient tab	113
		Graph tab	107
		Signals & Curves tab	111
		Username dialog	27
		V	
		Valid	344-345
		Valley	87, 174-175, 215
		Valley To Valley Slope	190
		ValleyToValley Slope	86, 214
		Value	85, 184, 187-188, 213
		Value Font	110, 259, 287, 336
		Variables	153
		Variance	233
		Vendor	13
		Verify	144
U			
Uncal	89, 208		
Uncal. Base	91		
Uncal. Response	91, 227		
Undo	306		
Unidentified Peaks	91		
Units ...15, 151, 162, 165, 209, 268, 272			
Units After Scaling	91		
Units Fon	110		
Units Font	259, 287		
Units Type	162, 164		
Unknown	248, 298		
Unprotected mode	361		
Unretained Peak Time	234		
Unretained Time	93		
Unselect	109		
Unselect All	68, 109		
Unzoom	156, 257, 285		
Up	81		
Update Calibration Response	273		
Update Retention Time	273		
Update the Station	35		
Use Default Font	106, 155		
Use ISTD	240		
Use Printer Setting	337		
Use Scale Factor	90, 209		

Verify Signatures	144
Version	46
VertexCountLimit	388
Vial No.	300
View	57, 88
menu	30, 103, 317, 323, 329
Visible	162, 164
Voltage Range	279

W

W05	224
WAITING	279
Warn before running already	105
measured sequence	
Warn when maximum zoom	105
reached	
Weight	272
Weighting Factor	272
Weighting Method	241, 248, 275
What's New tab	39
WHATSNEW.TXT	39
Window	
menu ...	116, 237, 278, 293, 317, 326
Window header	118
Window title bar	361
Windows Default	108
Without compressing	51, 53
Without Unknown Sample	248

Y

Y Multiplier	138
---------------------------	------------

Z

Zoom Button	106
Zoom In	156, 257, 285
Zoom Out	156, 257, 285